

Cell Sorter

Operator's Guide

Chapter 1 Overview	
Chapter 2 Preparation	
Chapter 3 Basic Operation	
Chapter 4 Configuring Experiments	
Chapter 5 Analysis	
Chapter 6 Sorting	
Chapter 7 Window Description	
Chapter 8 Maintenance	
Appendix A Operating Principles	
Appendix B Miscellaneous	
Index	

Table of Contents

Software Revision History	8
Using the PDF Manual	11
Model Name and Function Table	12
Sample Line Support	12
Components and Documentation	13

Chapter 1 Overview

Main Features	14
Cell Sorter Block Diagram	16
System Configuration	17
Name and Function of Parts	18
Front Panel	18
Internal View (Front)	19
Internal View (Side)	23
Rear Panel	24
Fluidics Cart	25
Sorting Chip	27
Main Window	28
Cell Sorter Software Operating Environment	29
LCD Monitor	30

Chapter 2 Preparation

Checking Device Connections	31
Filling the DI Water Tank	32
Filling the Sheath Fluid Tank	32
Filling the Ethanol Tank	33
Preparing a Sample	33
Loading a Sample Tube in a Sample Tube Holder	33
Preparing Automatic Setup Beads	34
Sample Temperature Control	34
Preparing Collection Tubes	35
Loading 8-well Strips (SH800S, SH800Z, and SH800)	37
Attaching the Splash Guard (for use with multi-well plates) (SH800SP and SH800ZP)	38
Loading a Multi-well Plate (SH800SP and SH800ZP)	38
Loading a 384-well PCR Plate (SH800SP and SH800ZP)	39
Removing and Attaching the Plate Holder Adapter (SH800ZP and SH800SP)	39
Configuring Users	40
Adding Users	40
Editing User Settings	41
Changing Passwords	41

Chapter 3 Basic Operation

Workflow	43
System Startup	44
Logging In	45
Initial Instrument Setup	46
Using the Sorting Chip	46
De-bubbling the Sheath Filter	48
Cleaning the Sample Line	49
Automatic Calibration	50
Creating an Experiment	52
Creating a New Experiment	52
Acquiring Data	53
[Acquisition] Tab Main Functions	54
Controlling Data Acquisition and Recording	55
Customizing the Worksheet	55
Analyzing Data	56
Sorting	58
Supported Sorting Methods for Sorting Chip Nozzle	
Diameter	58
Two-way Tube Sorting	59
Sorting onto 8-well Strips (SH800S, SH800Z, and SH800)	61
Sorting into Multi-well Plates (SH800SP and SH800ZP)	64
Performing Fluorescence Compensation	67
Flushing the Fluidics System	70
Shutdown	71
Specifying User Preferences	72
Exchanging the Sorting Chip	73
Inheriting Auto Calibration Results Between Users	74
Backing Up/Restoring the Database	76
Backing up the Database	76
Restoring a Database	76

Chapter 4 Configuring Experiments

Creating a Customized Experiment	77
Configuring an Experiment	80
Editing an Experiment	81
Editing a Sample Group	81
Editing a Sample Tube	81
Changing Component Settings	82
Exporting/Importing an Experiment	82
Exporting an Experiment	82
Importing an Experiment	83

Sharing Experiments	83
Searching Experiments	84
Saving Tube Data as FCS Files	84
Saving an Experiment as a Template	85
Configuring Detector Settings	86
Basic Configuration Method	86
Loading Reference Data Method	87
Adjusting Fluorescence Compensation	88
Using a Target Gate in the Gate Hierarchy	88
Adjusting Fluorescence Compensation Manually	89

Chapter 5 Analysis

Adding a Plot	91
Editing Plots	92
Changing Plot Type	93
Changing the Histogram Display View	93
Duplicating a Plot	94
Removing a Plot	94
Changing Axis Scale Type	94
Changing the Axis Parameters	95
Adjusting Scale Ranges	95
Adding Gates	97
Adding a Physical Gate	97
Copying a Physical Gate	98
Adding a Boolean gate to the Gate Hierarchy	98
Editing Gates	99
Displaying Multiple Plots Overlaid	101
Displaying Plots Overlaid (Overlay Display)	101
Customizing the Overlay Display	103
Deleting Overlay Layers	103
Deleting Overlay Gates	104
Displaying Statistics	104

Chapter 6 Sorting

Sorting Mode	106
Overview	106
Single Cell Mode	108
Normal Mode	109
Semi-Purity Mode	110
Purity Mode	110
Ultra Purity Mode	111
Semi-Yield Mode	111

Yield Mode	112
Ultra Yield Mode	112
Adjusting the Sort Position (SH800SP and SH800ZP)	113
Index Sorting	114
Configuring Index Sorting	114
Analyzing Index Sort Data	115
Adjusting Sort Parameters Manually (Experienced Users)	117
Configuring Custom Sorting Mode	117

Chapter 7 Window Description

[File] Window	119
[Information] window	119
[Print] Window	124
[Database] Window	125
[Help] Window	126
[Log] Window	126
[Create Experiment] Window	128
Main Window	130
[Experiment] Tab (ribbon)	130
[Cytometer] Tab (ribbon)	134
[Compensation] Tab (ribbon)	141
[Worksheet Tools] Tab (ribbon)	143
[Plot Tools] Tab (ribbon)	146
[Gate Tools] Tab (ribbon)	148
[Gates and Statistics] Table	149
[Acquisition] Tab (control pane)	149
[Experiments] Tab (control pane)	150
[Sort Control] Pane	153
Context Menu	157
Experiment Explorer Context Menu	157
Worksheet Context Menu	159

Chapter 8 Maintenance

Maintenance Schedule	161
Preparations for the Day Before Measurement	161
Maintenance for Each Startup Operation	161
Maintenance for Each Shutdown Operation	161
Monthly Maintenance	162
Periodic Maintenance	162
Miscellaneous	162
Fluidics Cart	163

Refilling the Sheath Tank	164
Emptying/Changing the Waste Tank	166
Refilling the DI Water Tank	167
Refilling the Ethanol Tank	169
Cleaning the Sample Fluidics System using Bleach	170
Cleaning the Collection Tube Holder with Embedded Cooling Agent (5 ml)	171
General Cleaning	172
Cleaning the Deflection Plates	174
Cleaning the Waste Catcher	176
Running Maintenance Mode	177
When a Cleaning Reminder is Displayed	180
Changing the Sample Loader O-Ring	181
Cleaning the Sheath Tank and Sheath Line	183
Changing Tank Air Filters	186
Cleaning the DI Water Tank and DI Water Line	187
Changing the PEEK Sample Line (PEEK sample line compatible models)	189
Changing the Sample Line and Probe (Non-PEEK sample line compatible models)	192
Changing the Probe Adapter (PEEK sample line compatible models)	195
Changing the Sheath Filter	196
Changing the DI Water Filter	198
Changing the Sintered Sheath Line Filter	200
Releasing Air in the Sheath Filter and Sorting Chip	202
Releasing Air in the DI Water Filter	203
Autoclaving the Sheath Filter and DI Water Filter	204
Ejecting the Sorting Chip Manually	206
Cleaning and Handling of Optical Filters	207
Disconnecting and Reconnecting the Fluidics Cart	208
Cleaning the Internal Sheath Line and DI Water Line using Ethanol	209
Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))	212

Appendix A Operating Principles

Fluid Flow	219
Sample Fluid	220
Sheath Fluid	220
Hydrodynamic Focusing	221
Droplet Sorting	222
Optical System	224
Light and Signal Generation	224
Laser Interrogation	225
Detection	225
Fluorescence Compensation	227

Appendix B Miscellaneous

Series Lineup	229
Optical Filter Patterns	230
Optical Filter Pattern 1	230
Optical Filter Pattern 2	231
Fluorescent Protein Optical Filter Pattern (Option)	231
Filter Pattern for BV/PE/PI	231
Fluorochrome Detection Matrix	232
Relationship between Sample Pressure and Sample Flow/ Event Rate	240
Troubleshooting	241
Error Messages	244
Optional Accessories	257
Software License	258
Environmental Notices	258
Index	259

The supplied installation disc includes the Operator's Guide for this unit in PDF format.
For more details, see "Using the PDF Manual" (page 11).

- Reproduction or duplication, in whole or part, of the software or operator's guide supplied with the unit, as well as renting or leasing of the software without the authorization of the right holder is prohibited under copyright law.
- Sony assumes no responsibility for damages, loss of income, or any claims from a third party arising out of use of the unit or supplied software.
- For complete terms and conditions of the warranty for the unit, refer to the warranty card included in the package.
- The software supplied with the unit cannot be used with any other units.
- Note that the specifications of the unit and supplied software are subject to change for improvement without prior notice.

Software Revision History

Functions Added in Version 2.1

The following functions have been added in Cell Sorter Software version 2.1.¹⁾

1) Name has changed from SH800 Software.

Multi-well plate sorting support using 70 µm sorting chip

Multi-well plate sorting, including 96-well plates, is now supported when using a 70 µm sorting chip.

Supported models

SH800SP / SH800ZP / SH800 (96-well plate models)

For details, see “Supported Sorting Methods for Sorting Chip Nozzle Diameter” (page 58).

Custom sorting mode

Custom sorting mode allows users to adjust the percentage of the droplet that will be used for making the sort decision of target events.

For details, see “Configuring Custom Sorting Mode” (page 117).

Plot overlay display function

Multiple plots can now be displayed overlaid.

For details, see “Displaying Multiple Plots Overlaid” (page 101).

Optical filter pattern selection screen

An optical filter pattern selection screen has been added to [Initial Instrument Setup], allowing you to perform auto calibration with the optical filter pattern used for measurement.

For details, see “Initial Instrument Setup” (page 46).

Support for 100 µm sorting chip (LE-C3210)

SH800 series running on software Version 2.1.0 or higher, support the use of 100 µm sorting chips from the LE-C32 series.

Sorting chip can be setup using either “Standard” setting or “Targeted” auto calibration settings. “Standard” setting can be used for sorting of small cells and “Targeted” setting can be used for sorting into multi-well plates and for sorting large cells.

New product

- LE-C3210
100 µm sorting chip (40/box)

For details, see “Using LE-C32 series sorting chips” (page 51).

Cleaning reminder

If ethanol cleaning has not been implemented for a set time, a cleaning reminder prompting you to perform ethanol cleaning is displayed on the login screen.

For details, see “Setting a Cleaning Reminder” (page 42) and “When a Cleaning Reminder is Displayed” (page 180).

Switchable views of histogram plots

Contour line view of histogram plots is not supported, allowing you to switch between outline view and filled area view.

For details, see “Changing the Histogram Display View” (page 93).

Troubleshooting wizard

If an error occurs during fluidics system check or during auto calibration, a wizard is displayed to help resolve the issue.

For details, see “If fluid flow does not start” (page 48) and “If an error occurs during auto calibration” (page 51).

Improved functionality

The following functions have also been added or expanded.

- Data for gates can be exported individually (page 100)
- [Detector & Threshold] has been added to [User Preference] (page 122)
- The auto calibration result image can be viewed on the [Chip Alignment Log] screen and [Sort Calibration Log] screen (page 126)
- Satellite droplet detection (page 140)
- Added function for configuring an instrument to obtain equivalent measurement results as another instrument by specifying data acquired from the other instrument as reference data (page 87)
- Added ethanol cleaning log to logs (page 127)

Functions Added in Version 1.8

The following functions have been added in SH800 Software version 1.8.

70 µm sorting chip support

Added support for 70 µm sorting chip for 2-way sorting.

Supported models

Standard feature: SH800S / SH800SP

Upgrade support: SH800Z / SH800ZP / SH800

For details, see “Supported Sorting Methods for Sorting Chip Nozzle Diameter” (page 58).

Consumables

- LE-C3107
70 µm sorting chip (40/box)

Ethanol cleaning supported for whole fluidics system

Cleaning wizard supports ethanol cleaning all flow channels in the fluidics system, including the sample line and the sheath line from the sheath tank to the main unit.

Supported models

SH800S / SH800SP / SH800Z / SH800ZP / SH800 (96-well plate models)

For details, see “Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))” (page 212).

Added inheritance function for auto calibration results between users

The auto calibration results can now be shared between different users using the same sorting chip.

For details, see “Inheriting Auto Calibration Results Between Users” (page 74).

Multi-well plate sorting support using 130 µm sorting chip

Multi-well plate sorting, including 96-well plates, is now supported when using a 130 µm sorting chip.

Supported models

SH800SP / SH800ZP / SH800 (96-well plate models)

For details, see “Supported Sorting Methods for Sorting Chip Nozzle Diameter” (page 58).

SH800 Software executable by all Windows accounts

SH800 Software can now be run from any Windows account.

Improved functionality

The following functions have also been added or expanded.

- Added single-cell mode for 3-droplet sorting using 100 µm/70 µm sorting chip (pages 108, 154)
- Added number of events within a gate to recording stop conditions (page 54).

- The display range of the Y axis (number of events) on histogram plots can be changed manually (page 96).
- Added sheath filter de-bubble function to initial setup when sorting chip is inserted (page 48).
- Added function to start recording in conjunction with 2-way sorting when sorting is started (pages 60, 121).
- Default axis type for fluorescence plots is configurable in [User Preference] (page 123).
- Added [Total Event], [Target Ratio], [Sort Rate], [Sort Efficiency], [Abort Count], and [Abort Rate] to the [Sorting Result] display in the [Tube Results] dialog (page 133).
- Added toggle function to the multi-well plate selection button (page 65).
- Added a function to restrict the sort calibration performed when the chip alignment is not completed successfully (page 135).
- Added function for switching the worksheet display to the assigned sample tube at the start of data acquisition (page 55).
- Added display function of the laser usage time (page 139).

Functions Added in Version 1.7

The following functions have been added in SH800 Software version 1.7.

Support for PEEK sample line

Support added for PEEK sample line that can be detached easily and safely.

A sample probe is also attached to the PEEK sample line.

For details, see “Sample Line Support” (page 12).

Support for 384-well plates (option)

Sample fluid can now be sorted into 384-well plates using this option.

For details, see “Preparing Collection Tubes” (page 35) and “Supported Sorting Methods for Sorting Chip Nozzle Diameter” (page 58).

Added sample line exchange to maintenance mode

The sample line can be replaced using a wizard when the unit is in maintenance mode.

For details, see “Running Maintenance Mode” (page 177).

Support for sample line cleaning at startup

Sample line cleaning can be performed during initial setup when a sorting chip is inserted.

For details, see “Initial Instrument Setup” (page 46).

Improved functionality

The following functions have also been added or expanded.

- Display of remaining sorting time (*page 154 and page 155*)
- Improved accuracy in sorting position adjustment for multi-well plate sorting (four corners of plates) (*page 113*)
- Additional test modes for multi-well plate sorting (*page 113*)
- Added option to skip rinsing of the outside of the sample probe during cleaning using DI water (*page 121*)
- Added multi-well pattern selection buttons for when configuring 8-well strip and multi-well plate sorting (*page 62 and page 65*)

Functions Added in Version 1.6

The following functions have been added in SH800 Software version 1.6.

Support for 130 µm sorting chips on all models

Added support for 130 µm sorting chip, even for 2-way sorting on 96-well plate models.

For details, see “Sorting” (page 58).

[Large Cell] can be selected in all sorting methods

[Large Cell] can now be selected in 8-well strip, slide glass, and 96-well plate sorting methods, in addition to the 2-way sorting method.

Functions Added in Version 1.5.1

The following functions have been added in SH800 Software version 1.5.1.

Supported by the SH800

The following functions were added.

- Index sorting
- Specifying a gate as calculation target for fluorescence compensation
- Data source indication
- Sample empty detection function
- Maintenance mode

Functions Added in Version 1.5

The following functions have been added in SH800 Software version 1.5.

Multi-well plate support (SH800 (96-well plate model))

Sample fluid can be sorted into 6, 12, 24, 48, and 96-well plates.

For details, see “Preparing Collection Tubes” (page 35) and “Sorting” (page 58).

Index sorting

An index sorting function has been added that links the data for sorted events to the individual well in which the events were sorted.

For details, see “Index Sorting” (page 114).

Specifying a gate as calculation target for fluorescence compensation

This function allows you to specify the event population for a gate as a negative or a positive control used for calculating fluorescence compensation.

For details, see “Using a Target Gate in the Gate Hierarchy” (page 88).

Data source indication

A data source component has been added to the Experiment Explorer that indicates the type of sorting collection device for saved data.

For details, see “Data sources” (page 159).

Sample empty detection function

This function detects when the level in the sample tube is too low, and stops data acquisition automatically.

For details about each item, see “[Sample Empty Detection]” (page 140) in the [Advanced Settings] dialog.

Maintenance mode

This function allows an operator to run ethanol cleaning or bleach cleaning without needing to log in to SH800 Software.

For details, see “Running Maintenance Mode” (page 177).

Using the PDF Manual

The Operator's Guide can be read on a computer with Adobe Reader installed.

You can download Adobe Reader free from the Adobe website.

- 1** Open the "Operator's Guide" folder in the installation disc.
- 2** Select and click on the Operator's Guide that you want to read.

If you have lost or damaged the installation disc, you can purchase a new one from your Sony distributor.

Model Name and Function Table

The following model names are used to refer to the various unit models in this document. The following functions are supported for each model.

Name	Model	2-way sorting	Multi-well plate sorting ¹⁾	Sample line	70 µm sorting chip support
SH800S	LE-SH800SA, SB, SC, SD, SE, SF, SG, SH	Yes	8-well strips (option) Slide glass (option)	PEEK sample line	Yes
SH800SP	LE-SH800SAP, SBP, SCP, SDP, SEP, SFP, SGP, SHP	Yes	96-well plates (6 to 96) 384-well plates (option)	PEEK sample line	Yes
SH800Z	LE-SH800ZA, ZB, ZC, ZD, ZE, ZF, ZG, ZH	Yes	8-well strips (option) Slide glass (option)	Sample line + sample probe (metallic) ²⁾ or PEEK sample line	No ⁴⁾
SH800ZP	LE-SH800ZAP, ZBP, ZCP, ZDP, ZEP, ZFP, ZGP, ZHP	Yes	96-well plates (6 to 96) 384-well plates (option)	Sample line + sample probe (metallic) ²⁾ or PEEK sample line	No ⁴⁾
SH800	LE-SH800AC, BC, CC, DC, EC, FC	Yes	8-well strips (option) Slide glass (option)	Sample line + sample probe (metallic) ²⁾ or PEEK sample line	No ⁴⁾
SH800 ³⁾ (96-well plate upgrade models)		Yes	96-well plates (6 to 96) 384-well plates (option)	Sample line + sample probe (metallic) ²⁾ or PEEK sample line	No ⁴⁾


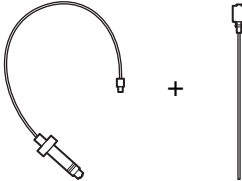
1) In this document 6, 12, 24, 48, 96, and 384-well plates are collectively referred to as “multi-well plates” or “96-well plates.”

2) Upgradeable to PEEK sample line.

3) SH800 models upgraded to support 96-well plate sorting are functionally equivalent to SH800ZP models.

4) Upgradeable to support 70 µm sorting chip.

Sample Line Support

	PEEK sample line (with probe)	Sample line + sample probe (metallic)
		
PEEK sample line compatible models	Yes	No
Non-PEEK sample line compatible models	No	Yes

Components and Documentation

System Components

The SH800 system comprises the following items.

- Main unit (1)
- Fluidics cart (1)

- Safety Guide (1)
- Operator's Guide (this document) (1)
- Installation disc (Operator's Guide, software) (1)

- Power cord (1)
- PC connection cable (1)
- Sample line (1)
For details about sample lines supported by each model, see "*Model Name and Function Table*" (page 12).
- Sample line (1) and sample probe (1) combination (non-PEEK sample line compatible models)
- Sheath filter (1)
- Sheath filter bypass line (yellow) (1)
- DI water filter bypass line (green) (1)
- DI water filter (1)
- Drip tray (1)
- DI water tank (1)
- DI water tank probe (1)
- Filter for DI water tank (1)

- Sheath tank (1)
- Ethanol tank (1)
- Ethanol tank probe (1)
- Air filter for ethanol tank (1)
- Waste tank (1)
- Waste tank nozzle (1)
- Air filter for waste tank (1)

- Sheath air line (clear) for fluidics cart (1)
- Sheath line (blue) for fluidics cart (1)
- Ethanol line (yellow) for fluidics cart (1)
- Waste line (red) for fluidics cart (1)
- Cart connection cable (1)

- Sample tube holder (0.5 ml) (1)
- Sample tube holder (1.5 ml) (1)
- Sample tube holder (5 ml) (1)
- Sample tube holder (15 ml) (1)
- Cleaning tube holder (30 ml) (1)

- Collection tube holder (5 ml) (1) (SH800, SH800Z, SH800ZP models)
- Collection tube holder with cooling agent (5 ml) (1) (SH800S, SH800SP models)
- Collection tube holder (15 ml) (1)

SH800SP and SH800ZP (standard equipment)

- 96-well plate holder (1)
- 96-well PCR plate holder (1)
- Plate holder adapter (1)
- Retaining screw (1)
- Splash guard (1)

SH800 and SH800ZP (options)

- 384-well plate holder (1)
- 384-well PCR adapter (1)

Documentation Structure

The system documentation comprises the following manuals.

- Safety Guide
Contains safety information, usage precautions, and specifications.
Read carefully before operating the unit.
- Operator's Guide
Describes the required preparation and basic operating procedures in order to use the instrument, adjustments and procedures for more advanced use of the instrument, and the configuration items in the Cell Sorter Software windows and dialogs.
"Chapter 3 Basic Operation" (page 43) provides an overview of the basic procedures, from setting up and calibrating the instrument through to acquiring data, analyzing data, and sorting.
Refer to the subsequent chapters for detailed information about operating procedures and instrument settings.

Main Features

The LE-SH800 series is a benchtop, high-speed, multilaser flow cytometer and cell sorter designed for research laboratory use. It employs a novel, replaceable, microfluidics, cell “sorting chip” that features high reliability and greatly simplifies cell sorting setup. It can be easily and quickly exchanged as needed for quick turnaround between measurements to minimize system downtime and to improve user work flow.

- Flow channels embedded within the chip for precision sheath fluid and sample fluid flow control
- High-uniformity droplet formation
- Automated chip loading and unloading system
- Easy optimization of system sorting and analysis conditions when changing sorting chip

The SH800 interrogates cells in samples using up to four lasers (405 nm, 488 nm, 561 nm, 638 nm), depending on the model. It measures forward scatter, back scatter, and six channels of fluorescent light. The wide range of excitation wavelengths supports the use of a wide variety of common fluorochromes for use in cell analysis.

The SH800 can measure intrinsic properties, such as size and internal complexity using forward scatter and back scatter analysis. It can simultaneously measure a wide range of extrinsic properties using fluorescent light emission analysis.

High-speed electronics allow the operator to specify multiple gates that isolate cells with defined characteristics for further processing. Gates can be applied to provide multi-parameter criteria for real-time identification of selected cell populations at speeds of 100,000 events per second (eps) for analysis and up to 10,000 eps for sorting. (100 µm sorting chip) The SH800 supports high-speed, two-way sorting to isolate populations for further study or testing.

An array of automated mechanisms is incorporated to enable easy use by non-specialist operators. These systems include:

- Sorting chip loading and alignment
- Sample tube loading
- Collection tube¹⁾ loading and alignment

- Fluorescence compensation
- Droplet formation, frequency, and calibration
- Sample line and sorting chip cleaning
- Temperature control of sample tube and collection tubes

1) In this document, sample tubes, 8-well strips, slide glass, and multi-well plates used for collecting sorted samples are collectively referred to as “collection tubes.”

Manual controls for fluorescence compensation and droplet formation are also supported in Cell Sorter Software.

The hardware user interface is designed to be as simple as possible to use. An LCD monitor on the front of the unit displays status information and notification of error messages.

The SH800 connects to standard 100 V to 240 V AC 50/60 Hz power supplies, and has a 17.5 °C to 27.5 °C (63.5 °F to 81.5 °F) recommended operating temperature range. It communicates with the host computer via the PC connection cable. The SH800 requires the connection of a laboratory compressed-air supply for operation. It can also be supplied using an external air compressor unit. An external fluidics cart, placed under or near the main unit, supplies the main unit with sheath fluid used during testing and ethanol used for cleaning, and collects the waste fluid from the main unit.

A single host computer running Cell Sorter Software controls all aspects of instrument operation, including cell analysis and sorting. The operator uses the software to create experiments for measuring data and to create plots and define statistics for analyzing data. The export functions are fully compliant with Flow Cytometry Standard version 3.0 or 3.1. The software provides quick access to all functions from the main window.

The SH800 has a compact footprint size. The benchtop unit has external dimensions (W × D × H) of 550 × 550 × 720 mm (21 3/4 × 21 3/4 × 28 3/8 in) and weighs 100 kg (220 lb).


Note

Exposing the main unit or the fluidics lines between the unit and the fluidics cart to drafts from a room air-conditioner may cause droplet formation to become unstable.

Tip

This document describes the basics of flow cytometry as it pertains to understanding how the SH800 cytometer operates. It is not intended as a reference on the subject of flow cytometry. There are many reference sources available, both in print and online, that you should consult for flow cytometry specifics, such as sample preparation, dye selection, analysis techniques, results publication methods, and other topics.


Cell Sorter Block Diagram

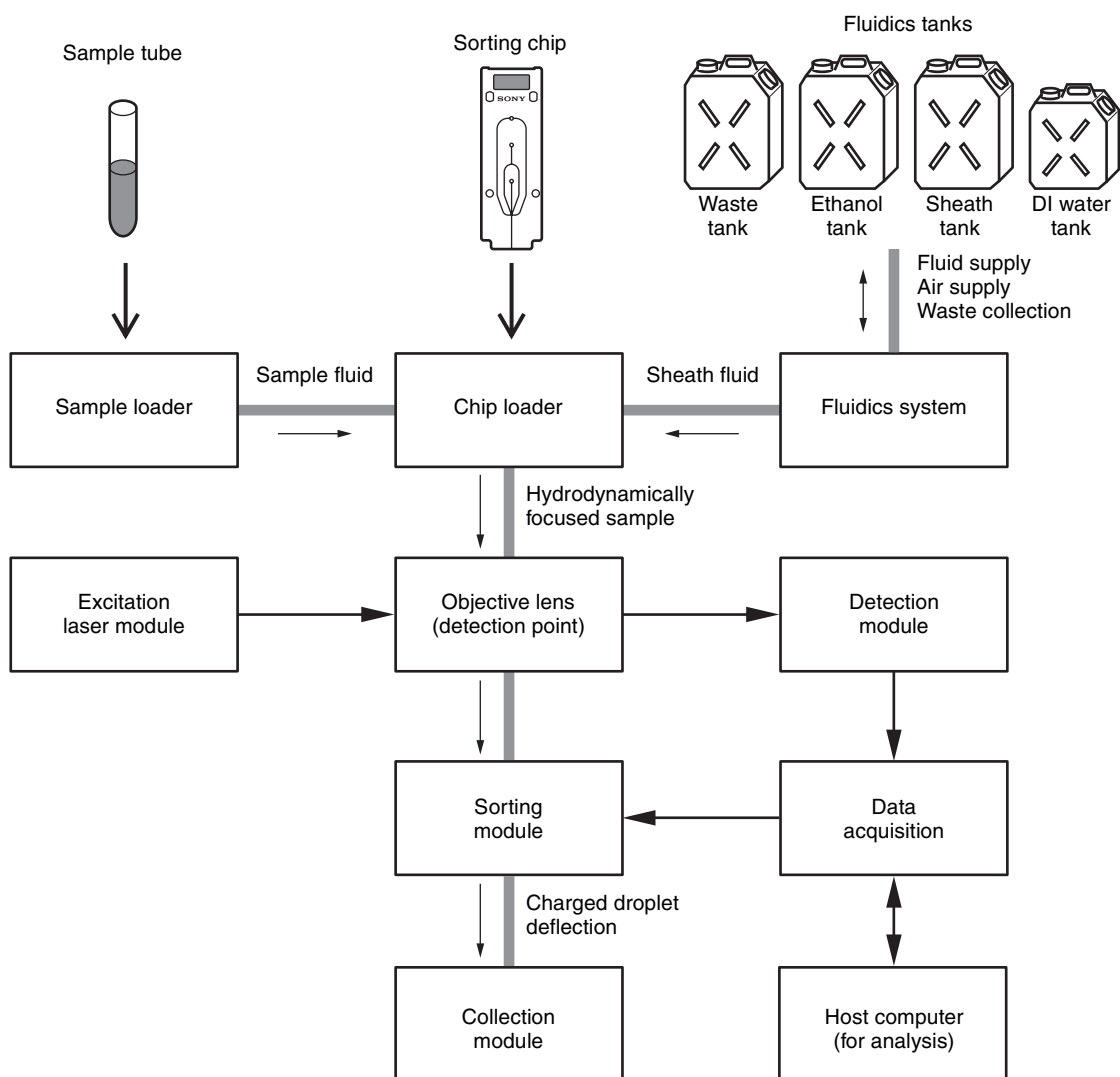

- The flow of waste fluid, ethanol, and DI water is not depicted in this diagram. All waste fluid is collected and pumped to the waste tank in the fluidics cart.


System Configuration

The SH800 system comprises the SH800 main unit, fluidics cart, host computer, and compressed air supply.

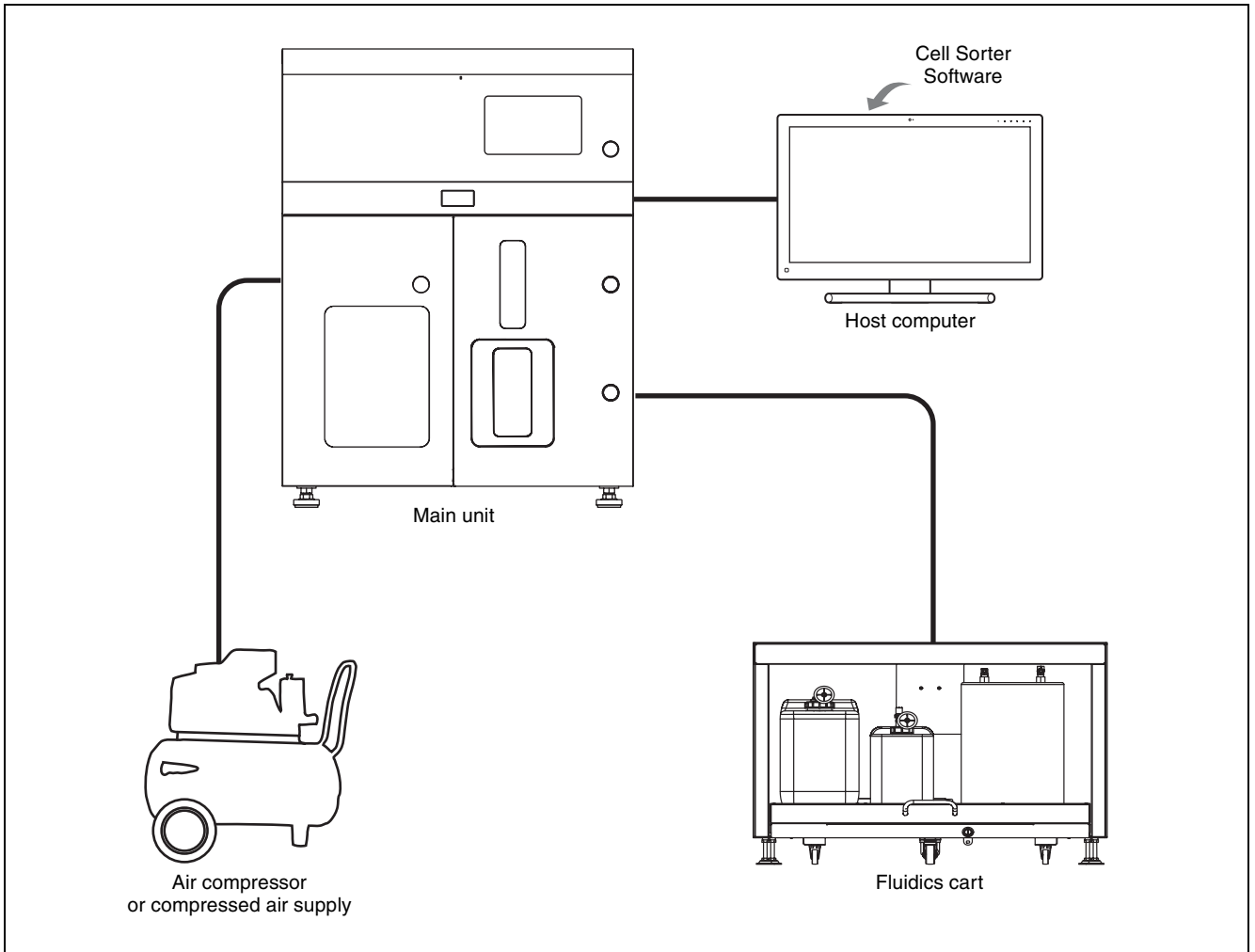

Name and Function of Parts

Front Panel

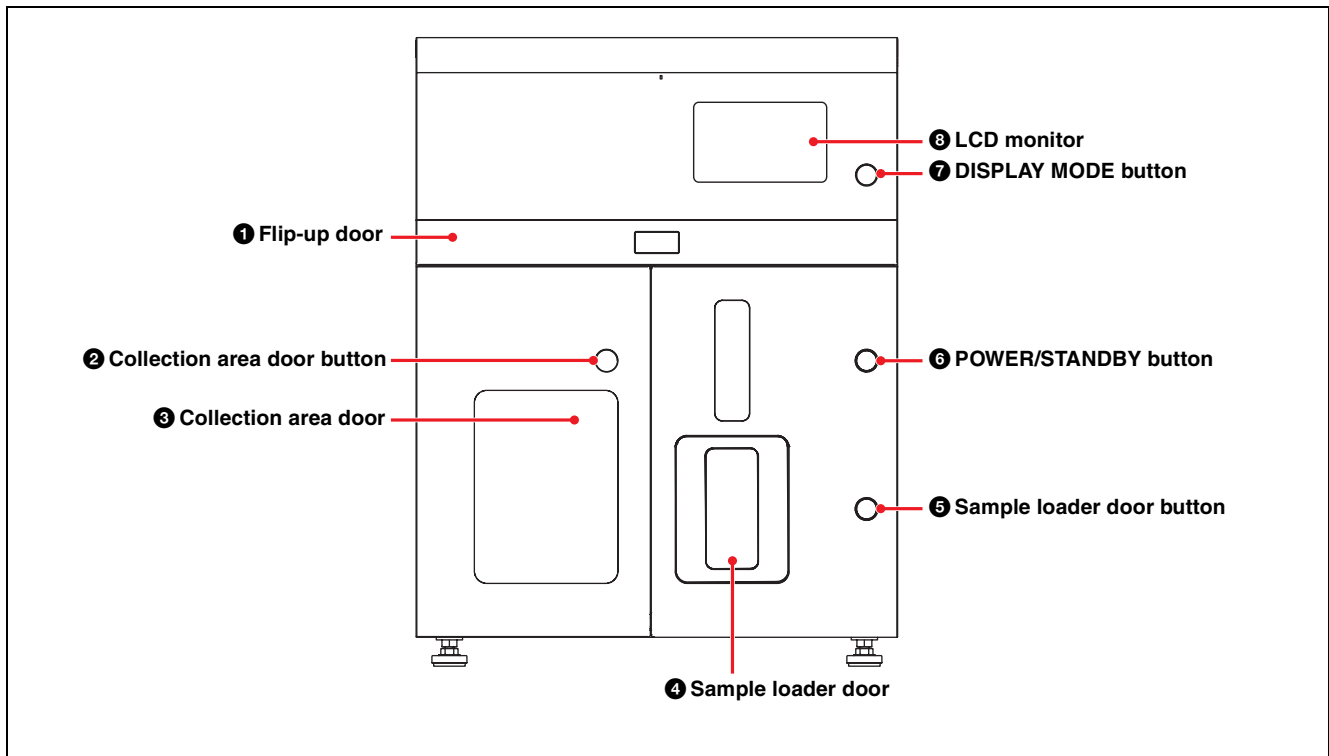

1 Flip-up door

Provides access to the chip loader for inserting and removing the sorting chip. The door also provides access to the optical filters, and the sample line and probe for routine maintenance.

Notes

- Make sure to close the flip-up door to run samples.
- The flip-up door opens and closes with assistance from the compressed air supply. Make sure the compressed air supply is turned on and supplying the rated air pressure when operating the door. For safety, a lock function operates to prevent the door from opening and closing when the air supply is turned off.

2 Collection area door button

Opens the collection area door.

3 Collection area door

Provides access to the collection area for loading and unloading collection tubes for sorting. It also provides access for cleaning and routine maintenance. A safety interlock switch disables operation while the door is open.

Note

A rubber packing seal is used to ensure the integrity of the collection area door. If the door does not open when the door button is pressed, the door may be sticking to the packing. In this case, open the door by simultaneously

pressing the button with one hand while pulling gently on the bottom edge of the door with your other hand.

4 Sample loader door

Provides access to the sample loader for loading and unloading samples for testing.

A safety interlock switch disables operation while the door is open.

5 Sample loader door button (OPEN/CLOSE)

Opens and closes the sample loader door. Once a sample has been lifted into the injection chamber within the sample loader and acquisition started, the button is disabled to prevent opening and closing of the sample loader door.

Note

Observe the following precautions when opening and closing the sample loader door.

- Turn on the power before operating the sample loader door button.
- When opening and closing the door while the power is turned off, make sure the compressed air supply is turned off. Failure to do so may result in damage or injuries.

6 Power button (POWER/STANDBY)

Switches the SH800 main unit power supply between ON and OFF. When power is ON, press and hold for about two seconds to switch Power OFF.

7 Display Mode button

Switches the LCD monitor display on/off.

8 LCD monitor

Displays instrument status information.


For details, see “LCD Monitor” (page 30).

CAUTION

- Do not place any objects in front of the doors.
- Be aware of the door edges when the flip-up door, collection area door, and fluidics maintenance door are open.


Internal View (Front)

The following illustration depicts the SH800SP.

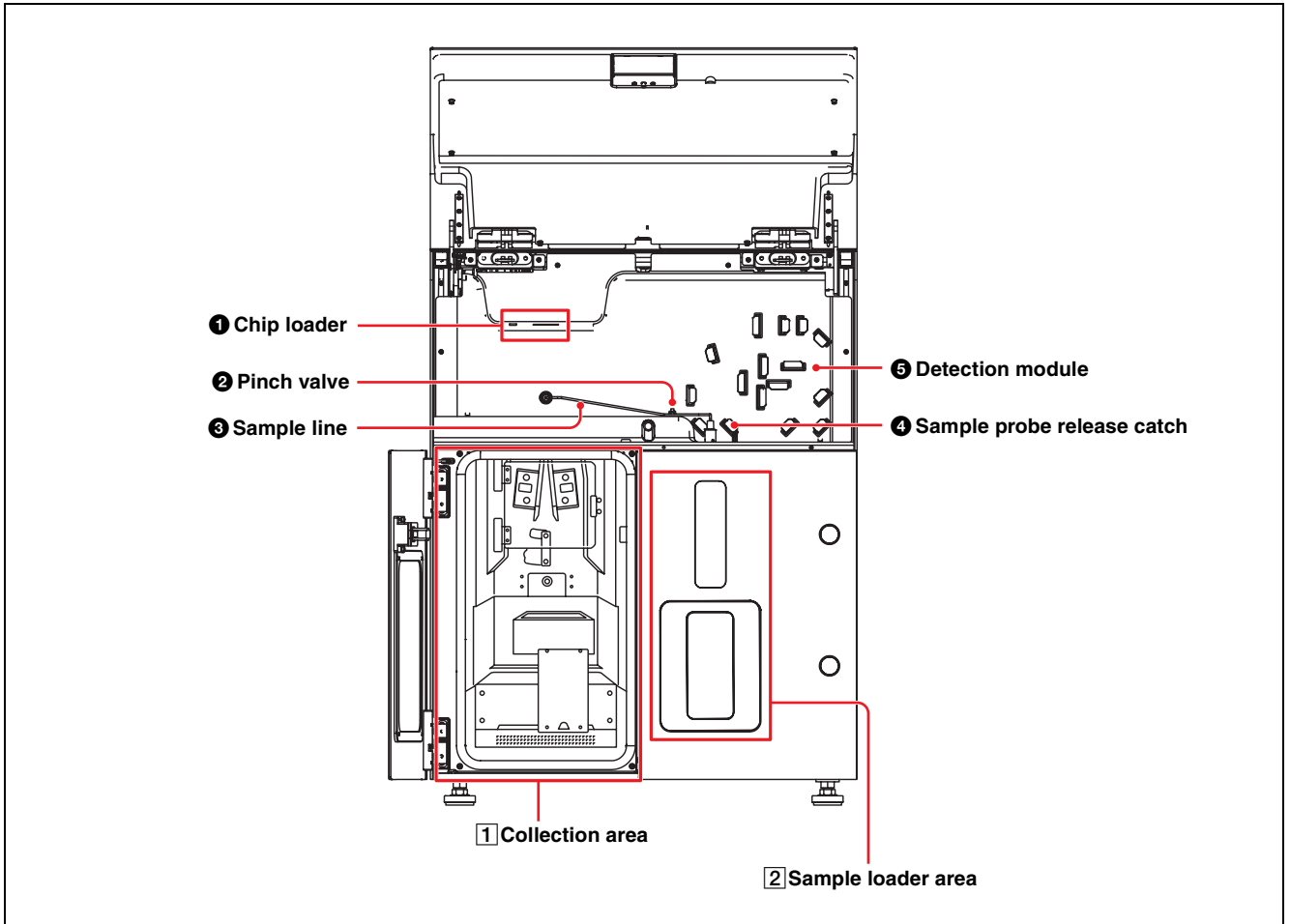

1 Chip loader

Loads and ejects the sorting chip. The chip is inserted in a top-loading slot.

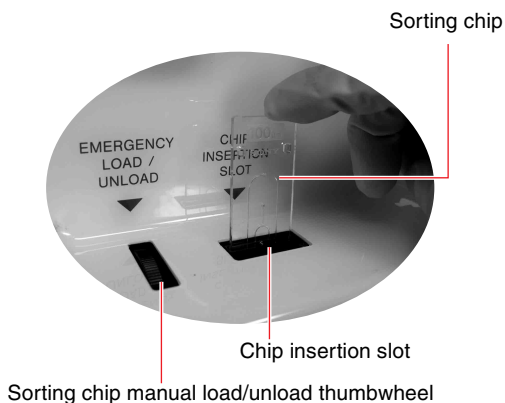

The chip unloading thumbwheel is used to eject the sorting chip if the chip will not unload automatically. For details, see “Ejecting the Sorting Chip Manually” (page 206).

2 Pinch valve

Controls the flow of sample fluid between the sample probe and the sorting chip.

3 Sample line

Transports the sample fluid from the sample loader tube to the sorting chip. The PEEK sample line combines the sample line and sample probe in a single component. The sample line and probe are maintenance items.

Notes


- Make sure the sample line is fully inserted into the pinch valve. Failure to do so may result in fluid leakage.
- Touching the sample line will directly affect the chip alignment. If you touch the sample line after the chip alignment has been adjusted, the chip must be realigned using the chip alignment function.
- Always leave the sample line connected when the unit is not in use for protection against dust.

For details, see “Changing the PEEK Sample Line (PEEK sample line compatible models)” (page 189) or “Changing the Sample Line and Probe (Non-PEEK sample line compatible models)” (page 192).

4 Sample probe release catch

Used when removing the sample probe or probe adapter.

For details, see “Changing the Sample Line and Probe (Non-PEEK sample line compatible models)” (page 192) or “Changing the Probe Adapter (PEEK sample line compatible models)” (page 195).

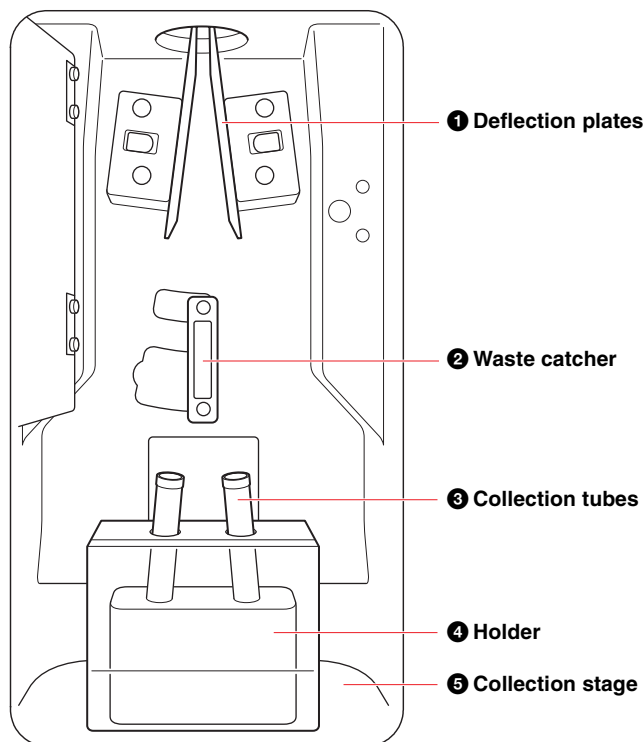
1 Collection area**SH800 model****5 Detection module**

Comprises an array of dichroic beam splitter longpass filters (LPF) to separate the collected fluorescent light into separate channels, and bandpass optical filters (BPF) to filter the light entering each detector. There are eight detectors: forward scatter (FSC), back scatter (BSC), and six fluorescent light (FL1 to FL6) channels for multi-color analysis. The optical filters can be changed to tailor the detection channels for specific measurements.


Notes

- The detection module uses components that are susceptible to static electricity, which may cause an error in operation. Do not touch the detection module during operation to prevent the transfer of static electricity.
- If you sense a discharge of static electricity from your body near the detection module during data acquisition, it is recommended that acquisition be paused and then resumed. This procedure restores normal gain settings, even in the event that static electricity affects the operation of the detection module.

SH800S and SH800Z models

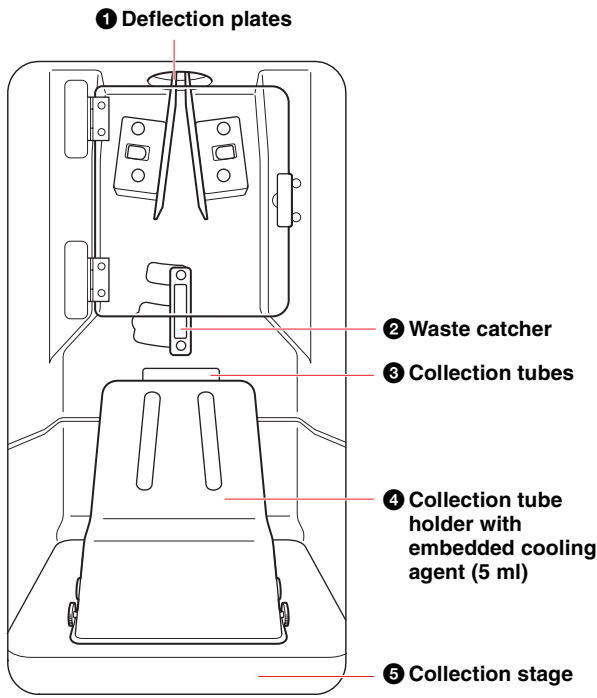

SH800SP and SH800ZP models

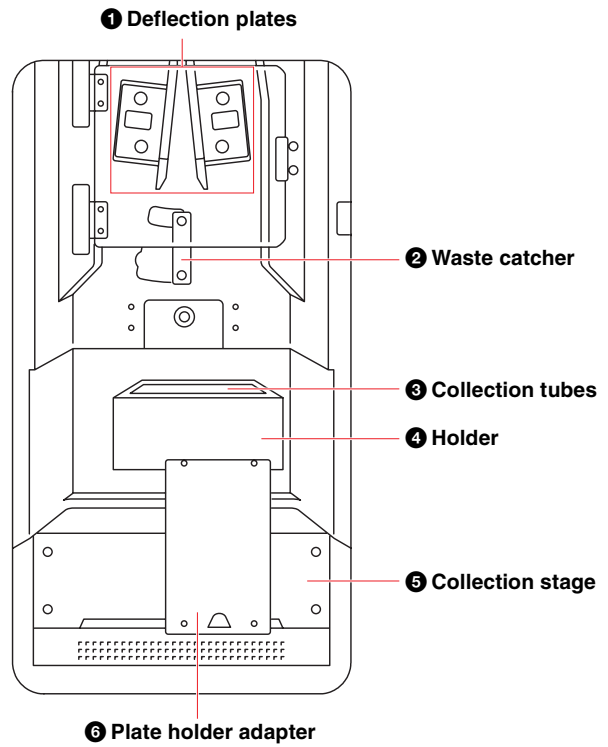

DANGER

Fluids may contain biological, chemical, or other agents. When handling collection tubes, always wear laboratory-use gloves, mask, protective goggles, and other protective clothing.

1 Deflection plates

Deflects electrostatically-charged droplets ejected from the nozzle of the sorting chip, sorting the droplets into the designated collection tubes.

Droplets that contain target cells for sorting are charged as they are ejected from the sorting chip nozzle, while other droplets containing non-target cells are uncharged and collected in the waste catcher. The target cells for sorting are specified using gates in Cell Sorter Software.

The trajectories of droplets that are sorted in the collection tubes are called side streams. For details, see “Side stream monitor” (page 154).

DANGER

The deflection plates are charged at an extremely high voltage. Contact with charged deflection plates can result in serious injury or electrical shock. Do not touch the deflection plates while power is connected to the main unit.

Note

Side stream calibration and data acquisition can be adversely affected by bright light shining directly on the side stream monitor. Do not place the front of the instrument near sources of bright light, such as lamps or unshaded windows.

Tip

Sorting may not occur correctly if there is any condensation on the deflection plates or if they are otherwise wet. Wipe off any condensation or water droplets.

2 Waste catcher

Collects the uncharged, non-target droplets. The collected droplets are passed to the waste tank in the fluidics cart.

3 Collection tubes

Collects the electrostatically-charged, target droplets. The SH800 supports the following collection tube supports. Each type of collection tube is used with a corresponding holder that mounts on the collection stage.


Collection tubes/ multi-well plates	Holder
15 ml conical tubes	Collection tube holder (15 ml)
5 ml round tubes	<ul style="list-style-type: none"> Collection tube holder (5 ml) Collection tube holder with embedded cooling agent (5 ml)
8-well strips	8-well strips holder ¹⁾
Slide glass	Slide glass holder ¹⁾
6-well plates	96-well plate holder
12-well plates	
24-well plates	
48-well plates	
96-well plates	

Collection tubes/ multi-well plates	Holder
8-well PCR strips	96-well PCR plate holder
96-well PCR plates	
384-well plates	384-well PCR plates ¹⁾
384-well PCR plates	384-well plate holder ¹⁾ + 384-well PCR adapter ¹⁾

1) Options.

For details about loading tubes and the types of collection tubes, see “Preparing Collection Tubes” (page 35).

2 Sample loader area

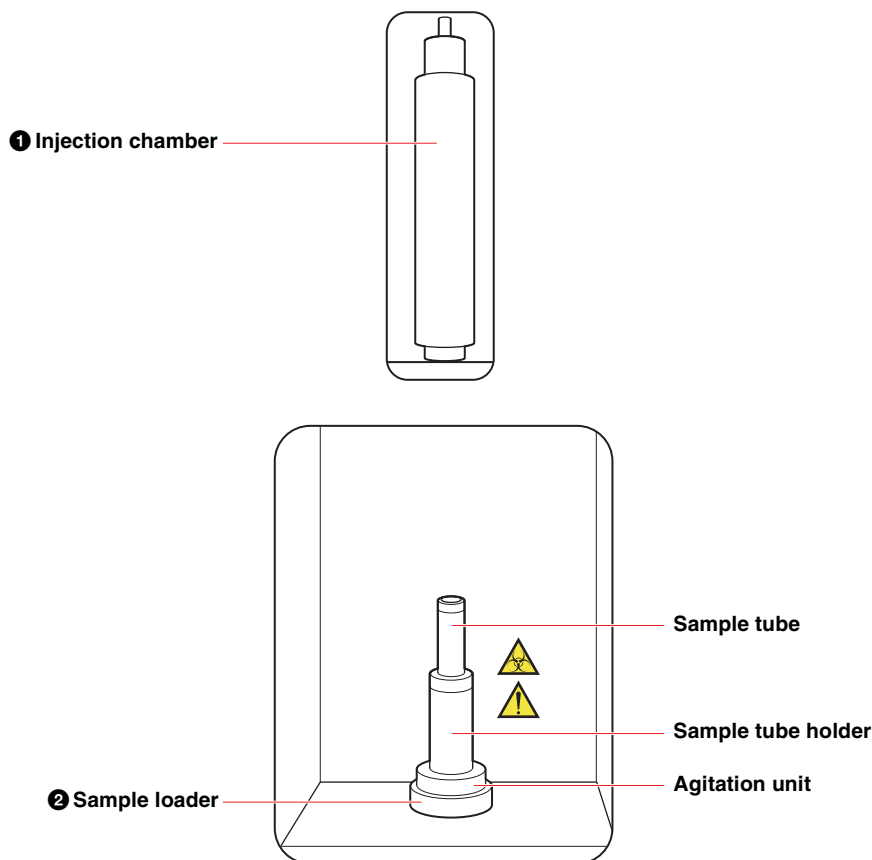

Notes

- Always wear laboratory-use gloves, mask, protective goggles, and other protective clothing when handling sample fluid to provide protection against biohazards.
- Clean the fluidics system regularly using a sodium hypochlorite liquid solution (bleach) to prevent clogging of the sample line and sample probe.
- Load the sample tube with its cap removed.

4 Holder

Holds the collection tube device.

5 Collection stage

Loads the collection tube device holder.

6 Plate holder adapter (multi-well plates)

Used for loading multi-well plate collection devices.

When running 2-way tube sorting, remove the adapter and load the corresponding collection tube holder.

1 Injection chamber

Extracts sample fluid from the sample tube, through the sample probe, and forces the fluid into the sorting chip.


2 Sample loader

Loads a sample tube holder containing a single sample tube containing sample fluid into the sample loader agitation unit.

Samples can be loaded in 0.5 ml or 1.5 ml micro tubes, 5 ml round tubes, or 15 ml conical tubes. A sample tube holder is provided for each size of sample tube.

For details, see “Loading a Sample Tube in a Sample Tube Holder” (page 33).

Internal View (Side)

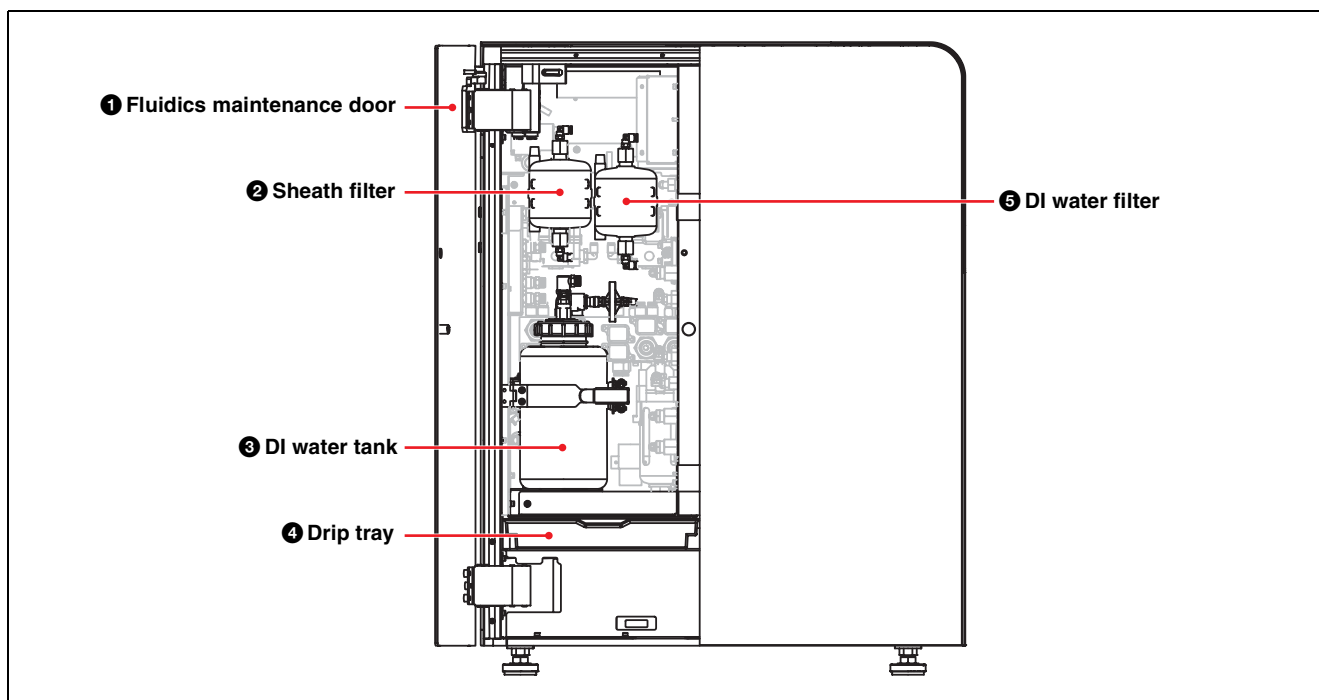

1 Fluidics maintenance door

Provides access to the fluidics system for performing maintenance.

CAUTION

Be aware of the door edges when the door is open.

2 Sheath filter

Filters the sheath fluid flowing between the sheath tank and the sorting chip. The sheath filter is a maintenance item.

3 DI water tank

Supplies deionized (DI) water used for flushing the fluidics system.

A sensor detects the level of DI water within the tank, and a message is displayed when the level drops below the minimum threshold.

The tank has a capacity of 1 liter (0.3 gallons US).

The DI water tank is a maintenance item.

4 Drip tray


Collects liquid spills, during routine maintenance, and leakages due to damaged tubing or improperly fitted connectors. The drip tray should be checked regularly for signs of leakage.

An alert is shown on the host computer if fluid is detected in the drip tray. If fluid is detected, resolve the source of the leak, clean the drip tray, and wipe clean all affected surfaces. The system must be restarted after a certain time has elapsed after the alert is raised.

5 DI water filter

Filters the DI water flowing from the DI water tank used for cleaning the sample probe. The DI water filter is a maintenance item.

Rear Panel

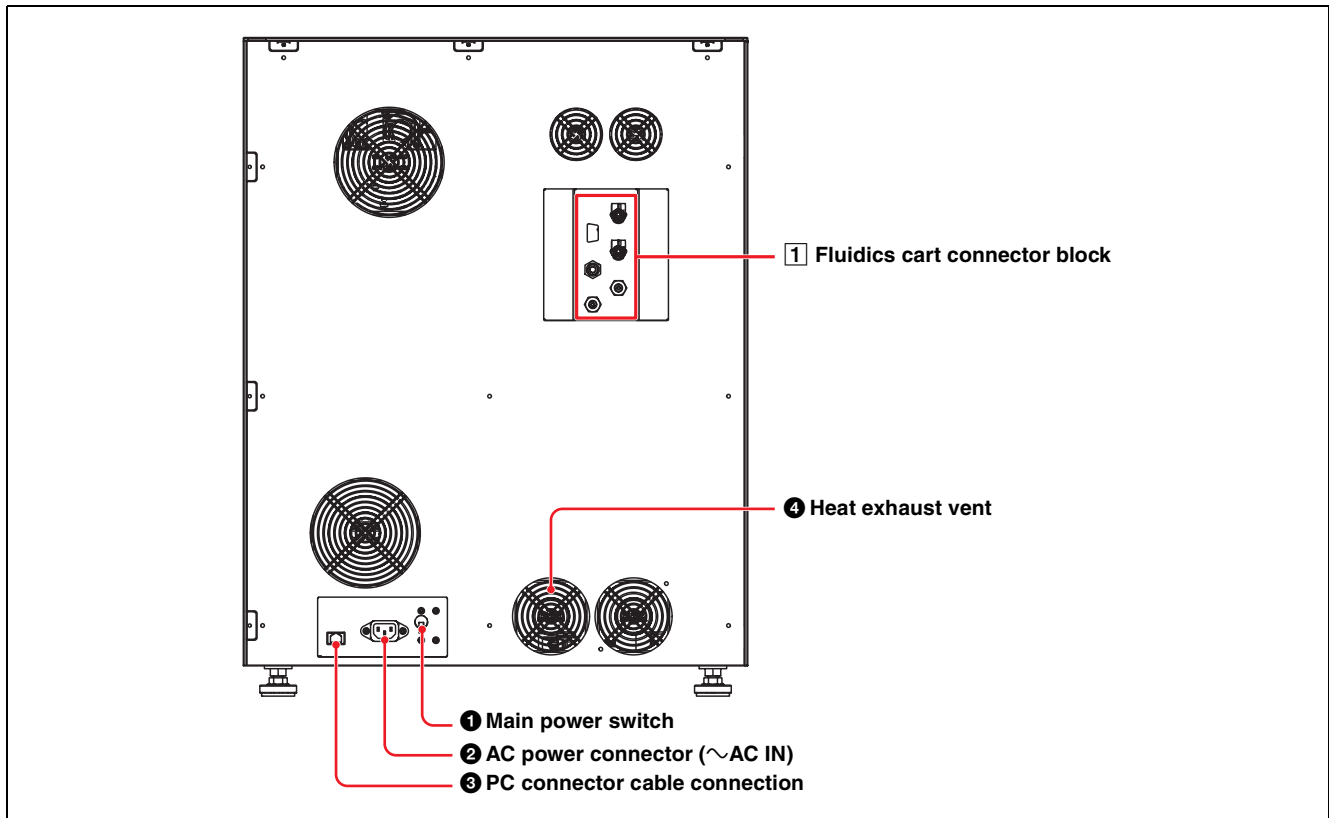

1 Main power switch (MAIN POWER)

Turns the power supply for the SH800 ON (up) and OFF (down). The breaker automatically disconnects the power supply if an electrical fault is detected.

2 AC power connector (~AC IN)

Connects to a standard AC power supply (100 V to 240 V AC, 50/60 Hz). Use only the supplied power cord to connect the SH800 main unit to a power supply.

Notes

- Use only within the specified supply voltage range.
- The power cord must be connected to a power outlet with an Earth connection.

3 PC connection cable connector (PC CONNECTION)

Connects directly to the host computer using the supplied PC connection cable.

Notes

- Do not disconnect the PC connection cable during operation. Data acquisition cannot be performed if the cable is disconnected.
- Do not connect any device other than the host computer using the PC connection cable. Device operation is not guaranteed if an external device is connected.

CAUTION

For safety, do not connect the connector to any peripheral device wiring that might apply a high voltage to this port.

Follow the instructions for this port.


4 Heat exhaust vent

Discharges heat generated from inside the unit.

Note

Maintain at least 10 cm (4 in.) space behind the main unit to allow for the dissipation of heat and to provide access to the power connector, main power switch, electrical connectors, fluidics cart connectors, and compressed air supply connectors.

1 Compressed air supply/fluidics cart connector block

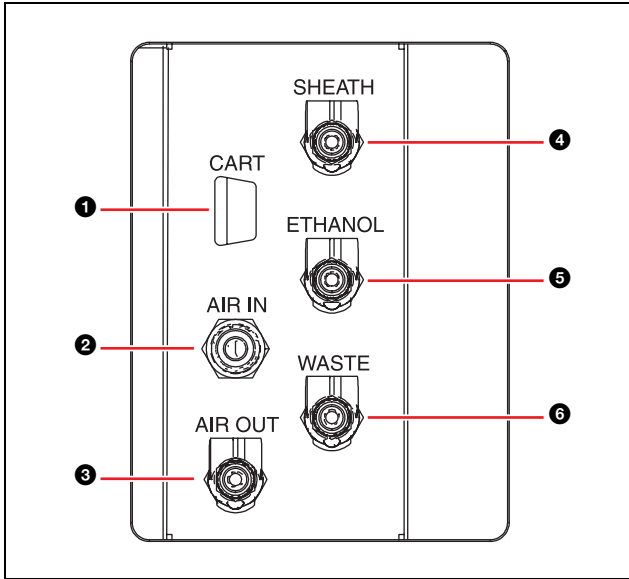

1 Cart connection cable connector (CART)
Connects to the cart connection cable connector on the fluidics cart. It is used to connect the sensors for the three fluid tanks in the fluidics cart.

Note

Use only the supplied cable to connect the main unit to the fluidics cart. Using third-party cables may cause incorrect operation or device failure.

2 Compressed air supply input connector (AIR IN)
Connects to the laboratory compressed air supply or a standalone air compressor.

The SH800 is designed to connect to 500 kPa to 650 kPa (73 psi to 94 psi) compressed air supply for stable output. Use polyurethane tubing with an outer diameter of 6 mm and an inner diameter of 4 mm.

3 Fluidics cart air line connector (AIR OUT)
Connects to the sheath air line connector on the fluidics cart (clear tubing). The air line supplies regulated compressed air for the sheath tank.

4 Fluidics cart sheath line (blue) connector
Connects to the SHEATH connector on the fluidics cart using the supplied sheath line (blue tubing).


5 Fluidics cart ethanol line (green) connector
Connects to the ETHANOL connector on the fluidics cart using the supplied ethanol line (green tubing).

6 Fluidics cart waste fluid line (red) connector
Connects to the WASTE connector on the fluidics cart using the supplied waste line (red tubing).

Note

The standard fluidics lines, sheath air line, and connection cable are about 2 m (6 ft) in length. It is recommended that the SH800 be placed on a table of height approximately 700 mm (28 in.) or greater, and that the fluidics cart be placed on the floor.

Fluidics Cart

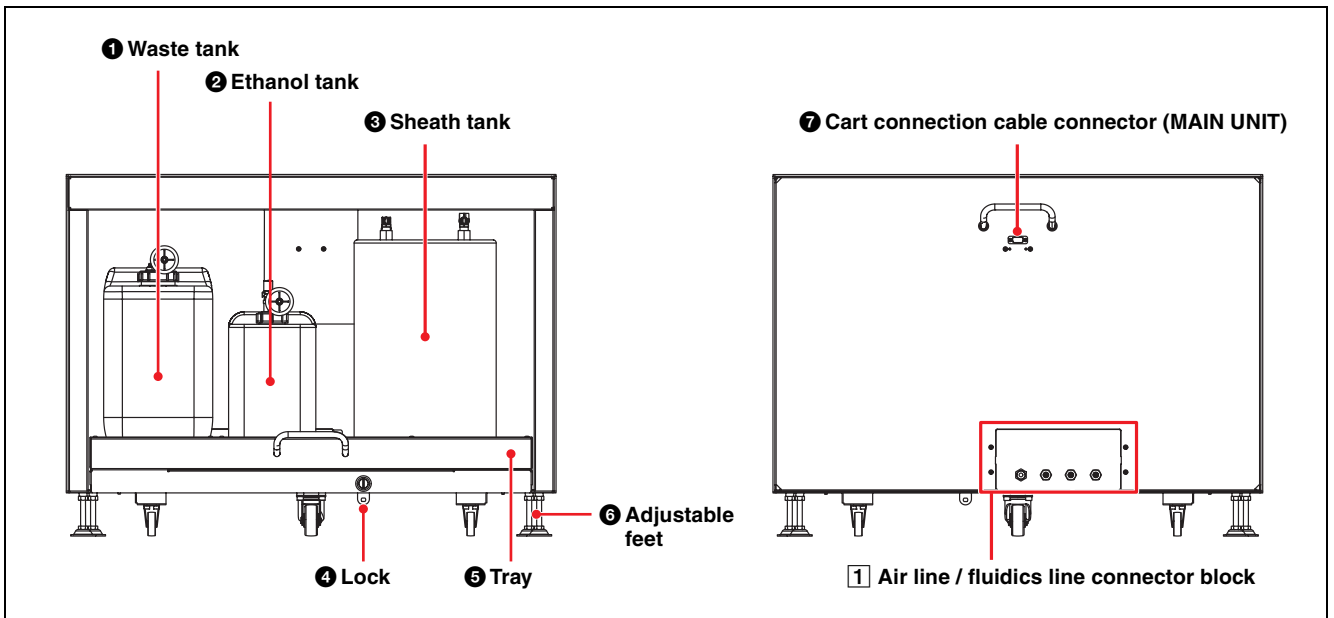

1 Waste tank
Collects waste fluid, comprising sample fluid, sheath fluid, and cleaning fluid, from various locations in the main unit. It also collects ethanol and DI water during cleaning. A sensor detects the waste fluid level for display on the LCD monitor and host computer. An

alarm is generated and SH800 operation automatically stops when the waste fluid rises above a set level. The tank has a capacity of 10 liters (2.6 gallons US).

DANGER

Waste fluid may contain biological, chemical, or other agents. Always wear gloves and other protective clothing, mask, and goggles, as required, when emptying or replacing the waste tank. The handling of waste fluid should be performed in accordance with biological hazard handling safety procedures.

2 Ethanol tank

Supplies the cleaning fluid (ethanol) used to clean the fluidics system during maintenance. A sensor detects the ethanol level for display on the LCD monitor and host computer. An alarm is generated when the ethanol falls below a set level. The tank has a capacity of 5 liters (1.3 gallons US).

3 Sheath tank

Supplies the sheath fluid used as the flow medium for transporting the sample fluid in a laminar flow through the sorting chip and into the collection tubes. The tank is pressurized using regulated compressed air supplied from the main unit.

A cart sensor detects the sheath fluid level for display on the LCD monitor and host computer. An alarm is generated when the sheath fluid falls below a set level. The tank has a capacity of 10 liters (2.6 gallons US).

Notes

- Never use the supplied sheath tank for anything other than the intended purpose.
- When opening the sheath tank cap, check that “Standby” is displayed on the LCD monitor or that the power to the main unit is turned off. If “Standby” is not displayed, place the unit in Standby mode using the following procedure.

① Click [Settings] on the [Cytometer] tab of the ribbon.

The [Cytometer Settings] dialog appears.

② Click [Advanced Settings].

The [Advanced Settings] dialog appears.

③ Click [Standby] on the [Pressure Options] tab.

④ Check that “Standby” is displayed on the LCD monitor.

When “Standby” is displayed, disconnect the fluidics cart air line (clear tubing) from the top of the sheath tank. For details, see “*Refilling the Sheath Tank*” (page 164).

- Do not force the lid down or place, place objects on the lid, or move the tank while the tank is pressurized.

Note

The sheath tank and sheath fluid line must not be disturbed to ensure stable, measurement conditions. Position the sheath fluid line where it will not be affected by drafts from air conditioning or vibration of the air compressor.

5 Lock

Locks the tray.

The lock prevents the tray from being withdrawn. To lock the tray, use a coin or similar object and turn to the LOCK position. To unlock, turn to the UNLOCK position.

6 Tray

Holds the sheath tank, DI water tank, and waste tank, and can be withdrawn when performing maintenance.

The tray is fitted with a lock to prevent it from being withdrawn accidentally.

See “*Maintenance*” (page 161).

7 Adjustable feet

Feet are adjustable to ensure the fluidics cart does not move during data acquisition.

8 Cart connection cable connector (MAIN UNIT)


Connects to the cart connection cable connector on the main unit. It is used to connect the sensors for the three fluid tanks in the fluidics cart.

CAUTION

Use only the supplied cable to connect the main unit to the fluidics cart. Using third-party cables may cause incorrect operation or device failure.

Note

The standard cable is about 2 m (6 ft) in length.

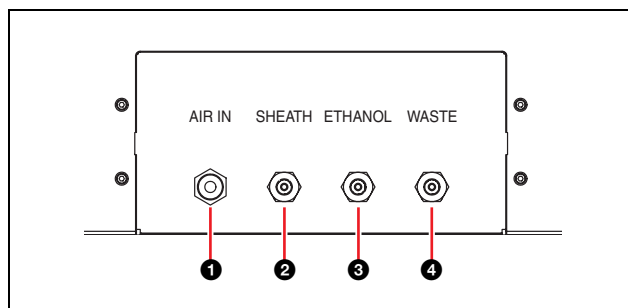1 Air line / fluidics line connector block**1 Fluidics cart air line connector (AIR IN)**

Connects to the fluidics cart air line connector (AIR OUT) on the SH800 (clear tubing). The air line supplies regulated compressed air for the sheath tank.

2 Fluidics cart sheath line (blue) connector

Connects to the SHEATH connector on the main unit using the supplied sheath line (blue tubing).

3 Fluidics cart ethanol line (green) connector

Connects to the ETHANOL connector on the main unit using the supplied ethanol line (green tubing).

4 Fluidics cart waste fluid line (red) connector


Connects to the WASTE connector on the main unit using the supplied waste line (red tubing).

Notes

- The standard fluidics lines and cable are about 2 m (6 ft) in length.
- Use only the supplied cable to connect the main unit to the fluidics cart. Using third-party cables may cause incorrect operation or device failure.

Sorting Chip

The SH800 uses a disposable sorting chip.

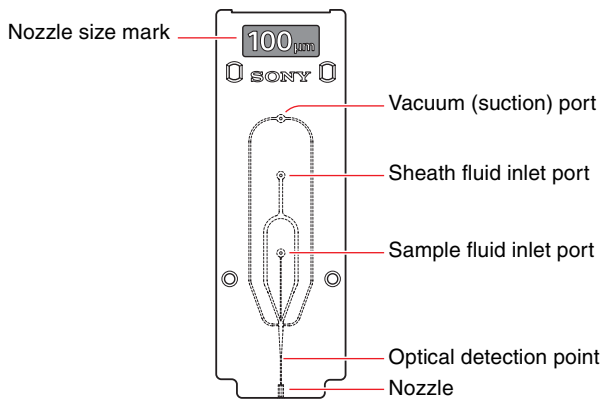

Sample fluid and sheath fluid enter the chip through inlet ports on the front surface of the chip. The sample fluid and sheath fluid are forced into the chip by air pressure from the sample loader and the sheath tank, respectively. The sample fluid and accompanying sheath fluid flows through the chip where the sample fluid is focused into a stream of single-file cells. The cells in the sample fluid then pass the optical detection point where they are interrogated by the lasers, before being ejected from the nozzle. A transducer vibrates the chip at an ultrasonic frequency to break the fluid stream into droplets for sorting.

A vacuum port provides negative pressure (suction) to remove bubbles and clogged samples from the fluidics channels in the sorting chip.

The nozzle size mark indicates the size of the nozzle located at the bottom of the chip.

The sorting methods that can be used vary depending on the nozzle diameter of the sorting chip and the size of cells. For details, see “Sorting” (page 58).

Note


Do not touch the sorting chip near the optical detection point. Fingerprints and smudges near the optical detection point can adversely affect data acquisition results.

Main Window

Notes

- The menus and buttons displayed vary depending on the user account privileges.
- The screenshots displayed in this document may vary from the actual software.

Operating window

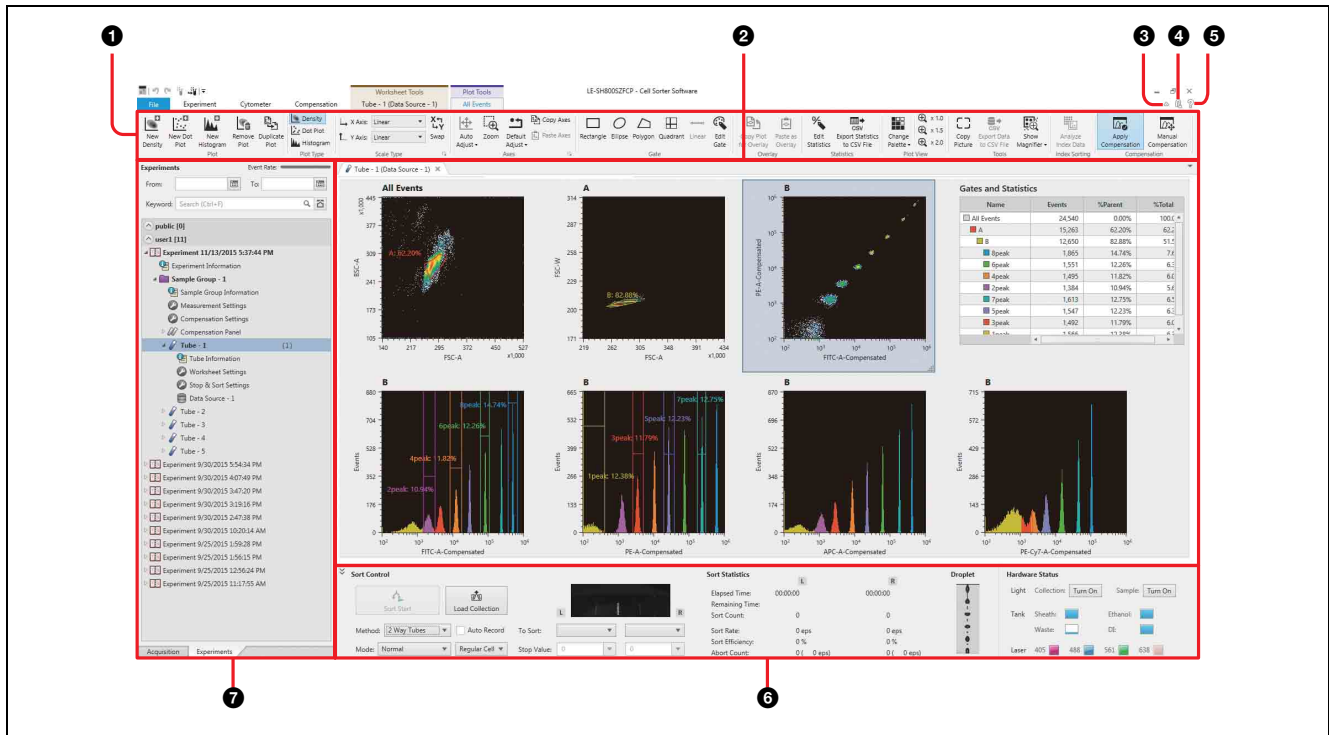

For details about the items displayed, see “Chapter 7 Window Description” (page 119).

1 Ribbon

Displays the tool buttons that can be used in the displayed window. The content of the ribbon varies depending on the selected tab.

2 Worksheet

Work area for displaying and editing the results from data acquisition and analysis.

3 Ribbon show/hide control (minimize/restore)

Clicking / alternately shows the ribbon or hides the ribbon.

4 Logout

Logs out the currently logged in user and displays the login window (page 45).

5 Help

Displays the Help window for accessing the Operator’s Guide.

6 Sorting control pane

Configures and controls droplet sorting. It also displays the current status of sorting, fluid tank levels, lights, and laser sources. The displayed items vary depending on the model and the sorting method. The sorting control pane is a collapsible window element. The panel can be shown or hidden as required.

Tip

When the unit is calibrated in analyzer mode, the sorting control pane is automatically hidden.

7 Experiment/Acquisition control area


Used to configure and control data acquisition and work with experiments.

[Acquisition] tab: Displays the controls used for acquiring and recording data.

[Experiments] tab: Displays experiments and the components in an experiment, used to acquire and analyze data, in a hierarchical list view.

[File] window

Example: [Information] tab

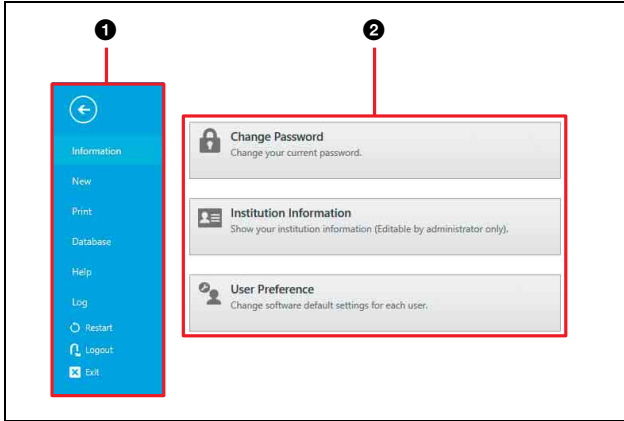

1 Menu items

Displays the following menu items.

[Information]: Displays commands used to configure user passwords and institution information, and to add/manage user accounts.

[New]: Creates a new experiment based on saved templates and/or recent experiments.

[Print]: Prints analysis results and other reports.

[Database]: Displays commands used to perform experiment database operations and to check hard disk free capacity.

[Help]: Displays commands used to display the Operator's Guide and to check software version information.

[Log]: Displays commands used to display the chip alignment log, sorting calibration log, and user log.

[Restart]: Restarts Cell Sorter Software.
When the software restarts, it loads the same conditions that existed at the time [Restart] was pressed, but without assigning any experiment as the active experiment.

Tip

The restart function should only be used when Cell Sorter Software is running very slowly or is otherwise unstable. It should not be used during normal operation.

[Logout]: Logs out the currently logged in user and displays the login window.

[Exit]: Exits Cell Sorter Software.

2 Operation area

Displays function buttons and information related to the menu item selected on the left side.

Cell Sorter Software Operating Environment

The host computer must satisfy the following requirements in order to run Cell Sorter Software.

Operating system (OS): Microsoft Windows 8/8.1/10 Operating System, Professional, 64 bit

CPU: Intel Core i7-3820QM 2.70 GHz or later

Cache: Level 3 cache memory 8 MB or greater

Memory: PC3-12800, 1600 MHz, 16 GB (8 GB × 2) or greater

Network interface: 1000BASE-T (1)

Main storage: 1 TB HDD (SATA 1 TB × 1) or greater

Monitor resolution: 1920 × 1080 (16:9 Full HD)

Graphics resolution: 1920 × 1080 (Full HD)

Camera: 1.3 megapixels or higher effective pixels

OS language setting: Symbol for the decimal point is “.” (period).

Note

If the decimal point is set to a character other than “.” (period), the following error message is displayed when starting Cell Sorter Software.

Invalid Language setting was detected.

Please change “Decimal symbol” setting to “.(period)” according to the Operator's Guide.

Press [OK] to close this application.

If this occurs, use the following procedure to change the decimal point to “.” (period), then restart Cell Sorter Software.

- 1 Open [Control Panel] > [Clock, Language, and Region] > [Region and Language].
- 2 Open [Additional settings...] on the [Formats] tab.
- 3 Set [Decimal symbol] to “.” (period) on the [Numbers] tab.

Tips

- For details about recommended antivirus software, contact your Sony distributor.
- It is the responsibility of the user to manage the Windows ID/password.

LCD Monitor

The LCD monitor on the front of the SH800 displays status information during operation.

The LCD monitor is turned on/off using the Display Mode button on the front panel.




Tip

If conflicting information is displayed in Cell Sorter Software display and on the LCD monitor, it does not necessarily indicate a failure or error.




Checking Device Connections

Check that the main unit, fluidics cart, host computer, and compressed air supply are connected correctly.

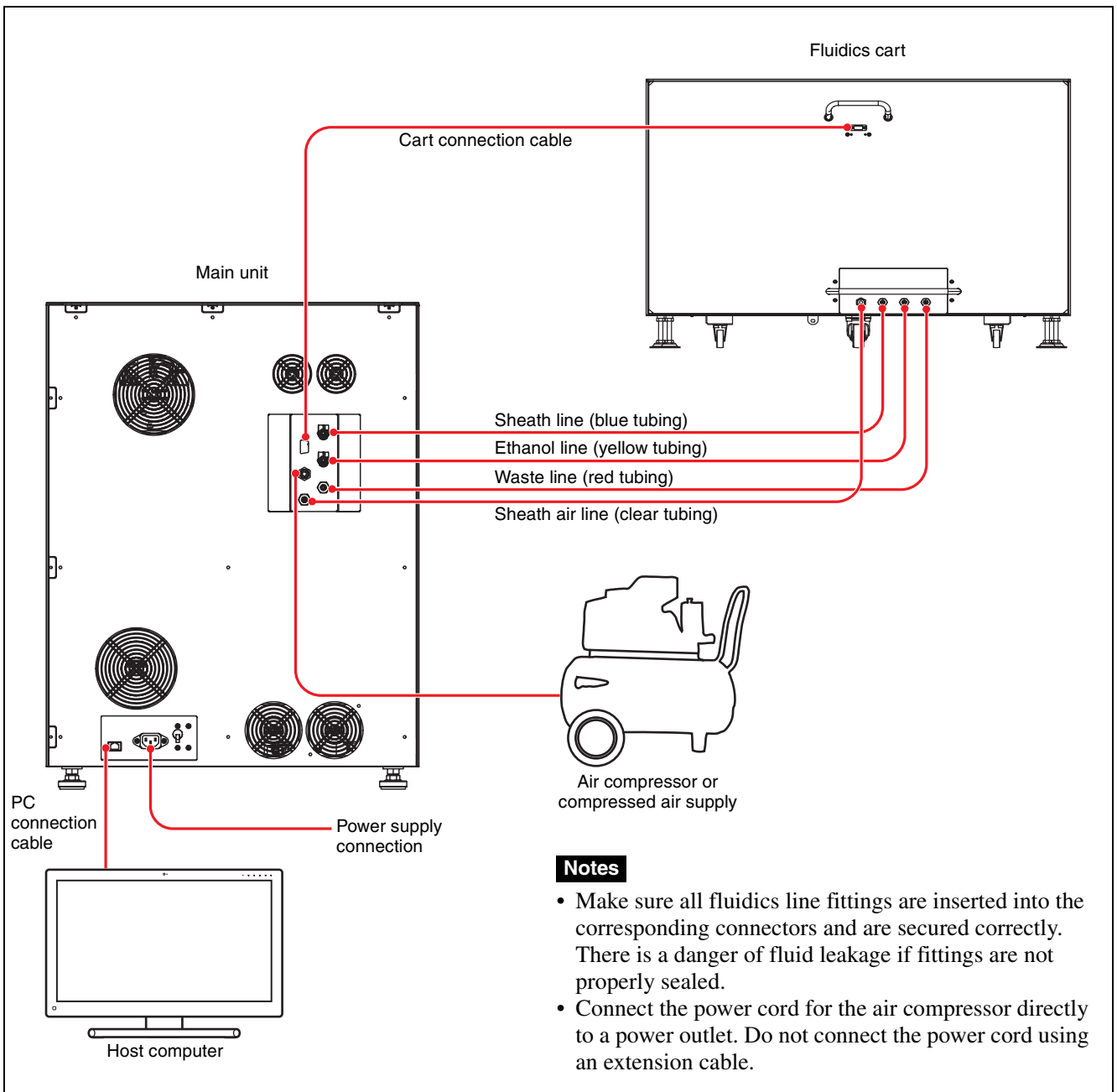

Filling the DI Water Tank

Fill the DI water tank in the main unit with DI water. DI water is used to flush sheath fluid, ethanol, and other cleaning agents from the fluidics system. Always check the level in the DI water tank and refill as required before starting/shutting down the SH800 and before running ethanol cleaning.

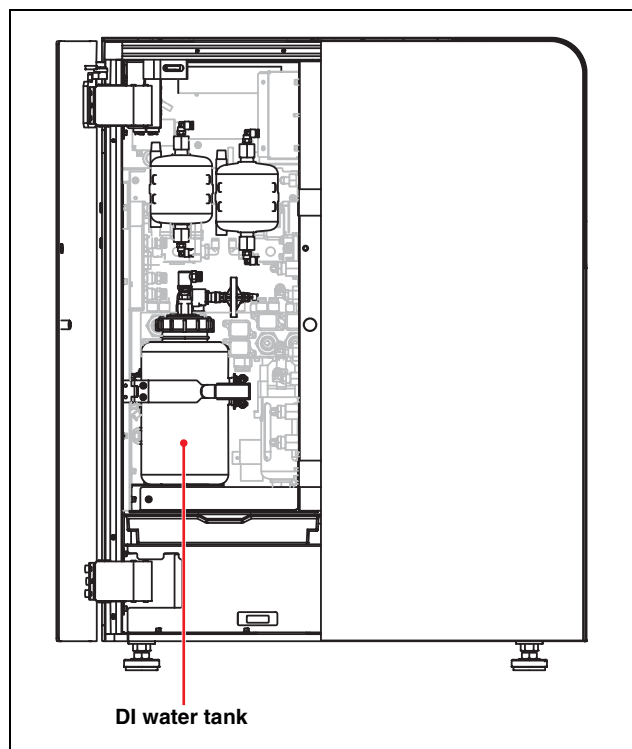

For details about refilling the DI water tank on the fluidics cart, see “Refilling the DI Water Tank” (page 167).

Tip


An alarm is raised if the level in the DI water tank drops too low.

Filling the Sheath Fluid Tank

Fill the sheath fluid tank on the fluidics cart with sheath fluid.

Sheath fluid is used as a transport media to deliver sample fluid to the sorting chip, and to clean the sample line.

Before starting and shutting down the SH800, make sure that there is sufficient sheath fluid in the sheath tank, and refill as required.

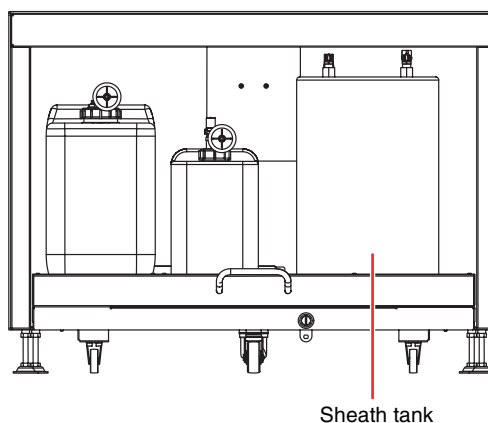

For details about refilling the sheath tank on the fluidics cart, see “Refilling the Sheath Tank” (page 164).

Notes

- Always place the sheath tank in the center of the tray area. The fluid level detector may not operate correctly if the tank is placed off-center.
- The sheath tank is not autoclavable.

Tip


An alarm is raised if the level in the sheath tank drops too low.

Filling the Ethanol Tank

Fill the ethanol tank on the fluidics cart with ethanol for ethanol cleaning.

Ethanol is used during ethanol cleaning to disinfect the fluidics system.

Always check the level in the sheath tank and refill as required before running ethanol cleaning.

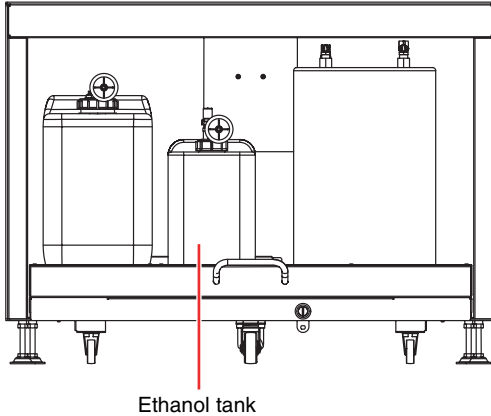

For details about refilling the ethanol tank on the fluidics cart, see “Cleaning the Internal Sheath Line and DI Water Line using Ethanol” (page 209).

Note

Always place the ethanol tank in the center of the tray area. The fluid level detector may not operate correctly if the tank is placed off-center.

Tip

An alarm is raised if the level in the ethanol tank drops too low.

Preparing a Sample

Loading a Sample Tube in a Sample Tube Holder

Insert the sample fluid for analysis or sorting in a sample tube, and place the tube in the corresponding sample tube holder.

The SH800 supports the following sample tubes.


Each sample tube is used with a corresponding holder that mounts in the sample loader.

Sample tube	Holder
0.5 ml micro tubes	Sample tube holder (0.5 ml)
1.5 ml micro tubes	Sample tube holder (1.5 ml)
5 ml round tubes	Sample tube holder (5 ml)
15 ml conical tubes	Sample tube holder (15 ml)


Load the sample tube into the corresponding size sample tube holder.

Note

Always remove the sample tube caps when performing measurements.

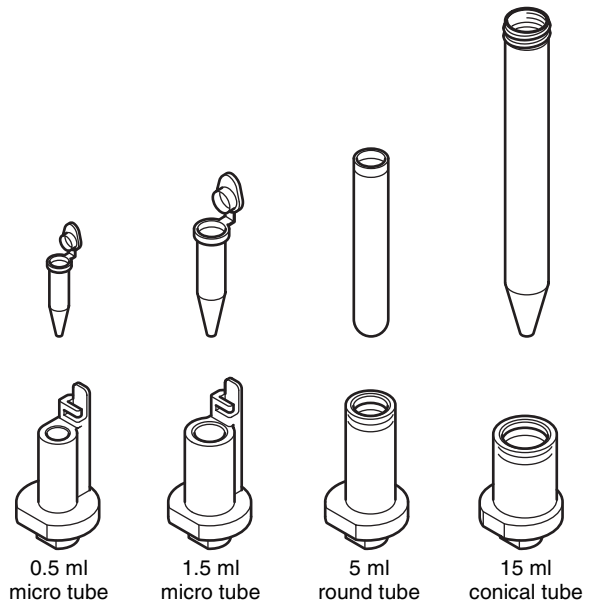

The tube holders for micro tubes have a recess at the top to hold the cap clear of the tube opening.

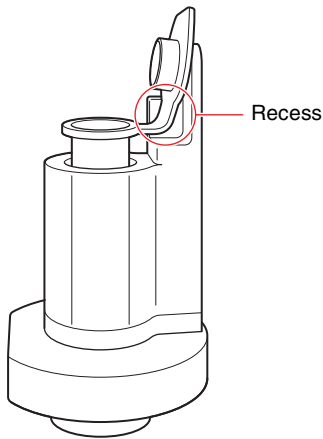

Notes

- Always wear laboratory-use gloves, mask, protective goggles, and other protective clothing when handling sample fluid to provide protection against biohazards. Also, take care when loading the sample tube in the tube holder to avoid sample fluid spills and breakage of the sample tube.
- Before using a tube, check that there is no dust or other foreign matter in the tube.
- Before place the sample tube in the sample tube holder, agitate sample to avoid trouble caused by aggregates.
- Before place the sample tube in the sample tube holder, check the volume of sample fluid to avoid introducing bubbles in fluidics line. The bubbles cause droplet formation to become unstable. Unstable droplet may cause fluid adhere to the surface of optical components, and stable performance cannot be maintained.

Preparing Automatic Setup Beads

Prepare automatic setup beads in advance if planning to run automatic calibration.

- 1 Filter the automatic setup beads using a 30 μm filter.
- 2 Insert 0.5 ml or more of undiluted, automatic setup beads in a sample tube, and place the tube in the corresponding sample tube holder.

Always use automatic setup beads in undiluted form.


Sample Temperature Control


The SH800 is fitted with thermal control units in the sample injection chamber and collection area to maintain sample fluid temperature control while running an experiment. The temperature control device can be set to 5 $^{\circ}\text{C}$ (41 $^{\circ}\text{F}$) or 37 $^{\circ}\text{C}$ (98.6 $^{\circ}\text{F}$) (collection stage: 5 $^{\circ}\text{C}$ (41 $^{\circ}\text{F}$) only). You configure settings in Cell Sorter Software.

Notes

- The temperature control function in Cell Sorter Software cannot be used to cool 8-well strip holders/slide glass holders, 96-well plate holders, 384-well plate holders, or 384-well PCR adapters. However, the holders contain an embedded cooling agent. You can cool a holder in a laboratory refrigerator (-20°C (-4°F)) beforehand, and then place the holder on the collection stage.
 - For 8-well strip holders/slide glass holders and 96-well plate holders:
 - Cooling time: 5 hours or longer
 - Temperature retention time: About 1 hour (varies with ambient temperature)
 - For 384-well plate holders and 384-well PCR adapters:
 - Cooling time: 8 hours or longer
 - Temperature retention time: About 1 hour (varies with ambient temperature)
- Note that holders/adapters may freeze if immediately placed on the collection stage when removed from the refrigerator.
- The collection tube holder with cooling agent (5 ml) contains an embedded cooling agent. Cool the holder in a laboratory refrigerator (-16°C (3 $^{\circ}\text{F}$) grade or cooler) beforehand for 8 hours or more, and then place the holder on the collection stage. When used with the temperature control function in Cell Sorter software, the holder provides about 2 hours cooling (will vary depending on ambient temperature).

For details about starting the SH800 instrument and Cell Sorter Software, see “System Startup” (page 44) and “Logging In” (page 45).

- 1 Click the  dialog launcher in the [Temperature Control] group on the [Cytometer] tab of the ribbon. The [Temperature Control Settings] dialog appears.
- 2 Select the temperature setting for the sample fluid, then click [Close].

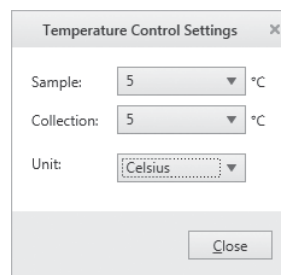

Sample: 5 $^{\circ}\text{C}$ (41 $^{\circ}\text{F}$) or 37 $^{\circ}\text{C}$ (98.6 $^{\circ}\text{F}$)
Collection: 5 $^{\circ}\text{C}$ (41 $^{\circ}\text{F}$) only

Note

When sorting with the temperature set to 5 $^{\circ}\text{C}$ (41 $^{\circ}\text{F}$), place the collection tube holder in a refrigerator to cool the holder before loading it on the collection stage.

- 3** To turn a temperature control device on/off, click [Sample] or [Collection] in the [Temperature Control] group on the [Cytometer] tab of the ribbon.

The buttons toggle between on and off.

Note

If the temperature control function is turned off, expected acquisition results may not be obtained, depending on the sample being analyzed.

Precautions when temperature control is set to 5 °C (41 °F)

- Cool the collection holder to approximately 5 °C (41 °F) in a refrigerator before loading it onto the collection stage (for the collection tube holder with embedded cooling agent (5 ml), cool to -16 °C (3 °F) grade or cooler). Cooling the collection holder beforehand significantly reduces the time required to cool the collection tubes.
- While the temperature is set to 5 °C (41 °F), do not remove the collection tube holder from the collection stage. Insert and remove the collection tubes directly from the collection tube holder.
- Loading a collection tube holder that is warm or that has become warm may result in a sudden change in temperature, and may prevent correct operation of the unit. When unloading and loading the collection tube holder, temporarily turn off the temperature control function, wait about 30 seconds, and then set the temperature to 5 °C (41 °F) again.
- The metallic parts of the 96-well plate holder become cold when the holder is cooled. Always load and unload by holding the plastic sides.

Preparing Collection Tubes

Prepare the collection tube device for collecting droplets containing target cells for sorting.


The SH800 supports the following collection tube devices. Each type of collection tube is used with a corresponding holder that mounts on the collection stage.

Collection tubes/ multi-well plates	Holder
15 ml conical tubes	Collection tube holder (15 ml)
5 ml round tubes	Collection tube holder (5 ml) Collection tube holder with cooling agent (5 ml)
8-well strips	8-well strips holder ¹⁾
Slide glass	Slide glass holder ¹⁾
6-well plates	96-well plate holder
12-well plates	
24-well plates	
48-well plates	
96-well plates	
8-well PCR strips	96-well PCR plate holder
96-well PCR plates	
384-well plates	384-well plate holder ¹⁾
384-well PCR plates	384-well plate holder ¹⁾ + 384-well PCR adapter ¹⁾


1) Option sale of this product has been discontinued (as of April 2019).

Common to all models


- Collection tube holder (15 ml)

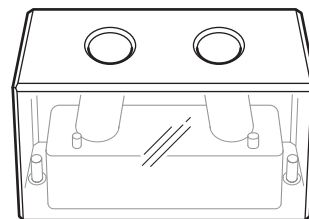

- Collection tube holder (5 ml)

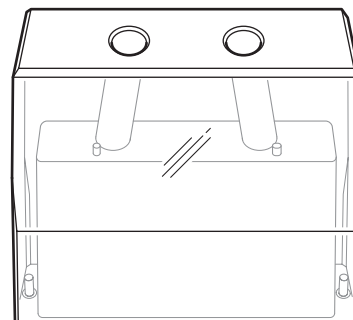

- Collection tube holder with cooling agent (5 ml)

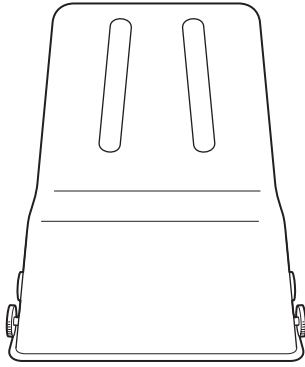

Use the collection tube holder with embedded cooling agent (5 ml) for more stable sorting if using a 70 μm sorting chip with the plain collection tube holder (5 ml).

When running 2-way tube sorting on the SH800SP and SH800ZP, the plate holder adapter must first be removed. See “Removing and Attaching the Plate Holder Adapter (SH800ZP and SH800SP)” (page 39).

SH800, SH800Z, and SH800S

- 8-well strips holder (with 8-well strips)




For details about loading 8-well strips, see “Loading 8-well Strips (SH800S, SH800Z, and SH800)” on page 37.

- Slide glass holder




SH800SP and SH800ZP


- 96-well plate holder (with 96-well plate)

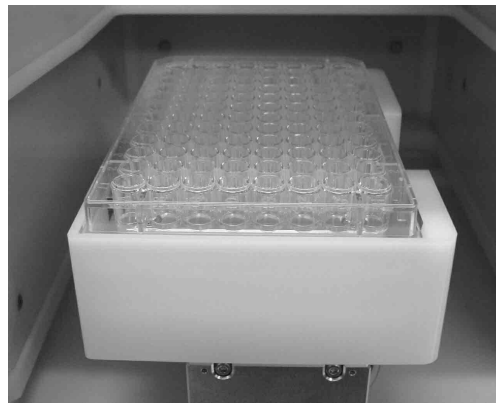

- 96-well PCR plate holder (with 96-well PCR plate)

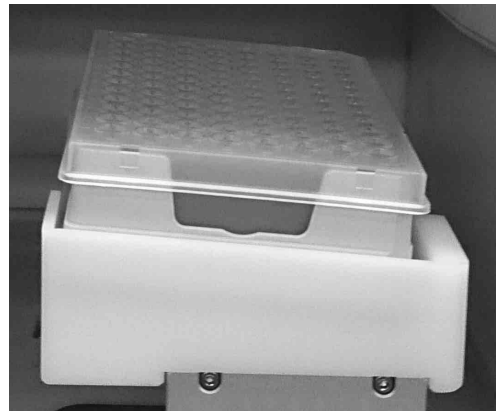

- 384-well plate holder (with 384-well plate)

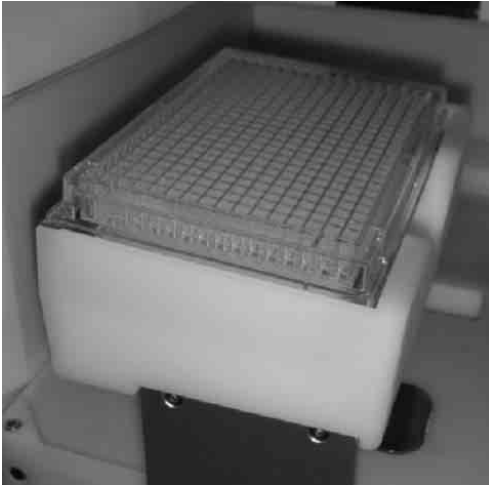

For details about loading tubes, see “Loading a Multi-well Plate (SH800SP and SH800ZP)” (page 38).

- 384-well plate holder (with 384-well PCR adapter and 384-well PCR plate)



For details about loading tubes, see “Loading a 384-well PCR Plate (SH800SP and SH800ZP)” (page 39).

When sorting into a 384-well plate or 384-well PCR plate, always adjust the sort position before starting. For details, see “Adjusting the Sort Position (SH800SP and SH800ZP)” (page 113).

DANGER

Fluids may contain biological, chemical, or other agents. When handling collection tubes, always wear laboratory-use gloves, mask, protective goggles, and other protective clothing.


Note

Before using a tube, check that there is no dust or other foreign matter in the tube.


Loading 8-well Strips (SH800S, SH800Z, and SH800)

Observe the following precautions when loading 8-well strips.

- 1 Insert the leading edge of the 8-well strip into the holder at an angle, and push the plate in all the way.

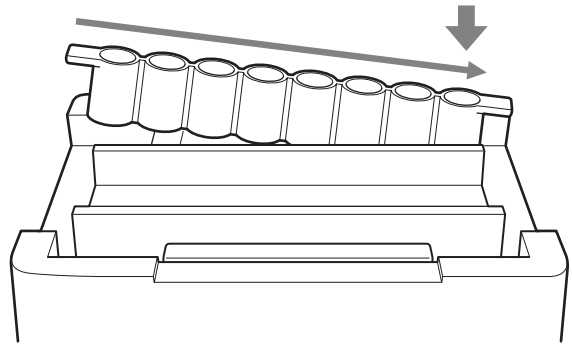


- 2 Orient the holder with the “FRONT” arrow towards you and place the holder on the collection stage.

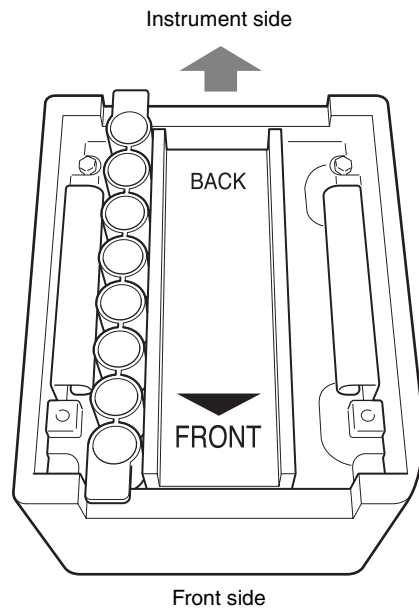




Attaching the Splash Guard (for use with multi-well plates)

(SH800SP and SH800ZP)


Attach the supplied splash guard when sorting into multi-well plates to prevent splashing of sample fluid.

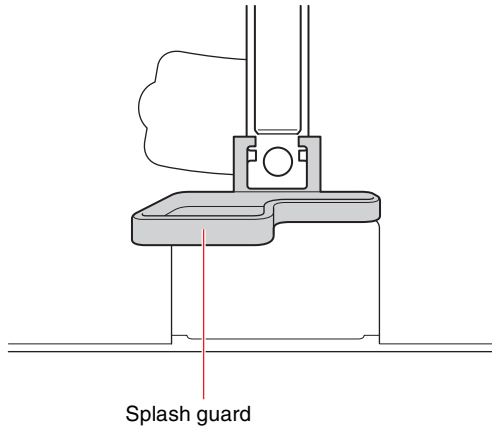


- 1 Hold the splash guard in the orientation shown in the diagram, and insert the protrusion on the splash guard into the notch on the waste catcher.

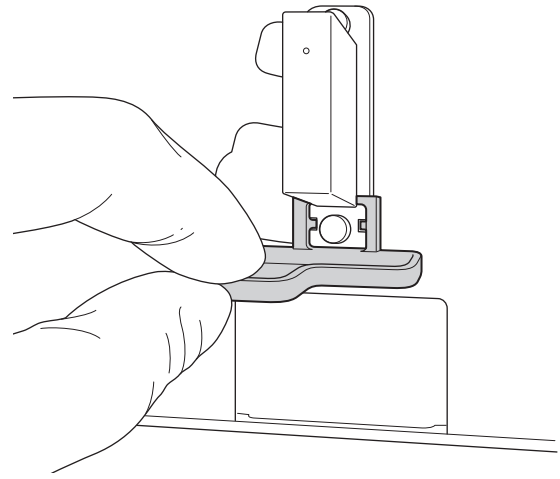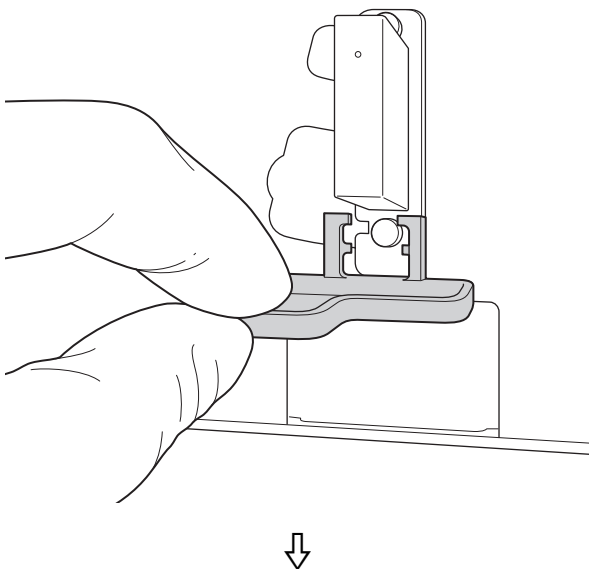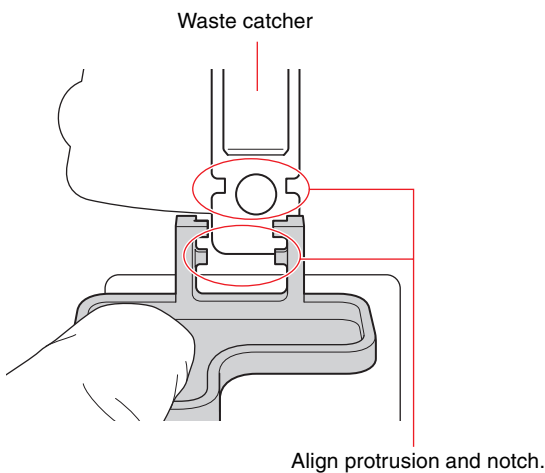

Insert the protrusions on both sides.

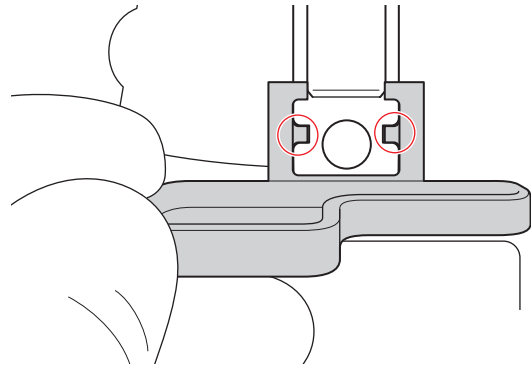

- 2 Push the splash guard in until it clicks into position.

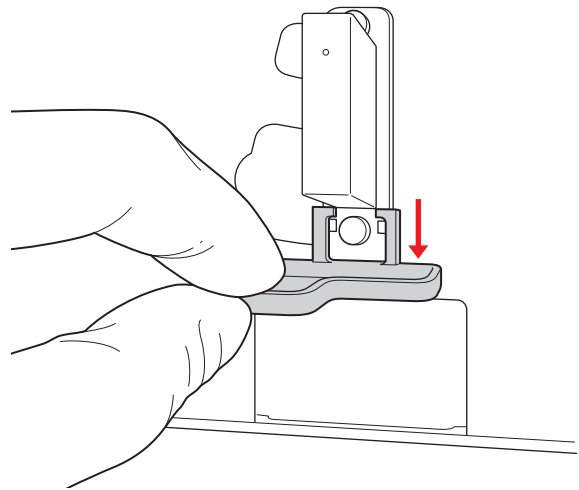

Tip


Remove the splash guard when unloading multi-well plates after sorting is finished.

Loading a Multi-well Plate

(SH800SP and SH800ZP)

Place the multi-well plate into the plate holder with the chamfered corners toward the rear of the instrument.


Rear of instrument

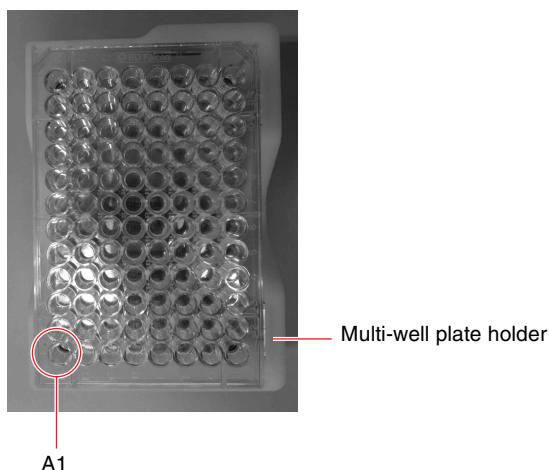**Notes**

- Always use a 384-well plate holder when using 384-well plates. Do not use a 96-well plate holder.
- Always use a 96-well PCR plate holder when using 96-well PCR plates. Do not use a 384-well plate holder.

Loading a 384-well PCR Plate (SH800SP and SH800ZP)

Place the 384-well PCR adapter in the 384-well plate holder, and place a 384-well PCR plate on the adapter. Place the 384-well PCR plate into the adapter with the A1 well at the front left of the instrument.

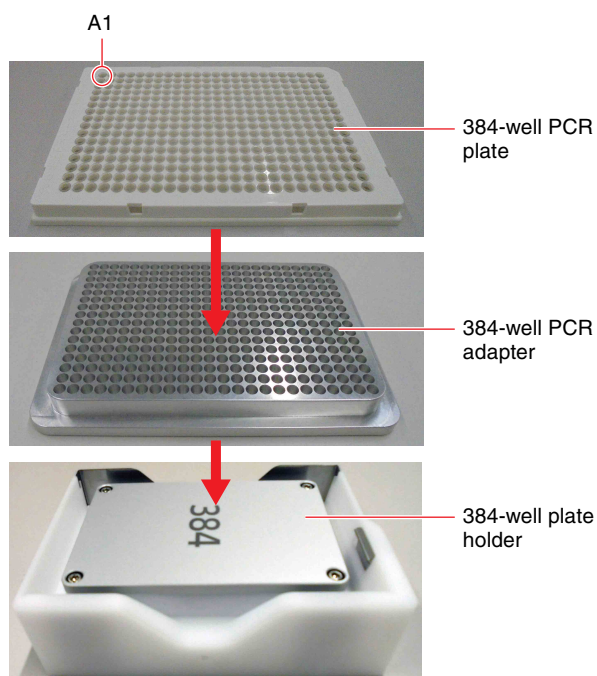**Note**


Always use a 384-well PCR plate holder and 384-well PCR adapter when using 384-well PCR plates. Do not use a 96-well PCR plate holder.

Removing and Attaching the Plate Holder Adapter (SH800ZP and SH800SP)

When running 2-way tube sorting, the plate holder adapter must first be removed.

Remove the screw securing the adapter using a coin or screwdriver, and remove the adapter.

For multi-well plate sorting, re-attach the adapter.

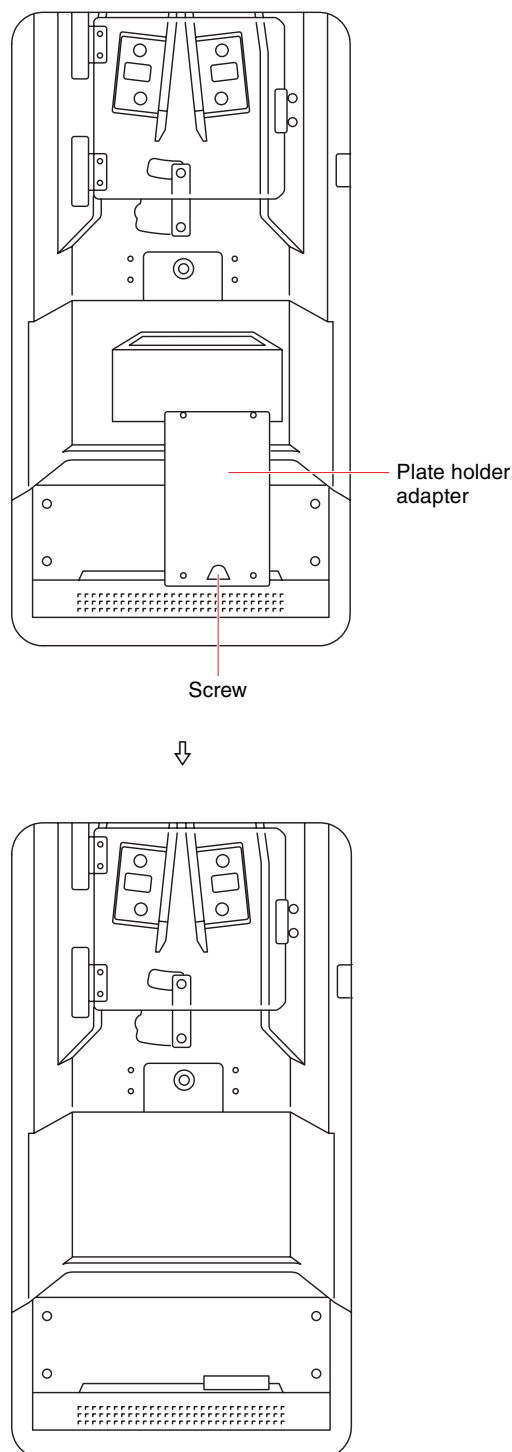

Configuring Users

This section describes how to manage user accounts in Cell Sorter Software.


For details about starting the SH800 instrument and Cell Sorter Software, see “System Startup” (page 44) and “Logging In” (page 45).

About the administrator account

The administrator account user name is set to “administrator” by default. The administrator account cannot be used to operate the SH800.

If the administrator account password has not been configured, the change password screen appears for the administrator account when launching Cell Sorter for the first time.

Specify each item and click [OK] to set the password.

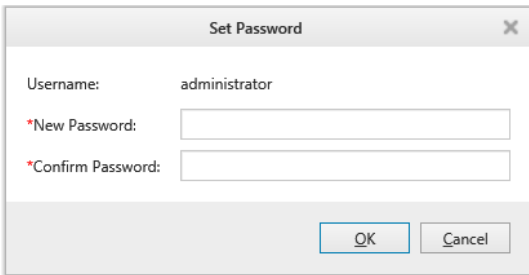


[New Password]: Enter a new password, comprising 8 to 20 alphanumeric characters (required item).

[Confirm Password]: Enter the same password again for confirmation (required item).

Adding Users


Only administrators can add new users.

- 1 Click the [File] tab, then click [Information] in the menu on the left.
- 2 Click [Account Settings].

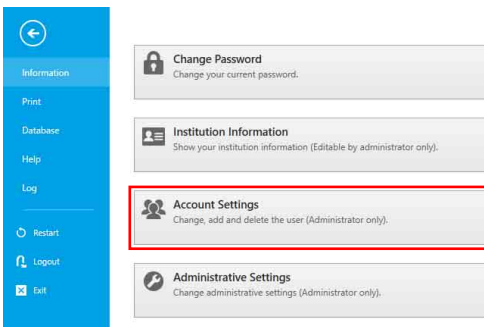


The [Account Settings] dialog appears.

- 3 Click [Add].

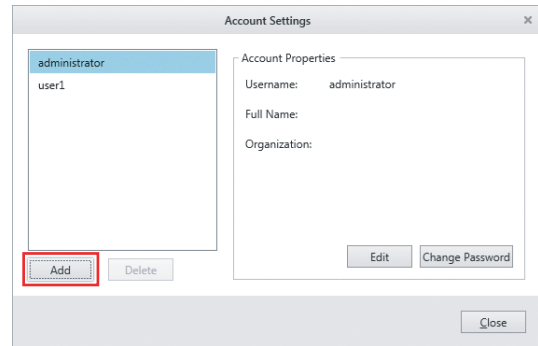

The [Add Account] dialog appears.

- 4 Set each field, then click [OK].

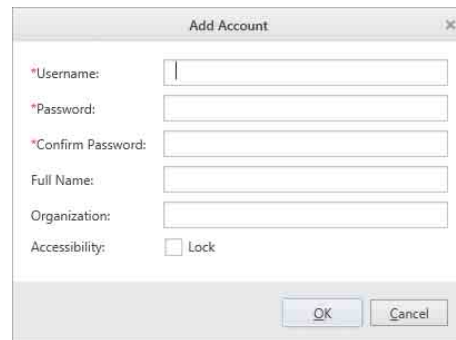

[Username]: Enter the user name, of up to 32 alphanumeric characters, used when logging in to Cell Sorter Software (required item).

[Password]: Enter the password, comprising 8 to 20 alphanumeric characters, used when logging in to Cell Sorter Software (required item).


[Confirm Password]: Enter the same password again for confirmation (required item).

[Full Name]: Enter the full name of the user.

[Organization]: Enter the name of the organization.

[Accessibility]: Place a check mark in the checkbox if you want to temporarily disable a general user account.

The user is registered in the user list in the [Account Settings] dialog.

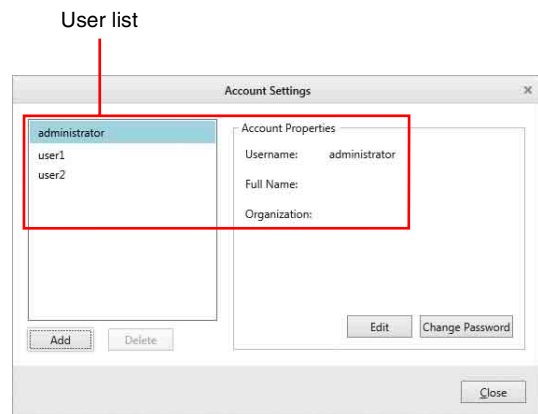

Tip

A warning dialog appears if the password is less than eight characters in length.

Editing User Settings

Only administrators can edit user settings.

- 1 Select the user you want to edit in the [Account Settings] dialog, then click [Edit].

Selecting a user in the list displays the settings for the user in the area on the right. The [Edit Account] dialog appears.

- 2 Edit the settings, then click [OK].

For details, see step 4 (page 40) in “Adding Users”.

The account settings are updated.

Changing Passwords

Individual users can change their password used to log in to Cell Sorter Software.

- 1 Click the [File] tab, then click [Information] in the menu on the left.
- 2 Click [Change Password].

The [Change Password] dialog appears.

- 3 Set each field, then click [OK].

[Old Password]: Enter the current password (required item).


[New Password]: Enter a new password, comprising 8 to 20 alphanumeric characters (required item).

[Confirm Password]: Enter the same password again for confirmation (required item).

The password is updated.

Setting a Cleaning Reminder

The instrument must be cleaned periodically using ethanol to control contamination and side stream instability resulting from poor system maintenance. You can set to display a message (cleaning reminder) on the login screen after a certain preset time has elapsed without performing ethanol cleaning.

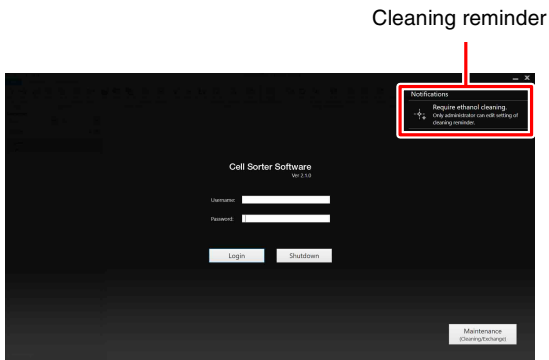


Tip

You can click the cleaning reminder to display the [Cleaning Reminder] screen and then start ethanol cleaning.

For details, see “When a Cleaning Reminder is Displayed” (page 180).


Use the following procedure to set the cleaning interval for sending cleaning reminder notifications. Only administrators can set the cleaning interval.

- 1 Click the [File] tab, then click [Information] in the menu on the left.
- 2 Click [Administrative Settings].

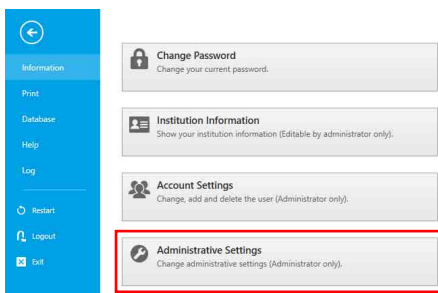

The [Administrative Settings] dialog appears.

- 3 Select [Cleaning Reminder], set each item, and then click [OK].

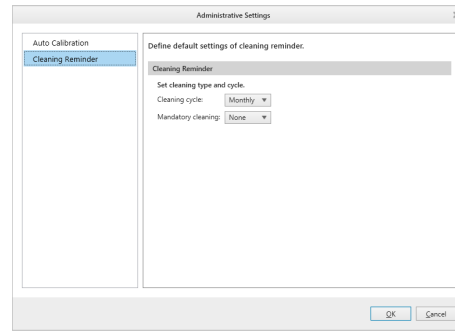

[Cleaning cycle]: Select the ethanol cleaning interval.

- None (no notifications)
- Daily
- Weekly
- Monthly (default value)

[Mandatory cleaning]: Select the mandatory ethanol cleaning interval.

If the interval specified here is exceeded, login is disabled until ethanol cleaning has been performed.


- None (no notifications) (default value)
- 1 week
- 2 weeks
- 1 month (30 days)

The ethanol cleaning intervals for sending cleaning reminder notifications are set.

You can check the cleaning history on the [Ethanol Cleaning Log] screen. For details, see “Ethanol cleaning log” (page 127).

This chapter describes the basic operating procedure for running a simple experiment, acquiring and analyzing data, and sorting selected populations in the sample.

Workflow

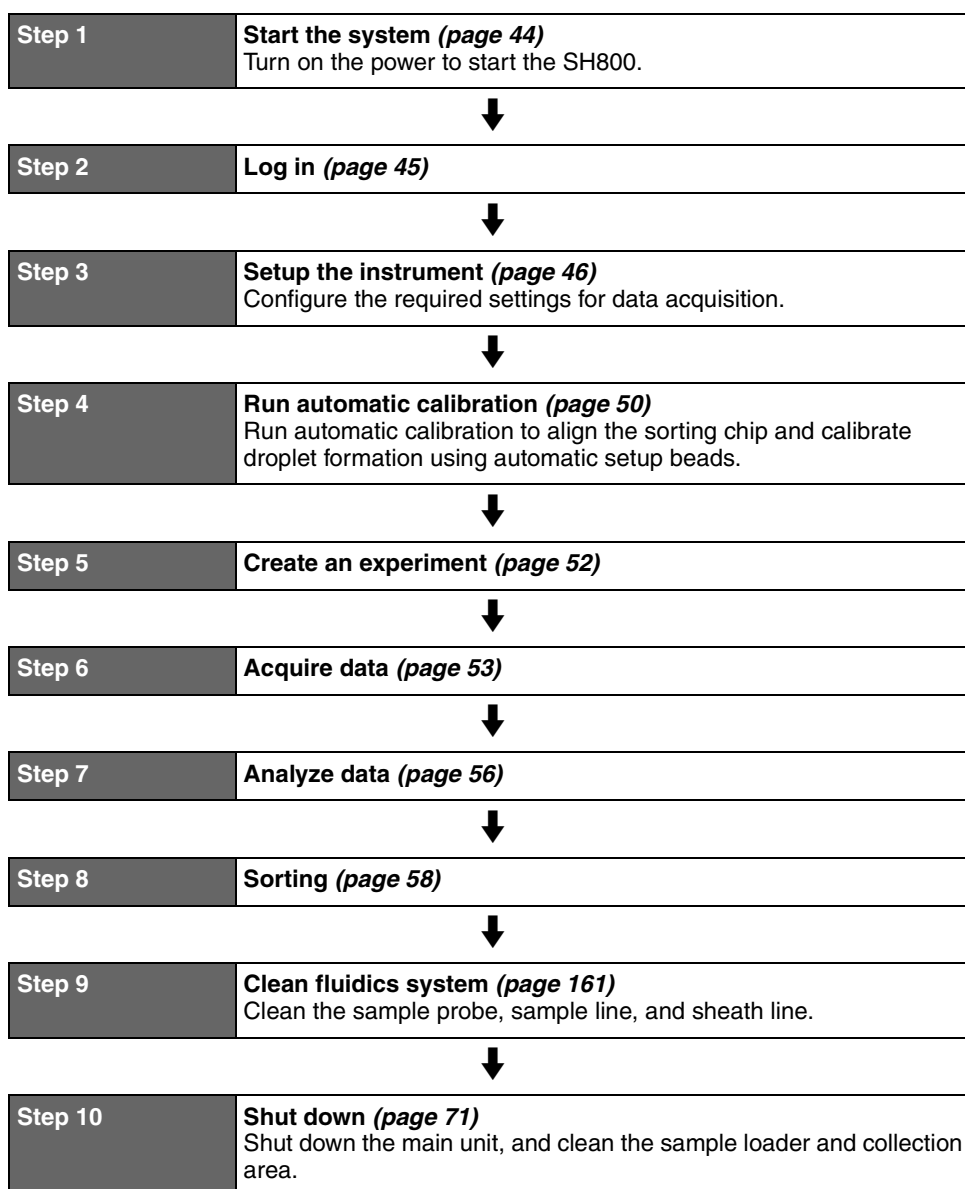

System Startup

Turn on the power supply and start the SH800 using the following procedure.


Note

Do not open the flip-up door at any time before turning on the compressed air supply. Doing so may damage the unit.

- 1 Check that the compressed air supply is connected and is supplying the rated pressure.
- 2 Check the level in the tanks on the fluidics cart, and refill or empty the tanks as required.

Also, check that the fluidics lines, sheath air line, and cart connection cable are connected correctly.

- 3 Turn on the MAIN POWER switch on the rear panel of the main unit.

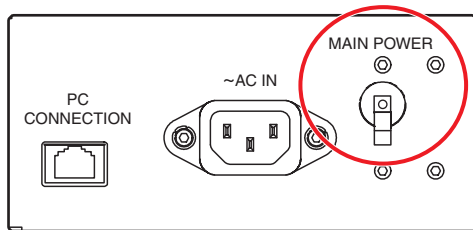


Note

During normal operation, the MAIN POWER switch can be left turned on all the time. Use the POWER/STANDBY button to turn power on/off on a daily basis.

- 4 Turn on the power using the POWER/STANDBY button on the front panel of the main unit.

The system goes through an initialization process during which self-diagnostics are run to test the proper working condition of all subsystems.

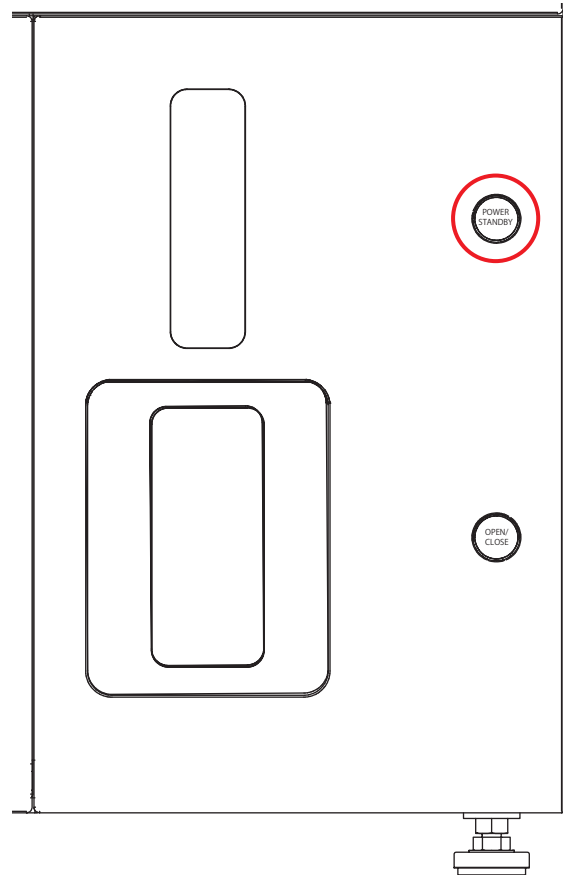

The SH800 status appears on the LCD monitor.




Logging In

Start Cell Sorter Software and log in to the SH800.

- 1 Check that “Standby” status is displayed on the LCD monitor.
- 2 Select Cell Sorter Software on the Windows Start screen to start the software.
The login window appears.
- 3 Enter your user name and password, then click [Login].



If login is successful, the [Initial Instrument Setup] wizard appears.

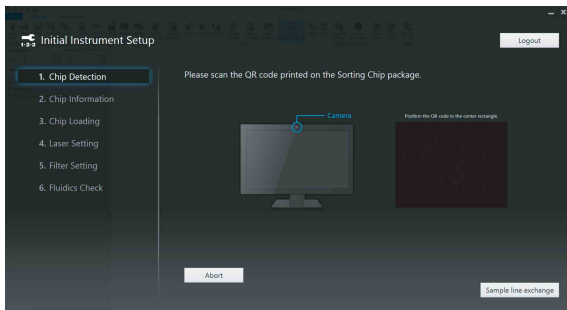


Tip

To change the sample line before initial instrument setup, click [Sample line exchange] and follow the instructions in the on-screen wizard.

For details, see “Changing the PEEK Sample Line (PEEK sample line compatible models)” (page 189) or “Changing the Sample Line and Probe (Non-PEEK sample line compatible models)” (page 192).

If periodic ethanol cleaning has not been performed

A cleaning reminder notification appears at the top right of the screen. Click the cleaning reminder to display the [Cleaning Reminder] screen and then start ethanol cleaning.

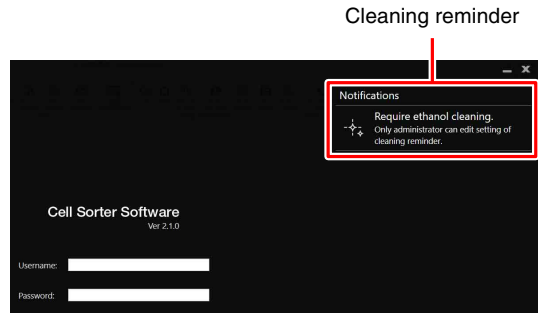

For details, see “When a Cleaning Reminder is Displayed” (page 180).

Initial Instrument Setup

Load a sorting chip, select the lasers to be used, and insert the optical filters in the required optical filter pattern by following the on-screen instructions in the [Initial Instrument Setup] wizard.

Using the Sorting Chip

- 1 Hold the QR code on the sorting chip packaging in front of the camera on the host computer.

QR code



The camera automatically scans the QR code.


Tip

You can click [Abort] to bypass instrument configuration and proceed directly to the main window. In this mode, Cell Sorter Software can be used only for analysis of existing recorded data in experiments.

When loading a sorting chip, the QR code on the chip packaging must be scanned using the built-in camera on the host computer to uniquely identify the chip and register the serial number and nozzle size of the chip with the SH800. This function maintains a history for the chip in order to ensure optimum performance.

When the QR code is successfully scanned, the product number and nozzle size are displayed, together with basic instrument settings. The serial number for the chip is registered internally.


- 2 Check that the displayed information is correct, then click [Next].

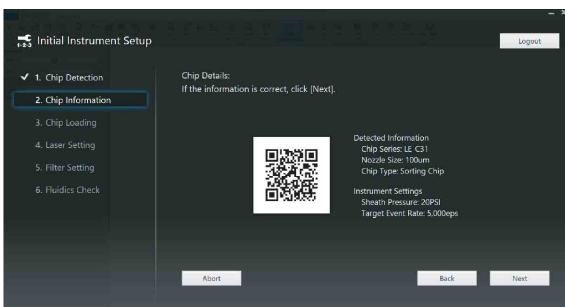


When [Next] is clicked, the currently loaded sorting chip is ejected from the chip loader.

Tip

If you wish to load a different chip, click [Back] to return to the previous screen, then scan the QR code for the different chip.

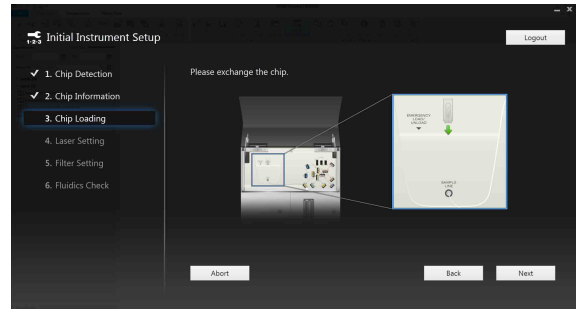

- 3 Open the flip-up door.
- 4 Remove the current sorting chip, if one is present, and insert the new chip into the chip slot at the top of the chip loader.

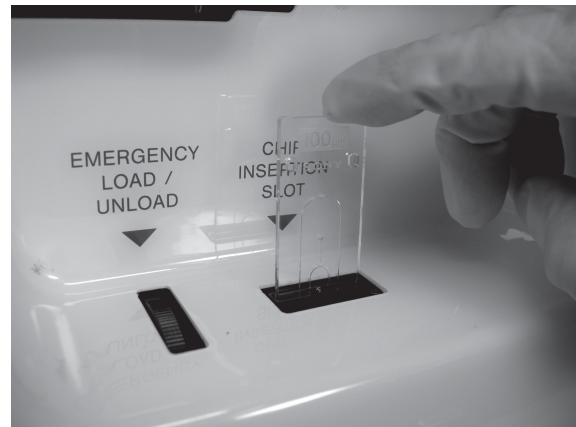

Insert the chip into the slot with the Sony logo and nozzle size mark at the top facing toward you. Take care not to insert the sorting chip facing backwards.

Place the previous chip in its original packaging. The chip and its corresponding QR code should be kept together at all times so that they are not mixed up with other chips.

Notes

- An error message appears if the sorting chip is inserted in the incorrect orientation.
- Do not insert any objects other than Sony sorting and cleaning chips designed for the SH800 into the chip insertion slot.
- The packaging for the loaded sorting chip should be stored in a safe location.
- Replace the sorting chip if it is discolored, dirty, or stained from previous use.
- Do not drop the sorting chip on a hard surface or bend the chip intentionally. If there is a problem


with the sorting chip, contact your Sony distributor.

- Always wear gloves and other protective clothing, mask, and goggles, when changing the sorting chip.
- If the sorting chip will not load correctly, run the instrument in maintenance mode as described in “*Running Maintenance Mode*” (page 177) to clean the instrument. If the problem persists, contact your Sony distributor for service.
- Replace the sorting if it is used for extended periods or after the usage period expires, and when it becomes clogged, discolored, or stained.

5 Click [Next].

The sorting chip is loaded and the [Laser Setting] screen appears.

6 Select the excitation lasers to be used for the experiment, then click [Next].

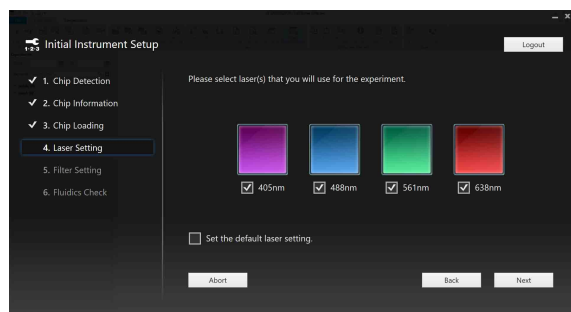

The excitation spectra of the fluorochrome labels you wish to detect largely determine the lasers needed for an experiment.

For a list of fluorochromes versus excitation laser wavelength for each model, see “*Fluorochrome Detection Matrix*” (page 232).

Tips

- The 488 nm excitation laser is used to measure forward scatter and back scatter light.
- It is recommended that lasers not required for an experiment be deselected.
- If the selected excitation lasers and [User Preference] settings do not match, [Set the default filter setting] is displayed. To make the currently selected excitation laser settings the default settings, place a check mark in [Set the default filter setting].

For details about [User Preference], see “*Specifying User Preferences*” (page 72) and “[Cytometer] tab” (page 121) of the [User Preference] dialog.


The [Filter Setting] screen appears.

7 Select the optical filter pattern setting, insert all the optical filters in the slots specified in “*Fluorochrome*

Detection Matrix” (page 232) and close the flip-up door, then click [Next].

The optical filter pattern settings that can be selected will vary depending on the lasers installed in the instrument.

Example: When 405 nm laser is installed

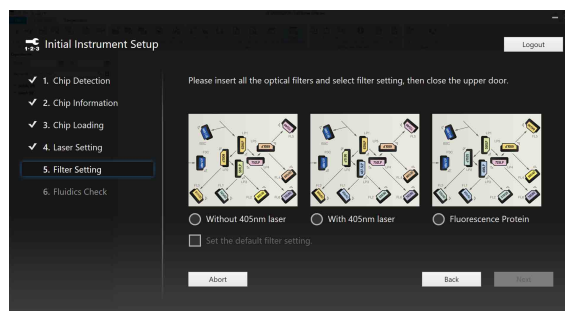

The image displayed shows the optical filter pattern to be installed.

Notes

- Make sure that the correct filter is inserted in each slot for the recommended filter pattern. Correct data cannot be obtained if a filter is inserted in the wrong slot.
- Exercise extreme care when handling the optical filters. The surfaces of each optical filter must be free of dust, smudges, fingerprints, and scratches for accurate data acquisition and cell sorting operation.

For details about handling optical filters, see “*Cleaning and Handling of Optical Filters*” (page 207).

Tip

If the selected optical filter pattern setting and [User Preference] settings do not match, [Set the default filter setting] is displayed. To make the currently selected optical filter pattern setting the default setting, place a check mark in [Set the default filter setting].

For details about [User Preference], see “*Specifying User Preferences*” (page 72) and “[Cytometer] tab” (page 121) of the [User Preference] dialog.

RESERVE slots

- On the following models, dummy filters are installed in the RESERVE slots.
SH800AC, SH800BC, SH800DC, SH800ZA, SH800ZB, SH800ZD, SH800ZH, SH800ZAP, SH800ZBP, SH800ZDP, SH800ZHP, SH800SA, SH800SB, SH800SD, SH800SH, SH800SAP, SH800SBP, SH800SDP, SH800SHP
- On the following models, filters not used in the current optical filter pattern should be stored in the RESERVE slots. Filter pattern 1 can also be used when the 405 nm laser is turned off to make efficient use of the six fluorescence detection


channels.

SH800CC, SH800EC, SH800FC, SH800ZC,
SH800ZE, SH800ZF, SH800ZG, SH800ZCP,
SH800ZEP, SH800ZFP, SH800ZGP, SH800SC,
SH800SE, SH800SF, SH800SG, SH800SCP,
SH800SEP, SH800SFP, SH800SGP

For details about filter pattern 1 and 2, see
“Optical Filter Patterns” (page 230).

The [Fluidics Check] screen appears.

8 Click [Next].

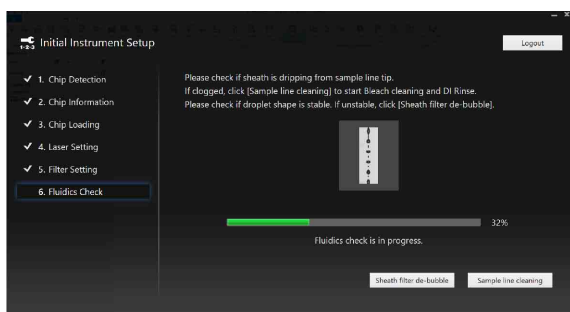


When fluidics check starts, sheath fluid droplets appear from the tip of the sample probe. If droplets do not appear, the sample line or the sample probe may be clogged. Click [Sample line cleaning] on the bottom right of the screen to clean the sample line.

During the fluidics check, the sorting chip de-bubble function is run periodically, making the droplets appear unsettled. If droplet formation is unstable, even when the sorting chip is de-bubbled, air may have accumulated in the sheath filter. Click [Sheath filter de-bubble] at the bottom of the screen to de-bubble the sheath filter.

When the fluidics check is completed, the [Auto Calibration] screen appears.

If fluid flow does not start

[Troubleshooting] is displayed at the bottom left of the screen. Click [Troubleshooting] to launch the troubleshooting wizard. Follow the on-screen instructions to resolve the issue.

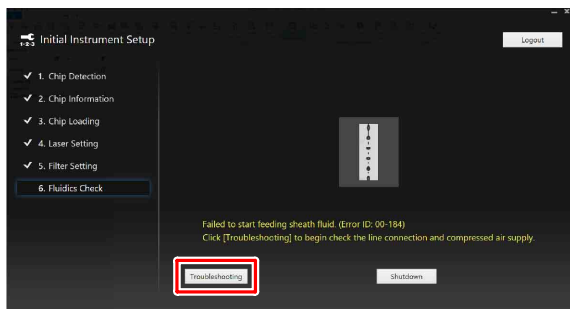


Tip

If you click the [Shutdown] button, the SH800 will shut down.

De-bubbling the Sheath Filter


You can de-bubble the sheath filter before running automatic calibration.

1 Click [Sheath filter de-bubble] on the [Fluidics Check] screen of the [Initial Instrument Setup] wizard.

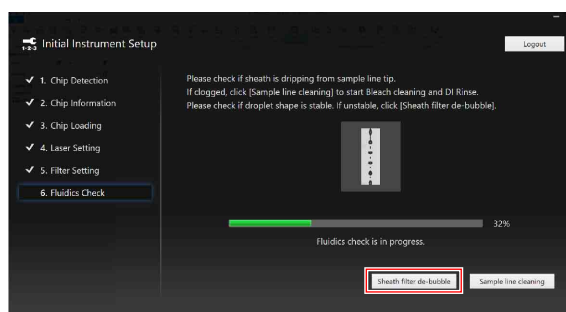

The [Sheath filter De-bubble] wizard appears.

2 Click [Start] to begin de-bubbling the sheath filter.

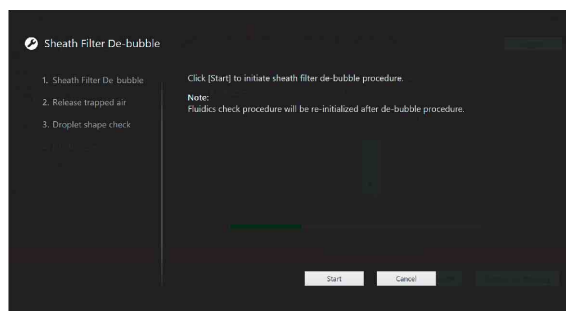

3 Click [Start] to begin de-bubbling the sheath filter. When de-bubble is completed, click [Next].

4 Follow the on-screen instructions and perform steps 4 to 8 of “To release air in the sheath filter manually” in “Releasing Air in the Sheath Filter and Sorting Chip” (page 202).

5 When all air has been released, click [Next].

6 Check that the droplet formation displayed on the screen is stable, then click [Next].

If unstable, click [De-bubble again] and repeat steps 3 to 5.

7 Click [OK] to exit the [Sheath filter De-bubble] wizard.

Tip


Running the sheath filter de-bubble function initializes the fluidics check in the Initial Instrument Setup.

When the [Sheath filter De-bubble] wizard finishes, the [Fluidics Check] screen appears. When the fluidics check is completed, the [Auto Calibration] screen appears.

Cleaning the Sample Line


You can clean the sample line before running automatic calibration.

- 1 Click [Sample line cleaning] on the [Fluidics Check] screen of the [Initial Instrument Setup] wizard.

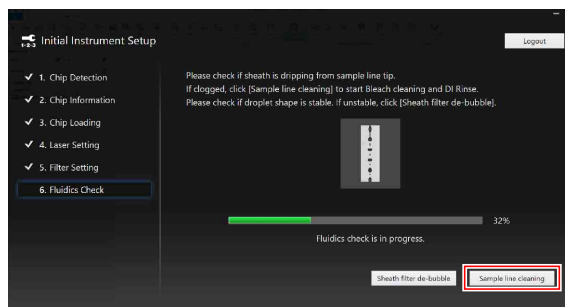

The [Sample Line Cleaning] wizard appears.

- 2 Click [Start] to begin cleaning the sample line.

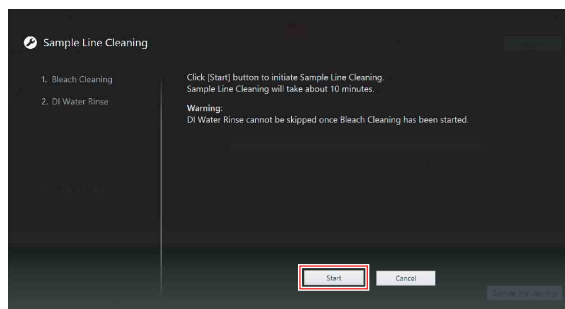

Follow the on-screen instructions.

In sample line cleaning, the fluidics system is cleaned using a sodium hypochlorite solution (bleach cleaning) and DI water, in that order.

Tips

- You can skip each cleaning step by clicking [Skip], but cleaning with DI water after cleaning with bleach cannot be skipped.
- You can also clean the system again after running the cleaning steps by clicking [Cleaning again].
- For details about cleaning using Sodium Hypochlorite solution, see “*Cleaning the Sample Fluidics System using Bleach*” (page 170).


When sample line cleaning finishes, the [Fluidics Check] screen appears. When the fluidics check is completed, the [Auto Calibration] screen appears.

Automatic Calibration

After configuring the excitation lasers and emission detection optical filters, the startup procedure performs sorting chip alignment, droplet calibration, sort delay calibration, and side stream calibration on the [Auto Calibration] screen.

Note

If the SH800 is installed in a BCC300 Biosafety Cabinet, wait 30 minutes after checking the fluidics lines before running auto calibration to allow the temperature within the cabinet to stabilize.

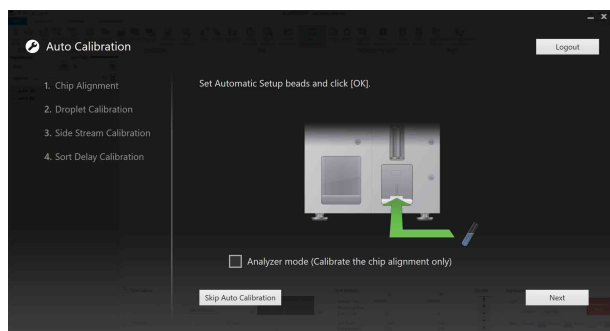


Tips

- You can click [Skip Auto Calibration] to bypass automatic calibration and proceed directly to the Cell Sorter Software main window. Bypassing automatic calibration is useful for quickly accessing the main window to perform analysis of existing data. Do not bypass automatic calibration if you plan to run any new samples.
- You can align the sorting chip and calibrate the sorting system later using [Chip Alignment] and [Sort Calibration], respectively, on the [Cytometer] tab of the ribbon from the main window.

1 Insert 0.5 ml or more of undiluted, automatic setup beads in a sample tube, and place the tube in the corresponding sample tube holder.

For details about sample tubes, sample tube holders, and automatic setup bead preparation, see “Preparing a Sample” (page 33).

2 Press the sample loader door button to open the sample loader door, and place the sample tube holder in the sample loader.

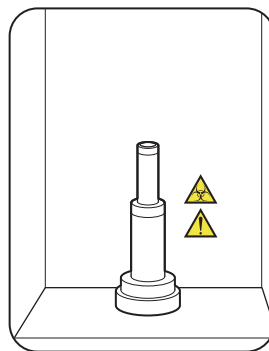

Align the D-shaped notch on the bottom of the sample tube holder with the slot in the top of the agitation unit when loading the sample tube holder.

Notes

- Always wear laboratory-use gloves, mask, protective goggles, and other protective clothing when handling sample fluid to provide protection against biohazards.
- Also, take care when loading the sample tube in the tube holder to avoid sample fluid spills and breakage of the sample tube.
- Before using a tube, check that there is no dust or other foreign matter in the tube.
- Do not subject the instrument to vibration during operation. Shock or vibration may adversely affect alignment and calibration.

3 Press the sample loader door button to close the sample loader door.

Tip

The sample loader door will close automatically, if you leave the door open, when you click [OK] in Cell Sorter Software in the next step to begin calibration.


4 Click [OK] on the [Auto Calibration] screen.

Tip


If you are planning to analyze samples without sorting them, you can select the [Analyzer mode (Calibrate the chip alignment only)] checkbox and click [OK] to bypass part of automatic calibration. This displays the main window in Analyzer mode.

Using LE-C32 series sorting chips

A screen for selecting auto calibration settings appears. Select [Standard] or [Targeted], and click [OK]. “Targeted” is a setting suitable for multi-well plate sorting and sorting of large cells.

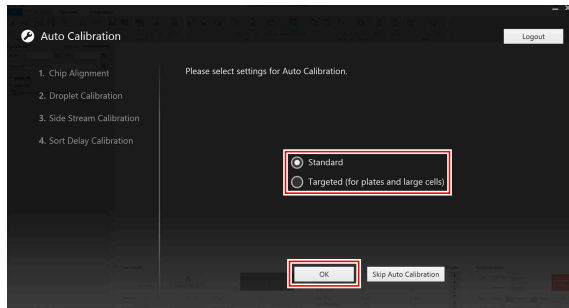

The automatic calibration sequence starts. A progress bar displays the status at each stage of the calibration process.

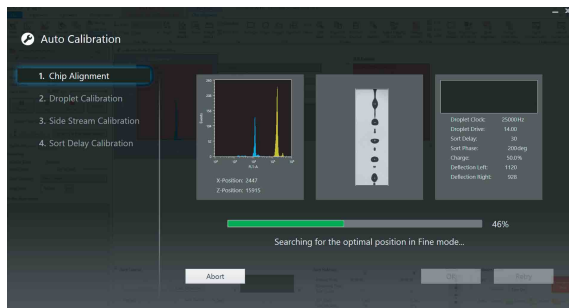

As automatic calibration proceeds, a histogram for the automatic setup beads, the droplet camera image and the side stream monitor image are displayed.

Tip

During automatic calibration, you can click [Abort] to return to the previous screen.

When automatic calibration finishes, [Abort] is disabled, and [Retry] and [OK] are enabled. Clicking [Retry] restarts the automatic calibration procedure.


5 When calibration finishes, click [OK].

The [Create Experiment] window appears for building a new experiment.


Note

During automatic calibration, precision measurement is performed using the side stream monitor. Accordingly, do not place any bright light sources (halogen lamps, for example) within the following angular ranges. Doing so may cause automatic calibration to malfunction.

Top view

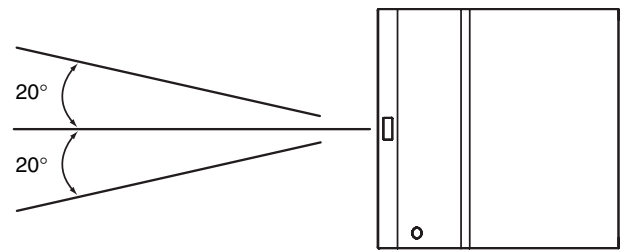

Side view

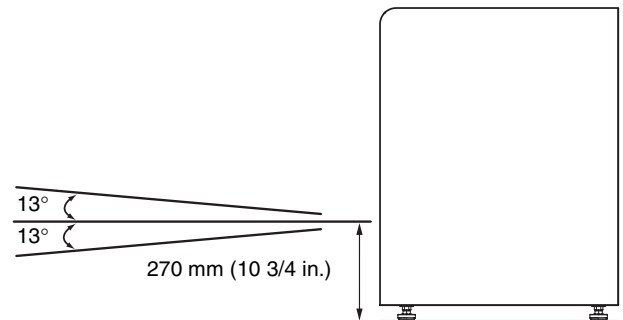


If an error occurs during auto calibration

If an error occurs during sorting chip alignment or side stream calibration, [Troubleshooting] appears at the bottom left of the screen.

Click [Troubleshooting] to launch the troubleshooting wizard. Follow the on-screen instructions to resolve the issue.

If [Troubleshooting] is not displayed when an error occurs, see “*Error Messages*” (page 244).

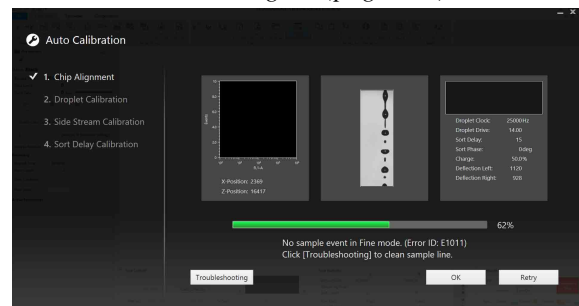

Tip

The content in the Troubleshooting wizard will vary depending on the type of error.


Creating an Experiment

When calibration finishes, you create an experiment that is used to acquire data from samples in the Create Experiment window.

Creating a New Experiment

This section describes how to create a new experiment using existing templates and/or recent experiments.

- 1 If the [Create Experiment] window is not displayed, click [New] on the [File] tab of the ribbon.
- 2 Create an experiment.

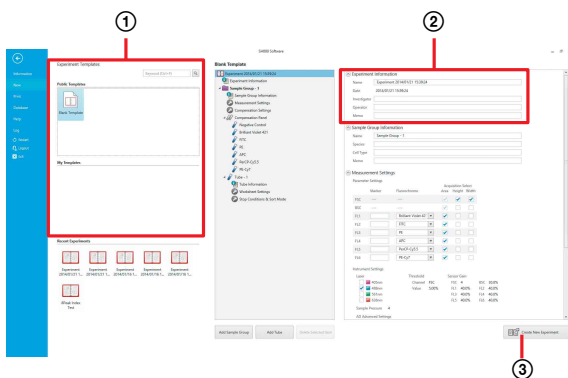

- 1 Select a template.
Clicking a template or an experiment on the left displays the structure of the selected experiment on the right side of the window.


For details about changing the structure of an experiment, see “Creating a Customized Experiment” (page 77).

- 2 Enter a name for the experiment in [Name].
You can configure sample groups, sample tubes, pulse parameters, and other settings for data acquisition.

For details about each item, see “[Create Experiment] Window” (page 128).

- 3 Click [Create New Experiment].
If “Blank Template” was selected, the [New Experiment Startup Procedure] dialog appears. Proceed to step 3.
If an existing template or experiment was selected, the main window appears. See “Acquiring Data” (page 53).

- 3 Select the action to perform, then click [OK].

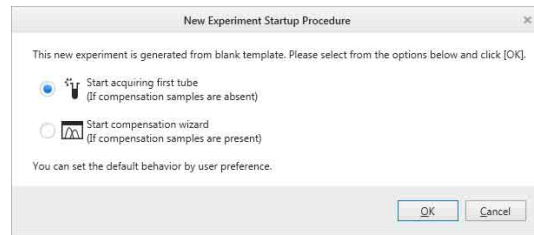


[Start acquiring first tube]: Displays the main window and assigns the first tube in preparation for data acquisition.

[Start compensation wizard]: Displays the [Compensation Wizard] to perform fluorescence compensation in preparation for measurement.

For details about the [Compensation Wizard], see “Performing Fluorescence Compensation” (page 67).

The experiment is created, and the main window appears.

The created experiment is automatically made the active experiment, and the first tube is automatically assigned as the active tube.

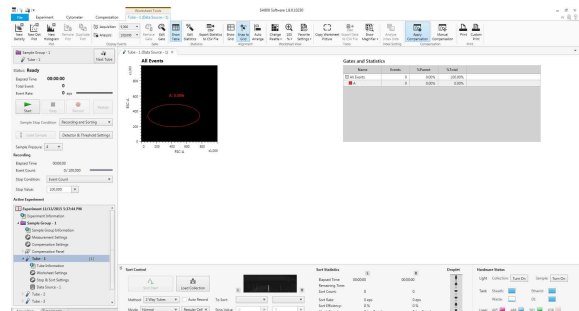

To create an experiment using an existing experiment

- 1 In step 2-1 of the procedure above, select an experiment you want to use from [Recent Experiments] instead of a template.

[Recent Experiments] displays the most recently used experiments in reverse chronological order.



- 2 Check the experiment details, and enter a name for the experiment in [Name] under [Experiment Information].

3 Click [Create New Experiment].

The experiment is created, and the main window appears.

To save an experiment as a template

You can save an experiment as a shared template (Public Templates) or a private template (My Templates).

For details, see “Saving an Experiment as a Template” (page 85).

Acquiring Data


After an experiment is created, you acquire and record data for samples in the experiment from the main window. Recorded data is displayed in real-time, and can also be exported for analysis at a later time.

Before starting acquisition

Check the following items:

- Check the fluid levels in each of the tanks.
- Check that the main unit and fluidics cart are connected correctly.

You control data acquisition and recording in the [Acquisition] tab of the Experiment/Acquisition control pane.

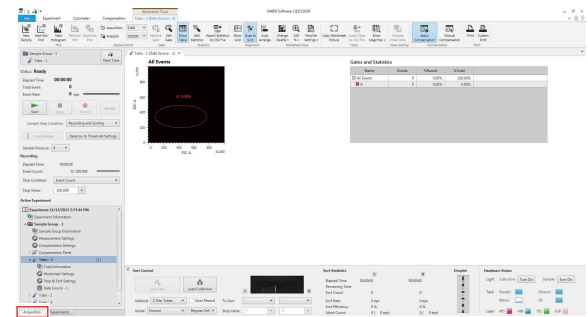


[Acquisition] tab

1 Insert sample fluid in a sample tube, and place the tube in the corresponding sample tube holder.

For details about sample tubes and sample tube holders, see “Preparing a Sample” (page 33).


2 Place the sample tube holder in the sample loader.

If the sample loader door is closed, press the sample loader door button to open the sample loader door.

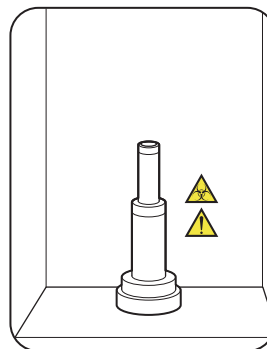


3 Press the sample loader door button to close the sample loader door.

Tip

The sample loader door will close automatically, if you leave the door open, when you click  (Start) in Cell Sorter Software to start acquisition.

4 Specify the measurement stop condition in [Sample Stop Condition].

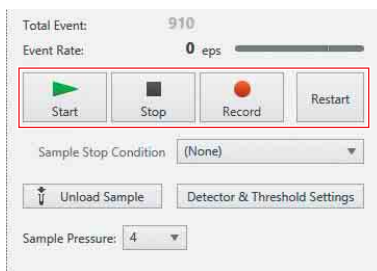

The following options are available.

- None (data acquisition will stop automatically when the number of detected events exceeds 100,000,000 or the elapsed time exceeds 12:00:00, whichever occurs first)
- Recording (stops data acquisition automatically when the recording stop condition is satisfied)
- Sorting (stops data acquisition automatically when the sorting stop condition is satisfied)
- Recording and Sorting (stops data acquisition automatically when both the recording and sorting stop conditions are satisfied)

Tip

[Recording] cannot be selected for index sorting. For details about index sorting, see “*Index Sorting*” (page 114).

For details, see “*Configuring Detector Settings*” (page 86).

5 Configure the detector settings.


Adjust the gain, threshold, and other settings.

6 Specify the recording stop condition under [Recording].


The following options are available.

- None
- Elapsed Time (total elapsed time)
- Event Count (total number of events)
- Gated Event Count (number of events for the specified gate)


You can also enter a value directly from the keyboard.

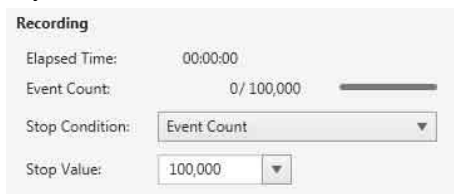

When [Gated Event Count] is selected, select the gate for the automatic stop condition and specify the number of events within the selected gate.

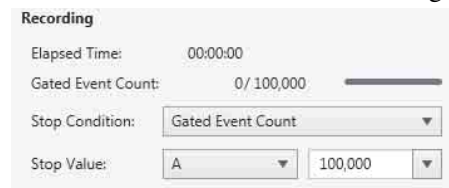

7 Click (Start).

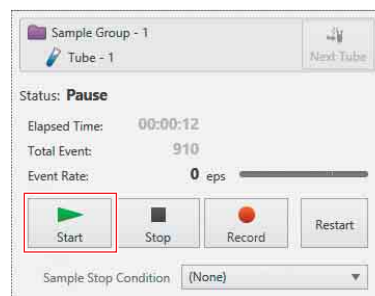


Sample fluid starts flowing from the sample tube into the sorting chip, and data acquisition starts.

8 Click (Record).

Data recording starts.

If an automatic stop condition was specified, data acquisition and recording stops automatically when the stop condition has been satisfied.

Tip

 (Record) cannot be selected for index sorting. For details about index sorting, see “*Index Sorting*” (page 114).


9 If an automatic stop condition was not specified (None), click (Stop) to stop data acquisition and recording.

The sample tube unloads automatically.

[Acquisition] Tab Main Functions

During data acquisition, the data is displayed on plots and in the [Gates and Statistics] table on the worksheet. You use the buttons on the [Acquisition] tab, as required, to control data acquisition while monitoring the data acquisition status on the worksheet. You can change the type of plots, edit gates on plots, and customize the content displayed in the [Gates and Statistics].


Controls data acquisition and recording.

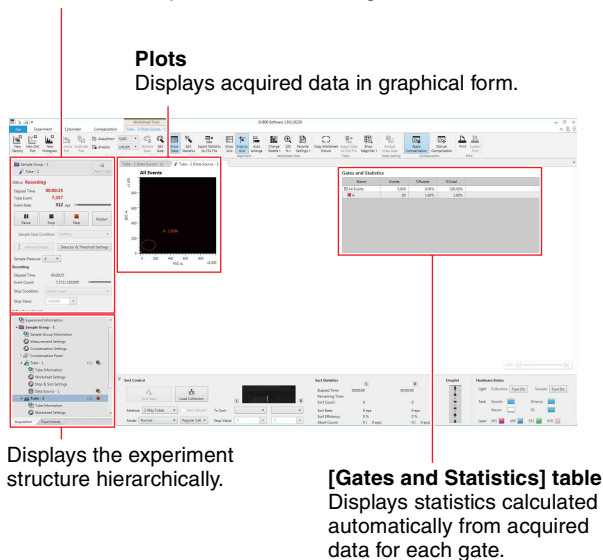
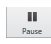

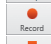
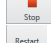



Controlling Data Acquisition and Recording


After data acquisition starts, the elapsed time since the start of data acquisition, the total number of events acquired, and the event rate measured in events/second (eps) are displayed on the [Acquisition] tab. You control acquisition, as required, while monitoring the acquisition status.

For details about the displayed items, see “[Acquisition] Tab (control pane)” (page 149).

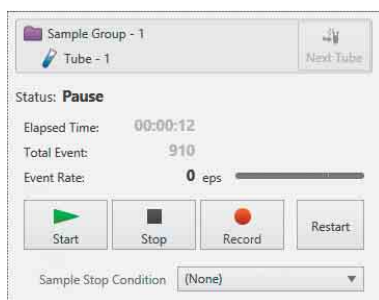

-  **(Start):** Starts/resumes data acquisition. This button does not start data recording.
-  **(Pause):** Pauses data acquisition. The elapsed time continues counting while acquisition is paused.
-  **(Stop):** Stops data acquisition manually.
-  **(Record):** Records the acquired data.
-  **(REC stop):** Stops data recording manually.
-  **(Restart):** Resets the total events and elapsed time counters to 0 and restarts data acquisition. Clicking this button while recording will delete all data recorded up to that point.


Tips

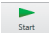- If an automatic stop condition has been specified, data acquisition and/or recording stops automatically when the corresponding stop condition has been satisfied.
- Clicking the  (Start) button to start data acquisition automatically switches the worksheet display to the assigned sample tube data, even if another sample tube data is displayed on the worksheet.

Customizing the Worksheet

You can change the type of plots, modify gates on plots, and customize the content in the [Gates and Statistics] table displayed on the worksheet during data acquisition. You customize the items on the worksheet using the tabs on the ribbon.

Example: [Worksheet Tools] tab

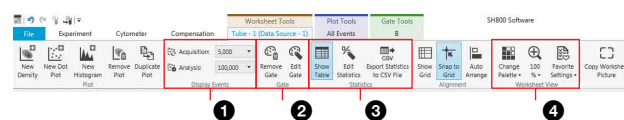

- 1 [Display Events]**
Sets the number of events displayed on plots.
 - 2 [Gate]**
Tools for editing gates.
 - 3 [Statistics]**
Shows/hides the [Gates and Statistics] table, customizes the content displayed in the table, and exports the table as a CSV file.
 - 4 [Worksheet View]**
Changes the view and scale of plots on the worksheet.
- For details, see “[Worksheet Tools] Tab (ribbon)” (page 143).

Analyzing Data


After data acquisition finishes, you analyze the acquired data events. The most recent data is displayed on the worksheet immediately after acquisition finishes. To analyze data for a past experiment, search for and select the target experiment on the [Experiments] tab of the Experiment/Acquisition control pane.

Plots

Displays acquired data in graphical form. Plots can be added/edited.

Gates

Groups events with common characteristics, isolating them from all other events.

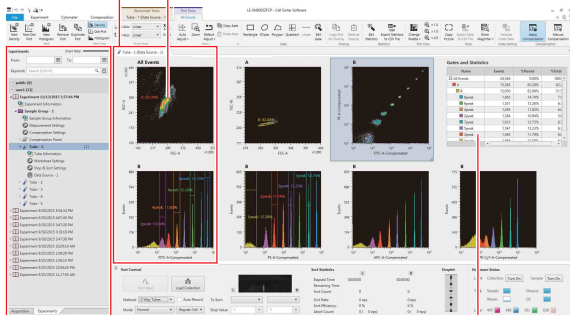

[Experiments] tab

Displays experiments that can be selected for analysis.

[Gates and Statistics] table


Displays statistics calculated automatically from acquired data for each gate. You can customize the statistics displayed in the table.

This section describes how to add plots to the worksheet, how to add gates to a plot, and how to display statistics for the gates.


For details, see “Analysis” (page 91).

1 Select the [All Events] plot on the worksheet.

The [All Events] plot is the top level in the hierarchy. All plots for gates are derived from [All Events]. If necessary, adjust the axes scales using the tool buttons in the [Axes] group on the [Plot Tools] tab of the ribbon.

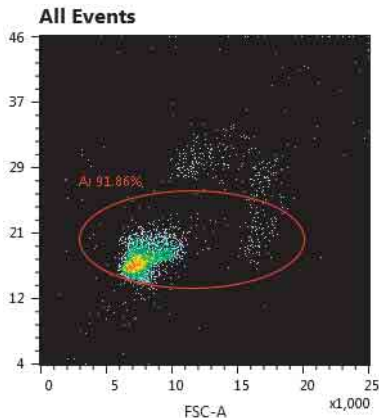

2 Adjust the gate size and position to surround the target population.

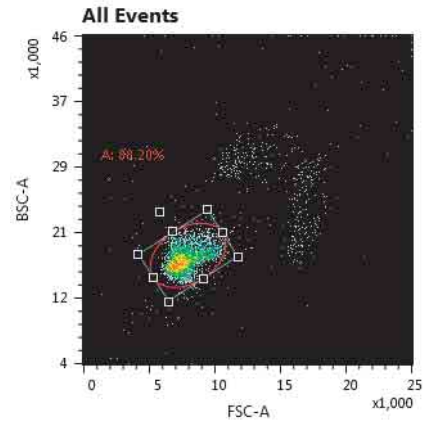

3 Double-click the gate in the plot to add a child plot for that gate.


You can also add a child plot from the context menu by right-clicking the gate in the [Gates and Statistics] table.

Tips

- You can add a new plot using the tool buttons in the [Plot] group (page 143) on the [Plot Tools] tab or the [Worksheet Tools] tab of the ribbon.
- You can change the type of the plot using the tool buttons in the [Plot Type] group (page 146) on the [Plot Tools] tab of the ribbon.


4 Click the axis labels and select the parameters to be displayed from the popup menu.

Adjust the axes scales to display the events on the plot, as required.

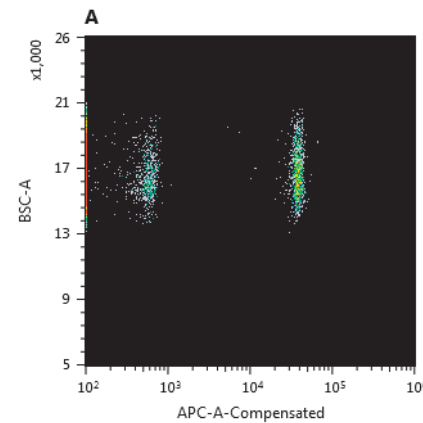

5 Select a gate type in the [Gate] group of the [Plot Tools] tab, and draw gates around both of the populations in the child plot.


The gates are automatically named in alphabetical order. You can change the name and color of gates in the [Property Window] dialog by right-clicking a gate and selecting [Properties] from the context menu.

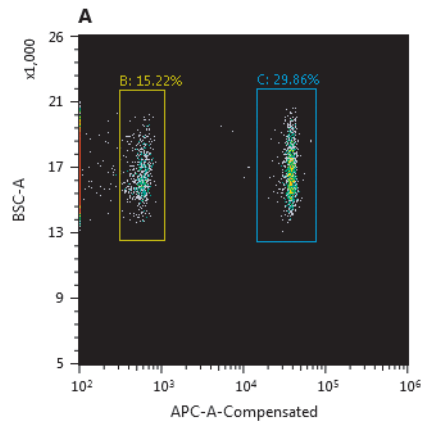

- 6 Click [Edit Statistics] in the [Statistics] group on the [Worksheet Tools] tab or [Plot Tools] tab.

The [Statistics Editor] dialog appears.

- 7 Select the statistics to display in the [Gates and Statistics] table.

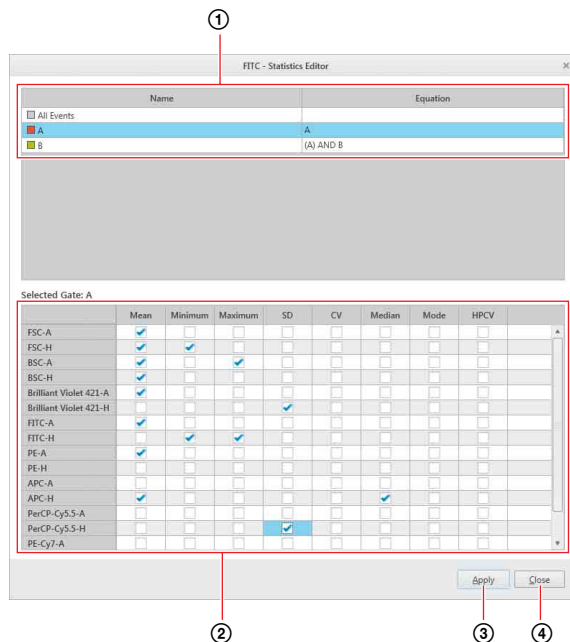

- ① Select a gate.
- ② Place check marks in the checkboxes for the statistics to display.
- ③ Click [Apply] to update the [Gate and Statistics] table.
- ④ Click [Close] to close the dialog.
The selected statistics are displayed in the [Gates and Statistics] table.

Name	Events	%Parent	%Total	Mean	Maximum	CV
All Events	5,000	0.00%	100.00%			
A	4,477	89.54%	89.54%			
FSC-A				8,052		12.27%
FSC-H				77,516		13.32%
BSC-A				17,539		
BSC-H				141,886		7.67%
CD56: Brilliant...						-151.91%
CD45: FITC...					7,981	
CD45: FITC...					71,213	
C	1,484	33.15%	29.68%			

Sorting

Once events have been isolated on plots using gates, the events can be sorted into different collection tubes for further analysis. Sorting enables the targeted cells to be recovered from a sample to increase the concentration of the cells or to isolate cells for increased purity.

Supported Sorting Methods for Sorting Chip Nozzle Diameter

The sorting methods that can be used vary depending on the nozzle diameter of the sorting chip and the size of cells.

SH800S, SH800Z, and SH800

Sorting method		2 Way Tubes		8 Well Strips		Slides	
Cell size settings		Regular Cell	Large Cell ²⁾	Regular Cell	Large Cell ²⁾	Regular Cell	Large Cell ²⁾
Sorting chip nozzle	100 µm	Yes	Yes	Yes	Yes	Yes	Yes
	130 µm	Yes	–	–	–	–	–
	70 µm ¹⁾	Yes	–	–	–	–	–

1) A 70 µm sorting chip is supported in SH800S, SH800, and SH800Z models that have been upgraded to support the 70 µm sorting chip.

2) [Large Cell] is not supported for LE-C32 series sorting chips.

SH800SP and SH800ZP

Sorting method		2 Way Tubes		96 Well Plate		384 Well Plate	
Cell size settings		Regular Cell	Large Cell ²⁾	Regular Cell	Large Cell ²⁾	Regular Cell	Large Cell ²⁾
Sorting chip nozzle	100 µm	Yes	Yes	Yes	Yes	Yes	Yes
	130 µm	Yes	–	Yes	–	Yes	–
	70 µm ¹⁾	Yes	–	Yes	–	Yes	–

1) A 70 µm sorting chip is supported in SH800SP and SH800ZP models that have been upgraded to support the 70 µm sorting chip.

2) [Large Cell] is not supported for LE-C32 series sorting chips.

Notes

- [Large Cell] is an assistive mode for the sorting of large-sized cells when using the 100 µm sorting chip. Use it if you are having trouble sorting large cells. However, this function does not guarantee complete accuracy in sorting large cells. A rough guide for the size of large cells for which to use the [Large Cell] setting is 15 µm or larger.
- When sorting large cells using a 100 µm sorting chip (LE-C32 series), use the [Targeted] setting in automatic calibration.

For details, see “Using LE-C32 series sorting chips” (page 51).

Recommended conditions for Regular Cell sorting (96/384-well single cell mode)


In 96/384-well single cell mode sorting, the following sorting conditions are recommended.

Event rate	800 eps to 1,200 eps
Cell size	15 µm or smaller (100 µm, 130 µm sorting chip) 10 µm or smaller (70 µm sorting chip)
Sorting mode	Single Cell
Sample pressure (sample fluid pressure)	6 or lower

Two-way Tube Sorting

- 1 Check that data acquisition and recording are stopped or paused.

If data acquisition is in progress, click the (Stop) to stop data acquisition/recording or (Pause) to pause data acquisition/recording.

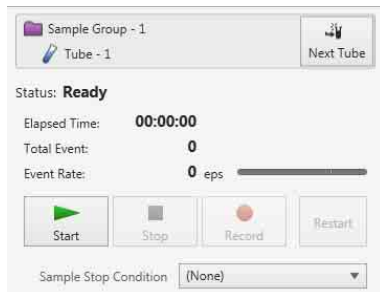


- 2 Place the collection tubes into the corresponding holder.


For details, see “Preparing Collection Tubes” (page 35).

- 3 Open the collection area door, place the holder on the collection stage, and close the door.


SH800, SH800Z, and SH800ZP

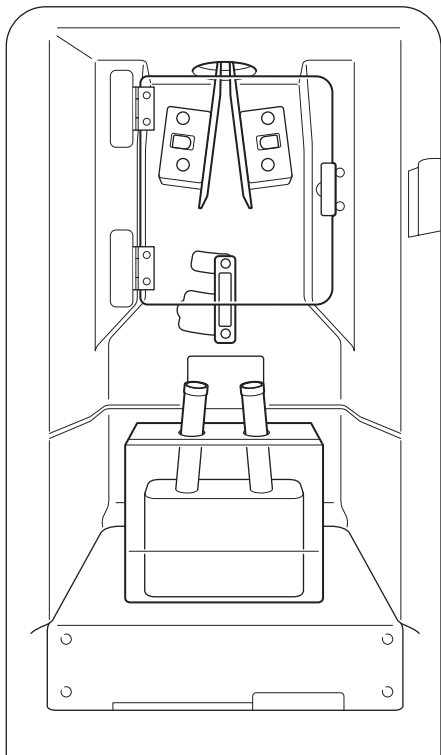

SH800S and SH800SP

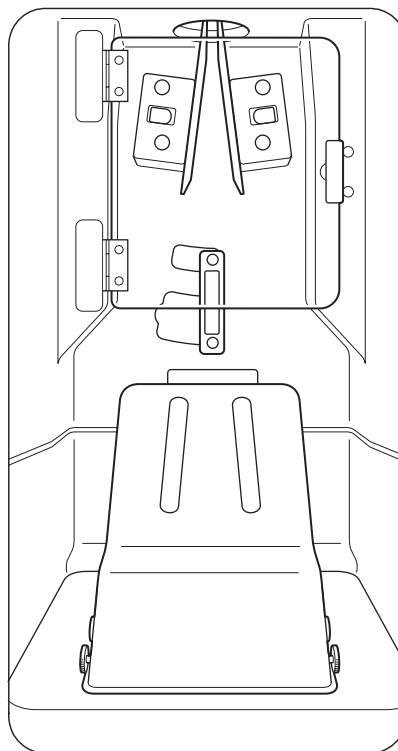

- 4 Click [Load Collection] in the [Sort Control] pane.

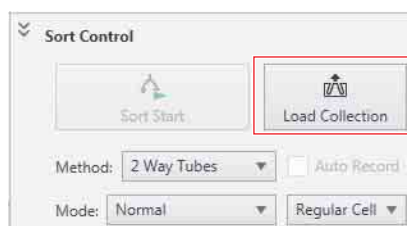


The collection tube device is loaded.

Note

Do not open the collection area door during sorting.


- 5 Select [2 Way Tubes] in [Method].

Always select the sorting method matching the collection tube device loaded on the collection stage.

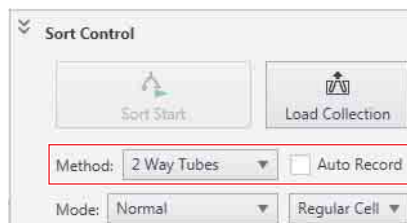

The sorting parameters vary depending on the selected sorting method.

To start recording at the same time as sorting, place a check mark in the [Auto Record] checkbox. [Sort Start] changes to [Sort & Record Start].

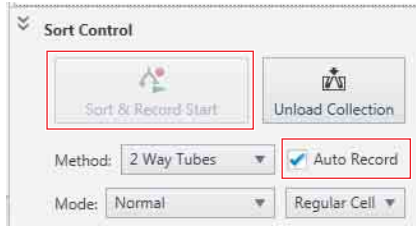

6 Set each parameter.

For details about each item, see “[Sort Control] Pane” (page 153).



[Mode]: Selects the sorting mode and cell size.

When sorting larger cells using a 100 μm /70 μm sorting chip, selecting [Single Cell (3 drops)] may improve the sorting performance. The cell size is fixed to [Large Cell] when [Single Cell (3 drops)] is selected.

To adjust the position for sorting of target cells within a droplet, you can adjust the settings using [Edit setting].

For details, see “[Sort Mode Setting] dialog” (page 156).


[To Sort]: Selects the gates for the target events to sort into the left and right collection tubes.

[Stop Value]: Specifies the event count conditions to stop recording automatically for each of the left and right collection tubes.


You can select a value or enter a value from the keyboard.

Entering a value of “0” disables the stop condition.

7 Click [Start] or [Resume] in the data acquisition control pane to begin data acquisition.

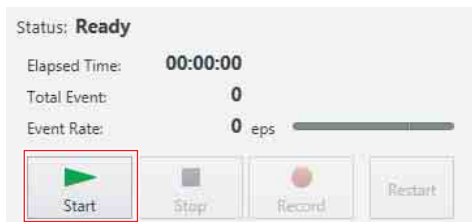


8 Click [Sort Start] or [Sort & Record Start] in the [Sort Control] pane to begin sorting.

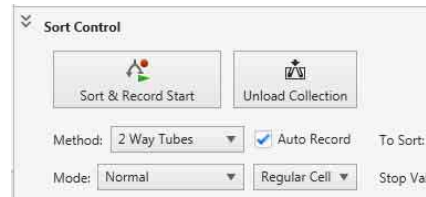

Basic sorting status is displayed, together with progress bars if an automatic stop condition is specified.

Sort Statistics	L	R
Elapsed Time:	00:00:00	00:00:00
Remaining Time:		
Sort Count:	0	0
Sort Rate:	0 eps	0 eps
Sort Efficiency:	0 %	0 %
Abort Count:	0 (0 eps)	0 (0 eps)


You can monitor the side streams and droplet flow in the main window.

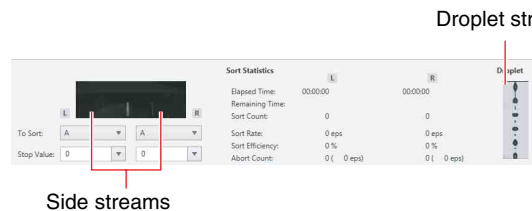

Notes

- If the SH800S and SH88SP are upgraded to the 70 μm sorting chip, the side stream laser turns off by default. To turn the laser on, see [Side Stream Laser] in “[Cytometer Settings] dialog” (page 138).
- When using the collection tube holder with embedded cooling agent (5 ml), turn the side stream laser off.

Double-clicking the image displays the side streams in a separate window for monitoring droplets.

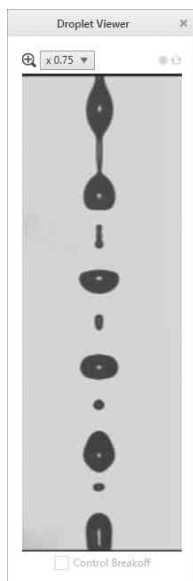

If an automatic stop condition is configured, a stop message appears when the stop condition is satisfied and sorting stops.

- 9 Click [OK] to close the confirmation message dialog.

Tip

If an event is detected that falls within both gates selected as the target for the left and right side streams, the left side stream takes precedence.



Note


Do not sort sample fluid above the permitted volume into the tube and do not overfill the tube with buffer solution to prevent sample fluid overflow from the tube.

For details about cleaning if an overflow occurs, see “Cleaning the Collection Tube Holder with Embedded Cooling Agent (5 ml)” (page 171).

Sorting onto 8-well Strips (SH800S, SH800Z, and SH800)

- 1 Check that data acquisition and recording are stopped or paused.

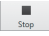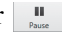If data acquisition is in progress, click the  (Stop) to stop data acquisition/recording or  (Pause) to pause data acquisition/recording.

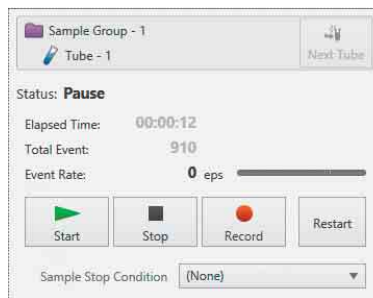

- 2 Place the 8-well strips or glass slides onto the corresponding holder.

For details, see “Preparing Collection Tubes” (page 35).

- 3 Open the collection area door, place the holder on the collection stage, and close the door.

When 8-well strip holder is loaded



When slide glass holder is loaded




Notes

- Do not sort sample fluid above the permitted volume into the wells and do not overfill the wells with buffer solution to prevent sample fluid overflow from the wells.
- Take care not to spill samples when removing the 8-well strips or glass slides from the multi-well plate holder.
- Always load 8-well strips or slide glass according to the method selected in [Method] on the [Sort Control] pane.
- Do not open the collection area door during sorting.

- Do not attempt to remove the 8-well strips or slide glass while the collection stage is moving.


4 Select the appropriate plate sorting method in [Method].

Always select the sorting method matching the collection tube device loaded on the collection stage.

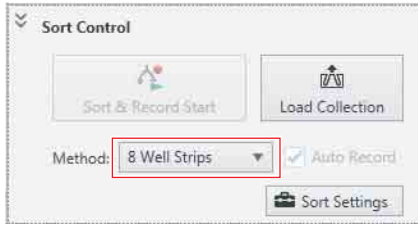

The sorting parameters vary depending on the selected sorting method.


5 Click [Sort Settings].

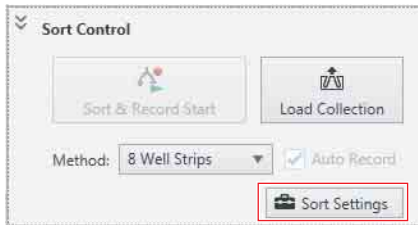

The [Sort Settings] dialog appears.

6 Configure the settings for sorting into wells on the [Plate Sort Settings] tab.

The description in this section shows an 8-well strip as an example. The operation is the same for glass slides.

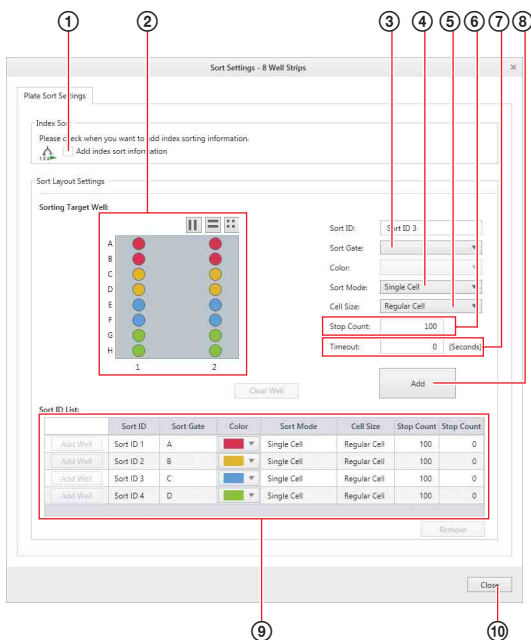

① If using index sorting, enable the [Add index sort information] checkbox.

In index sorting, you can associate the cells sorted in each well with the events for those cells on plots.

For details, see “Index Sorting” (page 114).

② Select the target well to set.

You can select the well using the following buttons or by clicking the well on the palette.

☐ : Select wells in alternate columns.

☐ : Select wells in alternate rows.

☐ : Select wells in groups of four.

You can select multiple wells using the Shift and Ctrl keys.

The ID is displayed in [Sort ID].

You can change the value in [Sort ID] manually.

③ Select the gate whose population you want to sort into the well in [Sort Gate].

You can change the color of the gate in [Color].

④ Select a sorting mode in [Sort Mode].

The default is “Single Cell.”

“Single Cell” mode is recommended for reliable sorting.

When sorting larger cells using a 100 μm sorting chip, selecting [Single Cell (3 drops)] may improve the sorting performance.

To adjust the position for sorting of target cells within a droplet, you can adjust the settings using [Edit setting].

For details, see “[Sort Mode Setting] dialog” (page 156).

⑤ Select the size of the cells to be sorted in [Cell Size].

As a rough guide, select [Regular Cell] for cells up to 15 μm in size, and [Large Cell] for larger cells.

The cell size is fixed to [Large Cell] when [Single Cell (3 drops)] is selected in [Sort Mode].

Note

When you select [Large Cell], a charge setting suited for sorting larger cells is used. However, this function does not guarantee complete accuracy in sorting large cells.

⑥ Enter the number of events to sort in [Stop Count].


Note

If a value exceeding the values below is entered in [Stop Count], a message appears warning you of the risk that sample fluid may overflow the well (operation does not stop).


Sorting chip	8-well strips	Slides
100 μm nozzle	50,000 events	300 events

- ⑦ Enter the timeout time in [Timeout] in units of seconds.
- ⑧ Click [Add].
The settings are added to [Sort ID List].
- ⑨ Repeat steps ② to ⑧ to set parameters for any other wells if necessary.
You can modify the parameters for a well by clicking the corresponding cells in [Sort ID List].
- ⑩ When finished, click [Close] to close the dialog.

- 7 Click [Start] or [Resume] in the data acquisition control pane to begin data acquisition.

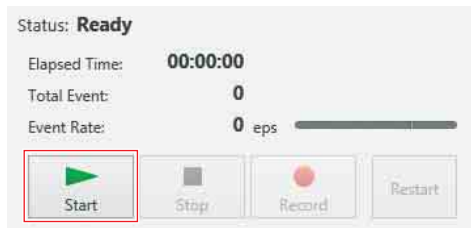


- 8 Click [Sort Start] or [Sort & Record Start] in the [Sort Control] pane to begin sorting.

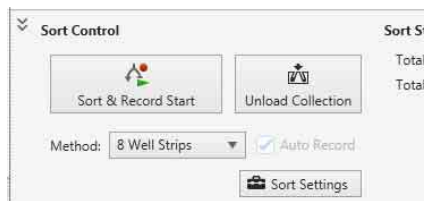


Tip

Starting sorting automatically starts recording.

Basic sorting statistics are displayed, together with progress bars if an automatic stop condition is specified.

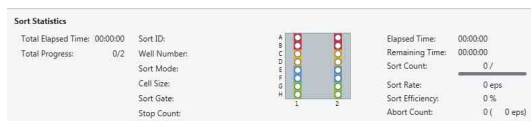


Double-clicking the image displays the side streams in a separate window for monitoring droplets.

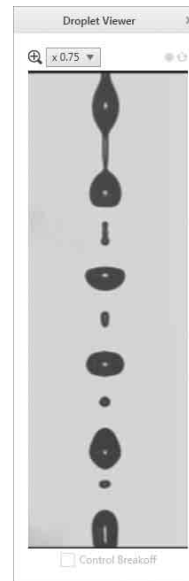

Double-clicking the plate image or right-clicking and selecting [Sorting Monitor] from the context menu displays the [Plate Sorting Monitor] window for monitoring progress. You can also display the window by right-clicking a tube in the tube list and selecting [Show Sorting Report] from the context menu.



After one set has been sorted, the following message appears.

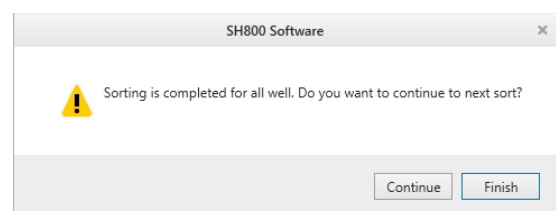

- 9 Click [Finish] to finish sorting.

To continue sorting, click [Continue].


When you click [Continue], the collection tube is unloaded, allowing you to load the next collection tube.

Sorting into Multi-well Plates (SH800SP and SH800ZP)

This section describes the procedure for sorting into a 96-well plate. The procedure is identical for sorting into 6, 12, 24, 48, and 384-well plates.

- 1 Check that data acquisition and recording are stopped or paused.

If data acquisition is in progress, click the (Stop) to stop data acquisition/recording or (Pause) to pause data acquisition/recording.

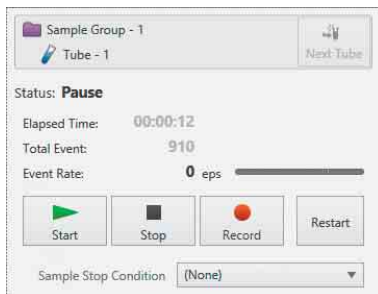

- 2 Attach the splash guard.

Always attach the splash guard when sorting into multi-well plates.


For details, see “Attaching the Splash Guard (for use with multi-well plates) (SH800SP and SH800ZP)” (page 38).

- 3 Place the 96-well plate into the corresponding holder.

For details, see “Preparing Collection Tubes” (page 35).

- 4 Open the collection area door, place the holder on the collection stage, and close the door.

96-well plate holder

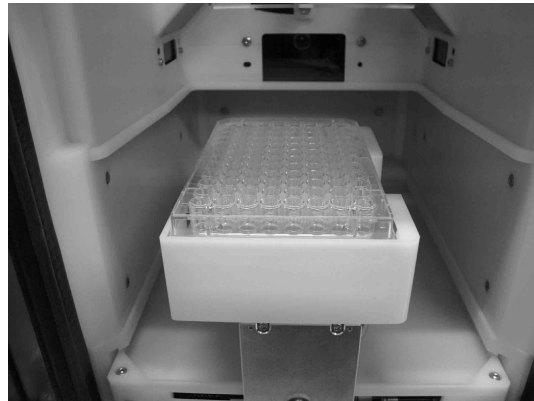

Notes


- If the multi-well plate is loaded in the incorrect orientation, correct sorting is not possible. Check the orientation when loading a multi-well plate.
 - Do not sort sample fluid above the permitted volume into the wells and do not overfill the wells with buffer solution to prevent sample fluid overflow from the wells.
 - Take care not to spill samples when removing the multi-well plate from the multi-well plate holder.
 - Securely mount the holder adapter to ensure successful sorting.
 - Always load a multi-well plate according to the method selected in [Method] on the [Sort Control pane].
 - Do not open the collection area door during sorting.
 - Do not attempt to remove a multi-well plate while the collection stage is moving.
- 5 Select the appropriate plate sorting method in [Method].

Always select the sorting method matching the collection tube device loaded on the collection stage.



The sorting parameters vary depending on the selected sorting method.


6 Click [Sort Settings].

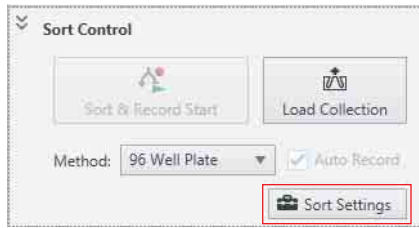

The [Sort Settings] dialog appears.

7 Configure the settings for sorting into wells on the [Plate Sort Settings] tab.

The description in this section shows a 96-well plate as an example. The operation is the same for all other multi-well plates.

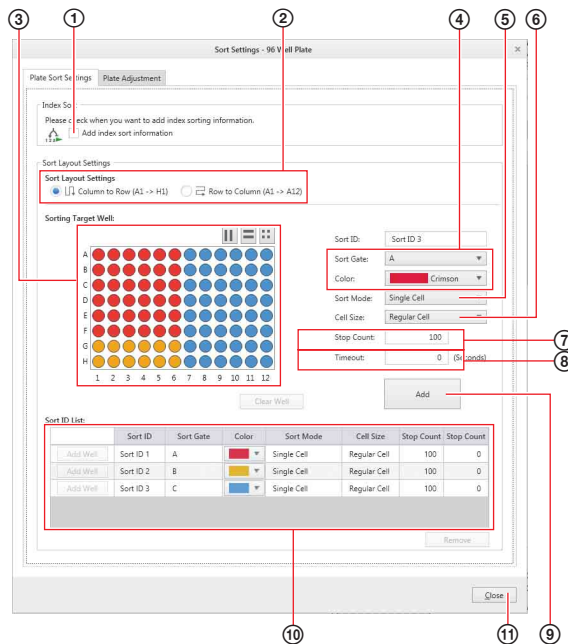

① If using index sorting, enable the [Add index sort information] checkbox.

In index sorting, you can associate the cells sorted in each well with the events for those cells on plots.




For details, see “Index Sorting” (page 114).

② Select the sorting sequence in [Sort Layout Settings].

- [Column To Row]: Sorting occurs in the sequence A1 → B1 → C1...H1 → A2 → B2...G12 → H12.
- [Row To Column]: Sorting occurs in the sequence A1 → A2 → A3...A12 → B1 → B2...H11 → H12.

③ Select the target well to set.

You can select the well using the following buttons or by clicking the well on the palette. You can select multiple wells using the Shift and Ctrl keys.

-  : Select wells in alternate columns.
The selection toggles between odd columns and even columns each time the button is clicked.
-  : Select wells in alternate rows.
The selection toggles between odd rows and even rows each time the button is clicked.
-  : Select wells in groups of four.
The selection toggles as follows each time the button is clicked.
Top left → Top right → Bottom left → Bottom right → Top left...

The ID is displayed in [Sort ID]. You can change the value in [Sort ID] manually.

④ Select the gate whose population you want to sort into the well in [Sort Gate]. You can change the color of the gate in [Color].

⑤ Select a sorting mode in [Sort Mode]. The default is “Single Cell.” “Single Cell” mode is recommended for reliable sorting.

When sorting larger cells using a 100 μm/70 μm sorting chip, selecting [Single Cell (3 drops)] may improve the sorting performance. To adjust the position for sorting of target cells within a droplet, you can adjust the settings using [Edit setting].

For details, see “[Sort Mode Setting] dialog” (page 156).

⑥ Select the size of the cells to be sorted in [Cell Size]. As a rough guide, select [Regular Cell] for cells up to 15 μm in size, and [Large Cell] for larger cells. The cell size is fixed to [Large Cell] when [Single Cell (3 drops)] is selected in [Sort Mode].

Note

When you select [Large Cell], a charge setting suited for sorting larger cells is used. However, this function does not guarantee complete accuracy in sorting large cells.

⑦ Enter the number of events to sort in [Stop Count].


Note

If a value exceeding the values below is entered in [Stop Count], a message appears warning you of the risk that sample fluid may overflow the well (operation does not stop).
96 Well Plate: 50,000 events

- ⑧ Enter the timeout time in [Timeout] in units of seconds.
- ⑨ Click [Add].
The settings are added to [Sort ID List].
- ⑩ Repeat steps ③ to ⑨ to set parameters for all wells. You can modify the parameters for a well by clicking the corresponding cells in [Sort ID List].
- ⑪ When finished, click [Close] to close the dialog.

For details about the [Plate Adjustment] tab, see “Adjusting the Sort Position (SH800SP and SH800ZP)” (page 113).

- 8 Click [Start] or [Resume] in the data acquisition control pane to begin data acquisition.

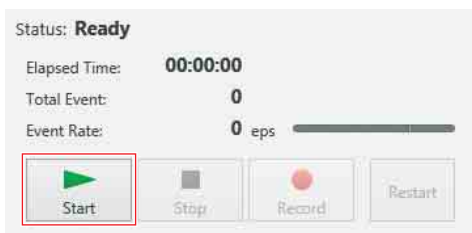


- 9 Click [Sort Start] in the [Sort Control] pane to begin sorting.




Tip

Starting sorting automatically starts recording.


Basic sorting statistics are displayed, together with progress bars if an automatic stop condition is specified.

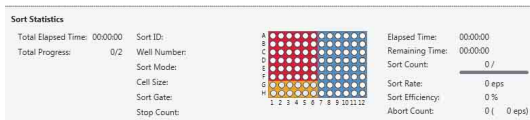


Double-clicking the image displays the side streams in a separate window for monitoring droplets.

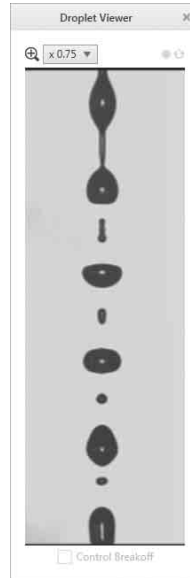

Double-clicking the plate image or right-clicking and selecting [Sorting Monitor] from the context menu displays the [Plate Sorting Monitor] window for monitoring progress. You can also display the window by right-clicking a tube in the tube list and selecting [Show Sorting Report] from the context menu.

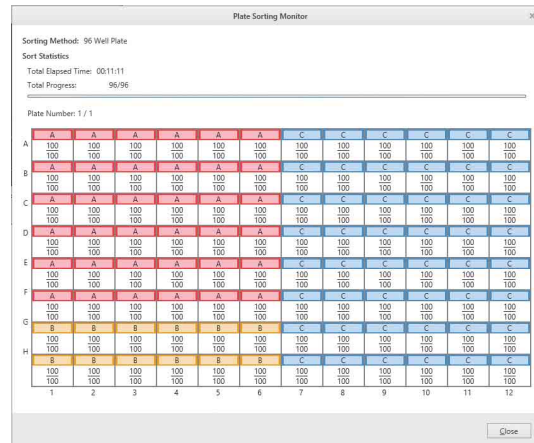

After one set has been sorted, the following message appears.

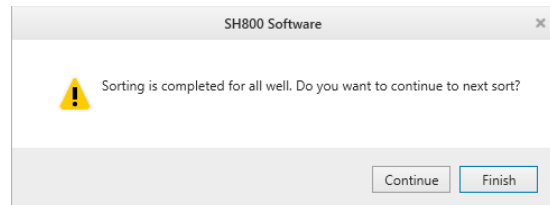

- 10 Click [Finish] to finish sorting.

To continue sorting, click [Continue].

When you click [Continue], the collection tube is unloaded, allowing you to load the next collection tube.

Performing Fluorescence Compensation

Fluorescence compensation is used to remove fluorescence spillover components from each channel to obtain the signal for the target fluorochrome only for each channel.

In fluorescence compensation, a compensation panel comprising an unstained negative control and a singly-stained, positive control samples for each fluorochrome are recorded.


This section describes the fluorescence compensation procedure using the [Compensation Wizard].

For details about using a gate for recording a control tube or adjusting fluorescence compensation manually, see “Adjusting Fluorescence Compensation Manually” (page 89).

Note


Do not change the gain of the channel detectors while recording the control tubes or samples for a compensation panel. The fluorescence compensation process is valid only for the instrument settings in place when all control tubes in a compensation panel are recorded. If the sensor gain of any detector is changed after recording the control tubes, the fluorescence compensation will no longer be valid and the control tubes will need to be loaded and recorded again.

- 1 Click [Compensation Wizard] on the [Compensation] tab of the ribbon.

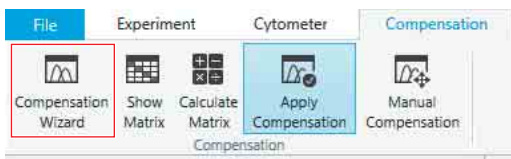

The [Compensation Wizard] launches.


- 2 Click [Next].

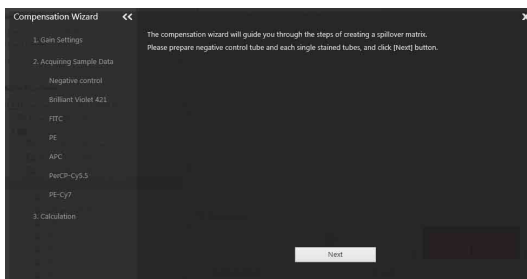

The following screen appears.


Follow the on-screen instructions to record control tubes.

The operating procedure is displayed.

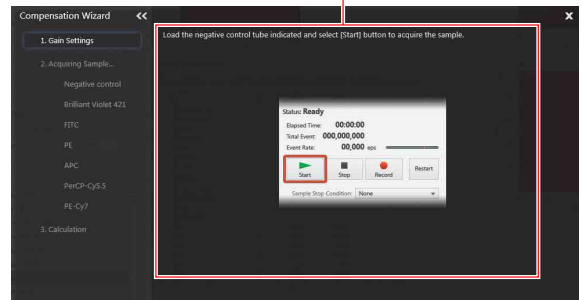


A control tube is automatically assigned as the active tube and the worksheet for the tube opens.

- 3 Load the control tube displayed in the wizard, then click [Start] on the data acquisition control panel.

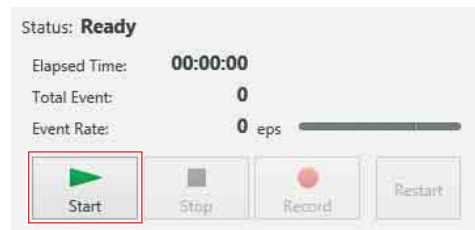


Tip

If, for any reason, you choose not to use a fluorochrome channel, you can click [Skip] to bypass the control tube.

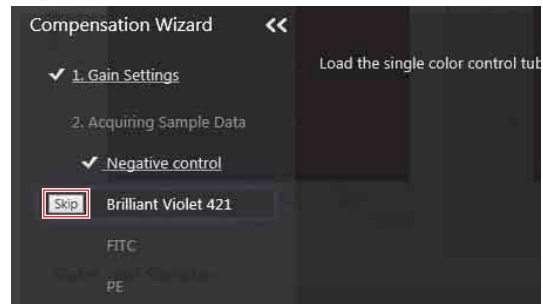


Data acquisition starts for the tube and data points are added to the events on the worksheet in real-time.

- 4 Click [Detector & Threshold Settings].

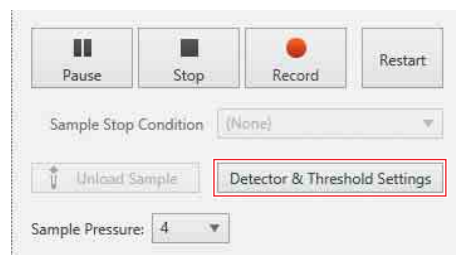


The [Detector & Threshold Settings] dialog appears.

- 5 Adjust the sensor gain of the FSC and BSC detectors to place the population on scale.

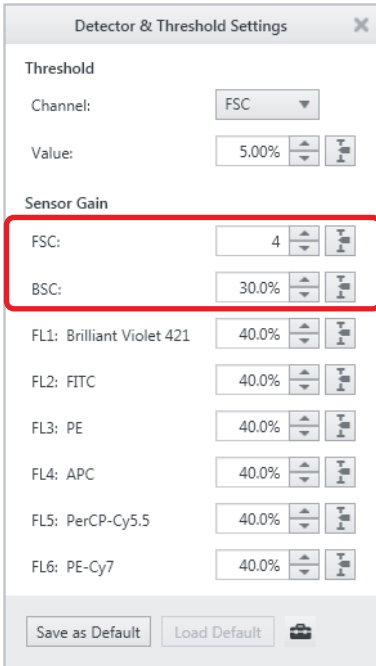


You adjust the gain while viewing the FSC vs. BSC density plot. You can also adjust the axis display scales to display a magnified portion of the plot.

- 6 Adjust the trigger threshold level to remove background noise without affecting the targeted population.

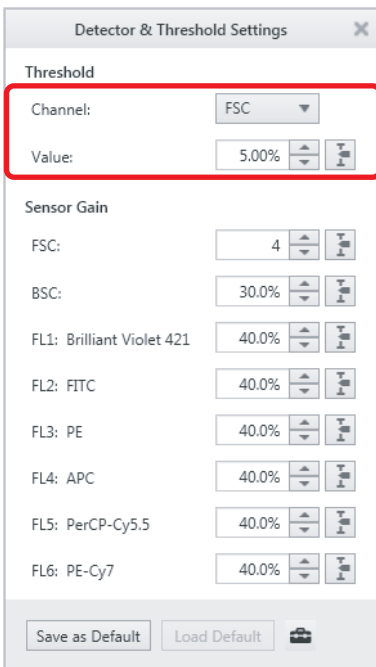


- 7 Adjust the size and position of the gate automatically added to the plot to encompass the desired population, then click [Next] in the [Compensation Wizard].

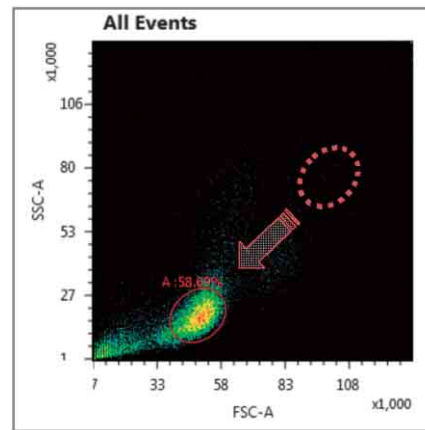


- 8 Adjust the gain levels of the FL1 to FL6 fluorescence channels to place the negative populations on scale in all channels.


You adjust the gain while viewing the histogram plots for each fluorescence channel.

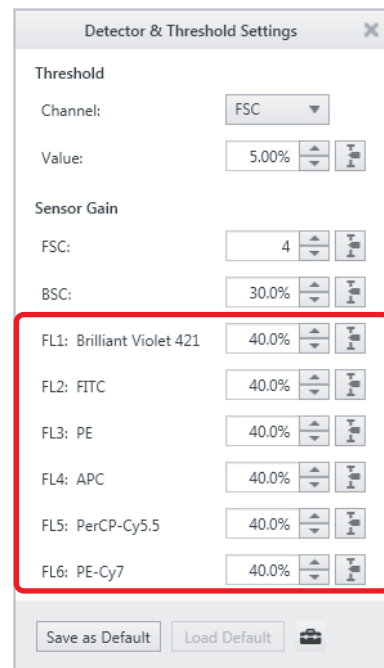


The negative peaks represent the signals detected due to cellular autofluorescence. The negative peaks should be on scale, and not wedged up against the axis in all channels.

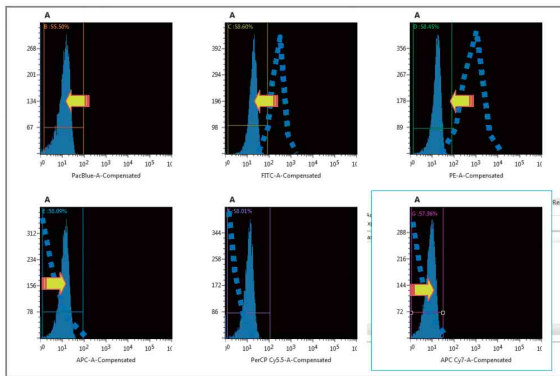

9 After adjusting the detector gain settings, click [Next].

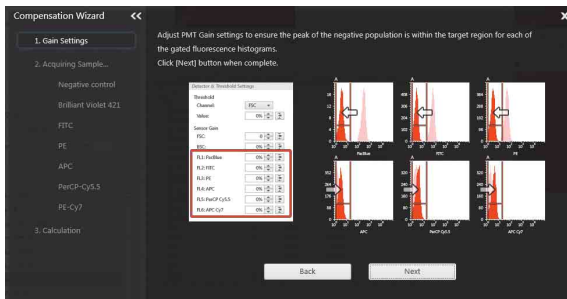

10 Click [Record] in the data acquisition control pane.

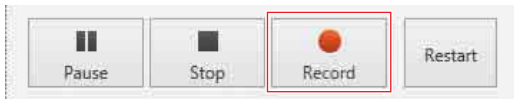


The SH800 starts recording the data acquired for the negative control tube. Recording stops automatically after satisfying the record stop condition when sufficient data has been recorded.

Tip


You can stop data acquisition and recording manually using [Stop] if a sufficient number of events have been recorded and the plot view has stabilized.

11 Click [Next].

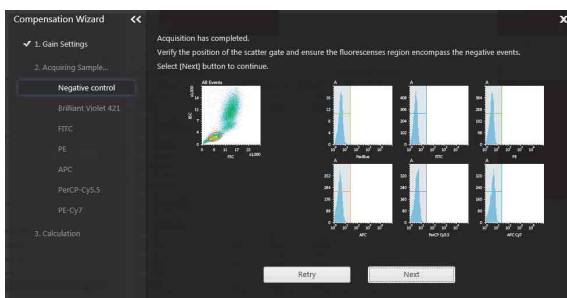

Data acquisition and recording for the negative control tube is completed. You are then prompted to load and record the positive control tubes.

12 Load the next control tube displayed in the wizard, then click [Start] on the data acquisition control pane.

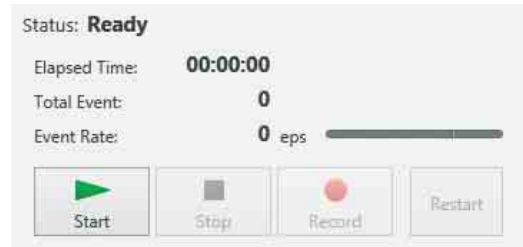

The positive controls must be loaded in the sequence displayed in the Experiment Explorer, as prompted by the [Compensation Wizard].

13 Check that data acquisition has started, then click [Record] in the data acquisition control pane.




Recording stops automatically after satisfying the record stop condition when sufficient data has been recorded. The positive population is displayed on the histogram plot.

Tip

You can stop data acquisition and recording manually using [Stop] if a sufficient number of events have been recorded and the plot view has stabilized.


14 Check the position of the gate on the density plot and the gate for the positive population on the histogram, then click [Next].

Adjust the gate, as required.

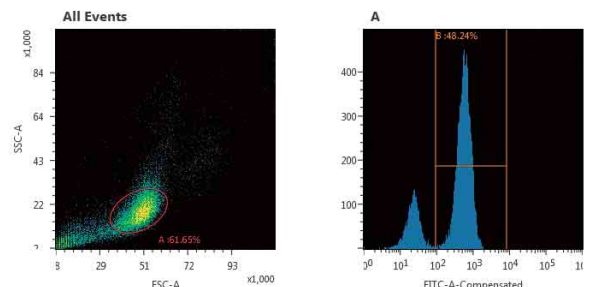


15 Repeat steps 11 to 14 to record each of the control tubes in sequence.

16 After recording all control tubes, click [Calculate Matrix] on the [Compensation] tab of the ribbon.

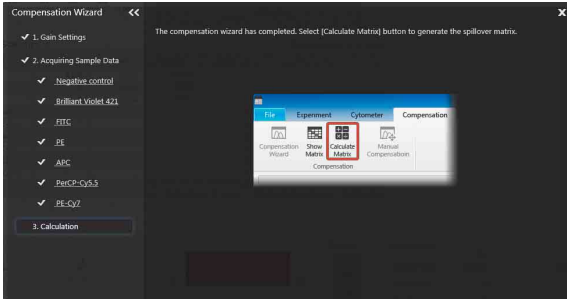

The [Calculate Compensation Settings] dialog appears.


17 Check the target compensation panel, then click [Calculate].

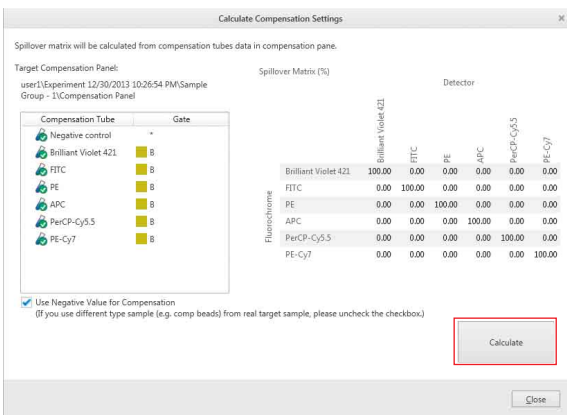

The spillover matrix and the negative fluorescence compensation values are calculated.

18 Check the fluorescence compensation values, then click [Close].

19 Check the compensation matrix values.

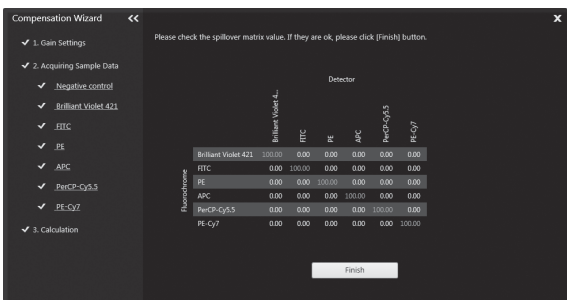

20 Click [Finish] to exit the [Compensation Wizard].

Flushing the Fluidics System

When measuring samples, the sorting chip, sample line, and sample probe should always be flushed with sheath fluid after unloading a sample tube and before loading the next sample tube to eliminate carryover contamination between samples.

1 Check that data acquisition and recording are stopped.

The fluidics system cannot be washed during data acquisition.

2 Click [Probe Wash] in the [Cleaning] group on the [Cytometer] tab of the ribbon.



The sorting chip, sample line, and sample probe are flushed with sheath fluid.

For details about cleaning manually using bleach, see “Cleaning the Sample Fluidics System using Bleach” (page 170).

Shutdown


Shut down the SH800 when finished measurement for the day using the following procedure. The fluidics system should be cleaned according to the instructions in the Shutdown Wizard before shutting down the SH800 main unit.

The Shutdown Wizard describes the bleach cleaning procedure. For details, see “Cleaning the Sample Fluidics System using Bleach” (page 170).

Note

Always leave the sorting chip in the chip loader after shutting down. The chip seals the inlet ports when the SH800 is not in use.

- 1 Click [Unload Collection] to unload the collection stage and remove the collection tubes.

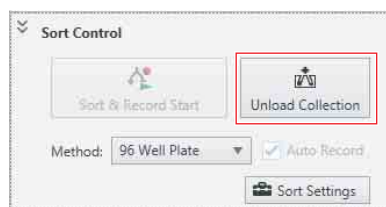


Tip

Remove the splash guard when unloading multi-well plates after sorting is finished.

- 2 Clean the sample loader, collection area, and the collection stage using ethanol to prevent salt precipitation within the unit and to prevent biological contamination.


For details about cleaning, see “General Cleaning” (page 172).

- 3 Click [Software and Hardware] in the [Shutdown] group on the [Cytometer] tab of the ribbon.

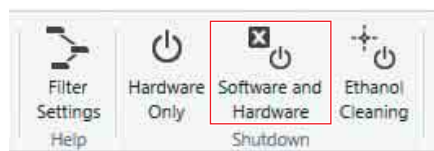

The Shutdown Wizard appears.


- 4 Click [Start].

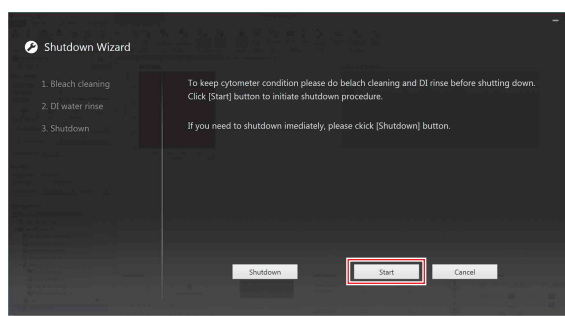

Tip

If you click [Shutdown], the SH800 shuts down without doing any cleaning.

- 5 Select the sodium hypochlorite cleaning sample tube size and click [Start].

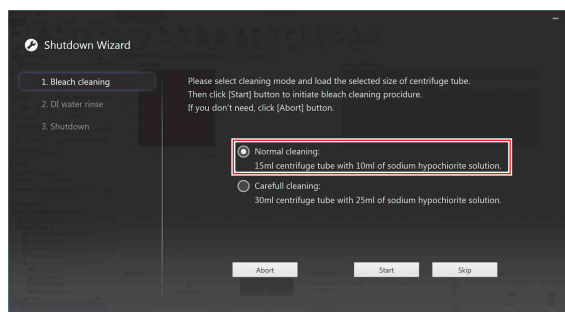


Bleach cleaning starts.

When bleach cleaning finishes, flush the system using DI water.

- 6 Select the DI water cleaning sample tube size and click [Start].

To shutdown the instrument with the sample probe filled with DI water, place a check mark in [Keep DI water inside sample probe].

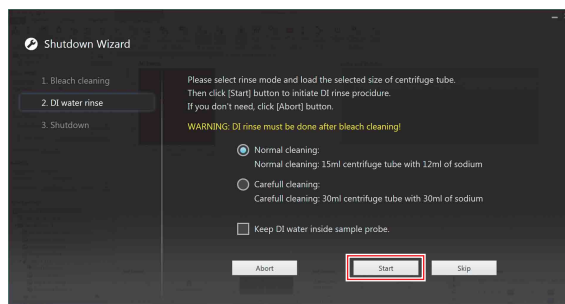


DI water cleaning starts.

When cleaning with DI water finishes, the following screen appears.

Note

After cleaning, the fluidics lines are filled with DI water to prevent the formation of salt deposits. Running data acquisition in this state may damage cells.

7 Click [OK] to shut down the SH800.

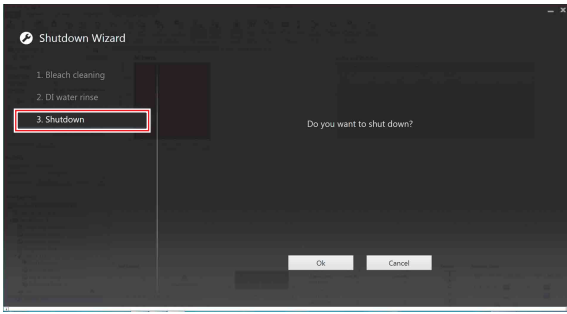


The fluidics system is depressurized and all fluidics lines are sealed.

Specifying User Preferences

Various configuration settings and values can be saved by each user in user preferences. User preferences are loaded automatically each time a user logs in, saving having to reconfigure the instrument each time you want to acquire or analyze data.


1 Click [Information] on the [File] tab of the ribbon, then click [User Preference].

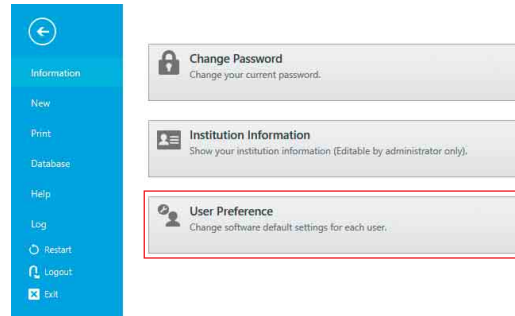

The [User Preference] dialog appears.

2 Set each item.

The values set here become the default values for each setting.

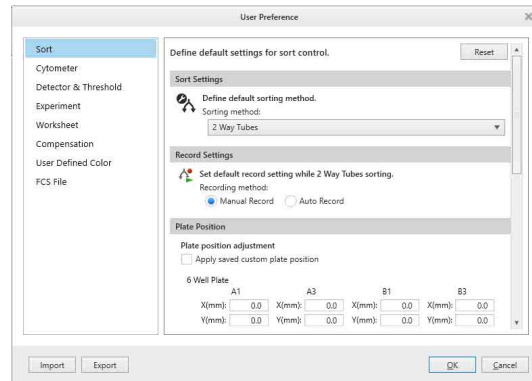

For details about each item, see “[User Preference] dialog” (page 120).

Exchanging the Sorting Chip

The sorting chip must be changed regularly to ensure an unobstructed flow path.


- 1 Click [Chip] in the [Exchange] group on the [Cytometer] tab of the ribbon.

A confirmation dialog appears.

- 2 Click [Yes].


The [Chip Exchange] wizard appears.

- 3 Scan the QR code on the packaging of the replacement chip.

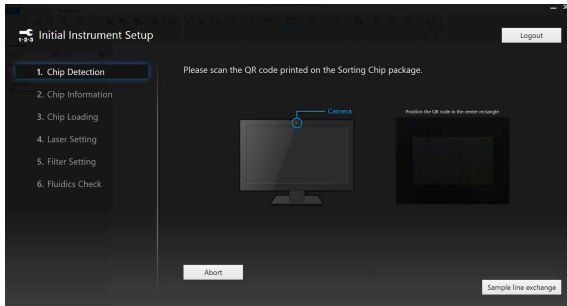


If the scan is successful, the chip information is displayed.

- 4 Click [Next].

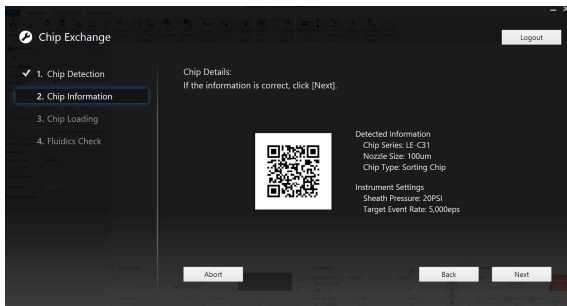


You are prompted to exchange the chip.

- 5 Open the flip-up door, and remove the current sorting chip.
- 6 Insert the replacement chip in the chip insertion slot, then close the flip-up door.

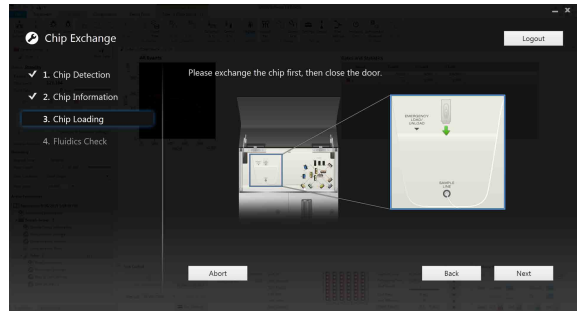

- 7 Click [Next].

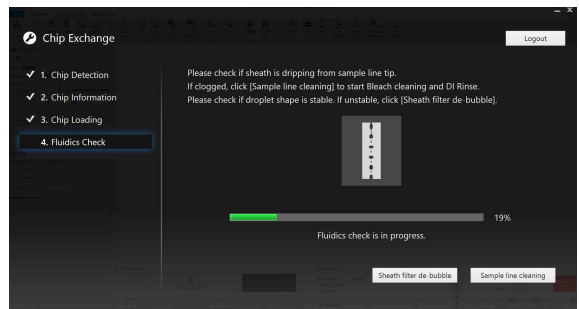


The fluidics check starts.

- 8 Click [Next].

The fluidics check and automatic calibration sequence begins. Follow the on-screen instructions to align the chip and calibrate the instrument.

Notes

- Do not insert any objects other than Sony sorting and cleaning chips designed for the SH800 into the chip insertion slot.
- The packaging for the loaded sorting chip should be stored in a safe location.
- Replace the sorting chip if it is discolored, dirty, or stained from previous use.
- Do not drop the sorting chip on a hard surface or bend the chip intentionally. If there is a problem with the sorting chip, contact your Sony distributor.
- Always wear gloves and other protective clothing, mask, and goggles, when changing the sorting chip.
- If the sorting chip will not load correctly, run the instrument in maintenance mode as described in “*Running Maintenance Mode*” (page 177) to clean the instrument. If the problem persists, contact your Sony distributor for service.
- The sorting chip must be changed regularly if subjected to prolonged exposure to irradiation by the lasers to prevent burn-in, especially by the 405 nm laser. You can monitor the length of time the chip has been exposed to the light from each laser in the [Chip Information] tab of the [Cytometer Settings] dialog, accessed by clicking [Settings] on the [Cytometer] tab of the ribbon.

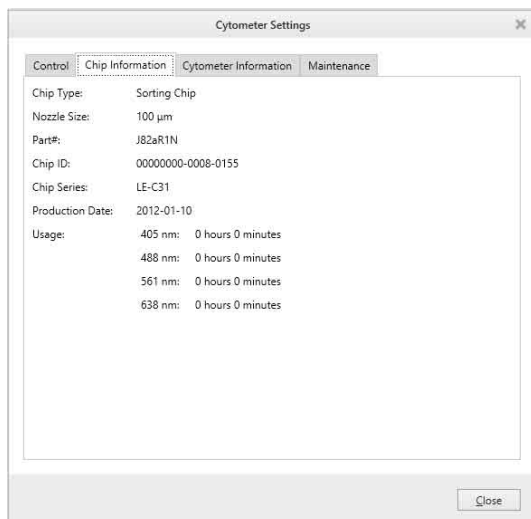

Inheriting Auto Calibration Results Between Users


When using the same sorting chip, the auto calibration results can be inherited between users. This reduces the time taken for auto calibration when another user logs in. Allowing auto calibration results to be passed between users is determined when logging out.

Tips

- The inherited auto calibration results varies depending on the implemented auto calibration sequence used. If chip position auto alignment only is performed, in Analyzer mode, only the chip position alignment result is inherited. If sorting calibration is also performed, then the chip position alignment result and sorting calibration result are inherited.
- An administrator can set whether the inherited auto calibration result is made the default settings when logging out.

For details, see [Administrative Settings] in “[Information] window” (page 119).

- 1 On the logout confirmation screen, place a check mark in the [Keep Auto Calibration result for next user.] checkbox and click [Yes].

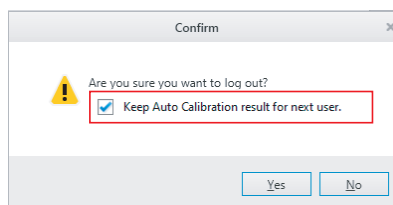


A confirmation message appears.

Notes


- If the level in the sheath tank is too low, droplet formation is unstable, the fluidics system stops due to an error, or other instrument status problem, the auto calibration result may not be inherited by another user, even if auto calibration was completed successfully.
- To enable the auto calibration result to be inherited by another user, run [Bleach Cleaning] and [DI Rinse] to clean the sample line before logging out.

- 2 Check the message details, then click [OK].

If the sample line has not been cleaned, click [Cancel] and then run [Bleach Cleaning] and [DI Rinse].

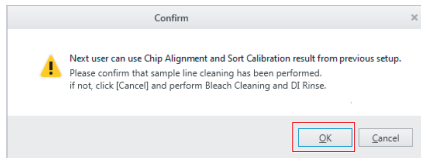

The login screen appears, displaying a timer showing when the auto calibration result will be cleared.

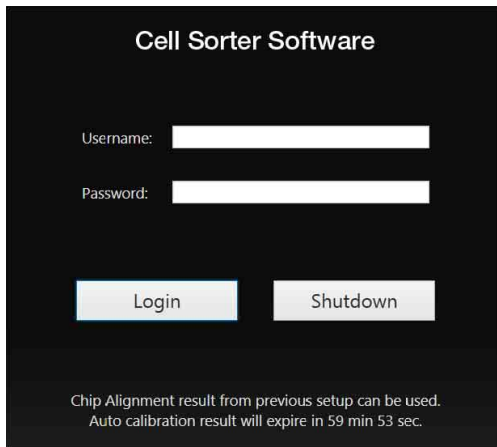


Tips

- When the auto calibration results are inherited, the sheath fluid flow continues.
- The auto calibration results are maintained for approximately 60 minutes.
- The auto calibration results may also be cleared, depending on the instrument status, even when the auto calibration results are inherited.
- If maintenance mode is run while the auto calibration results are inherited, the auto calibration results are cleared.

3 Log in.

The inherited auto calibration results are displayed.


4 Check the settings, then click [OK].

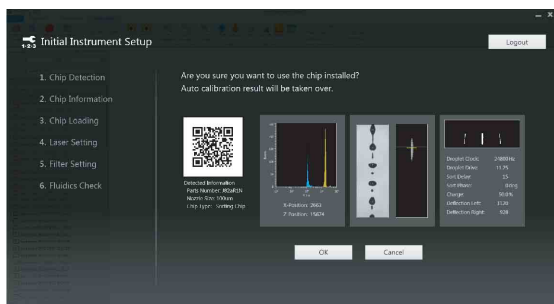

The [Create Experiment] window appears for building a new experiment.

Tips

- If only the chip position alignment result is inherited, the follow screen appears before the [Create Experiment] window appears.

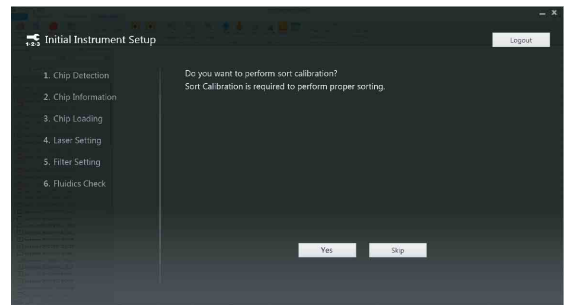

- To perform sorting, click [Yes] to run sorting calibration. After the sorting calibration is completed, the [Create Experiment] window appears.
- If [Skip] is clicked, the [Create Experiment] window appears immediately.

For details about sorting calibration, see “Automatic Calibration” (page 50).

Backing Up/Restoring the Database

You can back up all data in the Cell Sorter Software database. You can also restore all data from a backed up database, as required.


This operation is available only when logged in using an administrator account.

Tip

Backup/restore operations are supported on the NTFS-format file system. Other format file systems are not supported.


Backing up the Database

- 1 Click [Database] on the [File] tab of the ribbon, then click [Backup].

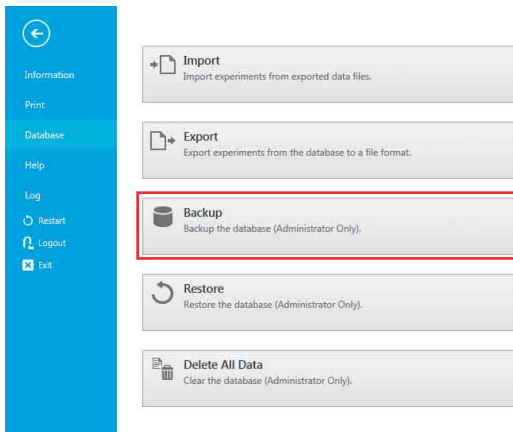

The [Backup Database] dialog appears.

- 2 Click (Browse) to specify the backup data destination, then click [Start].

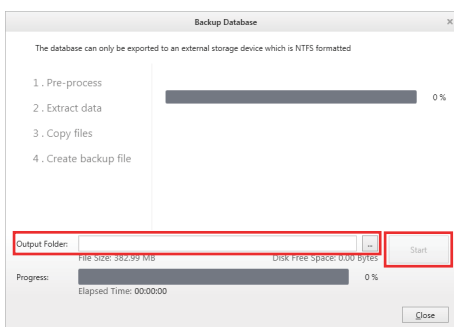

The database backup commences and a progress bar is displayed.

When the database is backed up, a confirmation message appears.


- 3 Click [OK].
- 4 Click [Close] to close the [Backup Database] dialog.

Restoring a Database

Tip


Restoring a previous database backup loads the user accounts, experiment data, templates, and other information in the backup, erasing and replacing all the information in the current database.

- 1 Click [Database] on the [File] tab of the ribbon, then click [Restore].

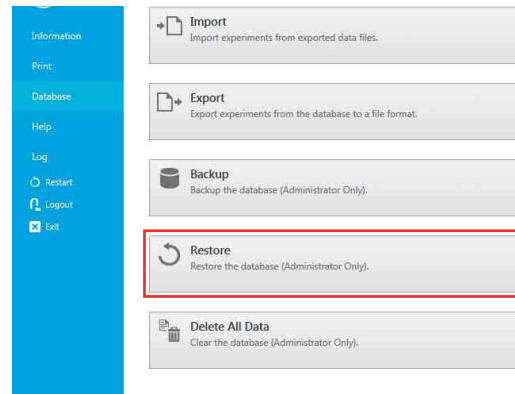

The [Restore Database] dialog appears.

- 2 Click (Browse) to specify the backup data to restore, then click [Restore].

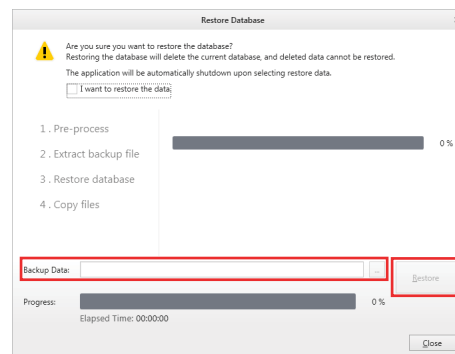

A confirmation message appears.

- 3 Click [Yes].
Cell Sorter Software launches and the database restore operation starts.
When the database is restored, a confirmation message appears.
- 4 Click [OK].
- 5 Click [Close] to close the [Restore Database] dialog.

This chapter describes how to customize experiments, and how to configure experiment components, detector settings, and fluorescence compensation.

For details creating an experiment quickly, see “Creating an Experiment” (page 52).

Creating a Customized Experiment

You can customize the structure of an experiment created using templates or recent experiments by adding sample groups and sample tubes to the experiment, and configuring the settings of each component.


You can also customize an experiment after it has been created. For details, see “Editing an Experiment” (page 81).

- 1 Click [New] on the [File] tab of the ribbon or click [New] in the [Experiment] group on the [Experiment] tab of the ribbon.

The [Create Experiment] window appears.

- 2 Select a template or a recent experiment.

Clicking a template or an experiment on the left displays the structure of the selected experiment on the right side of the window.

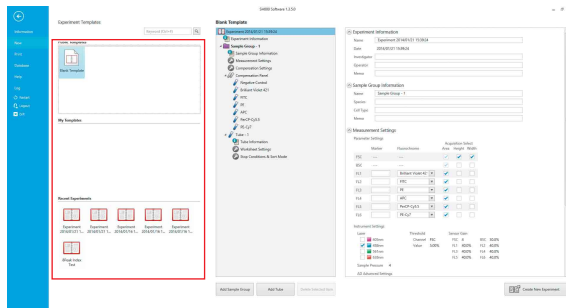

- 3 Modify the structure of the experiment, as desired.



For details about adding/deleting sample groups and sample tubes, see “To add a sample group” (page 78), “To add a sample tube” (page 79), and “To delete a sample group or tube” (page 80).

Tip

Changes made to the settings on the right side of the window are immediately reflected in the experiment structure.

4 Modify the experiment and sample group information, as desired.

Tip

The name of the experiment is supplied automatically, although it is recommended that you change it to a more appropriate name.


5 Select the detector channel pulse parameters for data acquisition in the [Measurement Settings] pane.

For details, see the [Measurement Settings] dialog description in “Configuring an Experiment” (page 80).

Tips


- The fluorochrome and marker names are optional, but recommended.
- Check that the fluorochrome and marker names assigned to detection channels correspond with the laser configuration and optical filter pattern. For details, see “Fluorochrome Detection Matrix” (page 232).

6 When finished building the experiment, click [Create New Experiment] in the bottom right-hand corner of the window.

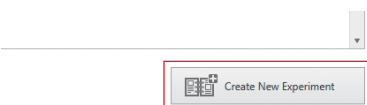

If “Blank Template” was selected, the [New Experiment Startup Procedure] dialog (page 52) appears.


If an existing template or experiment was selected, the main window appears. The created experiment is automatically made the active experiment, and the first tube is automatically assigned as the active tube.

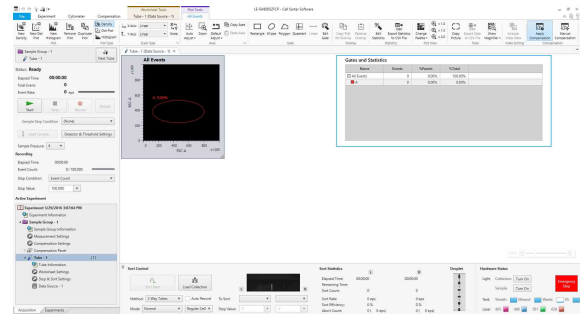


To add a sample group

Use the following procedure to add a sample group to an experiment.


- 1 Click [Add Sample Group] in the [Create Experiment] window.

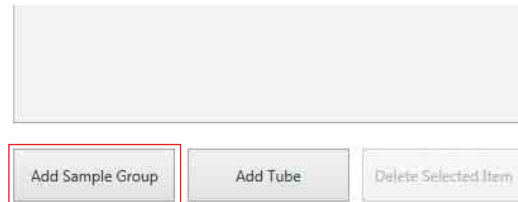

The [Add Sample Group] dialog appears.

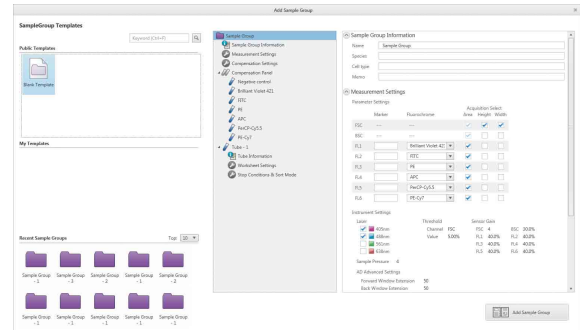

2 Select a sample group template or recent sample group in [Sample Group Templates].

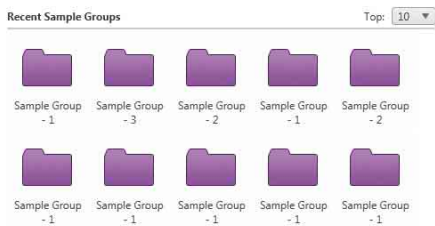


3 Enter a name for the sample group, and modify other settings as required.

The name of the sample group is a required item.



4 Select the detector channel pulse parameters for data acquisition in the [Measurement Settings] pane.

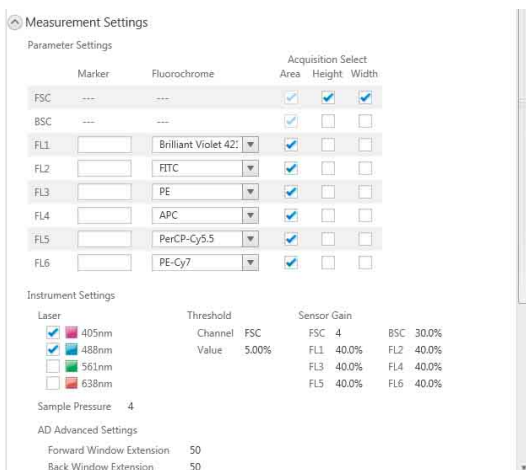

Only control tubes that have a selected detection channel pulse parameter are included in the compensation panel for the sample group in the experiment structure.

Tips

- The fluorochrome and marker names are optional, but recommended.
- Check that the fluorochrome and marker names assigned to detection channels correspond with the laser configuration and optical filter pattern. For details, see “*Fluorochrome Detection Matrix*” (page 232).

5 Modify the structure of the sample group, as required.

For details, see “*To add a sample tube*” (page 79) and “*To delete a sample group or tube*” (page 80).


6 When finished, click [Add Sample Group] in the bottom right-hand corner of the window.

The sample group is added to the experiment.

To add a sample tube

Use the following procedure to add a sample tube to the selected sample group.

1 Click [Add Tube] in the [Create Experiment] window.

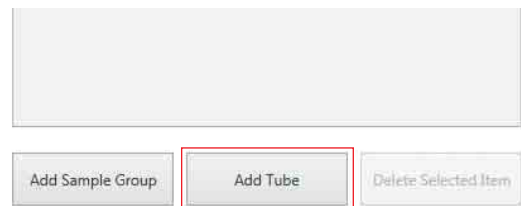

The [Add Tube] dialog appears.

- 2 Select a sample tube template or recent sample tube in [Tube Templates].



- 3 Enter a name for the sample tube, and modify other settings as required.


The name of the sample tube is a required item.



- 4 When finished, click [Add Tube] in the bottom right-hand corner of the window.

To delete a sample group or tube


Select the sample group or tube you want to delete in the structure tree in the [Create Experiment] window and click [Delete Selected Item].

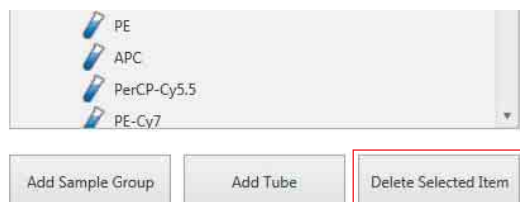

Configuring an Experiment


This describes how to configure each component in an experiment.

- 1 Expand the target experiment in the Experiment Explorer, and double-click [Measurement Settings] in the sample group you want to configure.

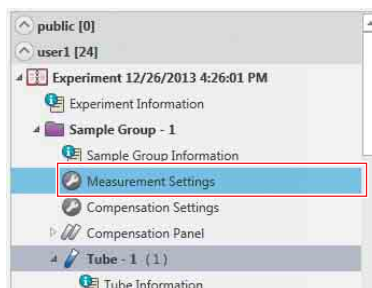

The [Measurement Settings] dialog appears.

- 2 Set each item.

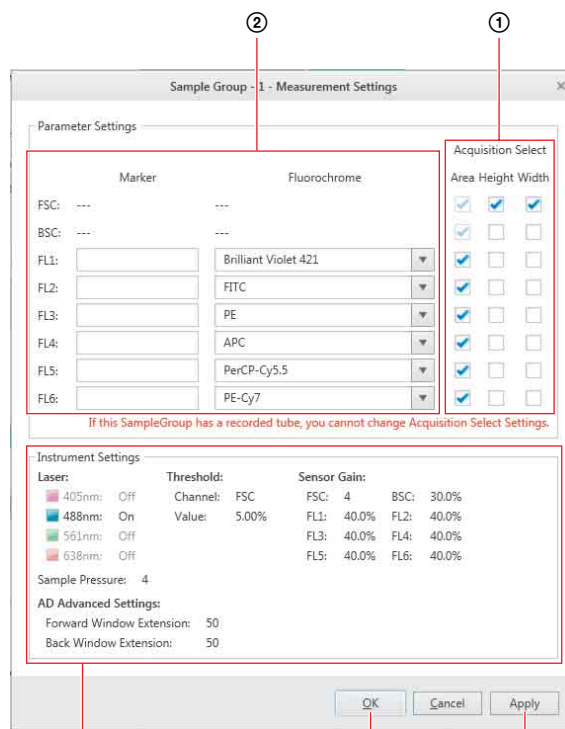

These parameters are configured in the [Detector & Threshold Settings] dialog ([page 86](#)).

- 1 Select the pulse parameters (Area, Height, Width) for each detection channel. The SH800 supports the acquisition of three pulse parameters (area, height, width) in the detector for each of the input channels (FSC, BSC, FL1 to FL6) for each tube in a sample

group. You select the detection channels for which to acquire data by selecting acquisition pulse parameters for those channels. Any combination of pulse parameters can be selected for each detector channel.

Tips

- The Area, Height, and Width pulse parameters are selected by default for the FSC detector channel. Only the Area pulse parameter is selected by default for all other detector channels.
 - The Area pulse parameters for the FSC and BSC detector channels are always enabled. Only detector channels with a pulse parameter selection will appear in the compensation panel.
- ② Specify a marker name and/or fluorochrome name for each fluorescence detector channel. You can also enter a fluorochrome name directly from the keyboard. Marker names and fluorochrome names, if specified, appear in the axis labels on plots, and fluorochrome names appear for control tubes in the compensation panel.

Tips

- Forward scatter (FSC) and back scatter (BSC) are measured using the incident laser light, not fluorescent light. Fluorochrome names and marker names cannot be entered for FSC and BSC.
 - Fluorochrome names and marker names can be modified before, during, or after data acquisition.
 - Marker names and fluorochrome names are optional, but recommended. If names are not specified, the fluorescence channel numbers (FL1 to FL6) will be displayed as the axis labels on plots by default, providing limited feedback as to the type of data displayed on plots.
- ③ Click [Apply] or [OK] to apply the measurement settings. Clicking [OK] applies the settings and closes the dialog.

Editing an Experiment

You can modify an existing experiment, as required. You can also edit an experiment using the [Experiment] tab of the ribbon.


For details about tools available on the [Experiment] tab, see “[Experiment] Tab (ribbon)” (page 130).

Editing a Sample Group

You can add, delete, and perform other edit functions on a sample group in an experiment.

Edit operations are performed using buttons in the [Sample Group] group on the [Experiment] tab of the ribbon.

You can also perform the same operations by right-clicking a sample group in the Experiment Explorer.

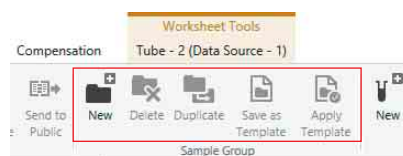


For details about each button, see “[Sample Group] group” (page 130) on the [Experiment] tab of the ribbon.

Editing a Sample Tube

You can add, delete, and perform other edit functions on a sample tube in an experiment.

Edit operations are performed using buttons in [Tube] on the [Experiment] tab of the ribbon.


You can also perform the same operations by right-clicking a sample tube in the Experiment Explorer.

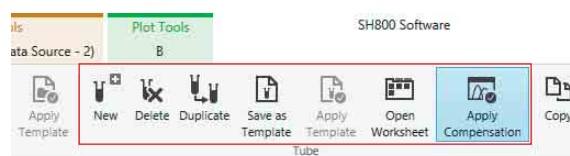

For details about each button, see “[Tube] group” (page 131) on the [Experiment] tab of the ribbon.

Changing Component Settings


- 1 Expand the target experiment in the Experiment Explorer, and double-click the component you want to configure.

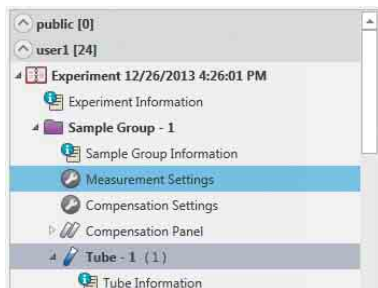

The corresponding dialog for the selected component appears.

- 2 Specify each parameter, then click [Apply] or [OK].

Example: If [Sample Group Information] is selected

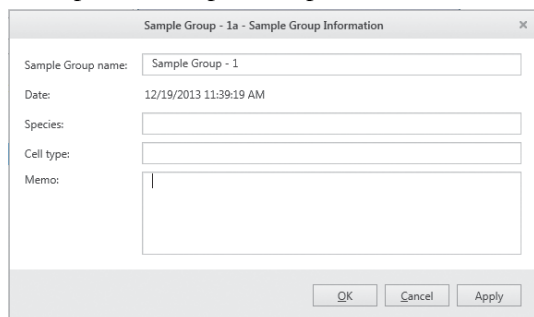

For details about items displayed in the dialogs, see “Experiment Explorer” (page 151).

For details about [Measurement Settings], see “Configuring an Experiment” (page 80).

Exporting/Importing an Experiment


You can import and export experiment data to back up the data, to reopen a previous experiment, or to review the experiment on another SH800 system.

Exporting an Experiment

When data acquisition for an experiment is finished, the data for the experiment can be exported to an external file.

- 1 Click [Database] on the [File] tab of the ribbon, then click [Export].

The [Export Experiment Data] dialog appears.

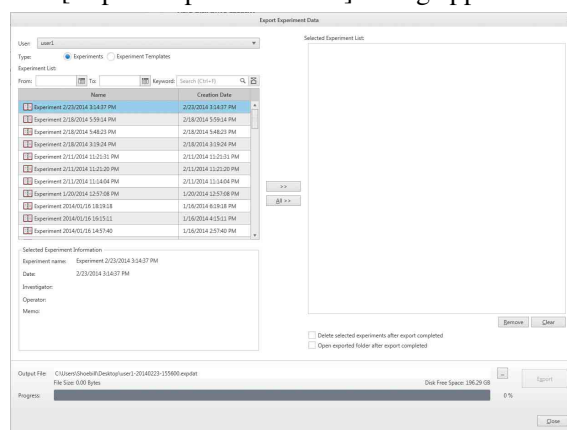

- 2 Select the owner of the data in [User], and select to export experiments or templates using the [Type] radio buttons.
- 3 Select an experiment you want to export in [Experiment List], and click [>>] to move it to [Selected Experiment List].

Clicking [All >>] moves all the experiments to [Selected Experiment List].

Tips

- To remove the exported experiments from the Experiment Explorer after the export is finished, place a check mark in [Delete selected experiments after export completed].
- To open the export destination folder after the export is finished, place a check mark in [Open exported folder after export completed].

- 4 Click (Browse) to specify the destination folder in which the experiment data is to be exported.

- 5 Click [Export].

The export status is displayed in the progress bar. A confirmation dialog appears when exporting is completed.

- 6 Click [OK].


- 7 Click [Close] to close the dialog.


Importing an Experiment

You can import previously exported experiment data into the database.

- 1 Click [Database] on the [File] tab of the ribbon, then click [Import].

The [Import Experiment Data] dialog appears.

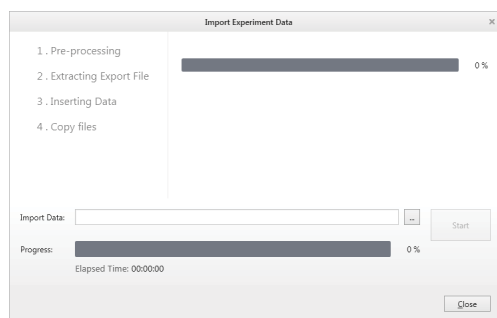

- 2 Click  (Browse) to specify the folder where the experiment data you want to import (.expdat file) is stored.

- 3 Click [Start].

The import status is displayed in the progress bar. A confirmation dialog appears when importing is completed.

- 4 Click [OK].

- 5 Click [Close] to close the dialog.


Sharing Experiments

You can share experiments with other users. Shared experiments allow multiple users to work with the experiment data on worksheets.


To share an experiment

- 1 Select the experiment you want to share in the Experiment Explorer, and click [Send to Public] on the [Experiment] tab of the ribbon.

Alternatively, you can right-click an experiment and select [Send to Public] from the context menu.

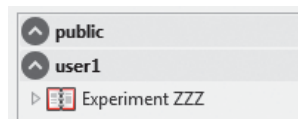

The shared experiment is duplicated in the [public] node in the Experiment Explorer.

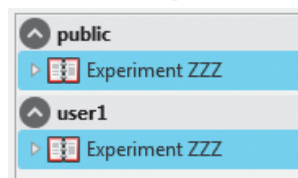


Tip

Other users can access the shared experiment under the [public] node. After an experiment is shared, the shared experiment and the original private experiment exist as independent experiments, with no interaction between the two.

Searching Experiments


You can search experiments by date and keyword.

- 1 Enter the search criteria in the Experiment Explorer of the [Experiments] tab on the left side of the main window.

Clicking  displays a calendar for specifying a date.



- 2 Click  to set optional addition search options.

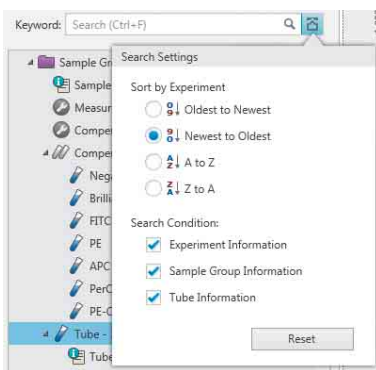

[Oldest to Newest]: Displays the search results in chronological order from oldest to newest.

[Newest to Oldest]: Displays the search results in chronological order from newest to oldest.

[A to Z]: Displays the search results in ascending alphabetical order.

[Z to A]: Displays the search results in descending alphabetical order.

[Search condition]: Selects the experiment components to search for the entered text keyword.

[Reset]: Resets the search options to their defaults.

Experiments matching the search criteria are displayed in the Experiment Explorer.


Saving Tube Data as FCS Files

You can export data recorded for tubes in flow cytometry industry standard FCS 3.0 or FCS 3.1 format for use in third-party applications.

Note

Cell Sorter Software does not support the importing of FCS format files.


- 1 Select the target tube in the Experiment Explorer on the [Experiments] tab on the left of the main screen.
- 2 Click [Export to FCS file] in the [Export] group on the [Experiment] tab of the ribbon.

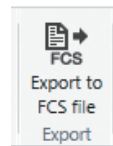

The [FCS File Export] dialog appears.

- 3 Place check marks in [Select target tubes] for tubes whose data is to be exported, specify the FCS format and export destination folder, then click [Export].

You can also select the required data source if there is more than one data source for a single tube.

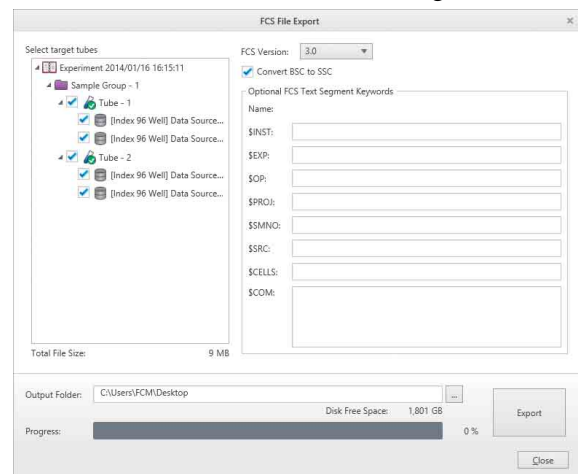

The tube data is exported.

The name of the exported file is the same as the name of the tube, with an “.fcs” filename extension.

Tips


- [Convert BSC to SSC]: Select to convert the data name of back scatter data from BSC (back scatter) to SSC (side scatter) when creating FCS files.
- The required FCS text segment keywords are exported automatically. The following optional FCS text segment keywords can also be specified.

Keyword	Description
\$INST	Institution at which data was acquired
\$EXP	Name of investigator initiating the experiment
\$OP	Name of flow cytometer operator
\$PROJ	Name of the experiment project
\$SMNO	Specimen (e.g., tube) label
\$SRC	Source of the specimen (patient name, cell type)
\$CELLS	Description of objects measured
\$COM	Comment

Saving an Experiment as a Template


You can save an experiment, sample group, or sample tube as a shared template (Public Templates) or a private template (My Templates.)

- 1 Select the target experiment in the Experiment Explorer on the [Experiments] tab on the left of the main screen.
- 2 Click [Save as Template] in the [Experiment] group on the [Experiment] tab of the ribbon.

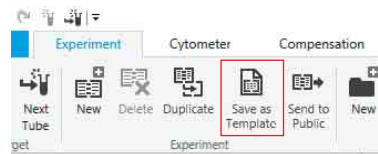

The [Save as Experiment Template] dialog appears.

- 3 Select the template type, enter a name, and then click [OK].

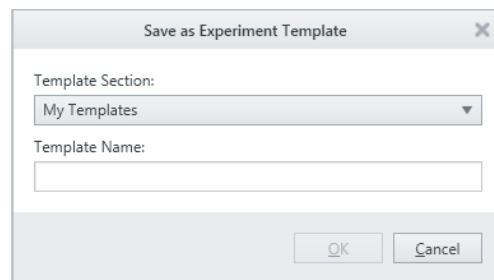

The experiment is saved as a template.

Tip

Overlay displays are not saved as templates.

Configuring Detector Settings

The following instrument settings must be optimized in an experiment to acquire accurate data.

- Trigger threshold channel
- Trigger threshold value
- Sensor gain of the forward scatter, back scatter, and fluorescence detectors

The instrument settings are saved with the sample group in an experiment. Duplicating an experiment or creating a new experiment using templates will load the saved instrument settings into the new experiment. Running an experiment using the same instrument settings as a previous experiment enables direct comparisons to be made between the results obtained, even if the experiments were conducted at different times.

Note

The gain parameter values directly affect the fluorescence compensation calculation of the spillover matrix. Do not change the gain of the channel detectors after or while recording the control tubes in a compensation panel. The fluorescence compensation process is valid only for the instrument settings in place when all control tubes in a compensation panel are recorded. If the gain of a detector for any channel is changed for any reason after recording the control tubes, the fluorescence compensation will no longer be valid and the control tubes will need to be recorded again.


For details about recording control tubes in a compensation panel, see “Performing Fluorescence Compensation” (page 67).

Basic Configuration Method

You can also configure many of the same parameters using the [Compensation Wizard] and on the [Control] tab of the [Cytometer Settings] dialog.


For details, see “Performing Fluorescence Compensation” (page 67) and “[Cytometer Settings] dialog” (page 138).

- 1 Click [Detector & Threshold Settings] in the [Acquisition] tab of the Experiment/Acquisition control pane.

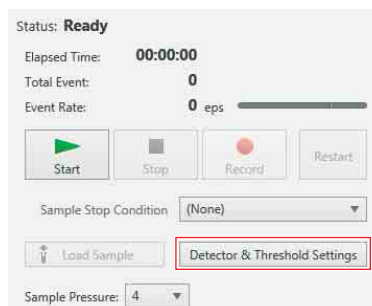

The [Detector & Threshold Settings] dialog appears.

- 2 Set each item.

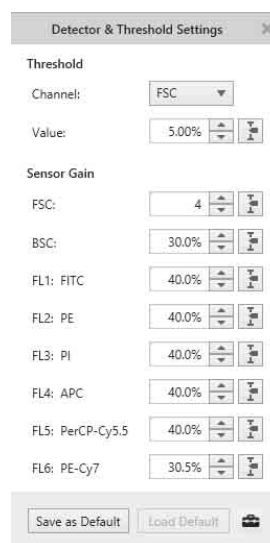

[Threshold]

Specifies the threshold for event detection.

[Channel]: Selects the channel (FSC, BSC, FL1 to FL6) used to trigger the detection of events. The default trigger channel is FSC.

[Value]: Sets the threshold value for the trigger channel as a percentage of the output signal of the detector for the trigger channel. The default value is 5.00%.

[Sensor Gain]

Specifies the sensor gain of the detectors.

[FSC]: Specifies the gain of the forward-scatter (FSC) detector. The gain setting has a 16-step range.

[BSC]: Specifies the gain of the back scatter (BSC) detector as a percentage of the output signal of the detector. The default value is 30.0%.


[FL1] to [FL6]: Specifies the gain of the corresponding fluorescence channel detector as a percentage of the output signal of the detector. The default value is 40.0%.

[Default Settings]

[Save as Default]: Saves the current settings in the [Detector & Threshold Settings] dialog as the default values.

The default values are used when creating an experiment or sample group using the Blank Template. The default values can also be configured in the [User Preference] dialog (page 120).

[Load Default]: Overwrites the current settings in the [Detector & Threshold Settings] dialog using the default values saved using the [User Preference] dialog. After loading, you can change the default values using [Save as Default]. [Load Default] is disabled if the default values have not been changed.

 **(Set):** Clicking this button displays the [User Preference] dialog (page 120) for setting user default values for the [Detector & Threshold Settings] dialog.

Tips

- Only the gain of fluorescence channels that have pulse parameters selected for acquisition can be adjusted. The gain settings of other channels are disabled.
- If the gain is too high, the detector output circuit may become saturated, providing incorrect signal output. In this case, reduce the gain and then readjust the instrument settings to find the optimum operating point.

3 Click [X] to close the dialog.

The settings are saved.

Loading Reference Data Method

You can load data acquired on another instrument as reference data in order to obtain equivalent measurement results.

Note

If the instrument is configured using this method, use the same kind of sample as the sample used to obtain the reference data. Note that because the detector characteristics will vary from instrument to instrument, data identical to the reference instrument data cannot be obtained.


Creating a gain adjustment template

You can create a gain adjustment template (template for instrument settings) for an experiment that includes data used as reference data. The experiment template contains a gain adjustment tube which includes the reference data.

1 In the Experiment Explorer on the [Experiments] tab, right-click a tube with acquired data to use as the reference, and select [Export Adjustment Template] from the context menu.

The [Export Adjustment Template] dialog appears.

2 Enter a name for the template in [Template Name], enter the folder path and file name in [Export File Path], and click [Export].

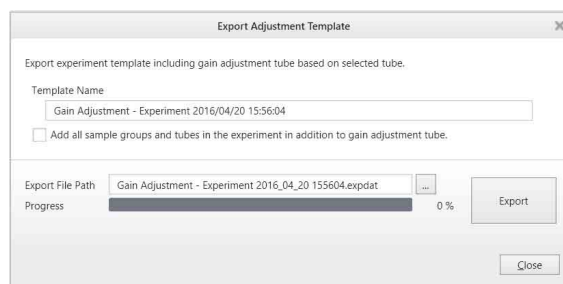

An experiment template file containing the reference data is exported.

Creating an experiment

You can create an experiment and apply instrument settings using an experiment template with gain adjustment tube.


Use the following procedure to configure the instrument.

1 Import an experiment adjustment template file containing reference data.

The imported template is added to [My Templates].

For details about importing a template, see “Importing an Experiment” (page 83).

2 Create a new experiment using the imported experiment adjustment template.

The gain adjustment tube is indicated by the  icon.

For details about creating an experiment, see “Creating an Experiment” (page 52).

3 Assign the gain adjustment tube as the active tube and perform data acquisition.

For details, see “Acquiring Data” (page 53).

4 Configure the detectors based on the reference data.

Configure the detector settings, including gain adjustment and threshold value adjustment. Two types of data, the reference data and acquired data are displayed on a plot for the gain adjustment tube.

Adjust the detectors until the acquired data matches the reference data.

For details about detector settings, see “Basic Configuration Method” (page 86).

5 Update the detector default values as required.

Click [Save] in the [Detector & Threshold Settings] dialog to save the current settings as the default values.

Tip


There are various other methods that can be used to optimize the detector gain settings for an experiment. For example, a more rigorous approach to detector gain optimization might be to measure the coefficient of variation (CV) for positive controls at various detector gain levels to determine the minimum gain levels for which the CV plateaus off in each detection channel sensor. A detector gain set at this minimum gain level will provide the optimal signal-to-noise ratio required for maximum sensitivity and maximum separation between negative and positive peaks.

Adjusting Fluorescence Compensation


Using a Target Gate in the Gate Hierarchy

You can specify a target gate within the gate hierarchy to be used as the negative control region or as a singly-stained positive control region for calculating the compensation matrix.


- 1 Right-click a gate on a plot for a control tube in the compensation panel to set the gate as the calculation target for the compensation matrix, and select [Set As Calculation Target] from the context menu.

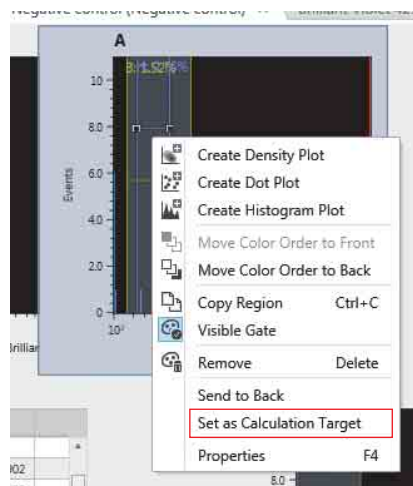

The selected gate is highlighted.

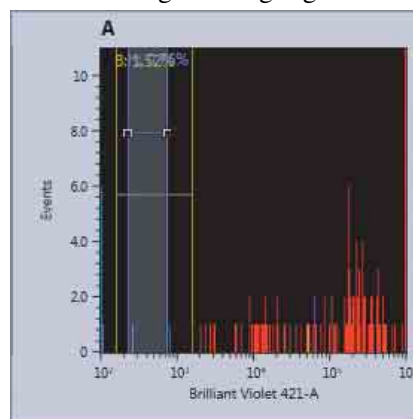


- Click [Calculate Matrix] on the [Compensation] tab of the ribbon.

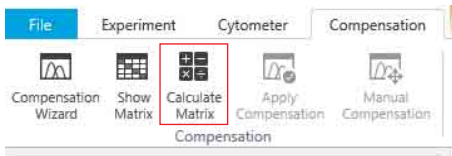

The [Calculate Compensation Settings] dialog appears.

- Check that the specified gate is the target of the calculation, and click [Calculate] to calculate the spillover matrix.

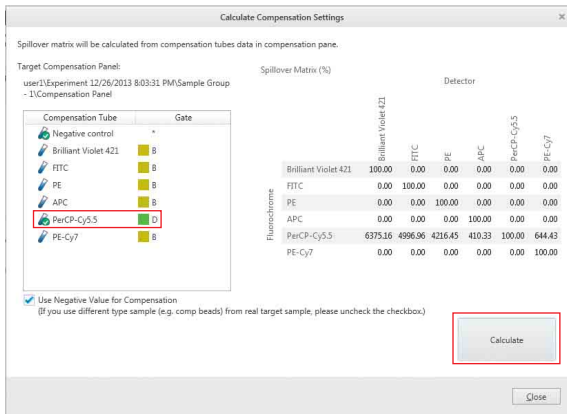

- Click [Close] to close the dialog.

Adjusting Fluorescence Compensation Manually

You can modify the compensation settings after using the [Compensation Wizard] or you can enter fluorescence compensation values manually to edit the spillover matrix directly.

Tips

- You can save the current fluorescence compensation settings using the [Compensation Settings] dialog.

For details about displaying the [Compensation Settings] dialog, see “To edit the spillover matrix directly” (page 90).


- Adjusting the position of events shown on a plot also adjusts the position of all events on other plots. Accordingly, you should add plots for all fluorescence parameters to the worksheet before adjusting compensation manually in order to monitor the affect of adjusting the compensation on each plot.

To adjust compensation graphically on plots

You can adjust fluorescence compensation graphically on the worksheet using a mouse.


- Display the worksheet for the tube whose compensation you want to modify.
- Enable [Apply Compensation] on the [Compensation] tab of the ribbon.
- Click [Manual Compensation] on the [Compensation] tab of the ribbon to activate manual compensation mode.

A diagonal line appears on plots that have fluorescence parameters on the X and Y axes.

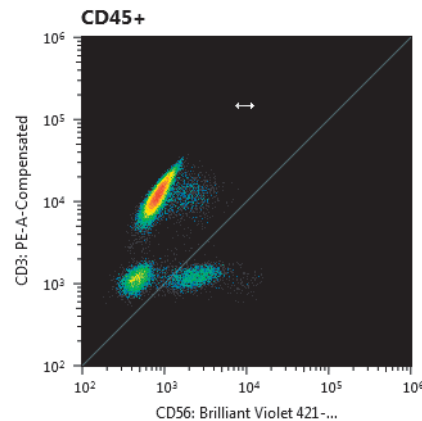

- Place the mouse cursor over the plot, and drag the plot to change compensation for the target parameter.

The mouse cursor changes to a horizontal double-ended arrow cursor in the left half of the plot, and a vertical cursor in the right half of the plot. The displayed events move on the plot to reflect the changed compensation.

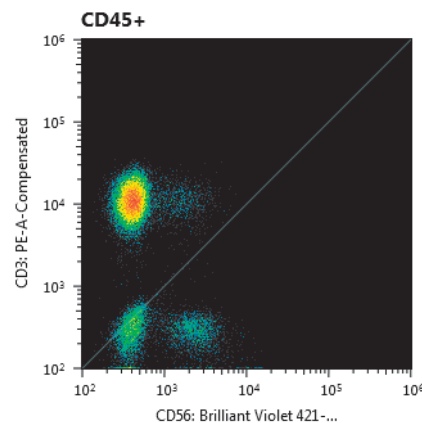

- When finished, click [Manual Compensation] again to deactivate manual compensation mode.

You can also deactivate manual compensation mode by selecting another worksheet to display.

Tips

- Adjusting the position of events shown on a plot also adjusts the position of all events on other plots. Accordingly, you should add plots for all fluorescence parameters to the worksheet before adjusting compensation manually in order to monitor the affect of adjusting the compensation on each plot.
- The actual compensation values in the spillover matrix can be monitored simultaneously in the [Compensation Settings] dialog. It is recommended that you save the compensation settings from the [Compensation Settings] dialog before adjusting compensation manually. You can also click [Reset] in the [Compensation Settings] dialog to reset the compensation settings.

To edit the spillover matrix directly

You can adjust the fluorescence compensation manually in the [Compensation Settings] dialog. You can adjust the values using visual feedback alone, or enter spillover matrix values directly to reproduce the compensation settings of another experiment.

Note

Editing the spillover matrix directly to adjust compensation is inherently difficult. Exercise care when editing the spillover matrix directly.


- 1 Display the worksheet for the tube whose compensation you want to modify.
- 2 Click [Apply Compensation] on the [Experiment] tab of the ribbon.

The fluorescence compensation obtained using the [Compensation Wizard] is applied to the plots on the worksheet.
- 3 Display the [Compensation Settings] dialog by doing one of the following.
 - Select [Compensation Settings] in the Experiment Explorer and click [Show Settings] on the [Experiment] tab of the ribbon.
 - Double-click [Compensation Settings] in the Experiment Explorer.
 - Select any component within the sample group and click [Show Matrix] on the [Compensation] tab of the ribbon.

The [Compensation Settings] dialog appears.

- 4 Click a cell in the matrix to select its value.

You can adjust a value using one of the following methods.

- Turn the mouse scroll wheel button.
- Enter a value directly in the spillover matrix.
- Click the  spin boxes.
- Drag the slider.



The plots on the worksheet are updated in real-time.

Tips

- The number of fluorochromes in the experiment determines the dimension of the spillover matrix. Spillover elements can be modified, but they cannot be added nor deleted.
- To save the current compensation settings, click [Save]. To load compensation settings for the sample group, click [Load].
- To reset the compensation settings, click [Reset].

- 5 Click [Close] to close the dialog.

Data is analyzed using various tools available on the worksheet.


This chapter describes gate settings used for analysis of the acquired data and for sorting, generating reports of data analysis, and related tasks.

For details about content displayed on the main window, see “Main Window” (page 130).

Adding a Plot

You can add plots to the worksheet.

- 1 Click a button for a new plot of the desired plot type to be added in the [Plot] group on the [Worksheet Tools] tab or the [Plot Tools] tab of the ribbon.

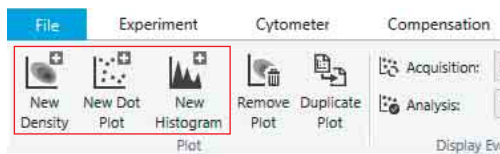

The plot is added to the worksheet. The new plot created represents the [All Events] population.

- 2 Change the layout of plots, as required.

You can change the position of plots by selecting the target plot and dragging it to a new location.



Edit the plots, add gates, and perform other actions on plots, as required.

See “Editing Plots” (page 92) and “Adding Gates” (page 97).

Editing Plots

Cell Sorter Software displays data on the following types of plots. You can change the plot type, change the axis scale, and modify other aspects of plots.


For details about the buttons on the [Plot Tools] tab of the ribbon, see “[Plot Tools] Tab (ribbon)” (page 146).

For details about adding gates to a plot, see “Adding Gates” (page 97).

Density plots

Density plots display a frequency distribution of fluorescence intensity for two acquisition parameters (bivariate distribution).



Density plots use color to show the distribution of the number of detected events having the same value. High-density regions are shown in red, low-density regions are shown in violet, and regions of intermediate density are shown using the colors of the visible spectrum in between. Density plots are most useful for displaying populations where the number of detected events is relatively large.

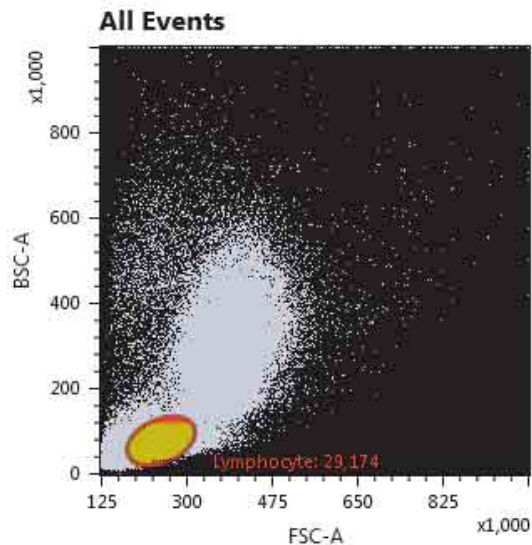

Histogram plots

Histogram plots display a frequency distribution of fluorescence intensity for a single acquisition parameter (univariate distribution).

Histogram plots are effective for identifying the peaks and troughs in measurement data.

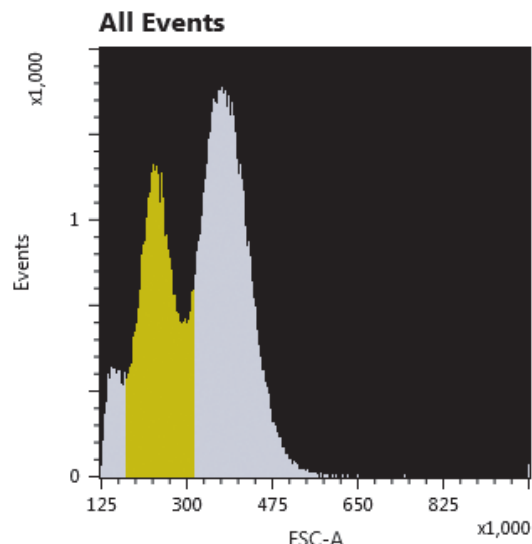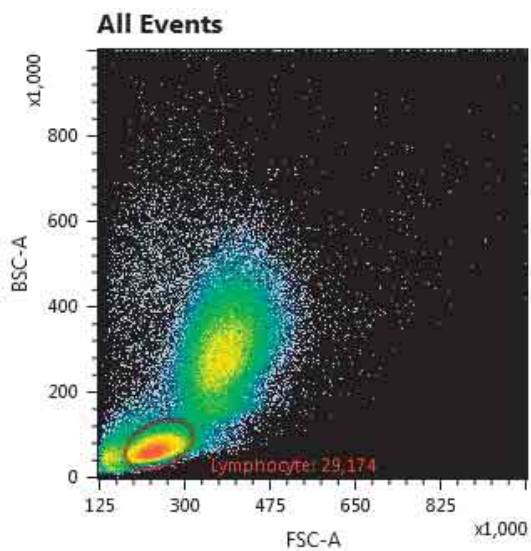

Dot plots

Dot plots display a frequency distribution of fluorescence intensity for two acquisition parameters (bivariate distribution).


Dot plots are most useful for displaying populations where the number of detected events is relatively small. Events on a dot plot are colored based on their classification (gate and population).

Tip

As the number of events on a dot plot increases, the plot may become saturated making it difficult to discern patterns. In this case, reduce the number of events displayed or use a density plot to view the data.

Changing Plot Type


- 1 Select a plot whose type you want to change on the worksheet.

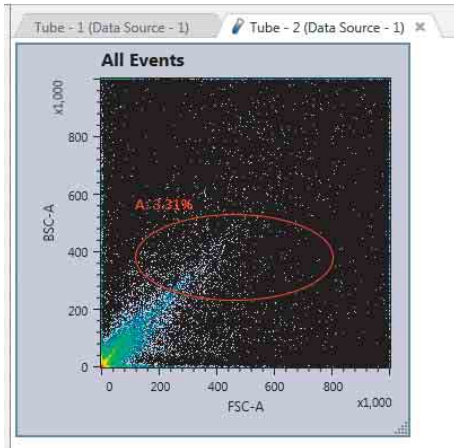


- 2 Click the target plot type in the [Plot Type] group on the [Plot Tools] tab of the ribbon.

You can also right-click a plot and select a plot type from the [Plot Type] submenu of the context menu.

For example, the diagram below shows a density plot being changed to a histogram plot.

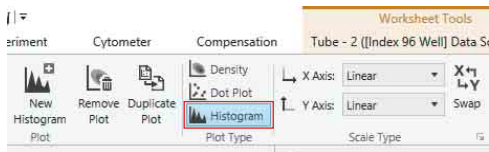

The type of the selected plot is changed.

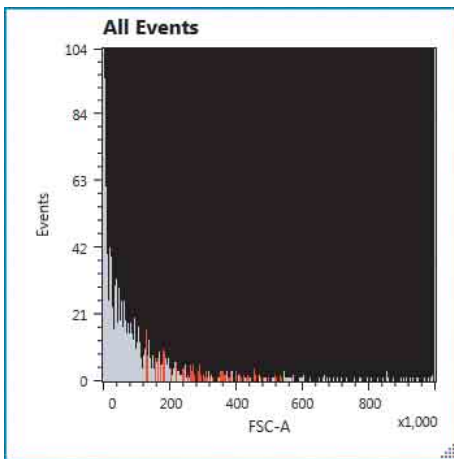

Tip


The plot type of an overlay display plot cannot be changed. To change the plot type, first remove the overlay.

For details about operation, see “Deleting Overlay Layers” (page 103).


Changing the Histogram Display View

You can select the following display views for histogram plots.


Filled area view

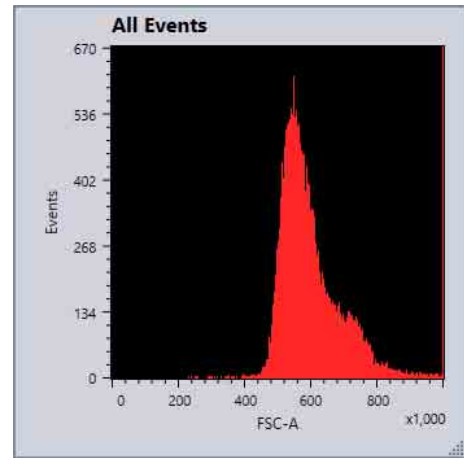

Outline view

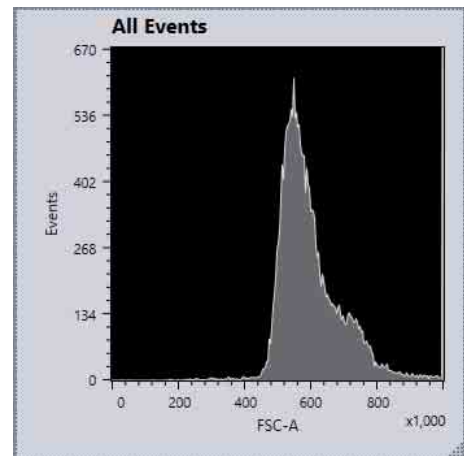


Select a histogram plot on the worksheet, and click [Histogram] repeatedly in the [Plot Type] group on the ribbon to toggle between filled area view and outline view.

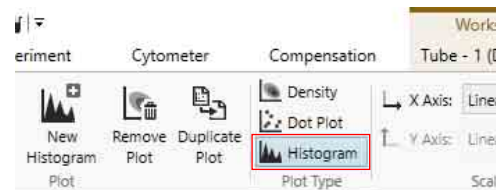

About color display of events


In filled area view

Events included within are gate are displayed in the specified colors.

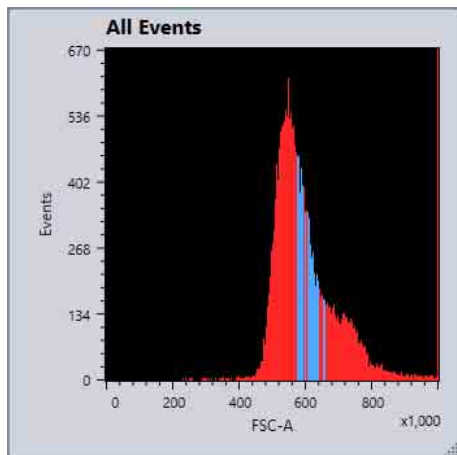

Outline view

The colors of each gate are not displayed.

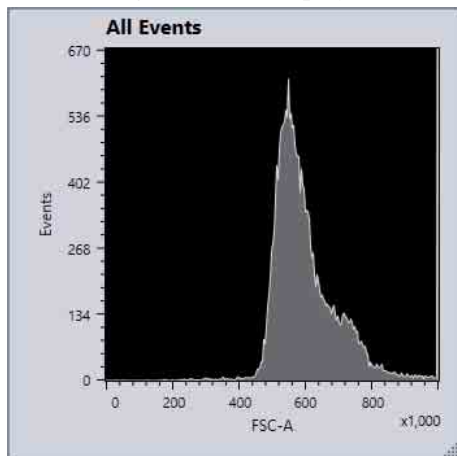


For details about each button, see “[Plot Tools] Tab (ribbon)” (page 146).

Removing a Plot


Select a plot to be deleted on the worksheet, then click [Remove Plot] in the [Plot] group on the [Plot Tools] tab of the ribbon.

Changing Axis Scale Type

Select a plot on the worksheet, then click the desired scale type for each axis in the [Scale Type] group on the [Plot Tools] tab of the ribbon.

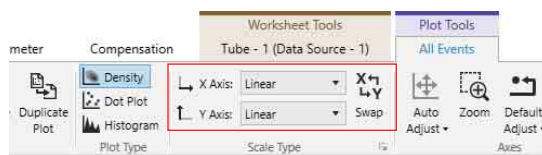


You can also change the scale by right-clicking the plot and selecting from the context menu.

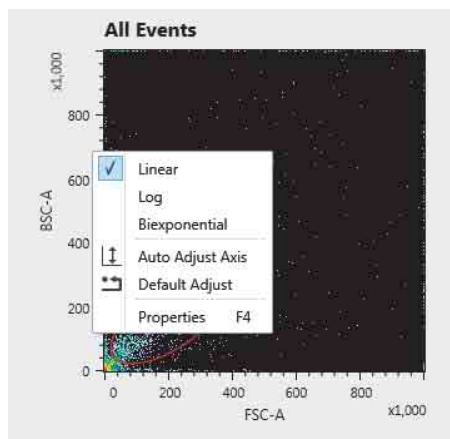

Duplicating a Plot

- 1 Select a plot to be duplicated on the worksheet, then click [Duplicate Plot] in the [Edit] group on the [Plot Tools] tab of the ribbon.

You can perform the same operation from the context menu by right-clicking the plot.

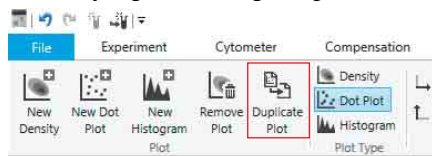

The selected plot is duplicated.

- 2 Edit the duplicated plot, as required, using the buttons on the [Plot Tools] tab of the ribbon.

You can perform the same operation from the context menu by right-clicking the plot.

For details about each item, see “[Scale Type] group” (page 146).


You can select the following scale types.

Tip

The axis scale type is applied only to the display of data on plots. Internally, all data is stored as linear data.


Linear

Linear scales are always used for the display of forward scatter and back scatter.

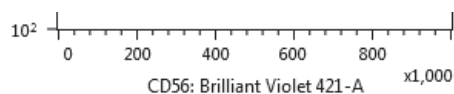


Log (logarithmic)

Logarithmic (log) scales are useful for displaying cells that exhibit fluorescence with intensities across a wide dynamic range. A logarithmic scale helps to make individual peaks easier to visualize. However, log scales cannot adequately represent data with a low mean and high variance.

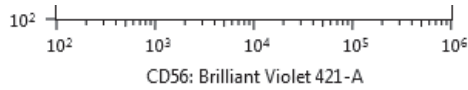

Biexponential


Biexponential scales combine a linear scale at the lower end of the axis (around zero) and a logarithmic scale at the higher end of the axis, with an algorithmic transition between the two scales. A biexponential scale makes it easier to analyze events close to zero or below zero that are difficult to observe on a logarithmic scale.

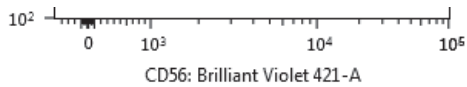

A common problem occurs when displaying fluorescence data on a logarithmic scale after applying fluorescence compensation. Fluorescence compensation introduces variations in measurement data due to variations in detector gain, background noise, and other factors. The result is that some data values can become close to zero or even negative. When values close to zero are plotted on a log scale, the values are wedged up against the axis making them difficult to discern. In addition, values below zero cannot be displayed at all. This can lead to these events, which are real events, being excluded from gates or other analysis. It can also lead to inaccurate fluorescence compensation when attempting to adjust compensation manually using visual cues. In these cases, a biexponential scale can be used to avoid some of the pitfalls of logarithmic scales, while maintaining the overall utility of logarithmic scales.

Changing the Axis Parameters

Click the axis title and select the desired parameter from the popup menu.

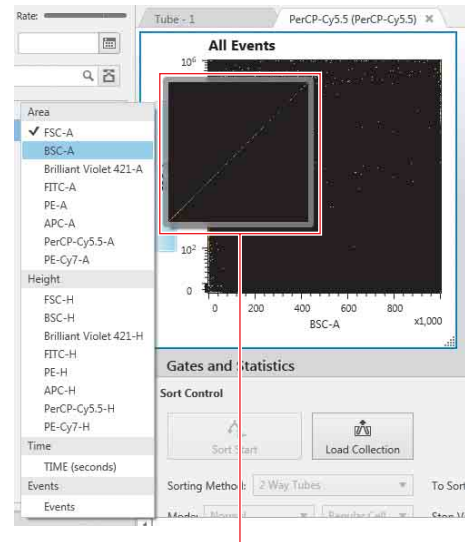


A preview of the plot appears as the mouse hovers over each

Adjusting Scale Ranges

To adjust a scale range automatically

Select a plot, then click [Auto Adjust] in the [Axes] group on the [Plot Tools] tab of the [Plot Tools] ribbon, and select the auto adjustment method.

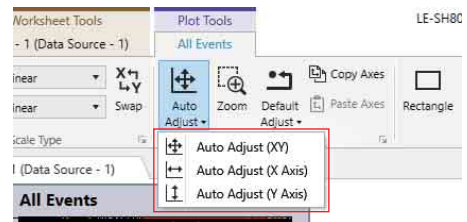


[Auto Adjust (XY)]: Automatically adjusts the scale of the X axis and Y axis.


[Auto Adjust (X Axis)]: Automatically adjusts the scale of the X axis only.

[Auto Adjust (Y Axis)]: Automatically adjusts the scale of the Y axis only.

To specify scale ranges manually


1 Select a plot, then click the  dialog launcher in the [Axes] group on the [Plot Tools] tab of the ribbon.

You can also right-click a plot, and select [Properties] in the context menu to display the [Property Window] dialog.

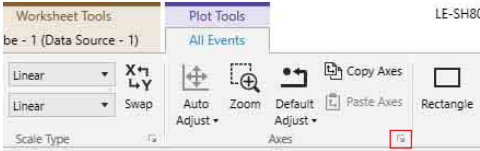

The [Property Window] dialog appears.


- 2 Enter values for the minimum and maximum values for the X axis and Y axis in [Axes].

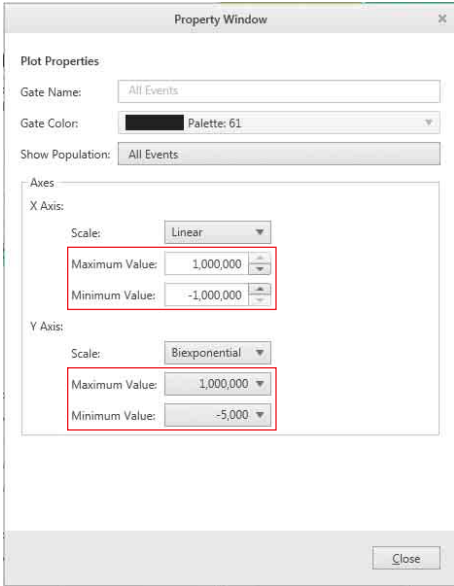

- 3 Click [Close] to close the dialog.

Tip


On histogram plots, by default the Y-axis range is adjusted automatically according to the data values, but the range can also be set manually. To set the range manually, clear the [Auto Adjust] checkbox, and enter the maximum value of the Y-axis manually. The minimum value is fixed at 0. Placing a check mark in the [Auto Adjust] checkbox selects automatic adjustment again.

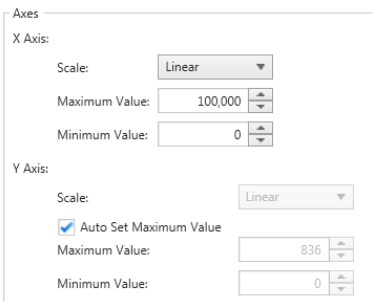


To zoom in using a mouse


You can enlarge a portion of the plot you want to display using the mouse.

- 1 Select a plot, then click [Zoom] in the [Axes] group on the [Plot Tools] tab of the ribbon.

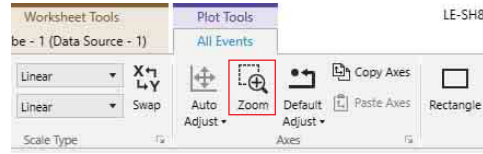


The mouse cursor changes to  when the mouse is placed over the selected plot.

- 2 Drag a selection rectangle over the portion of the plot to display then release the mouse.

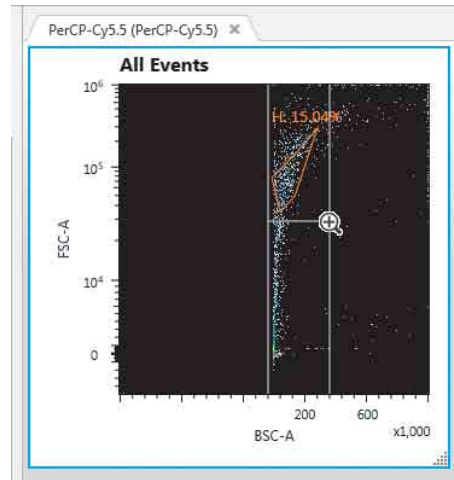

The selected range is scaled to fill the plot area.

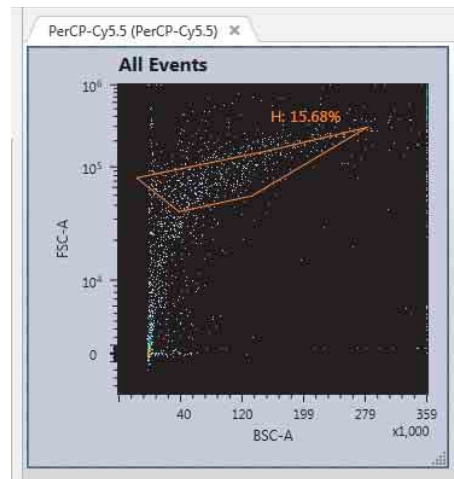

Tip

Gates change shape on plots to track the scale type and zoom factor, but may not be always reproduced accurately. In particular, calculation of polygon gates and ellipse gates are prone to error when a scale is changed from one type to another.

Adding Gates

You can isolate populations of events by drawing gates around the events on plots. Gates of different types (Rectangle, Ellipse, Polygon, Quadrant, Linear) are available which can be drawn to match the distribution of events. Adding a gate allows you to display a new plot representing the population of events that fall within the gate.

Gates are automatically added to the gate hierarchy in the [Gates and Statistics] table when they are added to plots.

You can also create Boolean gates which are defined by a logical combination of existing gates using Boolean logic. Boolean gates are defined not by drawing gates on plots, but by specifying an equation for the gate. The populations defined by Boolean gates can be displayed on plots, just like populations from physical gates. All statistics functions available for physical gates are also available for Boolean gates.

Physical gates and Boolean gates can be selected as the sorting criteria when sorting.


For details, see “Sorting” (page 106).

Adding a Physical Gate

You add a physical gate directly to plots.

- 1 Select a plot on which to add a gate.

Change the plot zoom, as required, in the [Worksheet View] group on the [Worksheet Tools] tab of the ribbon to make it easier to draw the gate.

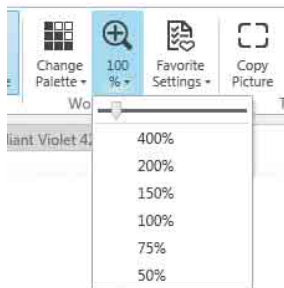

- 2 Click the desired gate type button in the [Gate] group on the [Plot Tools] tab of the ribbon, and draw a gate on the plot.

Alternatively, right-click the plot, then select [Create gate] and the desired shape of gate from the context menu.



[Rectangle]: Adds a rectangle gate.

[Ellipse]: Adds an ellipse gate.

[Polygon]: Adds a polygon gate. A polygon gate is added to a plot by clicking each vertex of the polygon using the mouse. Double-click to close the polygon.

[Quadrant]: Adds a quadrant gate. A quadrant gate is added to a plot by clicking the mouse at the intersection point of the four quadrants.


The four quadrants (“<gate_name>-Q1” to “<gate_name>-Q4”) are added to the plot.

Drag the intersection point to change the size of the quadrants.


[Linear]: Adds a linear gate. Linear gates can be added to histogram plots only.

The gate is added to the plot, together with a single statistic for the gate.

Example: Polygon gate

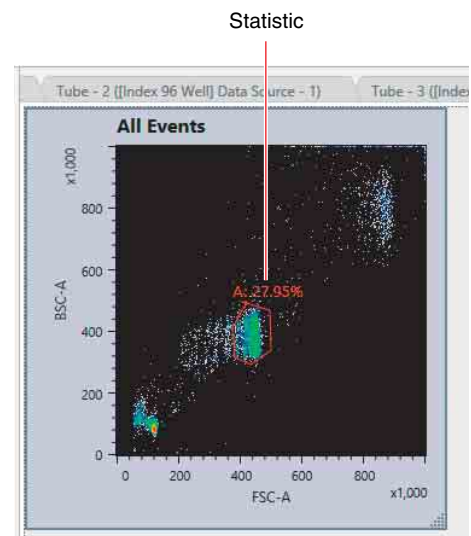


Example: Linear gate

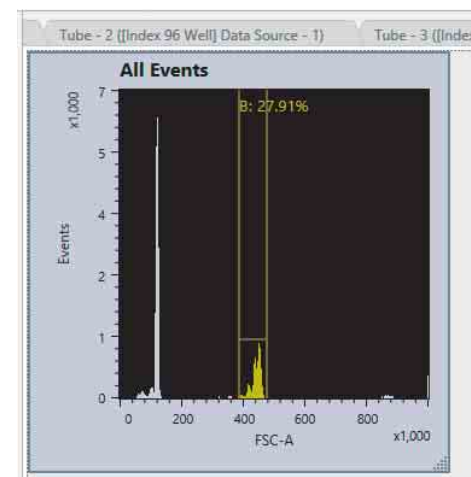

- 3 Edit the name of the gate, the displayed statistic, color, and other preferences, as required.

Select the gate, click [Edit Gate] in the [Gate] group on the [Plot Tools] tab of the ribbon to display the [Gate Editor] dialog, and then select the [Details] tab.

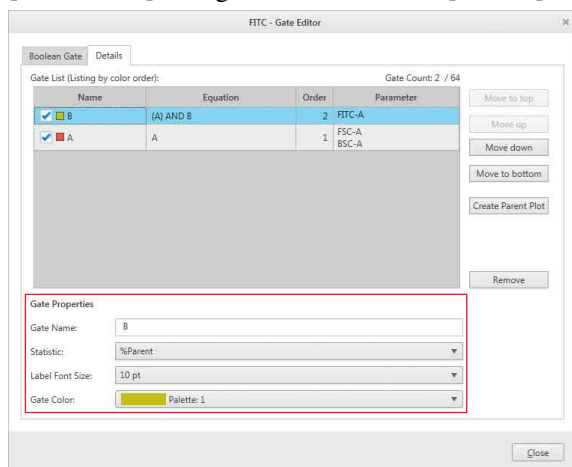

For details about each item, see “[Gate Editor] dialog” (page 145).

- Only linear gates can be pasted in a histogram. And linear gates cannot be pasted in a plot other than a histogram.

Adding a Boolean gate to the Gate Hierarchy


- 1 Select a plot to display the [Plot Tools] tab.
- 2 Click [Edit Gate] in the [Gate] group on the [Plot Tools] tab of the ribbon.



The [Gate Editor] dialog appears.

- 3 Enter the equation for the Boolean gate on the [Boolean Gate] tab.

Enter an equation as follows.

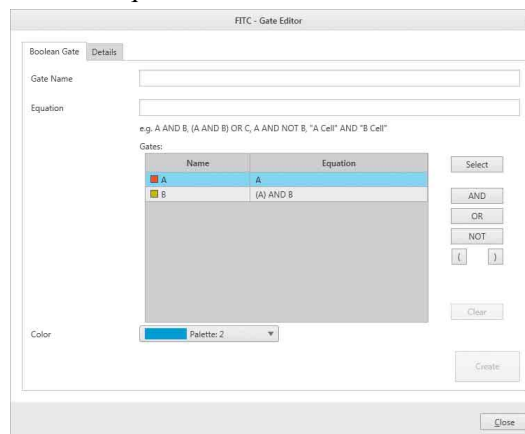

- ① Click the [Equation] text box to place the cursor in the box.
- ② Select a gate in the [Gates] list and click [Select], or double-click the name of the gate. The gate is added to the equation.
- ③ Click [AND], [OR], [NOT], [(, or]], as required. The logic operator is added to the equation. You can also enter the equation directly from the keyboard.
- ④ Repeat steps ② and ③ to build the equation for the gate.


Tip

The equations for all gates can be viewed on the [Details] tab.

Copying a Physical Gate

You can copy a gate drawn on a plot to another plot.

- 1 Right-click a gate to copy, then select [Copy Region] from the context menu.

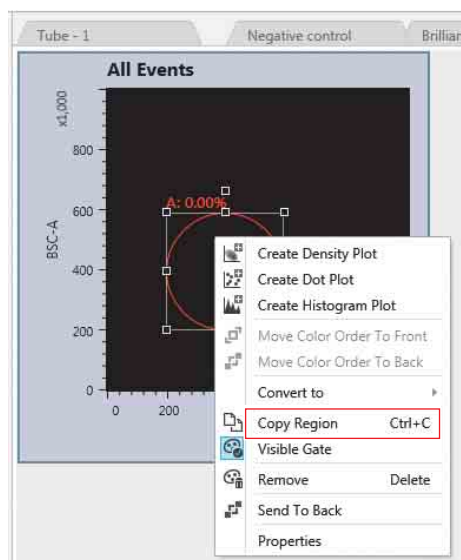

- 2 Paste the gate on the target plot.

Tips

- You can also right-click a gate in the [Gates and Statistics] table to copy a gate within the gate hierarchy using commands in the context menu.
- Adjustment to the scale of the copy destination is not performed. It is copied as a shape.

- ⑤ Change the name of the Boolean gate, as required.
Click in the [Gate Name] text box, and enter a new name.
The names of gates must be unique.
- ⑥ Select a color for the Boolean gate from the [Color] drop-down menu.
- ⑦ Click [Create].
If the equation specified for the gate is verified as a valid definition for a gate, the gate is added to the bottom of the gate hierarchy in the [Gates and Statistics] table.
- ⑧ Click [Close] to close the dialog.

4 Double-click the Boolean gate in the gate hierarchy.

A default FSC v. BSC density plot with linear X axis and Y axis is added to the worksheet.

5 Set the axis labels, scale types, and other properties for the new plot.

For details, see “Editing Plots” (page 92).

Editing Gates

You can change the shape of and settings for gates on plots.

To change the shape of a gate

Select the gate and drag the drawing handles to change the shape.

For quadrant gates, drag the intersection point of the quadrants to another position to change the gate.

To add/delete a vertex of a polygon gate, right-click the vertex point and select [Add Vertex] or [Remove Vertex] from the context menu.



To move a gate

Select the gate and drag the gate to the desired position.

To rotate a gate


You can rotate polygon and ellipse gates.

Select a polygon or ellipse gate and drag the rotation handle to rotate the gate.

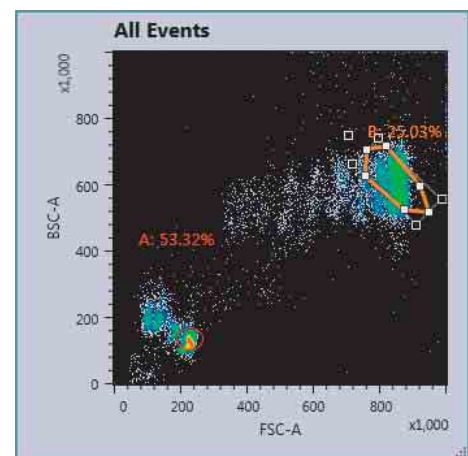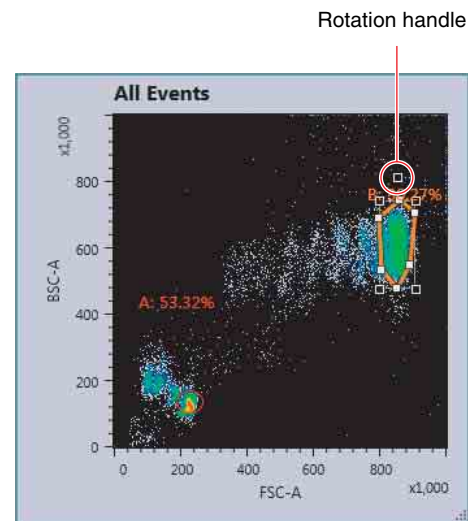

To change gate settings

Select the gate, click [Edit Gate] in the [Gate] group on the [Plot Tools] tab of the ribbon to display the [Gate Editor] dialog, and then select the [Details] tab.

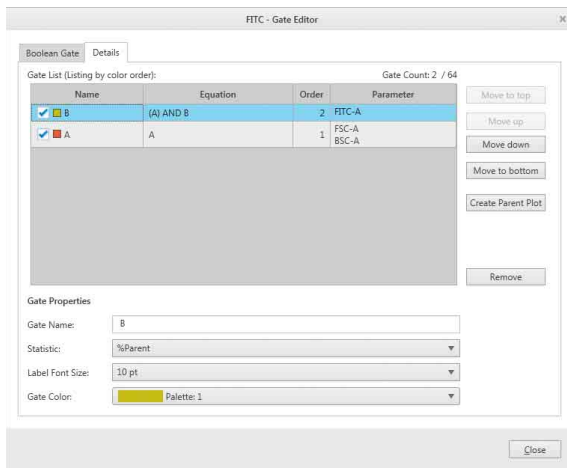

For details about each item, see “[Gate Editor] dialog” (page 145).

To hide a gate

Select a gate you want to hide, then click [Visible] in the [Style] group on the [Gate Tools] tab of the ribbon. Alternatively, right-click the gate you want to hide and select [Visible Gate] from the context menu.

To delete a gate

Select the gate to be deleted, then click [Remove Gate] in the [Gate] group on the [Gate Tools] tab of the ribbon. Alternatively, right-click the gate and select [Remove] from the context menu. You can also delete a gate using the Delete key on the keyboard.

Note

Deleting a gate will delete all data and statistics for the gate, including all child gates and plots derived from the gate.

Tip


There is an upper limit on the number of gates. When creating an overlay, overlay gates are created, reducing the number of gates that can be used on the overlay source worksheet. If the limit is exceeded, overlay display is not possible. Remove any gates that are no longer required.

For details about deleting overlay gates, see “Deleting Overlay Gates” (page 104).

To export the event data within a gate as a CSV file


You can export the event data within a gate as a CSV file on a gate-by-gate basis.

- 1 Select a gate whose event data you want to export, then click [Export Data to CSV File] in the [Tools] group on the [Gate Tools] tab of the ribbon.

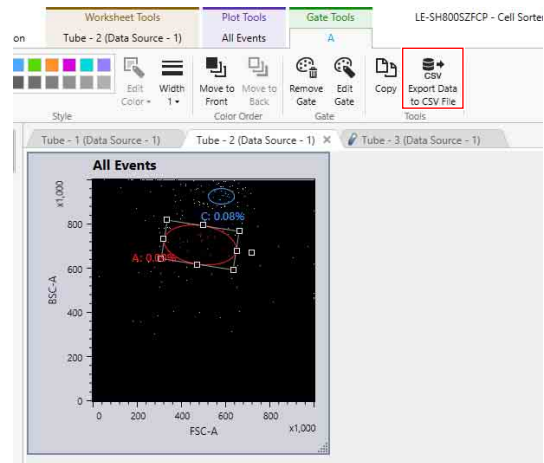

The [Exporting Data] dialog appears.

- 2 Click (Browse) to specify the folder to save the CSV file, then click [Export].

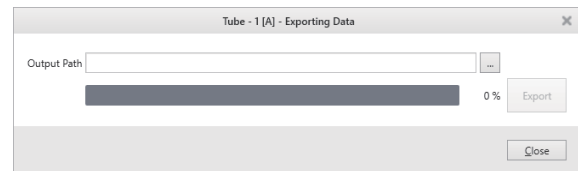

The CSV file is exported.


- 3 Click [Close] to close the dialog.

Displaying Multiple Plots Overlaid


You can display the data for up to eight plots overlaid so that you can compare data acquired in an experiment. Overlay display is supported for the following combinations of plots.

		Overlay source		
		Density plots	Dot plots	Histogram plots
Overlay destination	Density plots	Yes (displayed as dot plots)	Yes (displayed as dot plots)	No
	Dot plots	Yes (displayed as dot plots)	Yes	No
	Histogram plots	No	No	Yes

Example: Dot plot + dot plot

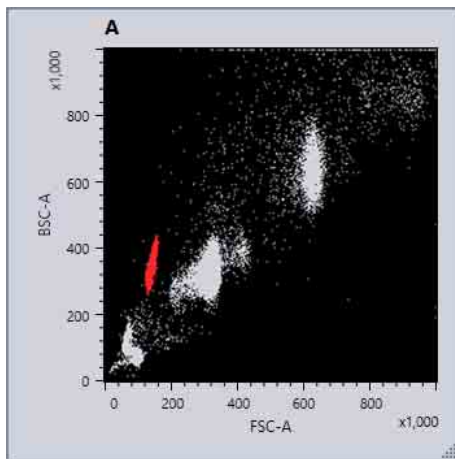

Example: Histogram plot + histogram plot

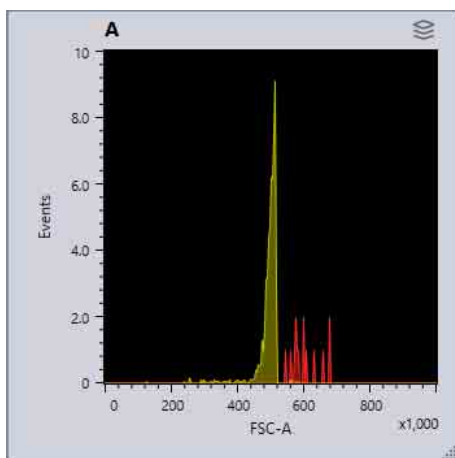



Tip

A plot of another tube can be displayed overlaid on the plot of the gain adjustment tube, but the plot of a gain adjustment tube cannot be overlaid on the plot of another tube.

Displaying Plots Overlaid (Overlay Display)


- 1 Open the worksheet of tube containing a plot that you want to display overlaid.

Open the worksheets for both the overlay source tube and overlay destination tube.
- 2 Select a plot from the overlay source, then click [Copy Plot for Overlay] in the [Overlay] group on the [Plot Tools] tab of the ribbon.

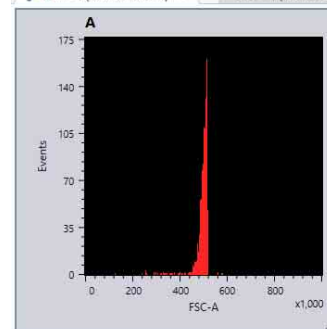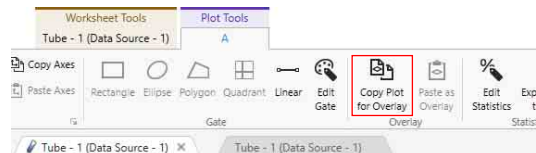


Tips

- [Copy Plot for Overlay] cannot be selected for a plot that has not been recorded. First acquire and record data. For details about operation, see *“Controlling Data Acquisition and Recording”* (page 55).
 - [Copy Plot for Overlay] cannot be selected for the All Events plot. First add a gate, and display a plot for the gate. For details about operation, see *“Adding Gates”* (page 97).
 - [Copy Plot for Overlay] cannot be selected for a plot that is displayed overlaid. First remove the overlay display. For details about operation, see *“Deleting Overlay Layers”* (page 103).
- 3 Select a plot from the overlay destination, then click [Paste as Overlay] in the [Overlay] group on the [Plot Tools] tab of the ribbon.

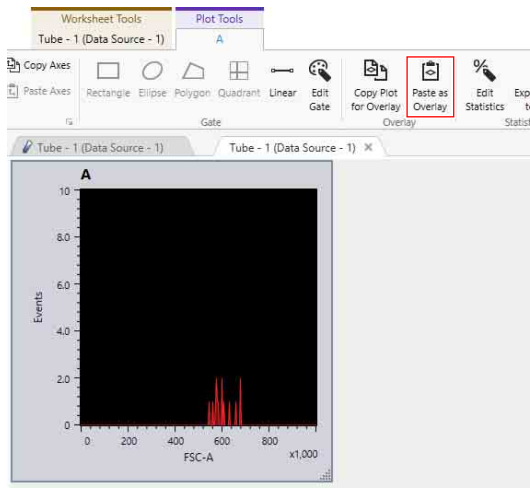

The plots are displayed overlaid. During overlay display, an overlay display icon is displayed at the top right of the plots. The scale and axes names correspond to the overlay destination.

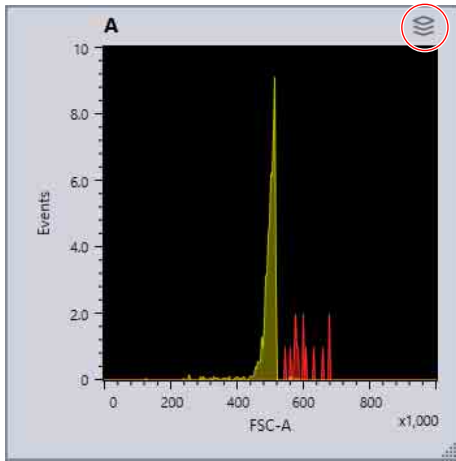


Tips

- Leave the overlay source experiment open; do not close it.
- Child gates in the overlay source are not displayed.
- When creating an overlay, overlay gates are created, reducing the number of gates that can be used on the overlay source worksheet. If the number of gates exceeds the limit, overlay display is not possible. In this case, remove any gates that are no longer required. For details about operation, see *“Deleting Overlay Gates”* (page 104).
- The layer information of the overlay source layer information is not exported when exporting the data of overlaid plots.
- The overlay source worksheet is not closed automatically when the overlay destination worksheet is closed. Close it manually.


To align the scales of histogram plots

On histogram plots, you can adjust the scales automatically or switch the scales to match the overlay destination scale.

- 1 Select a histogram plot, then click the  dialog launcher in the [Axes] group on the [Plot Tools] tab of the ribbon.


The [Property Window] dialog appears.

- 2 Select whether to adjust scales automatically using the [Auto Adjust] under [Maximum Value].

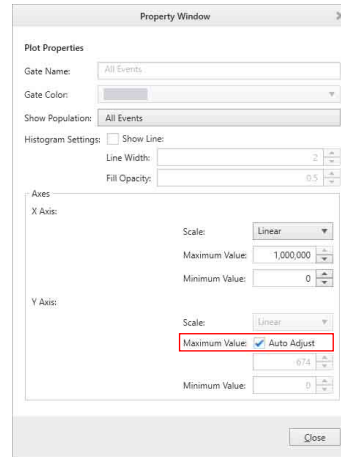


When on

The scales are adjusted automatically.

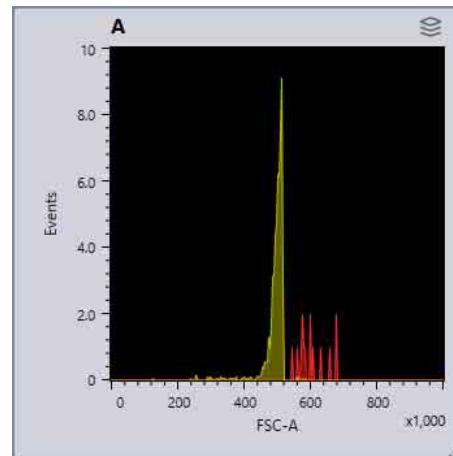


When off

The scales are aligned to the overlay destination.

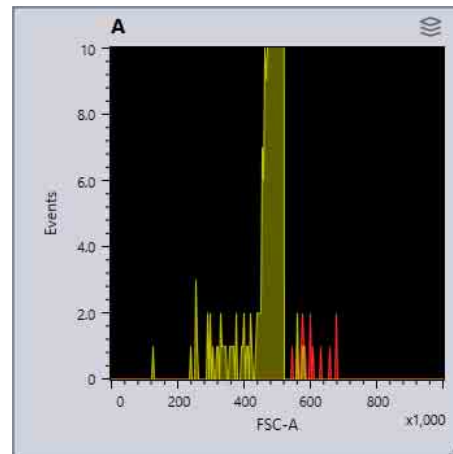


Customizing the Overlay Display

Click to display the overlay panel on the right side.

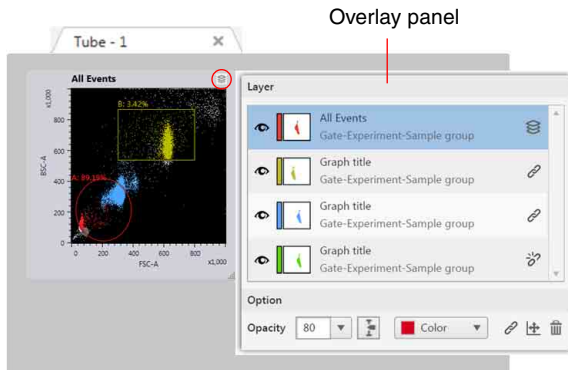

You can show/hide layers and change colors in the overlay panel.

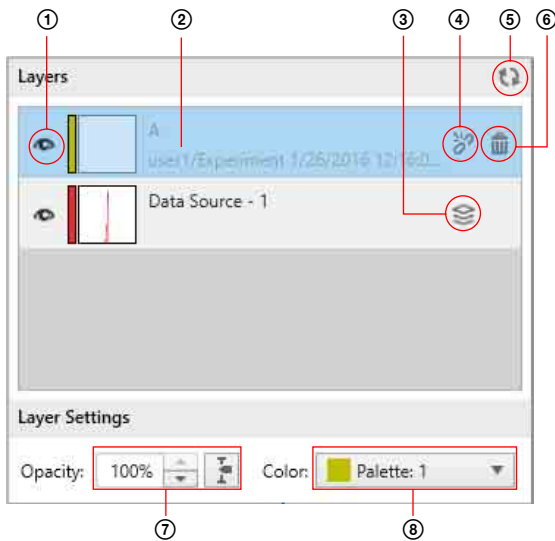

- ① **Layer show/hide selector**
Click repeatedly to toggle between show () and hide ().
- ② **Layer name**
Hover the mouse over the layer name to display user/experiment/sample group/sample tube/gate information.
- ③ **Base plot icon**
Indicates the base plot.
- ④ **Unload icon**
Indicates that overlay display is unavailable because the overlay source worksheet is not open.
- ⑤ **Refresh icon**
Updates the overlay display to the latest status.
- ⑥ **Delete icon**
Deletes the selected layer.
- ⑦ **[Opacity]**
Adjusts the opacity of the selected layer.

- ⑧ **[Color]**
Click to display a color palette in order to change the color of the selected layer.


Tip
If the overlay source worksheet is closed and the display is refreshed, the worksheet is reopened and the layer is shown. However, if the overlay gate of the overlay source is deleted, the display is not refreshed. Recreate the overlay display again.

Deleting Overlay Layers

You can delete a single overlaid layer or all layers.


To delete a single layer

Open the overlay panel, select the layer you want to delete, and click .

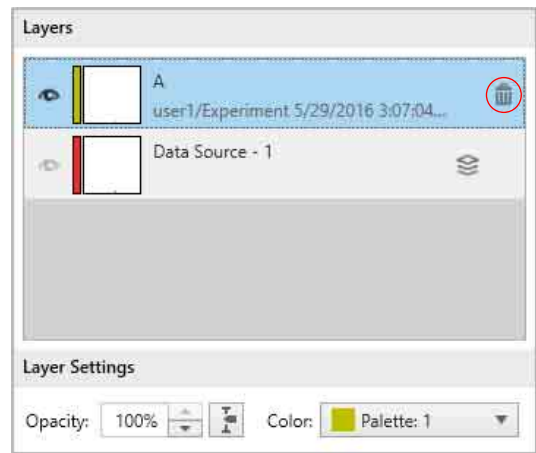

To delete all layers

Right-click the plot, and select [Clear overlay plots] from the context menu.

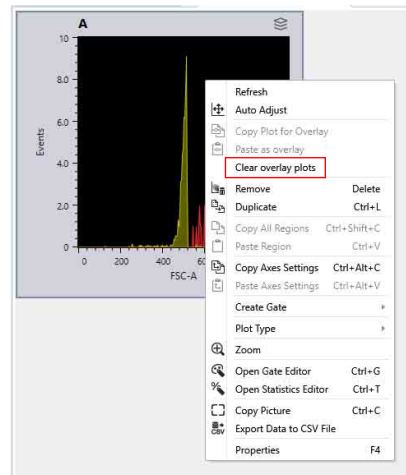


Deleting Overlay Gates

When creating an overlay, overlay gates are created, reducing the number of gates that can be used on the overlay source worksheet. If the number of gates exceeds the limit, overlay display is not possible. In this case, remove any gates that are no longer required.

- 1 Right-click the worksheet, then select [Overlay Gate List] from the context menu.

The [Overlay Gate List] dialog appears.

- 2 Select the gate to remove, and click [Remove].

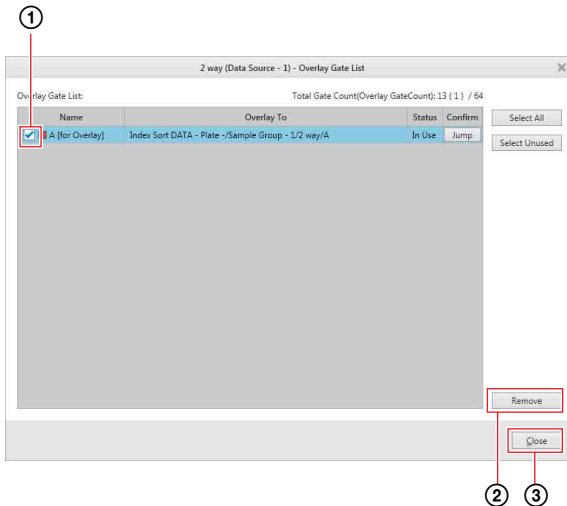

- 1 Place a check mark in the checkboxes for the gates to remove.
- 2 Click [Remove] to remove overlay gates that are no longer required.
- 3 Click [Close] to close the dialog.

Tip

The maximum number of gates is the total of physical gates and Boolean gates.

For details about deleting physical gates and Boolean gates, see “To delete a gate” (page 100).


Displaying Statistics

You can change the statistics items displayed for each gate in the [Gates and Statistics] table.

For details about statistics, see “[Gates and Statistics] Table” (page 149).


- 1 Click [Edit Statistics] in the [Statistics] group on the [Worksheet Tools] tab of the ribbon.

Alternatively, right-click the [Gates and Statistics] table and select [Open Statistics Editor] from the context menu.

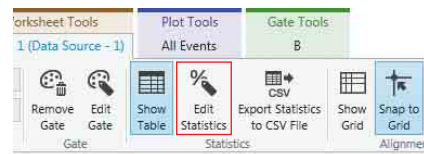

The [Statistics Editor] dialog appears.

- 2 Specify the items to be displayed in the [Gates and Statistics] table.

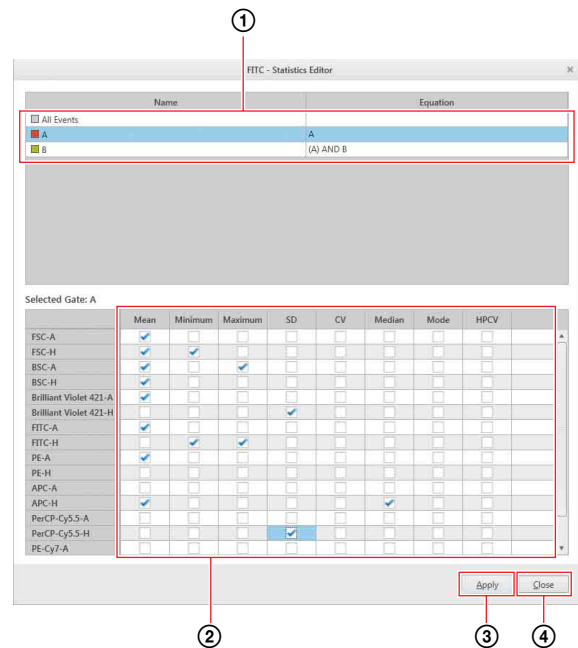

- 1 Select a gate.
The statistics values for the population in the selected gate are displayed in the lower area.
- 2 Place a check mark in the checkboxes for the statistics you want to display for each parameter.
- 3 Click [Apply] to update the statistics displayed in the [Gate and Statistics] table.
- 4 Click [Close] to close the dialog.

To change the width of columns

Drag the right edge of a table column to change the width of the column.

To change the display order of columns


Select the title of the column to move, and drag it to the desired location.

Tip

The [Name] column cannot be moved.

To export statistics as a CSV file

- 1 Click [Export Statistics to CSV File] in the [Statistics] group on the [Worksheet Tools] tab of the ribbon.

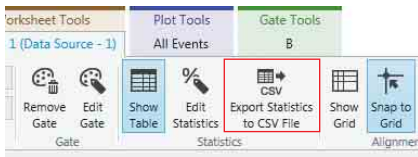

The [Exporting statistics table] dialog appears.

- 2 Click (Browse) to specify the folder to save the CSV file, then click [Export].



The CSV file is exported.

- 3 Click [Close] to close the dialog.

This chapter describes the various sorting modes and detailed sorting criteria.

The SH800 can sort cells contained in a selected subpopulation in the sample fluid at high speed. The instrument sorts droplets containing the selected target cells using gates drawn on plots on a worksheet. Sorted droplets can be collected in 5 ml round tubes, 15 ml conical tubes, and on multi-well plates. In addition, gates can be combined using Boolean logic to uniquely identify a population for sorting using multiple attributes, providing powerful sorting capabilities.

Sorting Mode

Overview

In an ideal sort, cells travel past the optical detection point in single file and are sorted one cell at a time in separate droplets. In practice, however, cells appear randomly and can occur in close proximity within the same droplet or in adjacent droplets (called coincidences).

When a cell is detected during sorting, it is generally not possible to know with great accuracy if the signal represents the passage of a large single cell, small multiple cells adhering to one another, or other non-cellular material. Accordingly, it is more common to refer to each signal detection not as a cell, but as an event.

The sorting mode sets the rules used when sorting coincidences. These rules determine the sorting decision for each event based on the type of event and the distance between adjacent events. For example, a decision might be made to sort an event regardless of an adjacent unwanted event if the goal is to isolate as many cells as possible from a small population. Conversely, a decision might be made to not sort an event if there is an adjacent unwanted event that could reduce the purity of the sorted population.

Tip

The trigger threshold setting in the [Detector & Threshold Settings] dialog (*page 86*) can be critical for effective sorting. If the threshold level is set too high, small cells and cell debris (noise) that fall below the threshold value may be unintentionally sorted together with events above the trigger threshold, depending on the sorting mode setting. To increase sorting purity, it is recommended that the trigger threshold value be set to a relatively low value.

About custom sorting mode

In addition to the preset sorting modes, each user can configure up to five custom sorting modes that allow you to adjust the position of target cells within a droplet for sorting.

For details about configuration, see “Configuring Custom Sorting Mode” (page 117).

The SH800 supports the following eight preset sorting modes for obtaining the desired sorting results. The modes provide different levels of purity¹⁾ and yield²⁾. Note that the droplets that are sorted in each sorting mode vary depending on the size of the nozzle of the sorting chip.

Single Cell ← Increasing purity

Increasing yield →

	Single Cell (page 108)	Ultra Purity (page 111)	Purity (page 110)	Semi-Purity (page 110)	Normal (page 109)	Semi-Yield (page 111)	Yield (page 112)	Ultra-Yield (page 112)
	3drops							
100 µm/ 70 µm sorting chip								
Number of droplets sorted	3	1	1	1	1	1	1	1
130 µm sorting chip								
Number of droplets sorted	3	1 to 3	1 or 2	1 or 2	1 or 2	1 or 2	2	3

△: Target cell, □: Non-targeted (conflicting) cell

: Sorted droplet

: Waste droplet

: Multi-drop inclusion

: Multi-drop exclusion

-----: Sorting regions

1) Purity measures the proportion of target cells collected as a percentage of all cells collected.

2) Yield measures the proportion of target cells collected as a percentage of target cells in the sample. It indicates the percentage of the target population that is collected.

A high-degree of purity of targeted cells can be obtained if the targeted cell events are separated from other

unwanted events. In this case, the droplets containing the targeted events are sorted, while all unwanted events and empty droplets are passed to waste disposal. In practice, however, there is a trade-off to be made between yield, recovery, and purity.

Sorting is a complex process. There are several factors that can affect the outcome. The position of a cell within a droplet can only be determined within a certain degree of accuracy, fluctuations can occur in the fluid flow due to the passage of large cells or variations in the sheath fluid pressure, and the droplet charging pulse timing may be slightly out of phase with droplet formation.

- If a targeted event appears close to the top or bottom edge of a droplet, the probability of correct recovery is reduced. The targeted event may actually be output in an adjacent droplet. One option is to sort successive droplets in order to catch the event to increase yield and recovery.

- If a targeted event appears close to the edge of a droplet and an unwanted event occurs in the adjacent droplet, or vice versa, the decision whether to sort the droplets or not depends on the goal: isolating a pure population or recovering as many targeted cells as possible. Sorting the targeted event will increase the yield but with a subsequent reduction in purity. Conversely, rejecting the targeted event will maintain high purity but with reduced yield. Likewise, a similar decision occurs if a targeted event and an unwanted event appear in the same droplet.
- If two, different targeted events for collection in separate collection tubes appear in the same droplet or close to one another in adjacent droplets, the sorting decision must determine which of the two targeted events is sorted correctly or whether to abort sorting of the two events.
- If two or more cells, including target and/or unwanted cells, appear at the optical detection point in very close proximity, a single event will be triggered if the detected signal does not fall below the threshold level between cells. In this case, it is impossible to distinguish the event as separate cells in a single-pass sort. If the event contains targeted and unwanted cells, it will exhibit some of the characteristics of both types of cell. If the sort criteria only define a limited subset of targeted cell characteristics, the event may satisfy that criteria and be sorted accordingly, with a corresponding reduction in purity due to the presence of the unwanted event. Sorting using multiple gates to isolate target cells for multiple characteristics can help isolate targeted events.

In addition, hardware-related factors can affect sorting. If two or more cells, including target and/or unwanted cells, appear at the optical detection point in reasonably close proximity, separate events will be triggered if the detected signal falls below the threshold level between cells. However, if the acquisition electronics are not yet ready to receive the second event, the physical presence of the second event is detected but the event is not analyzed. This time is called the dead time. These events are aborted automatically and cannot be sorted, reducing the potential yield.

From this discussion, it can be seen that sorting only a single targeted population where the targeted events occur in the center of droplets that contain no other unwanted events will always produce the best sort results in terms of purity. Sorting all events for a single targeted population will always produce the best results in terms of yield.

Single Cell Mode


In Single Cell mode, only droplets containing a single targeted cell within the center region of the cell are sorted (adjacent droplets are also sorted when using a 130 μm sorting chip) if the adjacent droplets are empty. If an adjacent droplet contains the same type of target cell, the droplet is not sorted.

100 μm /70 μm sorting chip

1-droplet sorting

A droplet is sorted if it contains a single targeted event located within the center region of the droplet and both of the adjacent droplets are empty.


One droplet is sorted for each event matching the sorting criteria.

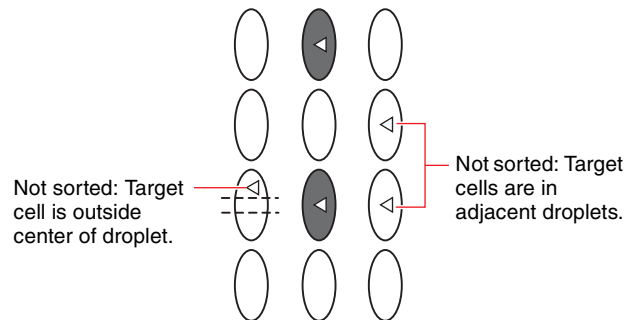

3-droplet sorting (3 drops)

A droplet is sorted if it contains a single targeted event and the adjacent and subsequent droplets on both sides are all empty.


Three droplets are sorted for each event matching the sorting criteria. In comparison with 1-droplet sorting, an improvement in sorting performance can be expected when sorting cells of 15 μm or larger, but this does not guarantee cells can be sorted.

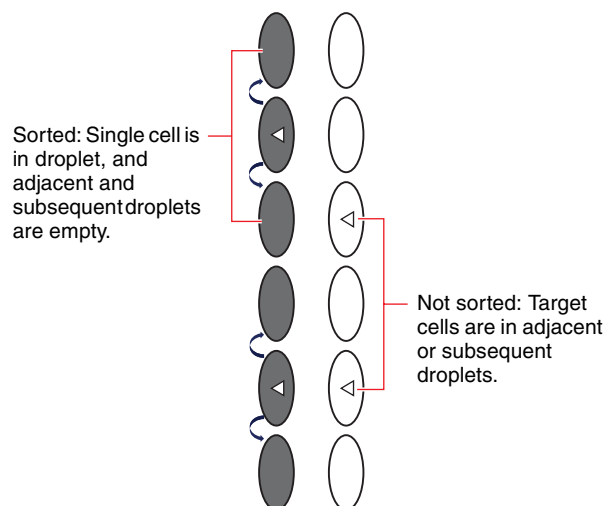

130 μm sorting chip

A droplet is sorted if it contains a single targeted event and the adjacent and subsequent droplets on both sides are all empty.

Three droplets are sorted for each event matching the sorting criteria.

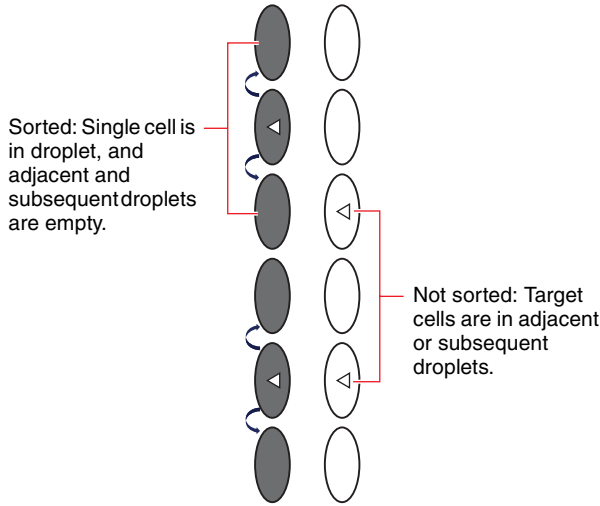


Normal Mode

Normal mode is the default sorting mode.

100 µm/70 µm sorting chip

A droplet is sorted if it contains one or more targeted events of a single type.

One droplet is sorted for each event matching the sorting criteria.

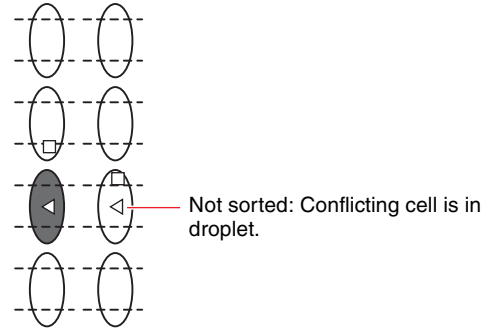


130 µm sorting chip

A droplet is sorted if it contains one or more targeted events of a single type. In addition, if the targeted event is near the edge of the droplet, the nearest adjacent droplet is also sorted if that droplet is empty or non-conflicting.

One or two droplets are sorted for each matching sorting criteria.

Sorted: Target cell is outside center of main droplet and nearest adjacent droplet is empty.

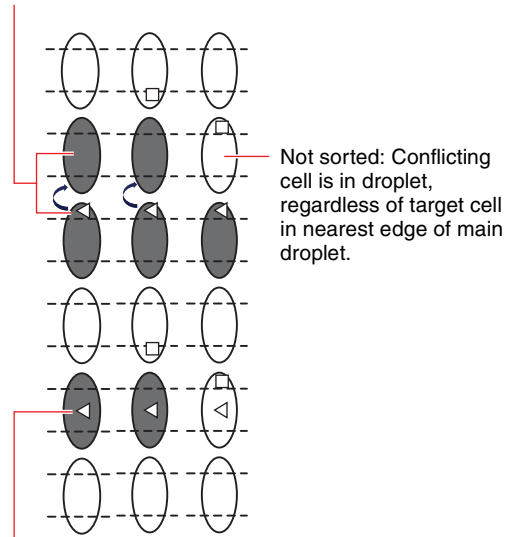


Sorted: Target cell is within center of droplet.

Semi-Purity Mode

In Semi-Purity mode, the purity achieved is higher than in Normal mode.


100 μm/70 μm sorting mode

A droplet is sorted if it contains one or more targeted events of a single type and the nearest edge region of both of the adjacent droplets is empty or non-conflicting. One droplet is sorted for each event matching the sorting criteria.

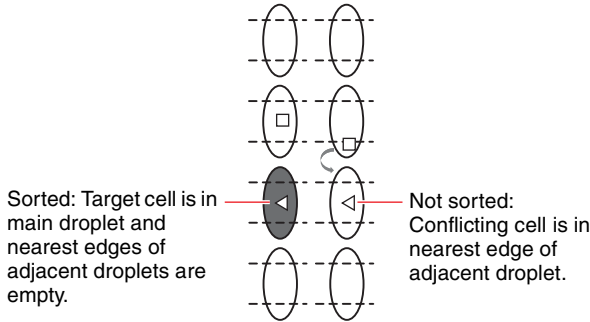

130 μm sorting mode

A droplet is sorted if it contains one or more targeted events of a single type and the nearest edge region of both of the adjacent droplets is empty or non-conflicting. In addition, if the targeted event is near the edge of the droplet, the nearest adjacent droplet is also sorted if the nearest edge region of the subsequent droplet is empty or non-conflicting. One or two droplets are sorted for each event matching the sorting criteria.

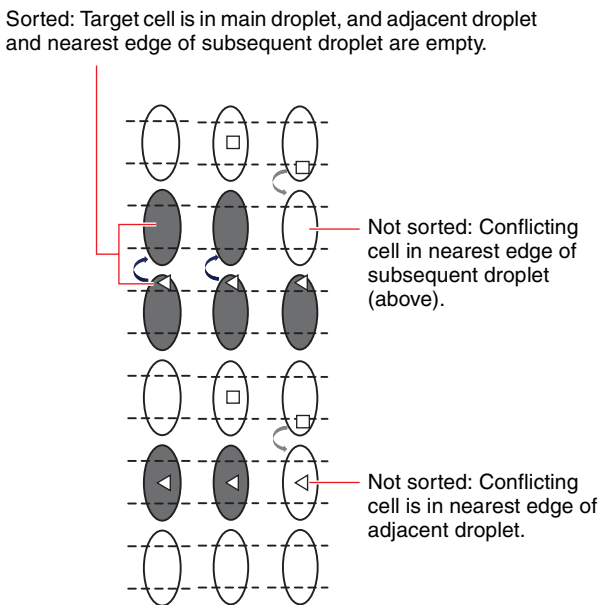


Purity Mode

In Purity mode, cells are sorted with high purity.

100 μm/70 μm sorting chip


A droplet is sorted if it contains one or more targeted events of a single type and the nearest edge region of both of the adjacent droplets is empty or non-conflicting. One droplet is sorted for each event matching the sorting criteria.

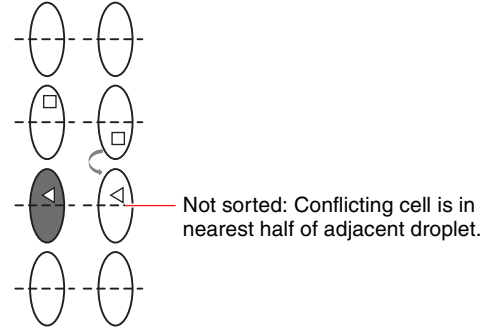

130 μm sorting chip

A droplet is sorted if it contains one or more targeted events of a single type and the nearest half of both of the adjacent droplets is empty or non-conflicting. In addition, if the targeted event is in one half of the droplet, the nearest adjacent droplet is also sorted if the nearest half of the subsequent droplet is empty or non-conflicting. One or two droplets are sorted for each event matching the sorting criteria.

Sorted: Target cell is in main droplet, and adjacent droplet and nearest half of subsequent droplet are empty.

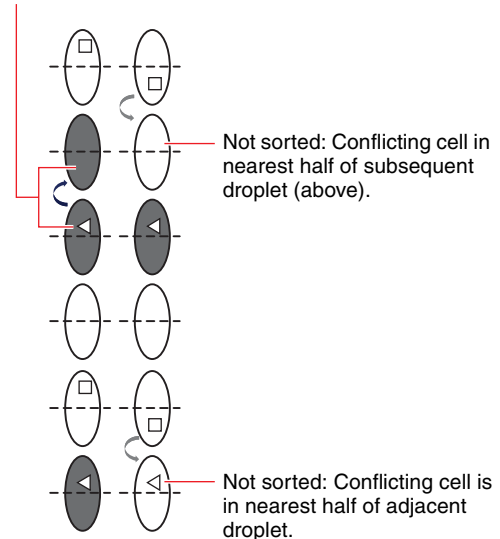


Ultra Purity Mode

In Ultra Purity mode, cells are sorted with the highest purity.

100 μm /70 μm sorting chip

A droplet is sorted if it contains one or more targeted events of a single type and the both of the adjacent droplets are empty or non-conflicting.


One droplet is sorted for each event matching the sorting criteria.

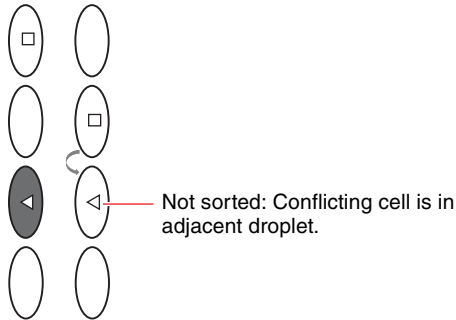

130 μm sorting chip

A droplet is sorted if it contains one or more targeted events of a single type and the both of the adjacent droplets are empty or non-conflicting. In addition, the adjacent droplet is also sorted if the subsequent droplet is empty or non-conflicting.


One, two, or three droplets are sorted for each event matching the sorting criteria.

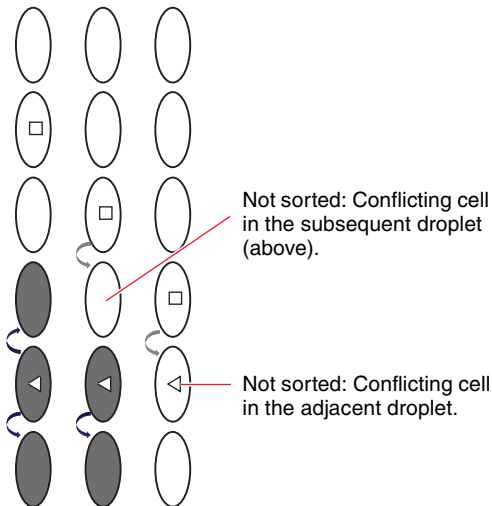

Semi-Yield Mode

In Semi-Yield mode, a droplet is sorted if it contains one or more targeted events, even if the droplet contains conflicting events. In addition, if a targeted cell is near the edge of the droplet, the adjacent droplet is also sorted, even if it too contains conflicting events.


One or two droplets are sorted for each event matching the sorting criteria.

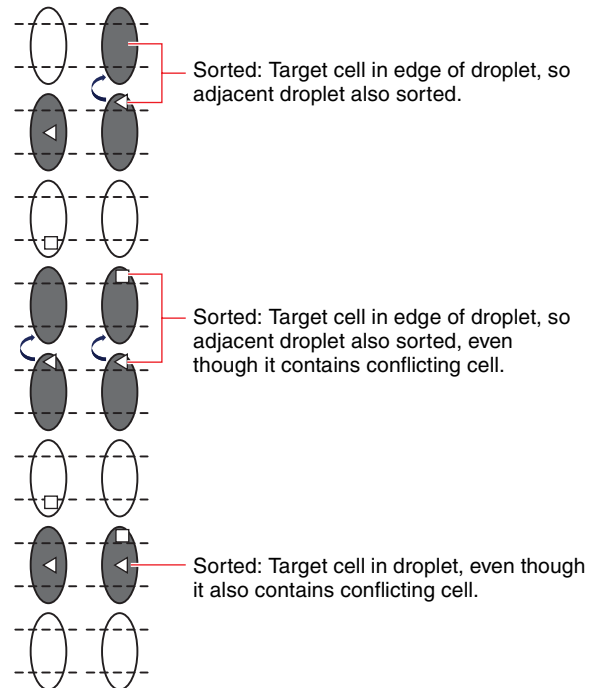

Yield Mode

In Yield mode, a droplet is sorted, together with the adjacent droplet nearest to the target event, if it contains one or more targeted events, even if the droplets contain conflicting events.

Two droplets are sorted for each event matching the sorting criteria.

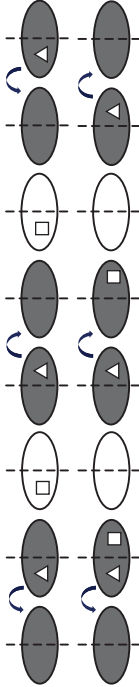

Ultra Yield Mode

In Ultra Yield mode, a droplet is sorted, together both adjacent droplets, if it contains one or more targeted events, even if the droplets contain conflicting events. Three droplets are sorted for each matching sorting criteria.




Adjusting the Sort Position (SH800SP and SH800ZP)

You can adjust the position where droplets fall within wells when sorting into a multi-well plate. The description in this section shows a 96-well plate as an example. The operation is the same for all other multi-well plates. When sorting into a 384-well plate or 384-well PCR plate, always adjust the sort position before starting.

Tip

The sort position cannot be adjusted when running 2-way tube sorting.


- 1 Place a cover over the multi-well plate so that you can see the position where droplets fall.
- 2 Select a multi-well plate sorting method in [Method] in the [Sort Control] pane, and click [Settings].

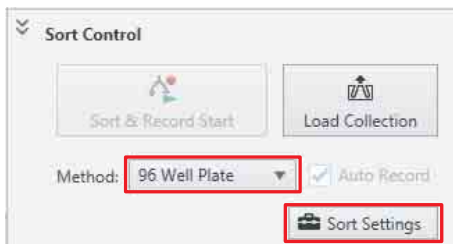

The [Sort Settings] dialog appears.

- 3 Adjust the sort position on the [Plate Adjustment] tab.

Adjust the position where droplets fall by visually monitoring the top of the multi-well plate.

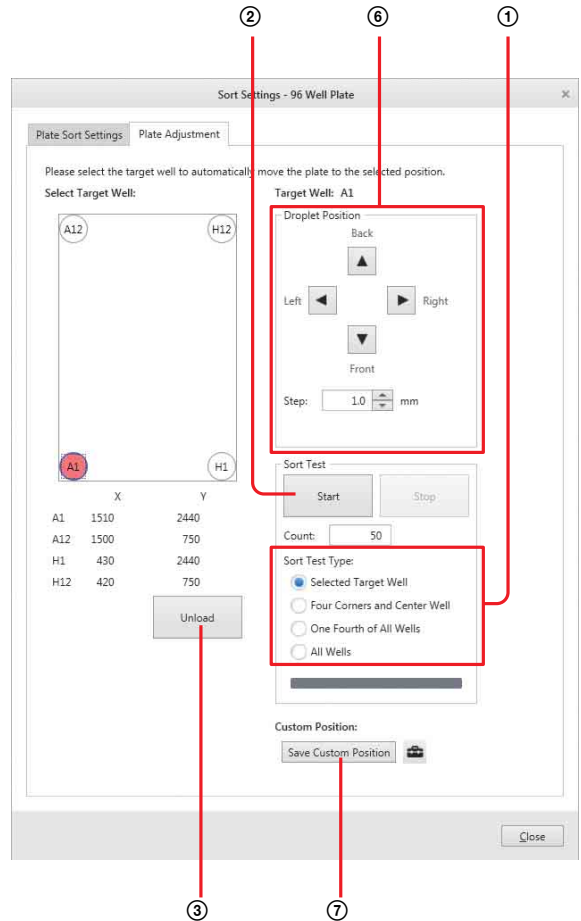


- ① Select the type of test sorting in [Sort Test] > [Sort Test Type].
 - [Selected Target Well]: Droplets fall in the well position selected in [Select Target Well]. Select this option to test sort into a single well at a time. When selected, select the well into which droplets should fall in [Select Target Well].
 - [Four Corners and Center Well]: Droplets fall into the four corner wells and the center well.
 - [One fourth of All Well]: Droplets fall into 1/4 of all wells.
 - [All Well]: Droplets fall into all wells.
- ② Click [Start] in [Sort Test]. Change the number of events to be sorted into each well in [Count], as required. Droplets start to fall in the position selected in step ①.
- ③ Click [Unload].
- ④ Visually check the position of the droplets on the multi-well plate cover.
- ⑤ Adjust the sort position of droplets using the arrow buttons in [Droplet Position]. For example, if the position for well A1 was 1 mm too far left and 1.5 mm too far up:
 - (1) Click well A1.
 - (2) Move the position 1 mm to the right using and 1.5 mm down using in [Droplet Position].

- (3) Click [Start] in [Sort Test].
- (4) Click [Unload].
- (5) Visually check the position of the droplets on the multi-well plate cover.
- (6) Repeat the procedure for wells A12 and H12, as required.

- ⑥ To set the adjusted position as the default, click [Save Custom Position].
- ⑦ Wipe off droplets after confirming their position.

Tip

Clicking  displays the [User Preference] dialog (page 120) for setting user default preferences.

- 4 When finished, click [Close] to close the dialog.

Index Sorting

You can associate the cells sorted in each well with the events for those cells on plots. Using the index sorting function allows you to select a specific well and view the data for cells collected in that well.


This section describes sorting using a 96-well plate, but the procedure is identical for sorting onto 8-well strips or 384-well plates.

Tip

Index sorting is not available for 2-way tube sorting.


Configuring Index Sorting

- 1 Select a sorting method in [Method] in the [Sort Control] pane, and click [Sort Settings].

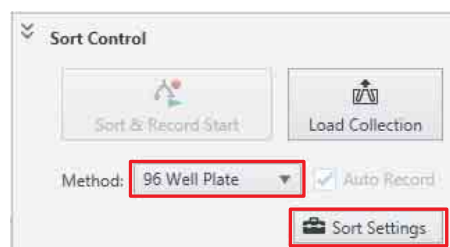


The [Sort Settings] dialog appears.

- 2 Place a check mark in [Add index sort information] on the [Plate Sort Settings] tab.

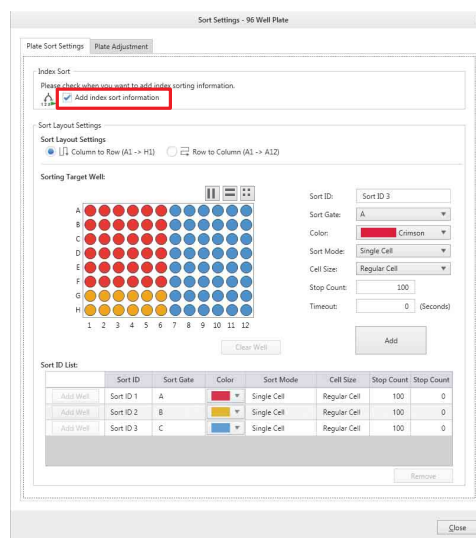


- 3 When finished configuring other items, click [Close] to close the dialog.

- Click [Start] on the data acquisition control pane to start acquiring data.

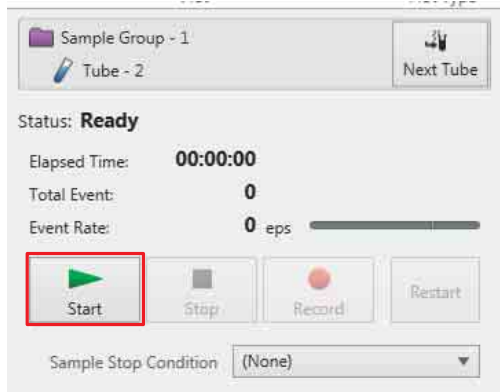

- Click [Index Sort Start] on the [Sort Control] pane to start sorting.

When configured for indexed sorting, “Index” is displayed on the button.

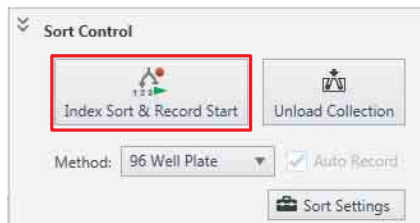

Tip

Starting sorting also automatically starts recording.




Analyzing Index Sort Data

This section describes how to analyze data acquired during index sorting.

This function allows you to examine the position on plots of cells sorted in a specific well.


- Click [Analyze Index Data] in the [Index Sorting] group on the [Worksheet Tools] tab of the ribbon.

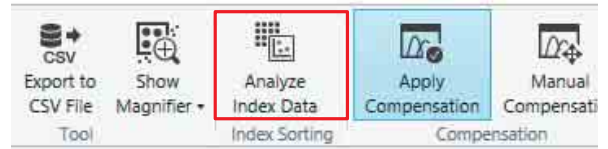

The [Index Analysis] dialog appears.

- Select a mode in [Analysis Mode].

The wells displayed vary depending on the selected mode.

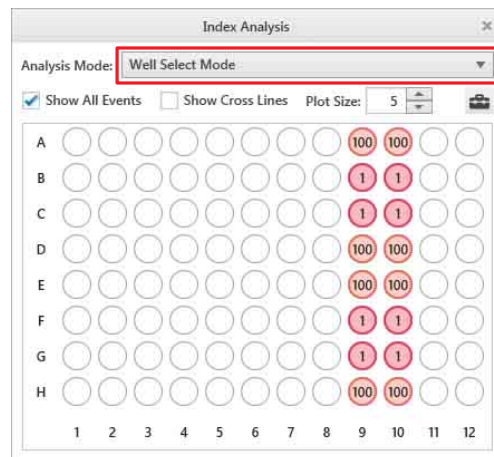

[Well Select Mode]: Select this mode when you want to display the position on plots of cells sorted in each well. You can select multiple wells at the same time.



[Gate Select Mode]: Select this mode to check which wells the cells within the selected gate are sorted.

If [Well Select Mode] is selected

Displays all wells in the multi-well plate. Clicking a well in the dialog displays the position on plots of the cells collected in that well.

Placing a check mark in [Show All Events] displays all data acquired during sorting (includes events that were sorted and events that were not sorted) on plots.


Specifies the size of the dot displayed on plots indicating the data position.

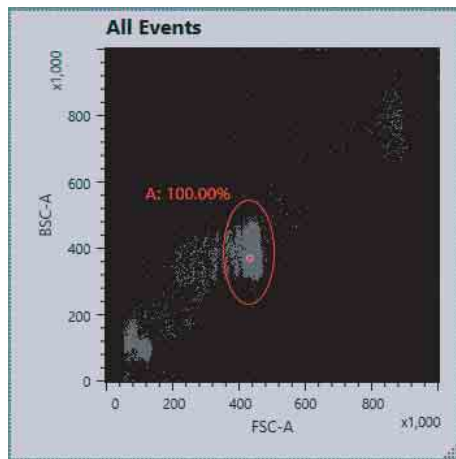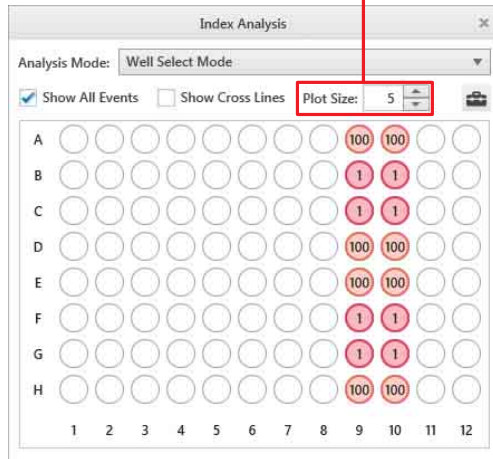

If [Gate Select Mode] is selected


Displays only those wells that contain events encompassed by the gate selected in [Selected Gate]. Placing a check mark in [Overlay Index Events] displays the wells into which the cells within the gate are sorted on plots.

Displays the wells into which the cells within the selected gate are sorted.

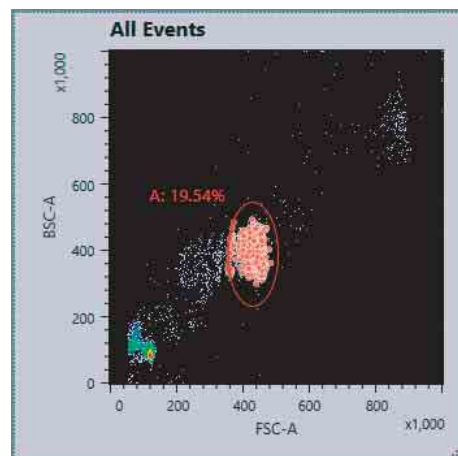


To view the dot position accurately


Placing a check mark in [Show Cross Lines] displays horizontal and vertical lines intersecting the dot position for more accurate positioning.

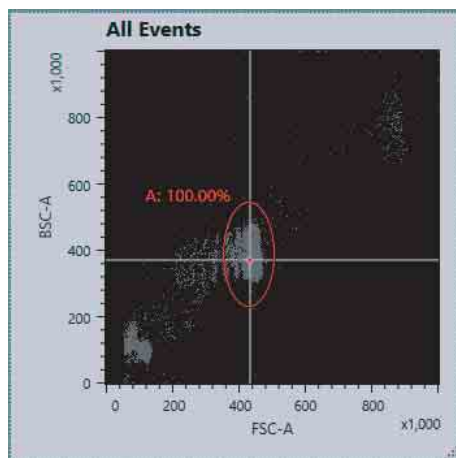

Tip

Clicking  displays the [User Preference] dialog (page 120) for setting user default preferences.

Tip


Clicking  displays the [User Preference] dialog (page 120) for setting user default preferences.

Adjusting Sort Parameters Manually (Experienced Users)


During startup, the SH800 automatically measures and calibrates the sorting parameters for sorting using automatic setup beads. However, there may be cases where you wish to override the automatic control system and adjust the sorting parameters manually.

These parameters are for experienced operators who are already familiar with manual sorting controls. The manual controls are located in the [Advanced Settings] window.

- 1 Click [Settings] on the [Cytometer] tab of the ribbon. The [Cytometer Settings] dialog appears.
- 2 Click [Advanced Settings].

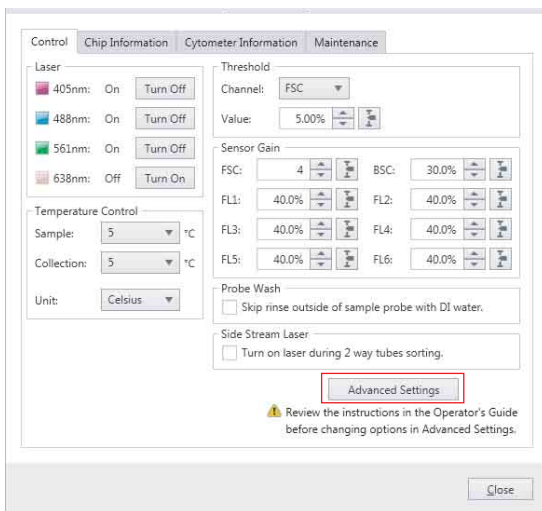

The [Advanced Settings] dialog appears.

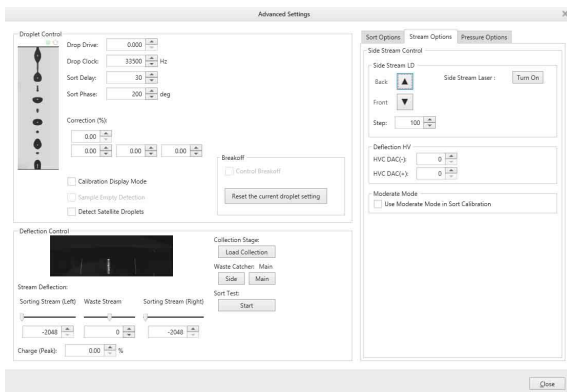

For details about each item, see “[Advanced Settings] dialog” (page 139).


Configuring Custom Sorting Mode

Each user can configure up to five custom sorting modes that allow you to adjust the position of target cells within a droplet for sorting.

Tip


Cells near the edge of droplet may cause unstable side streams. [Centering mask] in custom sorting mode allows you to adjust the balance between cell purity and yield.

- 1 Click [Information] on the [File] tab of the ribbon, then click [User Preference].

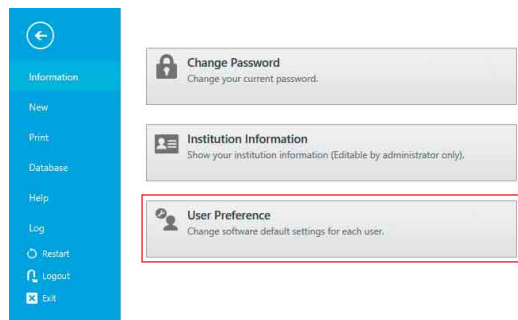

The [User Preference] dialog appears.


- 2 Select [Sort] from the menu on the left side, then click [Edit] in [Custom Sort Settings].

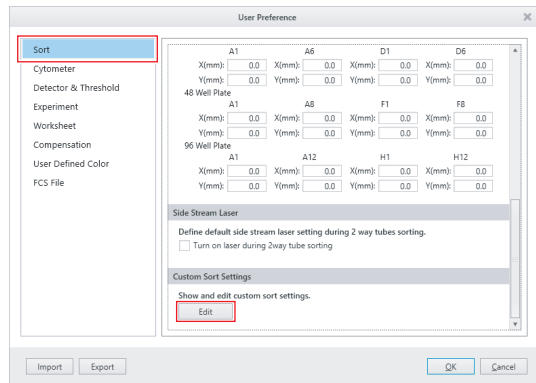

The [Custom Sort Settings] screen appears.

- 3 Set each item, then click [OK].

Select a base sorting mode in [Mode], then set a configuration name and the position range of cells within a droplet for sorting.

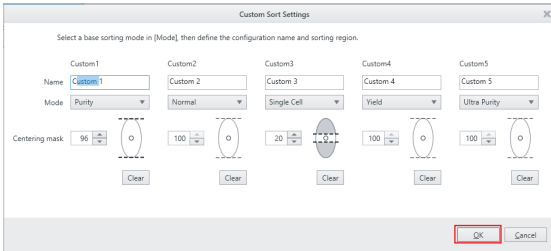

[Name]: Enter a name for the custom mode.

[Mode]: Select the base sorting mode.

When a sorting mode is selected, the default value for that mode is entered in [Centering mask].

[Centering mask]: Specify the position of target cells within a droplet for sorting.


Tips

- The [Name] field cannot be left blank. Always specify a name.
- Duplicated names cannot be specified.
- Specifying the same name as a preset sorting mode is not possible.

The screen returns to the [User Preference] dialog.

4 Click [OK].

The custom mode can now be selected for [Mode] on the [Sort Control] pane.

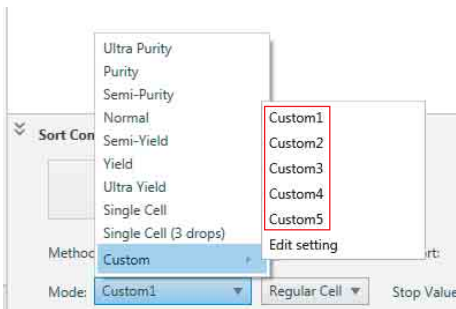

Tips

- The menus and buttons displayed vary depending on the user account privileges.
- The screenshots displayed in this document may vary from the actual software.

[File] Window


The [File] window (Backstage View) is displayed by clicking the [File] tab of the ribbon.

For details about [Restart], [Logout], and [Exit], see “[File] window” (page 29).


[Information] window

The [Information] window is displayed by clicking [Information] in the menu on the left side. The [Information] window is used to change login passwords, set institution information, and add/modify user accounts.

Example: Administrator account login

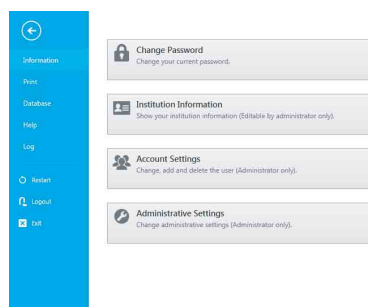

Example: User account login

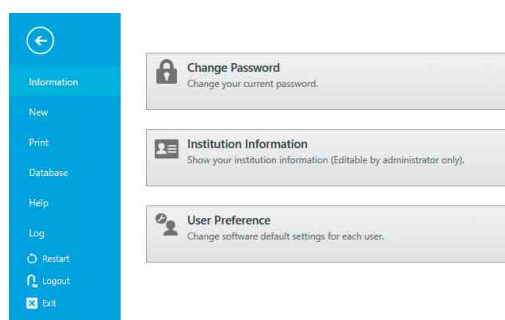

[Change Password]: Allows you to change the password used to log in to Cell Sorter Software. For details, see “*Changing Passwords*” (page 41).

[Institution Information]: Allows you to view information on the institution using the unit. You can also configure settings if you are logged in using an Administrator account. Settings cannot be configured if you are logged in using a general user account.

Click this button to display the [Institution Information] dialog (page 120).

[Account Settings]: Displays the [Account Settings] dialog (page 40) for managing user accounts for Cell Sorter Software. This button is displayed for Administrator accounts only.


[Administrative Settings]: Displays the “[Administrative Settings] dialog” (page 120) for configuring administrative settings on the instrument. This button is displayed only when logged in using an administrator account.

[User Preference]: Displays the [User Preference] dialog (page 120) for specifying software preferences, which can be specified independently by each user. This button is displayed for Administrator accounts only.

[Institution Information] dialog

The [Institute Information] dialog is displayed by clicking [Information] > [Institute Information] on the [File] tab of the ribbon.

The [Institution Information] dialog is used to specify information about the institution using the unit. Only administrators can modify this information.

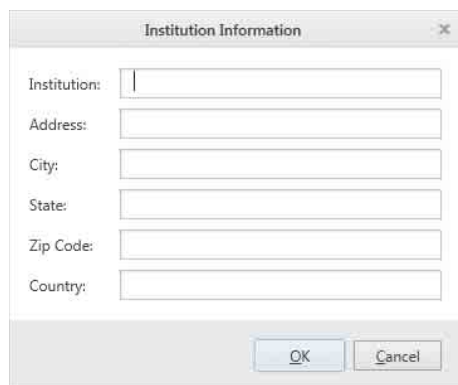

[Institution]: Enter the name of the institution.

[Address]: Enter the street address of the institution.

[City]: Enter the name of the city or town.

[State]: Enter the name of the state or province.

[Zip Code]: Enter the postal code number.

[Country]: Enter the name of the country or region.

[OK]: Applies the settings and closes the dialog.

[Cancel]: Cancels the settings and closes the dialog.


[Administrative Settings] dialog

The [Administrative Settings] dialog is displayed by clicking [Administrative Settings] in the [Information] window.

The [Administrative Settings] dialog is used to configure administrative settings on the instrument.

• [Auto Calibration]

Configures the inheritance of auto calibration results for other users.

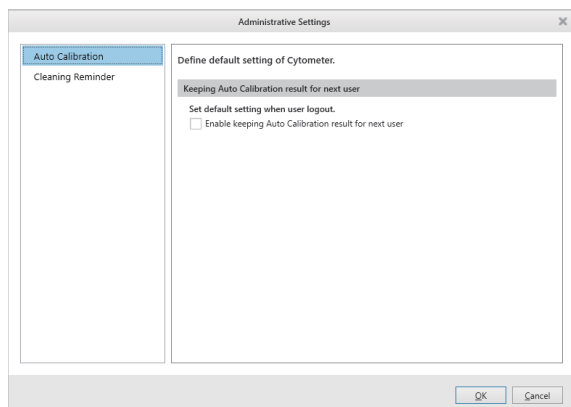


[Keeping Auto Calibration result for next user]

[Enable keeping Auto Calibration result for next user.]: Place a check mark in the checkbox to enable inheritance of auto calibration results.

• [Cleaning Reminder]

Sets the ethanol cleaning intervals for sending cleaning reminder notifications.

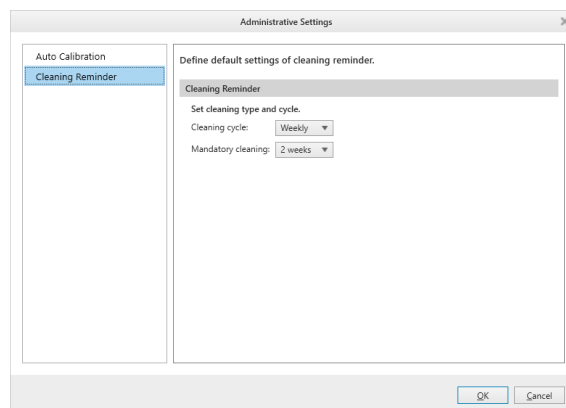

[Cleaning cycle]: Select the ethanol cleaning interval.

- None (no notifications)
- Daily
- Weekly
- Monthly (default value)

[Mandatory cleaning]: Select the mandatory ethanol cleaning interval.

If the interval specified here is exceeded, login is disabled until ethanol cleaning has been performed.

- None (no notifications) (default value)
- 1 week
- 2 weeks
- 1 month (30 days)

[User Preference] dialog

The [User Preference] dialog is displayed by clicking [Information] > [User Preference] on the [File] tab of the ribbon.


The [User Preference] dialog is used to specify software preferences, unique to each user. User preferences are loaded automatically when a user logs in, saving having to configure the instrument you want to perform acquisition, analysis, or sorting.

Tips

- You can click [Reset] to restore the default values for each setting.
- You can use the same [User Preference] settings as another user or another SH800 using [Import] / [Export].

[Sort] tab

Defines default settings related to sorting.

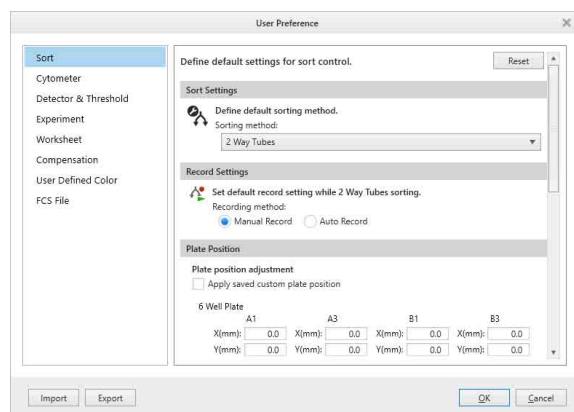

[Sort Settings]: Specifies the default sorting configuration.

[Sort method]: Selects the sorting method.

The available options vary depending on the model.

[Record Settings]

[Recording method]: Selects the recording method for 2-way sorting.

- Manual Record
- Auto Record (recording starts when sorting starts)

[Plate Position]: Adjusts the position of wells for each type of plate.

[Apply saved custom plate position]:

Applies saved plate position adjustment values.

[Side Stream Laser]

[Turn on laser during 2 way tube sorting]:

Selects whether to turn on the side-stream laser during 2-way sorting.

Tip

This item is displayed only on SH800S, SH800, SH800Z models that have been upgraded to support the 70 μ m chip.


[Custom Sort Settings]

[Edit]: Sets custom sorting modes that allow you to adjust the position of target cells within a droplet for sorting.

For details about configuration, see “Configuring Custom Sorting Mode” (page 117).

[Cytometer] tab

Defines default settings related to the instrument.

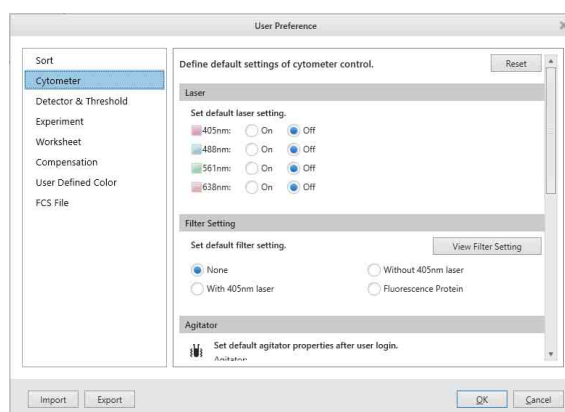

[Laser]

[Laser]: Turns the excitation lasers used for measurement on/off.

Tip

The laser settings are not inherited by the next user, even if inheritance of auto calibration results between users has been enabled. The settings follow the [User Preference] settings for those users. However, if all lasers are turned off, the laser settings are inherited.

[Filter Setting]: Selects the optical filter pattern setting.

Click [View Filter Setting] to display the selected optical filter pattern.

- None (no default setting)
- Without 405 laser (optical filter pattern 1)
- With 405 laser (optical filter pattern 2)
- Fluorescence protein

[Agitator]

[Agitator]: [Agitator]: Turns the agitation function on/off.

[Temperature Control]

[Sample Temperature Control]: Selects the temperature setting of the sample injection chamber.

- Off
- On (5 °C/41 °F)
- On (37 °C/98.6 °F)

[Collection Temperature Control]: Selects the temperature setting of the collection stage.

- Off
- On (5 °C/41 °F)

[Probe Wash]

[Skip rinse outside of sample probe with DI water.]: Sets whether to skip cleaning of the outer surface of the sample probe using DI water from the DI water tank when cleaning the probe. If a check mark is placed in this checkbox, you should load a sample

tube containing DI water, after cleaning the sample probe, to clean the outside of the probe.

Tip

When the probe is cleaned after auto calibration finishes, the outside of the probe is also cleaned, regardless of this setting.

[Area Light Control]


[Sample tube area light automatic control]: Specifies whether the sample loader light turns on automatically loading/unloading a sample.

- Off
- On

[Collection area light automatic control]: Specifies whether the collection area light turns on automatically when loading/unloading collection tubes.

[Detector & Threshold] tab

Defines default settings for the threshold value of the trigger channel and for the detectors.

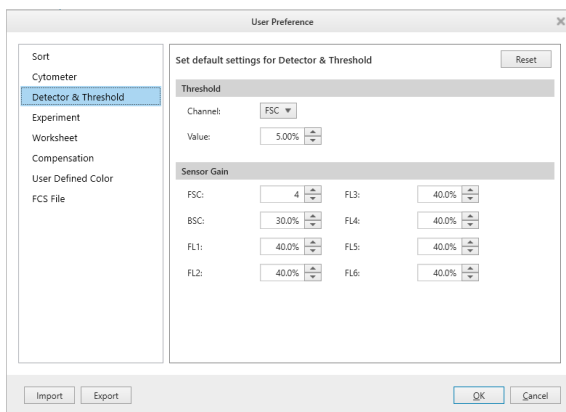

[Threshold]: Specifies the threshold for event detection.

[Channel]: Selects the channel (FSC, BSC, FL1 to FL6) used to trigger the detection of events.

[Value]: Specifies the threshold value for the trigger channel as a percentage of the output signal of the detector for the trigger channel.

[Sensor Gain]: Specifies the gain of each detector.


[FSC]: Specifies the gain of the forward-scatter (FSC) detector. The gain setting has a 16-step range.

[BSC]: Specifies the gain of the back scatter (BSC) detector as a percentage of the output signal of the detector.

[FL1] to [FL6]: Specifies the gain of the corresponding fluorescence channel detector as a percentage of the output signal of the detector.

[Experiment] tab

Defines default settings related to experiments.

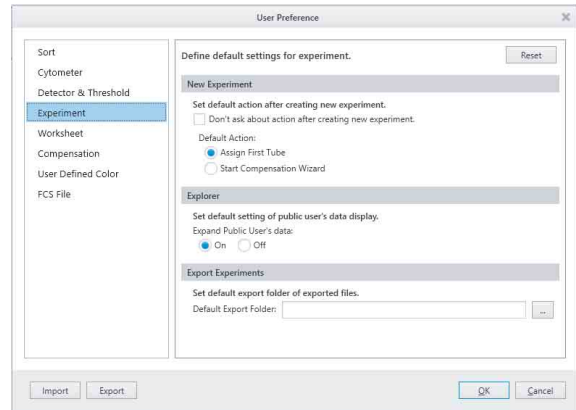

[New Experiment]

[Don't ask about action after creating new experiment.]: Sets whether to display a dialog prompting you to select the next action after creating an experiment using the Blank template.

[Default Action]: Selects the default action to take after creating an experiment.

- Assign First Tube
Assigns the first tube in preparation for data acquisition.
- Start Compensation Wizard
Select this to perform fluorescence compensation in preparation for measurements.

[Explorer]

[Expand Public User's data]: Selects whether to display the list of shared experiments.


- On
- Off

[Export Experiments]

[Default Export Folder]: Specifies the destination folder used when exporting experiment data.

[Worksheet] tab

Defines default settings related to the worksheet.

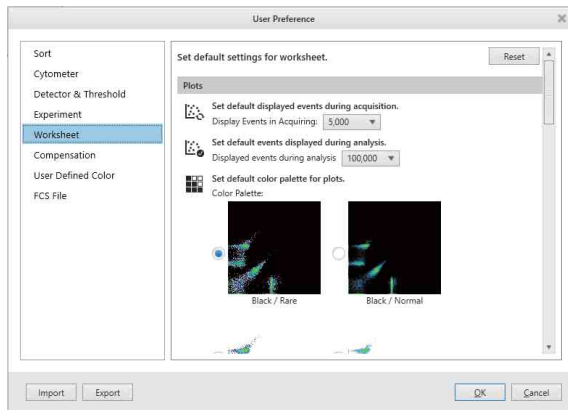

[Plots]

[Display Events in Acquiring]: Selects the number of events to display during data acquisition.

[Display events during analysis]: Selects the number of events to display when analyzing data.

[Color Palette]: Selects the color palette (background color and display of outlying (rare) events).

- Black/Rare
- Black/Normal
- White/Rare
- White/Normal

[Histogram Settings]: Makes settings for the display of histogram plots.

- For filled area view, select [No Line]. For outline view, select [Show Line]. When [Show Line] is selected, set the type of line in [Line Settings].
- [Line Settings]: Select the line thickness in [Width]. Set the opacity of the filled area in [Opacity]. You can check the type of line in [Preview].

[Gate]

[Label Font Size]: Selects the font size of text labels shown on gates. You can check the size in [Preview].

[Gate Line Width]: Selects the line width used to draw gates. You can check the line width in [Preview].

[Automatic zoom tracking function]:

Changes the shape and size of gates on plots to track the zoom size of plots and change of axis type.

[Default Scatter Plot]

[Plot Type]: Selects the default type of scatter plot added to a new worksheet.

- Density
- Dot Plot

[Axis Type]: Selects the scale type of an axis on scatter plots added to a new worksheet.

- Linear
- Log (logarithmic)
- Biexponential

[Gate type]: Selects the default gate type.

- Ellipse
- Polygon
- Rectangle

[Fluorescence Plots]

[Axis Type]: Selects the scale type of an axis on new plots displaying fluorescence data.

- Linear
- Log
- Biexponential

Tip

This setting is not applied to fluorescence plots for compensation panels.

[Index Analysis]

[Plot Size]: Specifies the size that events are displayed on plots (for emphasis) when using the index sorting function. You can check the size in [Preview].

[Alignment]

[Auto Arrange]: Selects whether to automatically align the plots and [Gates and Statistics] table to the grid of the worksheet.

- On
- Off

[Snap to Grid]: Selects whether to align the plots and [Gates and Statistics] table to the grid when moving or adding them on the worksheet.

- On
- Off

[Show Grid]: Selects whether to display the grid on the worksheet. The size of the grid is set in [Grid Size].


- On
- Off

[View]

[Expand]: Sets the zoom factor of the worksheet display.

[Compensation] tab

Defines default settings related to fluorescence compensation.

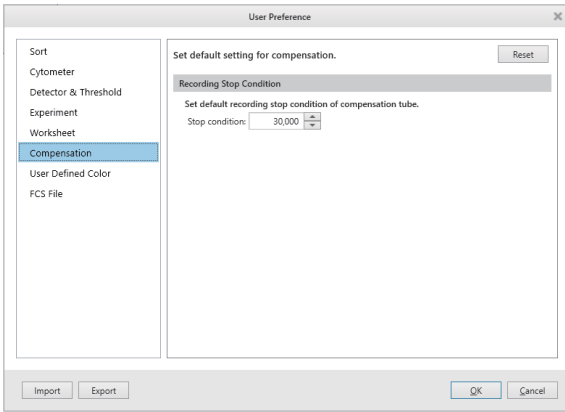


[Recording Stop Condition]

[Stop condition]: Selects the number of events before recording stops automatically.


[User Defined Color] tab

Specifies custom colors for gates.

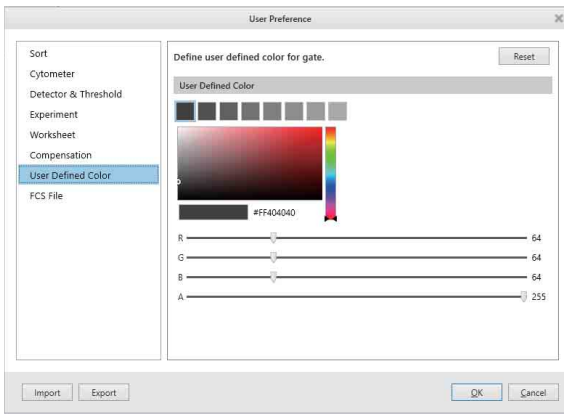

[FCS File] tab

Specifies default settings related to FCS files.

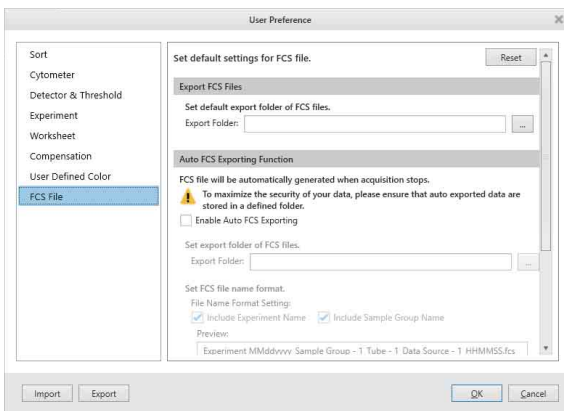

[Export FCS Files]

[Export Folder]: Specifies the destination folder used when exporting FCS files.

[Auto FCS Exporting Function]

[Enable Auto FCS Exporting]: Enables the automatic exporting of experiment data to FCS files.

Note that data security may be reduced if automatic export is enabled.

[Export Folder]: Specifies the destination folder used when exporting FCS files.

[File Name Format Setting]: Specifies the file naming format.

You can check the file name in [Preview].

- Include Experiment Name
- Include Sample Group Name

[FCS Version]: Selects the FCS file version.

- 3.0
- 3.1

[Convert BSC to SSC]: Selects whether to convert the data name of back scatter data from BSC (back scatter) to SSC (side scatter) when creating FCS files.

[Maximum Save Events]: Specifies the maximum number of events to save.


[Print] Window

Note

Printing cannot be executed during data acquisition and recording.

The [Print] window is displayed by opening the worksheet of the tube you want to print and clicking [Print] in the menu on the left side.

The [Print] window can print analysis results, statistics, and plots displayed on the worksheet.

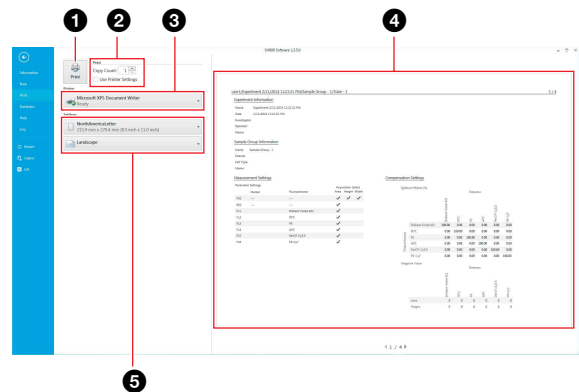

1 [Print] button

Sends the print job to the printer.

2 [Print]

Sets the number of copies to print.

Place a check mark in the [Use Printer Settings] checkbox to set advanced printer settings.

3 [Printer]

Selects the printer for printing.

4 Print preview area

Displays a preview image of the print job.

5 [Settings]


Sets the paper size and page orientation.

[Database] Window

The [Database] window is displayed by clicking [Database] in the menu on the left side.

The [Database] window is used to perform experiment database operations and to check hard disk free capacity.

Example: Administrator account login

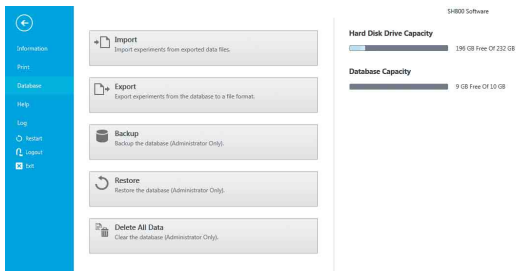

[Import]: Displays the [Import Experiment Data] dialog (page 125) for importing experiment data.

You can import experiment data created using another SH800 for use on this unit.

[Export]: Displays the [Export Experiment Data] dialog (page 125) for exporting experiment data.

You can export experiment data created on this unit for use by another SH800.

[Backup]: Displays the [Backup Database] dialog for backing up the experiment database. This button is displayed for administrator accounts only. Backing up the database allows the data to be restored at a later time, if necessary.

For details, see “Backing Up/Restoring the Database” (page 76).

[Restore]: Displays the [Restore Database] dialog for restoring a previously backed up experiment database. This button is displayed for administrator accounts only.

For details, see “Backing Up/Restoring the Database” (page 76).

[Delete All Data]: Displays the [Delete All Data] dialog (page 126) for resetting the experiment database. This button is displayed for administrator accounts only.

Note

Resetting the database removes all data, including user accounts, experiments, and analysis results.

[Hard Disk Drive Capacity]: Displays the available free space on the hard disk.

A warning will appear when the free space on the

hard disk falls below 50 GB. The assigning of sample tubes is disabled when the free space falls below 20 GB.


[Database Capacity]: Displays the available space on the database.

Up to 10 GB of data can be saved to the database in Cell Sorter Software. Various types of measurement conditions and configuration settings are saved in the database. Measurement data is not stored in the database, and has no effect on the available database capacity.

[Import Experiment Data] dialog

The [Import Experiment Data] dialog is displayed by clicking [Import] in the [Database] window.

The [Import Experiment Data] dialog is used to import data created on another SH800 into this instrument.

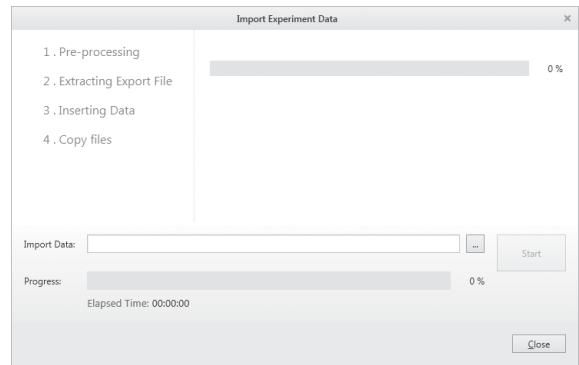

[Import Data]: Specifies the experiment data to import.

Click (Browse) to specify a file to import.

[Progress] bar: Displays the import progress.


[Start]: Starts importing the specified data.

[Close]: Closes the dialog.

[Export Experiment Data] dialog

The [Export Experiment Data] dialog is displayed by clicking [Export] in the [Database] window.

The [Export Experiment Data] dialog is used to export data created on this unit.

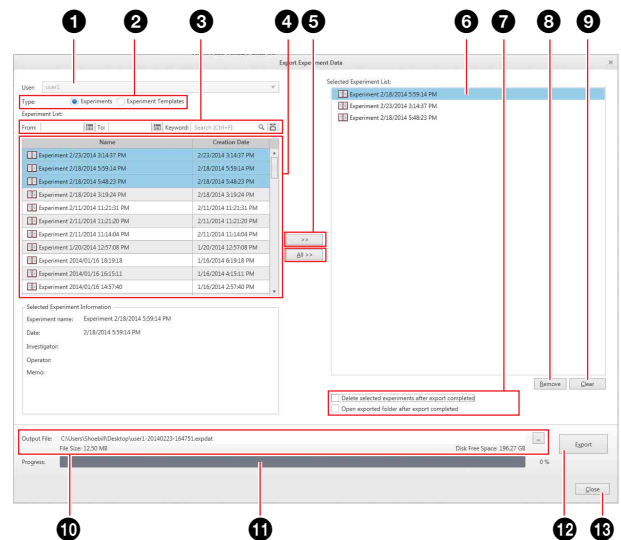

1 User

Selects the owner of the experiments to export.

2 [Type]

Selects whether to export experiments or experiment templates.

3 Search fields

Searches for experiments by date and keyword.

4 Experiment list

Displays a list of experiments that satisfy the specified search criteria.

Select the experiment(s) you want to export.

Selecting an experiment in the upper list displays information about the experiment in the lower list.

5 [>>] / [All >>]

Clicking [>>] adds the experiment data selected in the experiment data list to [Selected Experiment List].

Clicking [All >>] adds all the experiments in [Experiment List] to [Selected Experiment List].

6 [Selected Experiment List]

Lists the experiments to be exported.

7 [Delete selected experiments after export completed]

Deletes the exported experiments from the database after exporting has finished.

[Open exported folder after export completed]

Opens the export destination folder after exporting has finished.


8 [Remove]

Removes the experiment selected in [Selected Experiment List] from the list.

9 [Clear]

Clears the [Selected Experiment List].

10 [Output File]

Specifies the export destination file. Click  (Browse) to specify the file.

11 Progress

Displays the export progress.

12 [Export]

Starts exporting the experiment data.

13 [Close]

Closes the dialog.


[Delete All Data] dialog

The [Delete All Data] dialog is displayed by clicking [Delete All Data] in the [Database] window.

The [Delete All Data] dialog is used to remove all data, including user accounts and experiments.

Notes

- Resetting the database resets all data, including user accounts, experiments, and analysis results.
- After resetting the database, Cell Sorter Software automatically shuts down.

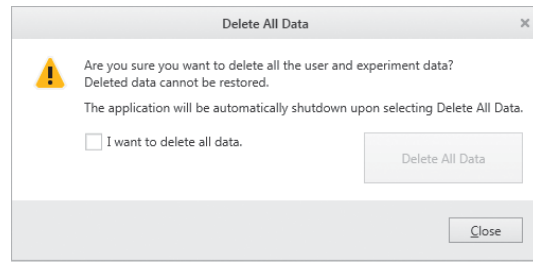

[I want to delete all data]: Place a check mark in the checkbox to confirm you want to reset the database.


[Delete All Data]: Resets the database, deleting all data.

[Close]: Closes the dialog.

[Help] Window

The [Help] window is displayed by clicking [Help] in the menu on the left side.

The [Help] window is used to display the Operator's Guide (this document) and to check software version information.

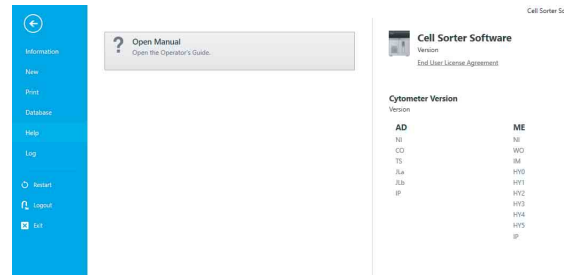

[Open Manual]: Displays the Operator's Guide.

[Cell Sorter Software]: Displays the Cell Sorter Software version information.


Clicking [End User License Agreement] displays the license agreement.

[Cytometer Version]: Displays version information for instrument components.

[Log] Window

The [Log] window is displayed by clicking [Log] in the menu on the left side.

The [Log] window is used to display the chip alignment log, sort calibration log, and user usage log.

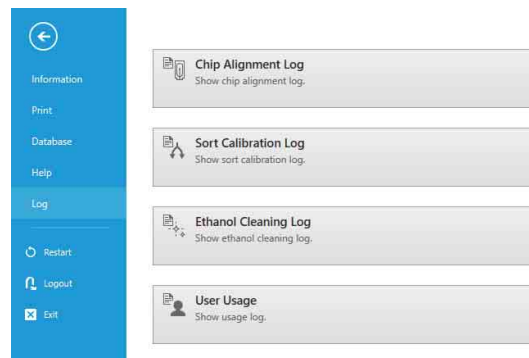

[Chip Alignment Log]: Displays the position alignment log for the sorting chip.

[Sort Calibration Log]: Displays the sort calibration log.

[Ethanol Cleaning Log]: Displays the ethanol cleaning history log.

[User Usage]: Displays the user usage log.

Chip alignment log

Displays the time and date when the sorting chip was aligned, together with other sorting chip information. Clicking a result in the [Result] column displays an image of the chip alignment result. Clicking [Export] exports the contents of the log to a CSV-format file.

Date/Time	Result	Action	User	Chip Type	Chip ID	Chip Status	Results Size	Chip Settings	Alignment Parameters
2017/04/04 09:38	Successful	Align	Administrator	Sorting Chip	20170404-0001-0001	LC-C31	70 µm	Without 405nm laser	X: 2415, Y: 4323, Z: 1000
2017/04/04 09:39	Successful	Align	Administrator	Sorting Chip	20170404-0002-0001	LC-C31	70 µm	Without 405nm laser	X: 2415, Y: 4323, Z: 1000

Ethanol cleaning log

Displays the time, date, name of user, and result for when ethanol cleaning was performed. This is displayed only when logged in using an administrator account.

Cleaning Type	Result	Start Time	End Time	Model Type	Serial Number	User
Ethanol cleaning of all fluidics lines	Successful	2017/04/04 10:31	2017/04/04 10:32	LC-S-000021C	999999	Administrator

User usage log

Displays the date and time users logged in to SH800 and other user information. Clicking [Export] exports the contents of the log to a CSV-format file.

User	Full Name	Organization	Login Date	Login Time	Usage time	Usage Sorting Chips	Usage Cleaning Chips
Administrator	Administrator	Organization	2017/04/04 10:31	2017/04/04 10:32	00:01:00	100 µm/000000-0000-0000	00:00:00
Administrator	Administrator	Organization	2017/04/04 10:31	2017/04/04 10:32	00:01:00	100 µm/000000-0000-0000	00:00:00

Sorting calibration Log

Displays the time and date when sorting was calibrated, together with other calibration information. Clicking a result in the [Result] column displays an image of the sort calibration result. Clicking [Export] exports the contents of the log to a CSV-format file.

Date/Time	Result	Action	User	Chip Type	Chip ID	Chip Status	Results Size	Chip Settings	Calibration Parameters
2017/04/04 09:38	Successful	Calibrate	Administrator	Sorting Chip	20170404-0001-0001	LC-C31	70 µm	Without 405nm laser	Droplet Clock: 240000Hz, Envelope Drive: 10.50V, Sort Delay: 33, Charge: 50.00pC, Deflection Left: 900µ, Deflection Right: 900µ

[Export Log] dialog


The [Export Log] dialog is displayed by clicking [Export] in the chip alignment log, sort calibration log, or user usage log dialogs. The [Export Log] dialog is used to export the contents of the displayed log as a CSV-format file.

[From] / [To]: Specifies the interval of the log to be exported. Click to enter a date.
[Output Folder]: Specifies the output destination folder. Click to browse for a destination folder.
Progress: Displays the export progress.
[Export]: Starts exporting the log.
[Close]: Closes the dialog.

[Create Experiment] Window

The [Create Experiment] window is displayed by clicking [New] on the [File] tab of the ribbon or [New] in the [Experiment] group on the [Experiment] tab of the ribbon.

The [Create Experiment] window is used to create a new experiment using templates.

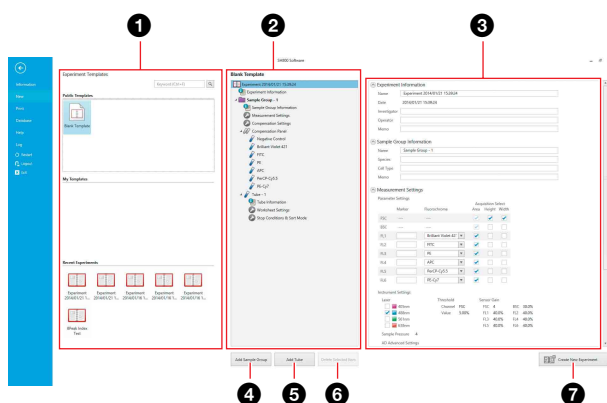

1 [Experiment Templates] list

Displays the available experiment templates and experiments that can be used as templates.

You can search the template list by keywords.

[Public Templates]: Displays experiment templates that were created as shared templates.

[Blank Template]: An empty template.

[My Templates]: Displays private experiment templates that you have created.

[Recent Experiments]: Displays recent experiments that can be used as templates.

2 Experiment structure

Displays the internal structure of the experiment template selected in the list.

You can add and delete sample groups and sample tubes, as required.

3 Experiment settings

Displays configurable information and instrument settings for the experiment.

You can also edit settings after an experiment has been created.

[Experiment Information]: Specifies the name and other basic information of the experiment. The date cannot be modified.

For details about each item, see “[Experiment Information] dialog” (page 152).

[Sample Group Information]: Specifies information relating to the sample groups.

For details about each item, see “[Sample Group Information] dialog” (page 152).

[Measurement Settings]: Specifies parameter settings and instrument settings.

For details about each item, see the “[Measurement Settings] dialog” (page 80).

[Tube Information]: Specifies information relating to the sample tubes.

For details about each item, see “[Tube Information] dialog” (page 152).

[Compensation Settings]: Displays compensation settings.

For details about each item, see “[Compensation Settings] dialog” (page 142).

[Sort & Stop Settings]: Displays the sorting and automatic stop conditions.

4 [Add Sample Group]

Displays the [Add Sample Group] dialog (page 129) for adding a sample group to the experiment structure.

5 [Add Tube]

Displays the [Add Tube] dialog (page 129) for adding a sample tube to the selected sample group in experiment structure.

6 [Delete Selected Item]

Deletes the selected item from the experiment structure.

7 [Create New Experiment]


Creates a new experiment with the specified settings.

The experiment is added to the Experiment Explorer in the main window. The created experiment is automatically made the active experiment.

[Add Sample Group] dialog

The [Add Sample Group] dialog is displayed by clicking [Add Sample Group] in the [Create Experiment] window.

The [Add Sample Group] dialog is used to add a sample group to the experiment structure.

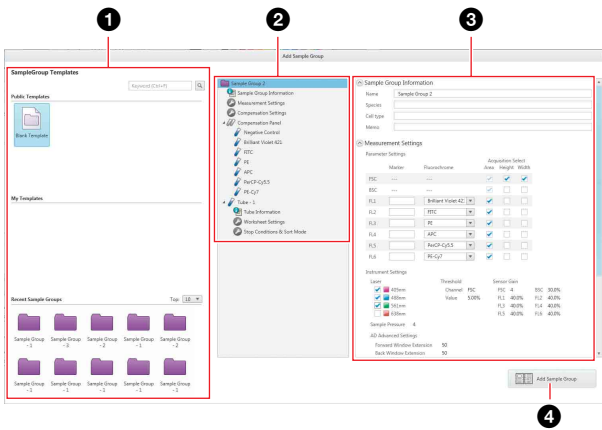

1 [Sample Group Templates] list

Displays the available sample group templates and recent sample groups that can be used as templates. You can search the template list by keywords.

[Public Templates]: Displays sample group templates that were created as shared templates.

[Blank Template]: An empty template.

[My Templates]: Displays private sample group templates that you have created.

[Recent Sample Groups]: Displays recent sample groups that can be used as templates.

2 Sample group structure

Displays the internal structure of the sample group template selected in the list.

3 Sample group settings

Displays configurable information and instrument settings for the sample group. You can also edit settings after an experiment has been created.

For details about configurations items, see “**3 Experiment settings**” (page 128) in the [Create Experiment] Window.


4 [Add Sample Group]

Adds a new sample group with the specified settings to the experiment structure.

[Add Tube] dialog

The [Add Tube] dialog is displayed by clicking [Add Tube] in the [Create Experiment] window.

The [Add Tube] dialog is used to add a sample tube to the selected sample group in the experiment structure.

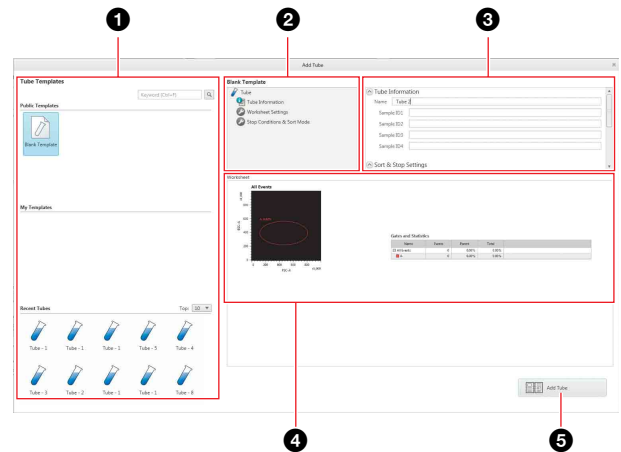

1 [Tube Templates] list

Displays the available tube templates and sample tubes that can be used as templates. You can search the template list by keywords.

[Public Templates]: Displays sample tube templates that were created as shared templates.

[Blank Template]: An empty template.

[My Templates]: Displays private tube templates that you have created.

[Recent Tubes]: Displays recent sample tubes that can be used as templates.

2 Sample tube structure

Displays the internal structure of the sample tube template selected in the list.

3 Sample tube settings

Displays configurable information and instrument settings for the sample tube. You can also edit settings after an experiment has been created.

For details about configurations items, see “**3 Experiment settings**” (page 128) in the [Create Experiment] Window.

4 Worksheet


Displays the plots and gate statistics configured for the selected template or sample tube.

5 [Add Tube]

Adds a new sample tube with the specified settings to the selected sample group in the experiment structure.

Main Window

The main window is where you acquire and record data, perform analysis, run cell sorting, and other functions. The main window has the following configuration.

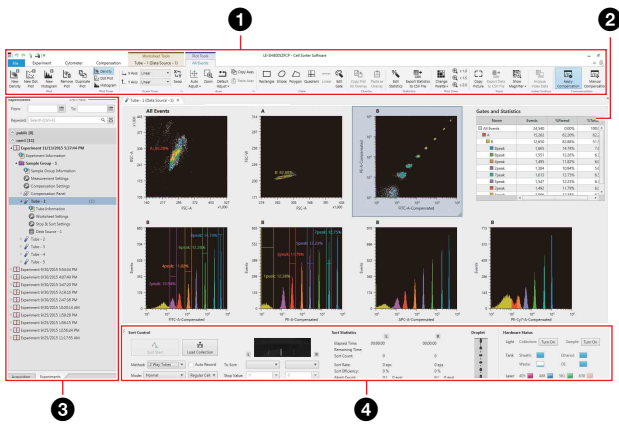

1 Ribbon

- [File] tab (page 119)
- [Experiment] tab (page 130)
- [Cytometer] tab (page 134)
- [Compensation] tab (page 141)
- [Worksheet Tools] tab (page 143)
- [Plot Tools] tab (page 146)
- [Gate Tools] tab (page 148)

2 Worksheet, [Gates and Statistics] table (page 149), context menu (page 157)

3 [Acquisition] tab (page 149) / [Experiments] tab (page 150), context menu (page 157)

4 [Sort Control] pane (page 153)

[Experiment] Tab (ribbon)

The [Experiment] tab of the ribbon has the following buttons.

[Target] group

Contains tool buttons for specifying the active sample tube used for data acquisition, recording, and sorting.




[Assign Tube]: Assigns the tube selected in the Experiment Explorer as the active tube for data acquisition, recording, and cell sorting.

[Next Tube]: Assigns the next tube. If there are no further tubes, it duplicates the current tube and assigns the new tube.

[Experiment] group

Contains tool buttons for creating, deleting, and performing other actions on experiments.

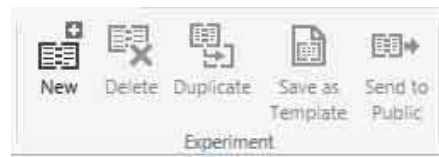

[New]: Displays the [Create Experiment] window (page 128) for creating a new experiment.

[Delete]: Deletes the experiment selected in the Experiment Explorer.


[Duplicate]: Duplicates the experiment selected in the Experiment Explorer.

[Save as Template]: Displays the [Save as Experiment Template] dialog (page 131) for saving the experiment selected in the Experiment Explorer as a template used when creating a new experiment.

[Send to Public]: Creates a duplicate of the selected experiment and places the duplicate under the [Public] node for sharing with other users.

[Sample Group] group

Contains tool buttons for creating, deleting, and performing other actions on sample groups.

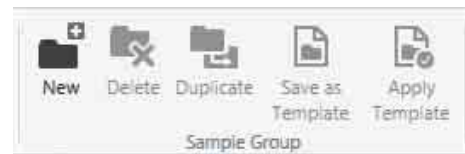

[New]: Creates a new, blank sample group in the experiment selected in the Experiment Explorer.

[Delete]: Deletes the sample group selected in the Experiment Explorer.


[Duplicate]: Duplicates the sample group selected in the Experiment Explorer.

[Save as Template]: Displays the [Save as Sample Group Template] dialog (page 131) for saving the sample group selected in the Experiment Explorer as a template used when creating a new experiment.

[Apply Template]: Displays the [Apply Template] dialog (page 132) for selecting a template for a new sample group selected in the Experiment Explorer.

[Tube] group

Contains tool buttons for creating, deleting, and performing other actions on sample tubes.

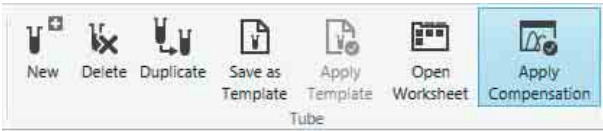

[New]: Creates a new, blank sample tube in the sample group selected in the Experiment Explorer.

[Delete]: Deletes the sample tube selected in the Experiment Explorer.

[Duplicate]: Duplicates the sample tube selected in the Experiment Explorer.

[Save as Template]: Displays the [Save as Tube Template] dialog (*page 131*) for saving the sample tube selected in the Experiment Explorer as a template used when adding a sample tube.

[Apply Template]: Displays the [Apply Template] dialog (*page 132*) for selecting a template for a new sample tube selected in the Experiment Explorer.

[Open Worksheet]: Opens the worksheet tab for the tube selected in the Experiment Explorer.

[Apply Compensation]: Applies fluorescence compensation to the plots on the worksheet for the tube selected in the Experiment Explorer.

[Settings and Information] group

Contains tool buttons for copying, pasting, and displaying settings and information.



[Copy] / [Paste]: Copies/Pastes the item selected in the Experiment Explorer.

[Apply to All Tubes]: Applies the worksheet settings for the tube selected in the Experiment Explorer to all sample tubes within the sample group.

Click [Worksheet Settings] for a tube in the Experiment Explorer and then click [Apply To All Tubes] to apply the same worksheet settings to the other tubes in the sample group for easy comparison during analysis.

[Show Information]: Displays the information dialog for the item ([Experiment Information], [Sample Group Information], etc.) selected in the Experiment Explorer.

You can also display the settings dialog by double-clicking [Experiment Information] and other items in the Experiment Explorer.

[Show Settings]: Displays the configuration dialog for the settings item ([Measurement Settings], [Compensation Settings], etc.) selected in the Experiment Explorer.


You can also display the settings dialog by double-clicking [Experiment Information] and other items in the Experiment Explorer.

[Show Results]: Displays the [Tube Results] dialog (*page 132*) for displaying the recording results and sorting results for the tube selected in the Experiment Explorer.

If [8 Well Strips], [Slides], [96 Well Plate], or [384 Well Plate] sorting was performed, the data source results that were recorded last are also displayed. If you click this button while a data source is selected, the [Data Source Results] dialog (*page 133*) appears, allowing you to view the results of each data source.

[Export] group

Contains a tool button for exporting recorded data for a sample tube.

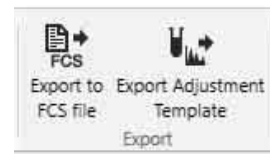

[Export to FCS Files]: Displays the [FCS File Export] dialog (*page 134*) for exporting the data recorded for the tubes in the experiment selected in the Experiment Explorer in FCS 3.0 or FCS 3.1 format.

Tip


The importing of FCS files is not supported in Cell Sorter Software.

[Export Adjustment Template]: Displays the [Export Adjustment Template] dialog (*page 87*) for creating a gain adjustment template (experiment template with reference data for instrument settings) from the experiment selected in the Experiment Explorer.

[Save as Experiment/Sample Group/Tube Template] dialog

The [Save as Experiment/Sample Group/Tube Template] dialog is displayed by selecting an experiment, sample group, or sample tube in the Experiment Explorer and clicking [Save as Template] on the [Experiment] tab of the ribbon or by right-clicking and selecting [Save as Experiment Template]/[Save as Sample Group Template]/[Save as Tube Template] from the context menu.

The [Save as Experiment/Sample Group/Tube Template] dialog is used to save the selected experiment, sample group, or sample tube as a template.

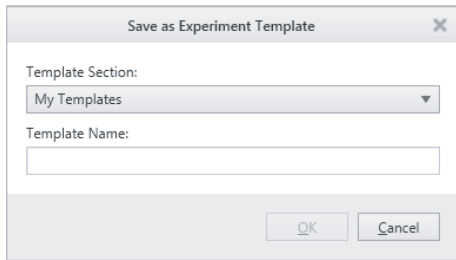


For details, see “Saving an Experiment as a Template” (page 85).

[Apply Template] dialog

The [Apply Template] dialog is displayed by selecting a new sample group or sample tube in the Experiment Explorer and clicking [Apply Template] on the [Experiment] tab of the ribbon.

The [Apply Template] dialog is used to select a template to apply to the sample group or sample tube.

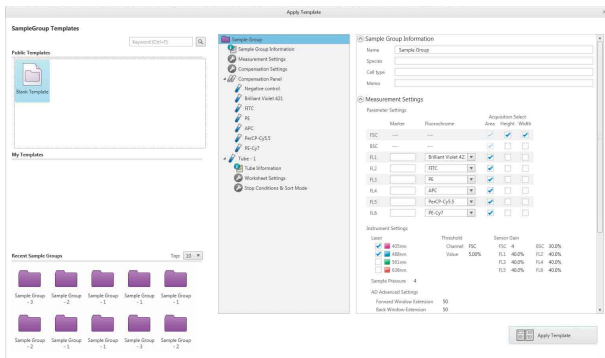

The configurable items are the same as the [Add Sample Group] dialog (page 129) or [Add Tube] dialog (page 129) for sample groups.


[Tube Results] dialog

The [Tube Results] dialog is displayed by selecting a sample tube in the Experiment Explorer and clicking [Show Results] in the [Settings and Information] group on the [Experiment] tab of the ribbon.

The [Tube Results] dialog is used to check the recording results and sorting results of the selected sample tube.

• [Basic Information] tab

Displays basic information identifying the sample tube.

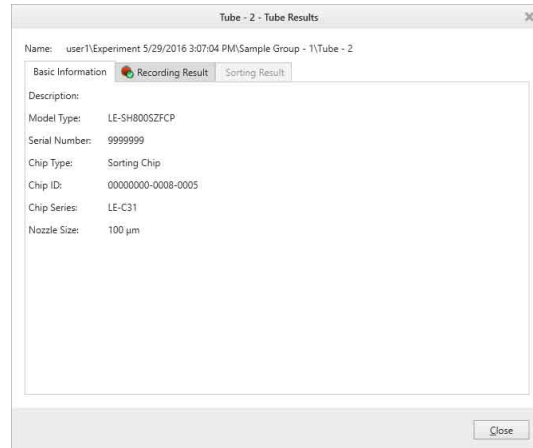


[Chip ID]: Displays the ID number of the sorting chip.

[Chip Series]: Displays the series name of the sorting chip.

[Nozzle Size]: Displays the nozzle size of the sorting chip.

• [Recording Result] tab

Displays recording results.

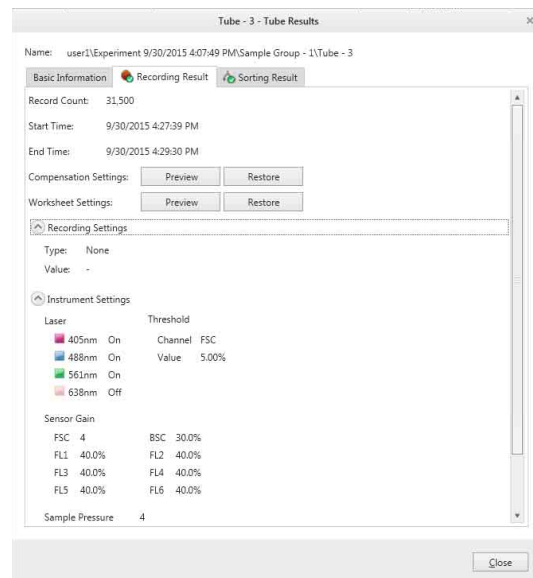

[Record Count]: Displays the number of events recorded.

[Start Time]: Displays the time when recording started.

[End Time]: Displays the time when recording ended.

[Compensation Settings]: Previews/restores fluorescence compensation settings, saved when data recording ends.

- **[Preview]:** Displays the fluorescence compensation settings.

- **[Restore]:** Restores the fluorescence compensation settings. Click [Yes] in the confirmation dialog to proceed.

[Worksheet Settings]: Previews/restores worksheet settings, saved when data recording ends.


- [Preview]: Displays the worksheet settings.
- [Restore]: Restores the worksheet settings. Click [Yes] in the confirmation dialog to proceed.

[Recording Settings]: Displays the recording settings.

[Instrument Settings]: Displays the instrument settings.

• **[Sorting Result] tab**

Displays the sorting results.

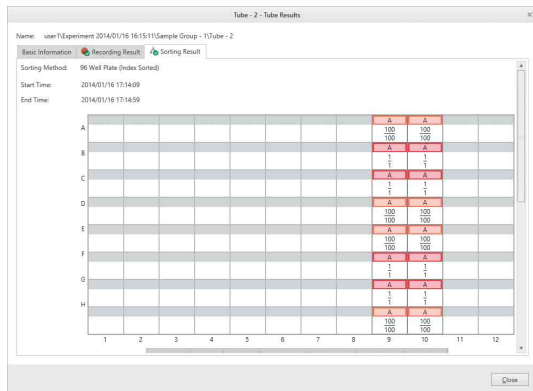

[Method]: Displays the sorting method.

[Start Time]: Displays the time when recording started.

[End Time]: Displays the time when recording ended.

[Sort Gate]: Displays the gate specified for sorting.

[Sort Mode]: Displays the sorting mode specified for sorting.

[Elapsed Time]: Displays the elapsed time since starting sorting.

[Total Events]: Displays the number of events detected during sorting.

[Target Ratio]: Displays the percentage of events detected during sorting that fall within the sort gate.

[Sorted Count]: Displays the number of events sorted.

[Sort Rate]: Displays the rate at which events are sorted in units of events per second (eps).

[Sort Efficiency]: Displays the number of sorts attempted as a percentage of targeted cells in the sample.

[Aborted Count]: Displays the number of target events in the sort gate that do not match the sort mode criteria and which are aborted during sorting.

[Abort Rate]: Displays the rate at which events are aborted during sorting in units of events per second (eps).

[Sort Result]: Indicates whether sorting is finished or not.

[Copy to Clipboard]: Copies the sorting results data to the clipboard as an image for pasting in other software, for example to produce reports.

[Save as CSV]: Saves the sorting results as CSV-format file.

[Compensation Settings]: Previews/restores fluorescence compensation settings, saved when sorting ends.

- [Preview]: Displays the fluorescence compensation settings.

- [Restore]: Restores the fluorescence compensation settings. Click [Yes] in the confirmation dialog to proceed.

[Worksheet Settings]: Previews/restores worksheet settings, saved when sorting ends.

- [Preview]: Displays the worksheet settings.

- [Restore]: Restores the worksheet settings. Click [Yes] in the confirmation dialog to proceed.


[Sort Settings]: Displays the sort settings.

[Instrument Setting]: Displays the instrument settings.

[Auto Parameters]: Displays automatically configured sorting parameters.

[Data Source Results] dialog

The [Data Source Results] dialog is displayed by selecting a data source in the Experiment Explorer and clicking [Show Results] in the [Settings and Information] group in the [Experiment] tab of the ribbon. The [Data Source Results] dialog allows you to view the results of each data source.

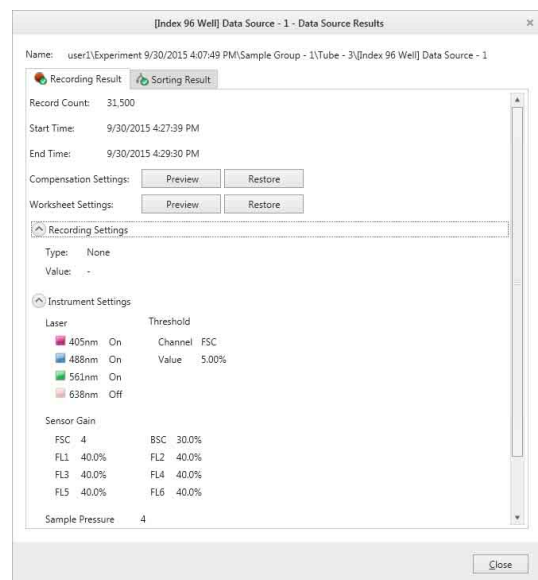


The [Data Source Results] dialog displays all the information from the [Tube Results] dialog, except the [Basic Information] tab.

For details, see “[Tube Results] dialog” (page 132).

[FCS File Export] dialog

The [FCS File Export] dialog is displayed by selecting an experiment in the Experiment Explorer and clicking [Export to FCS file] in the [Export] group on the [Experiment] tab of the ribbon or by right-clicking and selecting [Export FCS Files] from the context menu.

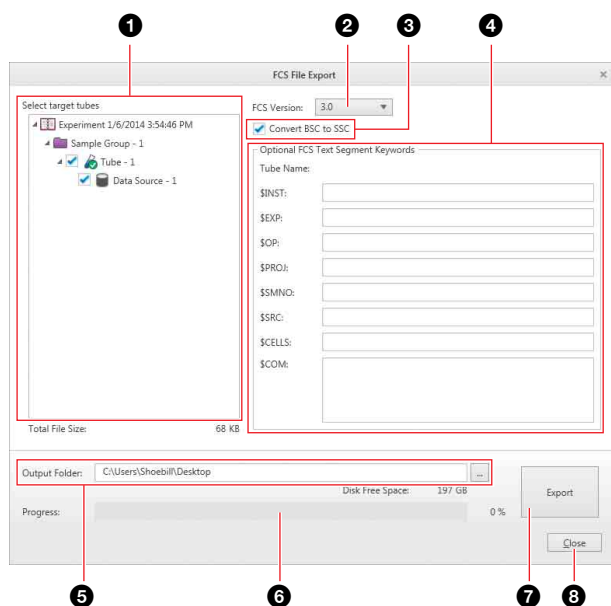

1 [Select target tubes]

Selects the sample tubes to export. Place a check mark in the checkbox for each target tube.

2 [FCS Version]

Selects the FCS file version (3.0 or 3.1).

3 [Convert BSC to SSC]

Selects whether to convert the data name of back scatter data from BSC (back scatter) to SSC (side scatter) when creating FCS files.


4 [Optional FCS Text Segment Keywords]

You can specify optional FCS text segment keywords, as required. Required FCS text segment keywords are exported automatically.

The following optional FCS text segment keywords can be specified.

- \$INST: Institution at which data was acquired
- \$EXP: Name of investigator initiating the experiment
- \$OP: Name of instrument operator
- \$PROJ: Name of the experiment project
- \$SMNO: Specimen (e.g., tube) label
- \$SRC: Source of the specimen (patient name, cell type)
- \$CELLS: Description of objects measured
- \$COM: Comment

5 [Output Folder]

Specifies the export destination folder. Click  (Browse) to specify the destination folder. The files are named in “<tube_name>.fcs” format.

6 Progress

Displays the export progress.

7 [Export]

Starts the export.

8 [Close]


Closes the dialog.

[Cytometer] Tab (ribbon)

The [Cytometer] tab of the ribbon has the following buttons.

[Load/Unload] group

Contains tool buttons for loading and unloading sample tubes and the collection tubes.

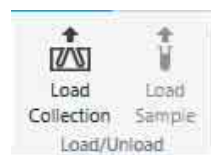


[Load Collection/Unload Collection]: Loads/unloads the collection tube in the collection area.

[Load Sample/Unload Sample]: Loads/unloads the sample tube in the sample loader.

[Light] group

Contains tool buttons for turning the collection stage and sample loader lights on/off.

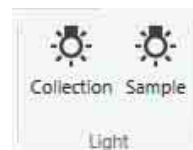

[Collection]: Turns the collection stage light on/off.

[Sample]: Turns the sample loader light on/off.

[Cleaning] group

Contains tool buttons for cleaning the fluidics system.



[Probe Wash]: Washes the sorting chip and the inner surface of the sample probe using sheath fluid, and washes the outer surface of the sample probe using DI water.

Tip

The outer surface of the sample probe is not washed when [Skip rinse outside of sample probe with DI water] is enabled.

[Bleach Cleaning]: Displays the [Bleach Cleaning] dialog (page 137) for cleaning the sorting chip, sample line, and sample probe with a Sodium Hypochlorite solution.

For details about cleaning using Sodium Hypochlorite solution, see “Cleaning the Sample Fluidics System using Bleach” (page 170).

[DI Rinse]: Displays the [DI Rinse] dialog (page 137) for flushing the sorting chip, sample line, and sample probe with DI water.

[Shutdown Rinse]: Displays the [Shutdown DI Rinse] dialog (page 137) for flushing the sorting chip, sample line, and sample probe with DI water before shutting down the instrument.

Note

After cleaning, the fluidics lines are filled with DI water to prevent the formation of salt deposits. Running data acquisition in this state may damage cells.

[De-bubble] group

Contains tool buttons for removing air bubbles from the sorting chip and sheath filter.



[Chip]: Removes air bubbles from the sorting chip.


[Sheath Filter]: Displays the [Sheath Filter De-bubble] dialog (page 137) for removing air bubbles from the sheath filter.

Tip

The sheath filter de-bubble function should be run at least three times after changing the sheath filter.

[Auto Calibration] group

Contains tool buttons for aligning the chip and calibrating sorting.

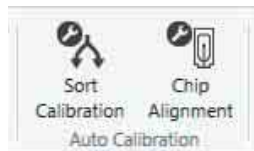

[Sort Calibration]: Displays the [Auto Calibration] wizard for calibrating sorting for performing droplet calibration, side stream calibration, and sort delay calibration automatically.

Run sorting calibration after inserting a sorting chip and the sorting chip alignment is successful.

For details, see “Automatic Calibration” (page 50).

Tip

Clicking the [Sort Calibration] button after inserting a sorting chip and while chip alignment is not completed successfully will display a warning message and sorting calibration will not be executed.

[Chip Alignment]: Displays the [Chip Alignment] screen of the [Auto Calibration] wizard for aligning the sorting chip.


Follow the on-screen instructions to align the sorting chip. During automatic alignment, the progress bar and status are displayed, allowing you to check the progress and current state.

[Temperature Control] group

Contains tool buttons for setting the temperature of the collection stage and sample injection chamber.


The following temperature control options are available.

- Collection stage: 5 °C (41 °F) only
- Sample injection chamber: 5 °C (41 °F) or 37 °C (98.6 °F)

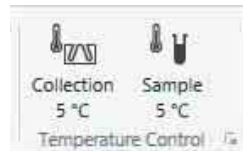

[Collection]: Turns the collection temperature control on/off and displays the current temperature setting.

[Sample]: Turns the sample injection chamber temperature control on/off and displays the current temperature setting.

Clicking the  dialog launcher displays the [Temperature Control Settings] dialog (page 137) for specifying the temperature settings.

[Agitate] group

Contains a tool button for controlling agitation of the sample fluid.



[Agitate]: Turns the sample loader agitation unit on/off.

[Safety] group

Contains a tool button for running a suction fan for safety.



[Vacuum Fan]: Displays the [Vacuum Fan Control] dialog (page 138) for setting a timer for operation of the vacuum fan.

The fan is used to extract aerosols (microparticles of biological and other matter in suspension created by sorting) from the collection area. Setting the duration to “0” runs the fan without the timer function.

[Exchange] group

Contains command buttons for changing the fluidics system components.



[Chip]: Ejects the sorting chip and displays the [Chip Exchange] wizard for inserting a different sorting chip.

For details, see “Exchanging the Sorting Chip” (page 73).

[Sample Line]: Displays the [Sample Line Exchange] wizard for changing the sample line.


For details, see “Changing the PEEK Sample Line (PEEK sample line compatible models)” (page 189) or “Changing the Sample Line and Probe (Non-PEEK sample line compatible models)” (page 192).

Tip

The sorting chip must be replaced or reinserted after changing the sample line.

[Settings] group

Contains tools button for configuring advanced settings related to the instrument.

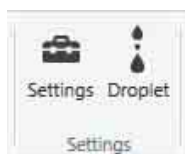


[Settings]: Displays the [Cytometer Settings] dialog (page 138) for configuring the instrument, adjusting sort parameters, and other instrument functions.

[Droplet]: Displays a detailed image of the droplets.

[Help] group

Contains a tool button for displaying the optical filter pattern.

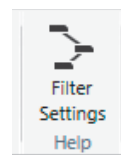

[Filter Settings]: Displays the configured optical filter pattern. This cannot be modified.

[Shutdown] group

Contains tool buttons used when shutting down the SH800.



[Hardware Only]: Displays the [Shutdown Wizard] screen for cleaning the SH800 and then shutting down the instrument only.

For details, see “Shutdown” (page 71).

[Software and Hardware]: Displays the [Shutdown Wizard] screen for cleaning the SH800 shutting down the instrument and Cell Sorter Software.

For details, see “Shutdown” (page 71).

[Ethanol Cleaning]: Displays the Ethanol Cleaning wizard for cleaning the internal sheath fluid line and DI water line with ethanol. The SH800 instrument and Cell Sorter Software automatically shut down after cleaning is finished.

For details, see “Cleaning the Internal Sheath Line and DI Water Line using Ethanol” (page 209).

Note

Use the following cleaning procedure to perform ethanol cleaning after sample measurement.


- ① Clean the sample line and sample probe using bleach solution (see “Cleaning the Sample Fluidics System using Bleach” (page 170)).
- ② Click [Shutdown Rinse] in the [Cleaning] group on the [Cytometer] tab of the ribbon to flush the system with DI water before shutdown (see “[Shutdown DI Rinse] dialog” (page 137)).
- ③ Click [Ethanol Cleaning] to clean the sheath line using ethanol (see above).

[Bleach Cleaning] dialog

The [Bleach Cleaning] dialog is displayed by clicking [Bleach Cleaning] in the [Cleaning] group on the [Cytometer] tab of the ribbon.

The [Bleach Cleaning] dialog is used to start cleaning of the sorting chip, sample line, and sample probe using a Sodium Hypochlorite (bleach) solution.

For details, see “Cleaning the Sample Fluidics System using Bleach” (page 170).

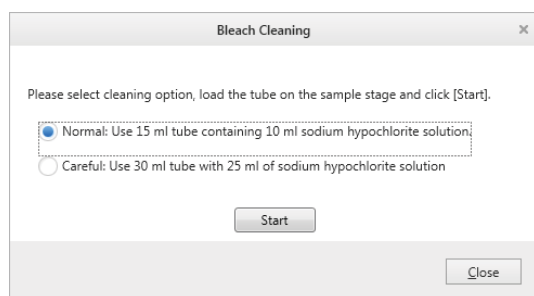


[DI Rinse] dialog

The [DI Rinse] dialog is displayed by clicking [DI Rinse] in the [Cleaning] group on the [Cytometer] tab of the ribbon.

The [DI Rinse] dialog is used to start flushing of the sorting chip, sample line, and sample probe using DI water.

Select the size of conical tube containing DI water to use, then click [Start] to load the tube and begin cleaning.

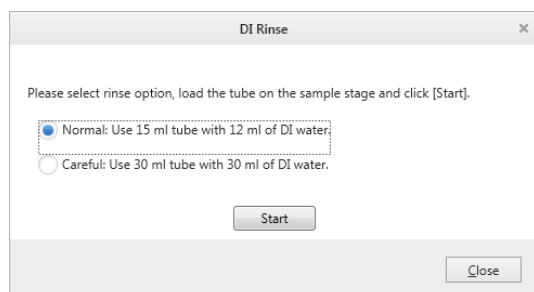


[Shutdown DI Rinse] dialog

The [Shutdown DI Rinse] dialog is displayed by clicking [Shutdown Rinse] in the [Cleaning] group on the [Cytometer] tab of the ribbon.

The [Shutdown DI Rinse] dialog is used to clean the sorting chip, sample line, and sample probe using DI water, and then automatically shutting down.


Select the size of conical tube containing DI water to use, then click [Start] to load the tube and begin cleaning.

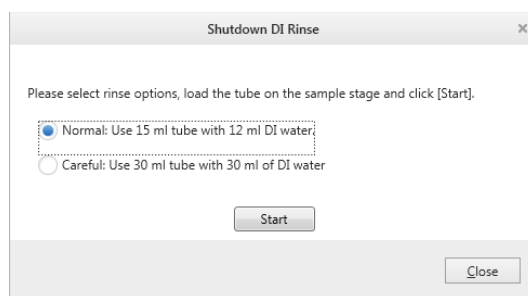

[Sheath Filter De-bubble] dialog

The [Sheath Filter De-bubble] dialog is displayed by clicking [Sheath Filter] in the [De-bubble] group on the [Cytometer] tab of the ribbon.


The [Sheath Filter De-bubble] dialog is used to remove air bubbles from the sheath filter. Click [Start] to begin the process.

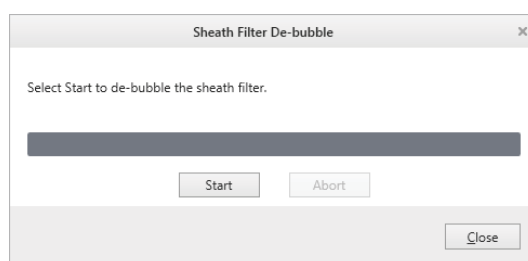


Tip

The sheath filter de-bubble function should be run at least three times after changing the sheath filter.

[Temperature Control Settings] dialog

The [Temperature Control Settings] dialog is displayed by clicking the  dialog launcher in the [Temperature Control] group on the [Cytometer] tab of the ribbon. The [Temperature Control Settings] dialog is used to set the temperature of the sample injection chamber and collection stage area.

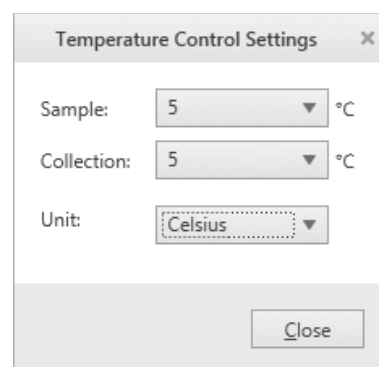

[Sample]: Sets the control temperature of the sample injection chamber to 5 °C (41 °F) or 37 °C (98.6 °F). The selected temperature values are displayed on the corresponding buttons in the [Temperature Control] group on the ribbon.

[Collection]: Selects the control temperature of the collection stage area. Only 5 °C (41 °F) is available.

[Unit]: Selects the units of temperature (Celsius or Fahrenheit).


[Close]: Closes the dialog.

[Vacuum Fan Control] dialog

The [Vacuum Fan Control] dialog is displayed by clicking [Vacuum Fan] in the [Safety] group on the [Cytometer] tab of the ribbon.

The [Vacuum Fan Control] dialog is used to set a timer for operation of the vacuum fan.

The fan is used to extract aerosols (microparticles of biological and other matter in suspension created by sorting) from the collection area.

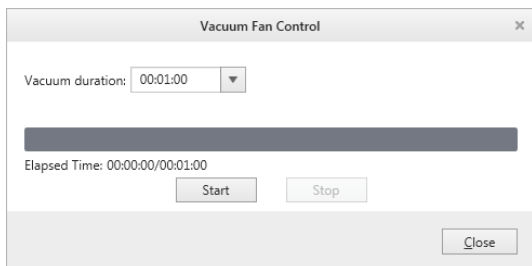

[Vacuum duration]: Set the duration for fan operation.

Setting the duration to “00:00:00” allows you to operate the fan for an unlimited duration. When you want to stop the fan in such cases, click [Stop].

[Elapsed Time]: Displays the elapsed time since starting the fan.

[Start]: Starts the fan.

[Stop]: Stops the fan.

[Close]: Closes the dialog.


[Cytometer Settings] dialog

The [Cytometer Status] dialog is displayed by clicking [Settings] in the [Settings] group on the [Cytometer] tab of the ribbon.

The [Cytometer Settings] dialog is used to configure instrument settings, adjust sort parameters, and other functions.

• [Control] tab

This tab is used to configure instrument settings.

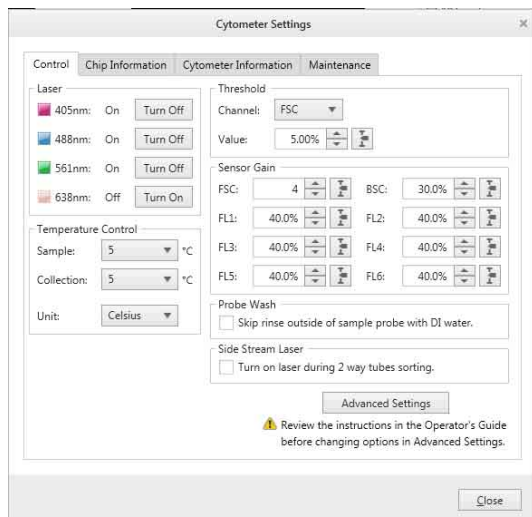

[Laser]: Turns each laser on/off.

[Temperature Control]: Sets the control temperature of the sample injection chamber and collection stage.

For details about each item, see “[Temperature Control Settings] dialog” (page 137).

[Threshold]: Sets the threshold trigger for detecting events.

- [Channel]: Selects the channel (FSC, BSC, FL1 to FL6) used to trigger the detection of events.
- [Value]: Sets the input threshold signal level of the trigger channel as a percentage of the maximum signal from the detection module.

[Sensor Gain]: Sets the gain of each detector.

- [FSC]: Selects the gain of the forward-scatter detector. The gain setting has a 16-step range.
- BSC, FL1 to FL6: Sets the gain of the corresponding back scatter and fluorescence detectors as a percentage of the maximum level.

[Probe Wash]: Sets probe wash options. To skip cleaning of the outer surface of the sample probe using DI water from the DI water tank when cleaning the probe, place a check mark in the [Skip rinse outside of sample probe with DI water.] checkbox. If a check mark is placed in this checkbox, you should load a sample tube containing DI water, after cleaning the sample probe, to clean the outside of the probe.

[Side Stream Laser]: Sets whether to turn the laser on during 2-way sorting.


Tip

This item is displayed only on SH800S, SH800, SH800Z models that have been upgraded to support the 70 µm chip.

[Advanced Settings]: Displays the [Advanced Settings] dialog (page 139) for manually adjusting sort parameters.


- **[Chip Information] tab**

Displays information about the loaded sorting chip. This is useful for checking the length of time the sorting chip has been in use.

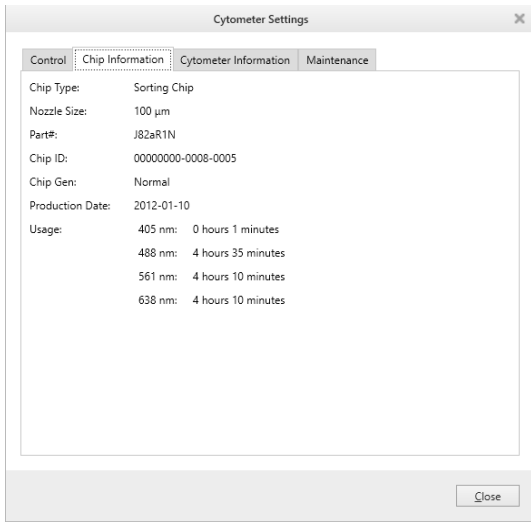


- **[Cytometer Information] tab**

Displays SH800 instrument information.

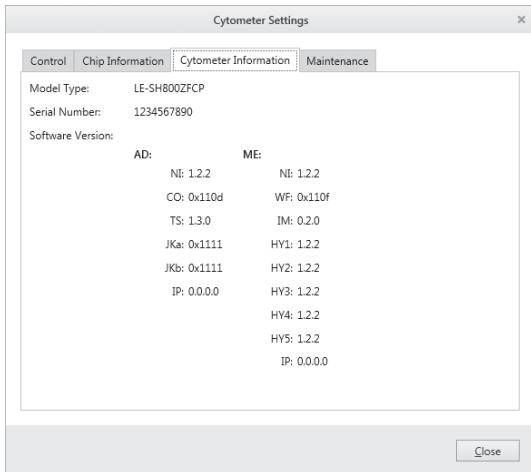

- **[Maintenance] tab**

Displays a function for de-bubbling the DI water filter and displays the laser usage time.

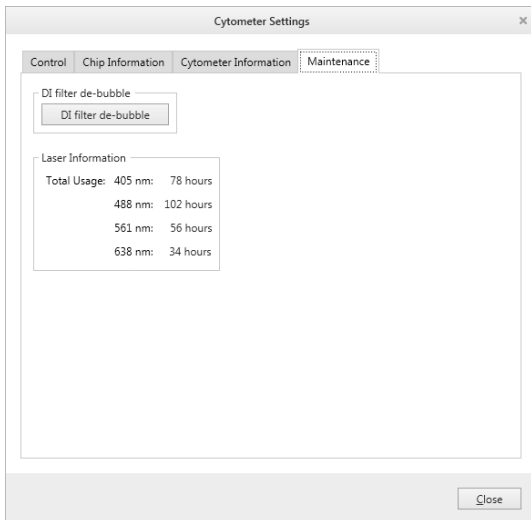

[DI filter de-bubble]: De-bubbles the DI water filter.

Click [DI Filter de-bubble] to display the [DI Filter De-bubble] dialog, and then click [Start].

Tip


The DI water filter de-bubble function should be run at least three times after changing the DI water filter.

For details, see “Changing the DI Water Filter” (page 198).

[Laser Information]: Displays the usage time of each laser.

- **[Advanced Settings] dialog**

The [Advanced Settings] dialog is used to adjust sort parameters manually.

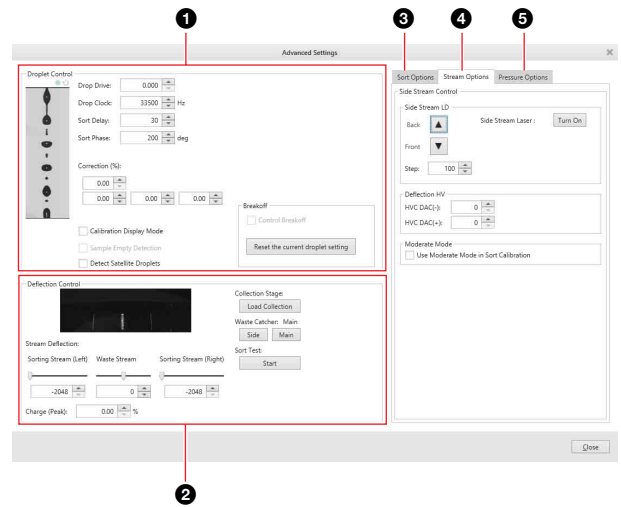

1 [Droplet Control]

Contains adjustments related to droplet formation.

[Drop Drive]

The drop drive setting adjusts the amplitude of the piezoelectric element drive voltage. The frequency sets the cycle time of vibration of the sorting chip. The vibration of the sorting chip causes the stream to break into droplets after the stream is emitted from the chip nozzle. The drop drive is optimized during auto calibration.

[Drop Clock]

The drop clock setting adjusts the frequency of the vibration of the piezoelectric element drive voltage. The frequency sets the cycle time of vibration of the sorting chip. The drop drive is optimized during auto calibration.

[Sort Delay]

The sort delay setting adjusts the number of vibrations applied to the sorting chip as cells pass between the optical detection point and the breakoff point. It is used to determine the timing of the electrostatic charging pulse applied to the sheath fluid.

The sort delay is optimized during auto calibration.

[Sort Phase]

The sort phase setting adjusts the phase difference of the charging pulse waveform as cells pass between the optical detection point and the breakoff point. It is used to determine the timing of the electrostatic charging pulse applied to the sheath fluid.

The sort phase is optimized during auto calibration.

[Correction (%)]

The correction setting adjusts the shape of the charging waveform applied to successive droplets.

- Top spinbox: Controls the overall charging waveform correction adjustment.
- Lower spinboxes: Fine-tunes the correction applied to the first droplet, second droplet, and third droplet, respectively.

When droplets are electrostatically charged, the next droplet receives an induced charge of the opposite polarity, which can affect the angle of deflection as the droplets pass through the deflection plates (called fanning). A defanning algorithm adjusts the charging waveform applied to successive droplets to reduce fanning.

[Calibration Display Mode]

The calibration display mode changes the image displayed by the droplet camera so that only droplets containing automatic setup beads are displayed.

This mode is used for fine manual control of the sort delay, sort phase, and related sort parameters.

[Reset the current droplet setting]

Saves the current droplet image as a reference image used by the Control Breakoff function. Clicking this button displays a confirmation message. Click [Yes] to save the current droplet image.

[Control Breakoff]

The control breakoff setting uses a feedback system to control the breakoff position automatically.

When this function is enabled, the drop drive is adjusted so as to maintain the optimized breakoff point and side streams obtained during auto calibration.

When auto calibration finishes, the Control Breakoff function is automatically enabled.

Tips

- The drop drive and drop clock cannot be adjusted manually while the Control Breakoff function is enabled.
- When breakoff feedback control is enabled, if the breakoff point moves and cannot be recovered within a predetermined time, sorting stops and an error message is displayed on the screen.

[Sample Empty Detection]

The empty detection function automatically stops data acquisition when the sample tube becomes empty. The function detects disturbances in the droplet stream as the sample fluid in the sample tube is used up, and then stops data acquisition.

Tips

- The sample empty detection function may stop data acquisition if cells of size 20 μm or larger are in the sample. If this occurs, disable the sample empty function to continue data acquisition.
- If a sample empty state is detected, click [Chip] in the [De-bubble] group on the [Cytometer] tab of the ribbon in Cell Sorter Software, and repeatedly de-bubble the sorting chip until the breakoff point returns to its original position.

[Detect Satellite Droplets]

The satellite droplet detection function detects irregular shaped droplets during sorting. When the function is enabled, disturbances in the droplets during sorting are detected, and sample flow is paused.

Tips

- In rare cases, the satellite droplet detection function may cause events to be aborted by mistake. If this occurs, disable the satellite droplet detection function to continue data acquisition.
- Satellite droplet detection is supported only under the following conditions.
 - When using a 70 μm sorting chip
 - When using an LE-C32 series chip with “Targeted” setting

2 [Deflection Control]

Contains controls for adjusting the side stream deflection angle, loading/unloading the collection stage, and running a sorting test.

[Stream Deflection]

Adjusts the angle of deflection of the two side streams.

[Charge (Peak)]

Sets the peak value of the charging voltage waveform.

[Collection Stage]

Loads/unloads the collection stage.

[Waste Catch]


Moves the collection area waste catcher to the side.

[Sort Test]

Runs a sorting test.

3 [Sort Options] tab

The sort option settings are used to adjust the position of the sorting chip, droplet image, and collection stage manually.

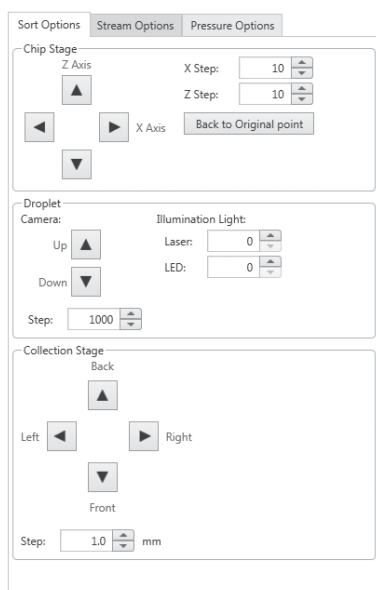


[Chip Stage]: Adjusts the positioning of the sorting chip within the chip loader.

[Droplet]: Adjusts the vertical positioning of the droplet camera, laser for droplet image fluorescence detection, and the brightness of the infrared LED for droplet imaging.

[Collection Stage]: Adjusts the horizontal position of the collection stage.

4 [Stream Options] tab

Adjusts side stream settings.

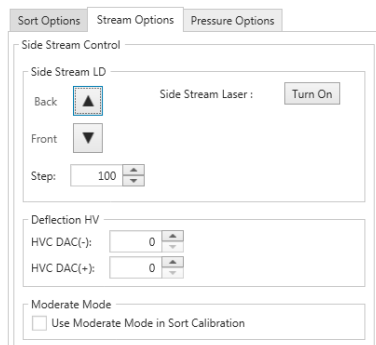


[Side Stream LD]: Adjusts the irradiation position of the side stream laser.

[Deflection HV]: Adjusts the voltage applied to each of the deflection plates.

[Moderate Mode]: Sets whether to use moderate mode for sorting calibration.

5 [Pressure Options] tab

Sets the sample fluid flow controls and the window extension settings.

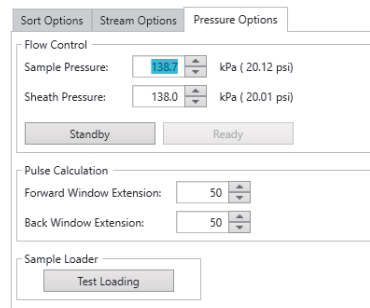

[Flow Control]: Adjusts the sample pressure and sheath pressure.

Clicking [Standby] stops the sheath and sample fluid flow. Clicking [Ready] starts sheath flow again in preparation for a sample.

[Pulse Calculation]: Sets the forward window extension (FWE) and back window extension (BWE) settings used to modify the pulse parameters for data acquisition.

Tip

The setting of the forward window extension and back window extension affect the data acquisition results for Area pulse parameters.

[Test Loading]: Loads the sample loader manually.


After loading, follow the instructions to unload the sample loader.

[Compensation] Tab (ribbon)

The [Compensation] tab of the ribbon has the following buttons.

[Compensation] group

Contains tool buttons related to fluorescence compensation.

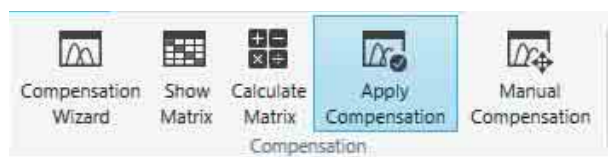

[Compensation Wizard]: Displays the [Compensation Wizard] for recording control tubes in a compensation panel.

The compensation settings should normally be configured using the [Compensation Wizard] to record a negative (unstained) control tube and up to six positive (stained) control tubes in a compensation panel. The [Compensation Wizard] acquires data for compensation samples and then automatically calculates the fluorescence compensation spillover matrix for the sample group.

For details, see “Performing Fluorescence Compensation” (page 67).

[Show Matrix]: Displays the [Compensation Settings] dialog (page 142) for viewing the spillover matrix, and loading/saving the spillover matrix.

[Calculate Matrix]: Displays the [Calculate Compensation Settings] dialog (page 142) for calculating the coefficients of the fluorescence compensation spillover matrix based on the compensation panel.


[Apply compensation]: Selects whether fluorescence compensation is applied to plots on the worksheet. Clicking this button toggles compensation on/off.

[Manual Compensation]: Toggles manual fluorescence compensation mode on/off for the plots on the worksheet.

[Compensation Settings] dialog

The [Compensation Settings] dialog is displayed by clicking [Show Matrix] on the [Compensation] tab of the ribbon.

The [Compensation Settings] dialog is used to view the spillover matrix, and load/save the spillover matrix.

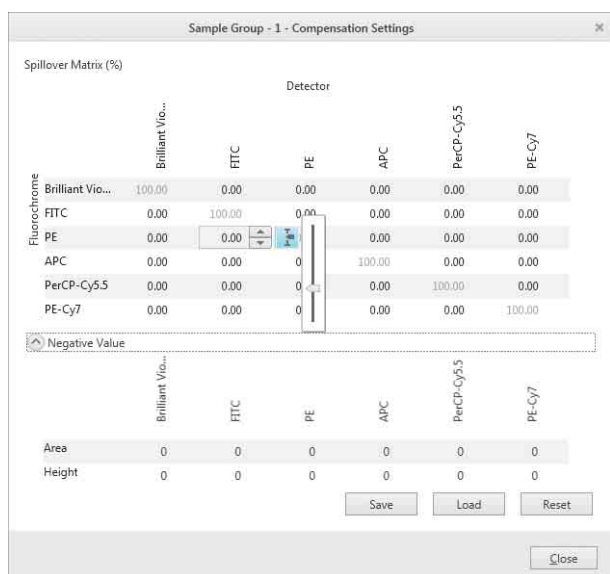

Tip

Fluorochromes cannot be added or deleted using the [Compensation Settings] dialog.

To make changes, see “[Measurement Settings] dialog” (page 80).

[Spillover Matrix (%)]: Displays the fluorescence compensation spillover matrix.

Spillover matrix values display the fluorescence intensity due to each fluorochrome as a percentage of the total intensity in each channel.

Spillover matrix coefficients can be modified manually.

Note

Changing compensation values directly is only recommended for experienced users.

[Negative Value]: Displays the spillover matrix for the negative control tube for both area and height pulse parameters.

[Save]: Saves the compensation settings as an XML file. You can load saved compensation settings into another sample group or another experiment.

[Load]: Loads previously saved compensation settings from an XML file.


[Reset]: Resets all compensation settings.

[Close]: Closes the dialog.

[Calculate Compensation Settings] dialog

The [Calculate Compensation Settings] dialog is displayed by clicking [Calculate Matrix] on the [Compensation] tab of the ribbon.

The [Calculate Compensation Settings] dialog is used to calculate the coefficients of the fluorescence compensation spillover matrix based on the compensation panel.

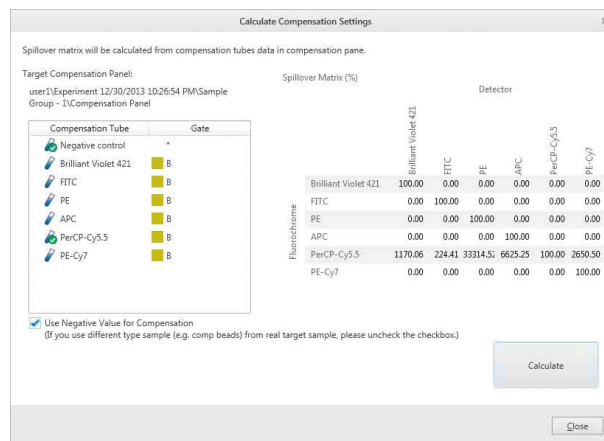

[Target Compensation Panel]: Displays the compensation panel for which to calculate fluorescence compensation.

[Use Negative Value for Compensation]: Enables calculation of the spillover matrix using a negative control. If an appropriate negative control is not available, deselect the checkbox to calculate the spillover matrix without using a negative control. The checkbox default setting is on (enabled).

[Spillover Matrix (%)]: Displays the spillover matrix. Spillover matrix values display the fluorescence intensity due to each fluorochrome as a percentage of the total intensity in each channel.

[Calculate]: Calculates the spillover matrix coefficients.

[Close]: Closes the dialog.

[Worksheet Tools] Tab (ribbon)

The [Worksheet Tools] tab of the ribbon has the following buttons.

[Plot] group

Contains tool buttons for adding, deleting, and duplicating plots.



For details, see “Adding a Plot” (page 91).

[New Density]: Adds a new density plot to the worksheet.

[New Dot Plot]: Adds a new dot plot to the worksheet.


[New Histogram]: Adds a new histogram plot to the worksheet.

[Remove Plot]: Removes the selected plot from the worksheet.

[Duplicates Plot]: Duplicates the selected plot on the worksheet.

[Display Events] group

Contains tool buttons for setting the number of events displayed on plots.

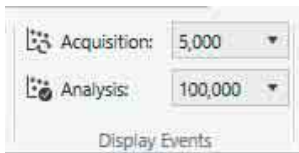

[Acquisition]: Sets the number of events to display on plot during data acquisition.

[Analysis]: Sets the number of events to display on plot when analyzing data.

[Gate] group

Contains tool buttons for editing gates.



For details, see “Editing Gates” (page 99).

[Remove Gate]: Deletes the selected gate.

Note

Deleting a gate will also delete all gates and plots that are derived from the selected gate.

[Edit Gate]: Displays the [Gate Editor] dialog (page 145) for editing gates.

[Statistics] group

Contains tool buttons for working with statistics.



For details, see “Displaying Statistics” (page 104).

[Show Table]: Shows/hides the [Gates and Statistics] table.


[Edit Statistics]: Displays the [Statistics Editor] dialog for configuring the attributes displayed in the [Gates and Statistics] table.

For details about the [Statistics Editor] dialog, see “Displaying Statistics” (page 104).

[Export Statistics to CSV File]: Displays the [Exporting statistics table] dialog (page 145) for exporting the data in the [Gates and Statistics] table as a CSV-format file.

[Alignment] group

Contains tool buttons for displaying a grid on the worksheet and for aligning displayed items with the grid.

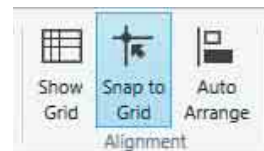


[Show Grid]: Displays the grid on the worksheet.

[Snap to Grid]: Aligns the plots and [Gates and Statistics] table on the worksheet with the grid when they are moved or added.


[Auto Arrange]: Automatically aligns the plots and [Gates and Statistics] table on the worksheet to the grid.

[Worksheet View] group

Contains tool buttons related to the display of items on the worksheet, and for saving the current settings.

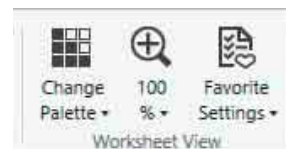

[Change Palette]: Displays a drop-down palette for changing the display style of the selected plot.

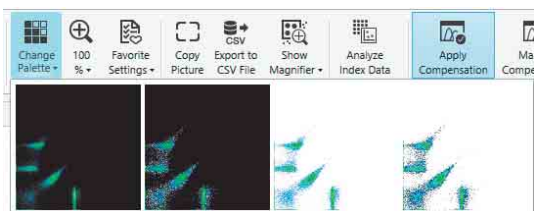

- Black background
- Black background with highlighted outlying events
- White background
- White background with highlighted outlying events


Highlighting outlying events is useful for identifying and analyzing rare events.

Zoom%: Displays a drop-down menu for changing the display zoom factor of the worksheet.

The current zoom factor is displayed on the button icon.

This setting is applied to all plots.

Click this button and select the zoom factor from the menu, or use the slider to set the zoom factor.

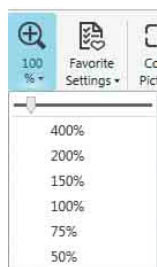

[Favorite Settings]: Allows you to store favorite worksheet settings.

- [Save as Favorite]: Saves the current worksheet state as the default settings.
- [Restore from Favorite]: Applies saved settings to the current worksheet.

[Tools] group

Contains tool buttons for copying, exporting, and displaying data on the worksheet.



[Copy Worksheet Picture]: Copies the selected worksheet to the clipboard for pasting in other software, for example to produce reports.

[Export Data to CSV File]: Exports the recorded data for the sample tube as a CSV file.


Clicking this button displays the [Exporting Data] dialog (*page 145*).

If conducting index sorting, you can export only the

index sorted data by selecting [Well Select Mode] in the [Index Analysis] dialog (*page 115*).

[Show Magnifier]: Shows/hides a worksheet magnifier.

You can also set the zoom factor.

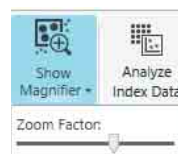

[Index Sorting] group

Contains a command button for analyzing data using the index sorting function.




[Analyze Index Data]: Displays the [Index Analysis] dialog for emphasizing the position on plots of cells sorted in each well in multi-well plates.

For details about the [Index Analysis] dialog, see “Analyzing Index Sort Data” (page 115).

[Compensation] group

Contains tool buttons related to fluorescence compensation.

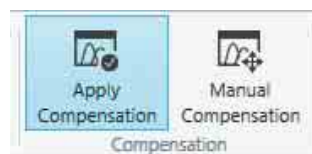


[Apply compensation]: Selects whether fluorescence compensation is applied to plots on the worksheet.

Clicking this button toggles compensation on/off.

[Manual Compensation]: Toggles manual fluorescence compensation mode on/off for the plots on the worksheet.

[Print] group

Contains tool buttons for printing statistics and plots displayed on the worksheet.

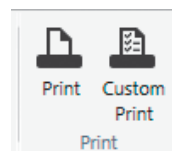

[Print]: Displays the [Print] window (*page 124*) for printing the worksheet.

[Custom Print]: Displays the [Custom Print] dialog (*page 146*) for selecting the items to print.


[Gate Editor] dialog

The [Gate Editor] dialog is displayed by clicking [Edit Gate] on the [Worksheet Tools] tab, [Plot Tools] tab, or [Gate Tools] tab of the ribbon.

The [Gate Editor] dialog is used to edit gates.

- **[Details] tab**

Edits attributes of gates.

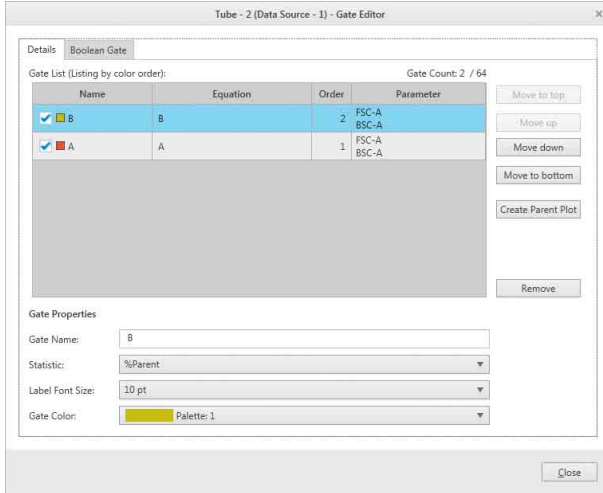

[Gate List]: Displays a list of the configured gates.

You can show or hide a gate on plots using the checkbox beside the gate name.

[Move to top]: Moves the gate with the selected color to the front on dot plots and histogram plots.

[Move up]: Moves the gate with the selected color one step toward the front on dot plots and histogram plots.

[Move down]: Moves the gate with the selected color one step toward the back on dot plots and histogram plots.

[Move to bottom]: Moves the gate with the selected color to the back on dot plots and histogram plots.

[Create Parent Plot]: Adds a parent plot to the worksheet.


[Remove]: Deletes the selected gate.

[Close]: Closes the dialog.

- **[Boolean Gate] tab**

Adds a Boolean gate to the gate hierarchy.

For details about the [Boolean Gate] tab, see “Adding a Boolean gate to the Gate Hierarchy” (page 98).

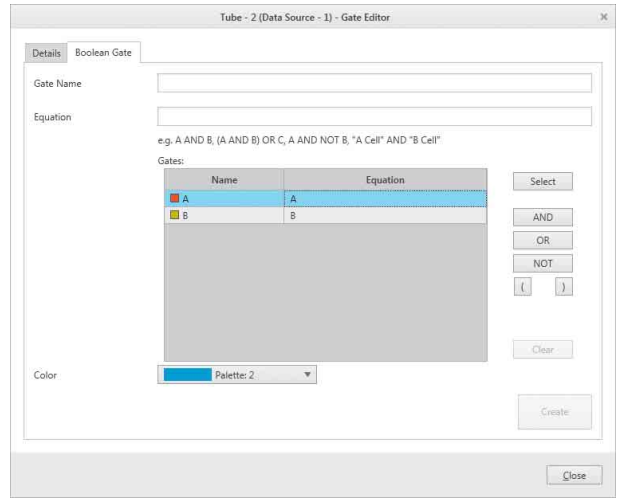

[Exporting statistics table] dialog

The [Exporting statistics table] dialog is displayed by clicking [Export Statistics to CSV File] on the [Worksheet] tab or [Plot Tools] tab of the ribbon.

The [Exporting statistics table] dialog is used to export the contents of the [Gates and Statistics] table as a CSV-format file.



[Output Path]: Specifies the export destination folder.

Click (Browse) to specify for a destination folder.

Progress bar: Displays the export progress.


[Export]: Starts the export.

[Close]: Closes the dialog.

[Exporting Data] dialog

The [Exporting Data] dialog is displayed by clicking [Export Data to CSV File] on the [Worksheet] tab of the ribbon.

The [Exporting Data] dialog is used to export recorded data for sample tubes as a CSV-format file.

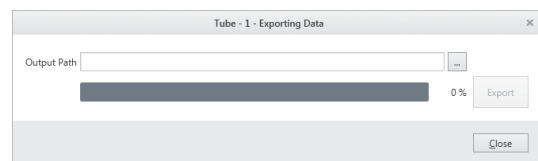

[Output Path]: Specifies the export destination folder.

Click (Browse) to specify for a destination folder.

Progress bar: Displays the export progress.

[Export]: Starts the export.

[Close]: Closes the dialog.

[Custom Print] dialog

The [Custom Print] dialog is displayed by clicking [Custom Print] in the [Print] group on the [Worksheet Tools] tab of the ribbon.

The [Custom Print] dialog is used to select items for printing.



[Custom Settings]: Selects the sample tubes for printing. Select the checkbox for a tube to print the worksheet for the tube.

[Print Experiment Information]: Selects whether to print the experiment information settings.

[Print Tube Information]: Selects whether to print the sample tube information settings.

[Print Sort Settings]: Selects print of the sort settings.

[Print Worksheets]: Selects printing of the worksheet settings.

[Print Statistics and Record Result]: Selects printing of the [Gates and Statistics] table and recorded results.

[Print Sort Result]: Selects print of the sorting results.

[Preview]: Displays a print preview image on the right side of the screen.

For details about other items, see “[Print] Window” (page 124).

[Plot Tools] Tab (ribbon)


The [Plot Tools] tab of the ribbon has the following buttons.

Refer to the [Worksheet Tools] tab for a description of the [Plot] group buttons (page 143), [Statistics] group buttons (page 143), [Tools] group buttons (page 144), and [Compensation] group buttons (page 144).

[Plot Type] group


Contains tool buttons for changing the type of the selected plot.

For histogram plots, select a histogram plot on the worksheet, and click [Histogram] repeatedly to toggle between filled area view and outline view.

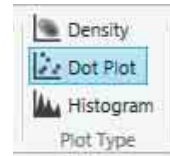

[Scale Type] group

Changes the type of scale displayed on axes on plots.

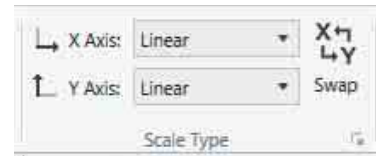

[X Axis] / [Y Axis]: Changes the type of scale displayed on the X axis and Y axis, respectively, of the selected plot.

For details about axis scale types, see “Changing Axis Scale Type” (page 94).

[Swap]: Swaps the position of the X axis and Y axis scales.

: Displays the [Property Window] dialog (page 147) for specifying axis types and scale ranges.

[Axes] group

Contains tool buttons for adjusting axis scales.



[Auto Adjust]: Automatically adjusts plot axes.

- [Auto Adjust (XY)]: Automatically adjusts the ranges of the X axis and Y axis of the selected plot to their minimum and maximum measured values.
- [Auto Adjust (X Axis)]: Automatically adjusts the ranges of the X axis of the selected plot to their minimum and maximum measured values.
- [Auto Adjust (Y Axis)]: Automatically adjusts the ranges of the Y axis of the selected plot to their minimum and maximum measured values.

[Zoom]: Zooms in on the area selected by dragging the mouse on a plot.

[Default Adjust]: Restores the default X axis and Y axis ranges. Click the down arrow to set the number of decades displayed on log scales.

[Copy Axes] / [Paste Axes]: Copies and pastes the plot axes into another plot.

: Displays the [Property Window] dialog (page 147) for specifying axis types and scale ranges.

[Gate] group

Contains tool buttons for drawing gates on plots and editing gate properties.



For details, see “Adding Gates” (page 97) and “Editing Gates” (page 99).

[Rectangle]: Creates a rectangular gate on a dot plot or density plot.

[Ellipse]: Creates an elliptical gate on a dot plot or density plot.

[Polygon]: Creates a polygonal gate on a dot plot or density plot.


[Quadrant]: Creates a quadrant gate on a dot plot or density plot.

[Linear]: Creates a linear gate on a histogram plot.

[Edit Gate]: Displays the [Gate Editor] dialog (page 145) for editing gates.

[Overlay] group

Contains tool buttons for overlay display of plots.

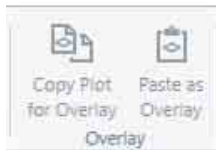

[Copy Plot for Overlay] / [Paste as Overlay]:

Copies and pastes an overlay source plate onto an overlay destination plot.

[Plot View] group

Contains tool buttons for changing the display of plots.




[Change Palette]: Displays a drop-down palette for changing the display style of the selected plot.


For details, see “[Change Palette]” (page 144).

x1.0, x1.5, x2.0: Displays the plot at 100% (default), 150%, and 200% size, respectively.

[Property Window] dialog

The [Property Window] dialog is displayed by clicking the  dialog launcher in the [Scale Type] group or [Axes] group on the [Plot Tools] tab of the ribbon. You can also display the dialog by right-clicking the worksheet, plot, gate, or a [Gates and Statistics] table cell, and selecting [Properties] from the context menu. The [Property Window] dialog is used to change the attributes of plots, and to specify the scale type and range of axes.

Example: Density plot / dot plot properties

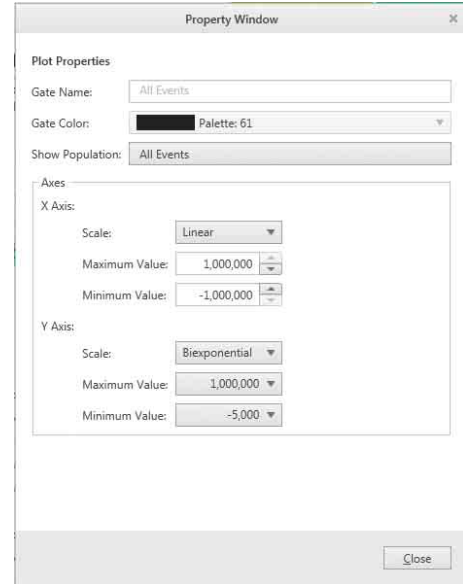

[Gate Name]: Displays the name of the gate (required). You can edit the name, as required. Note that the names of gates must be unique.

[Gate Color]: Selects the color of the gate on plots.


[Show Population]: Selects the population to display on plots.

[Axes]: Specifies the scale type and range of the X and Y axes.

- [Scale]: Selects the scale type.
- [Maximum Value]: Specifies the maximum value of the scale.
- [Minimum Value]: Specifies the minimum value of the scale.

[Close]: Closes the dialog.

Example: Histogram plots

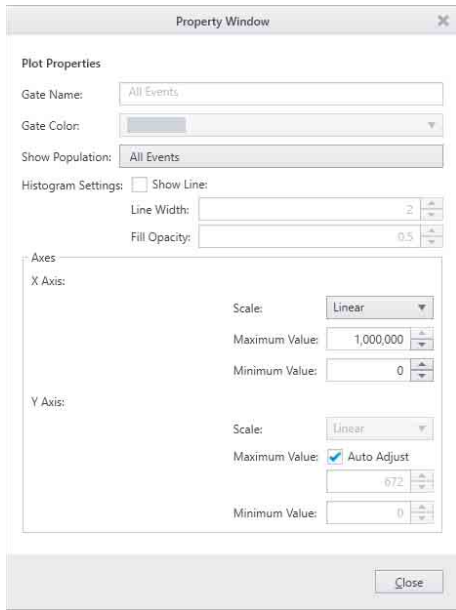

- [Gate Name]:** Displays the name of the gate.
- [Gate Color]:** Displays the color of the gate.
- [Show Population]:** Selects the population to display on plots.
- [Histogram Settings]:** Sets the display view of histogram plots.
 - **[Show Line]:** Toggles between outline view and filled area view.
 - **[Line Width]:** Selects the line thickness.
 - **[Fill Opacity]:** Selects the opacity of the filled area.
- [Axes]:** Specifies the scale type and range of the X and Y axes.
 - **[Scale]:** Selects the scale type.
 - **[Maximum Value]:** Specifies the maximum value of the scale.
 - **[Minimum Value]:** Specifies the minimum value of the scale.
 - Placing a check mark in **[Auto Adjust]** selects automatic adjustment of the Y axis.
- [Close]:** Closes the dialog.


[Gate Tools] Tab (ribbon)

The [Gate Tools] tab of the ribbon has the following buttons.

Refer to the [Worksheet Tools] tab for a description of the [Gate] group buttons (page 143).

[Label] group

Contains tool buttons for specifying gate labels.

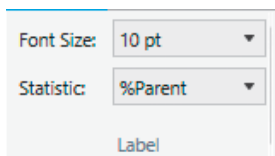


[Font Size]: Selects the font size of text labels.

[Statistic]: Selects the statistic that is displayed beside gates on plots.

[Type Convert] group

Contains tool buttons for converting the selected gate to a different gate type.

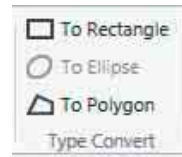


[Style] group

Contains tool buttons for changing the display style of gates.




- [Visible]:** Shows/hides the selected gate.
- Color palette:** Changes the color of the selected gate.
- [Edit Color]:** Allows you to edit the color in the lower color palette (user defined colors) to any color. Click a color in the color palette, click [Edit Color], then select a new color and click [Apply].

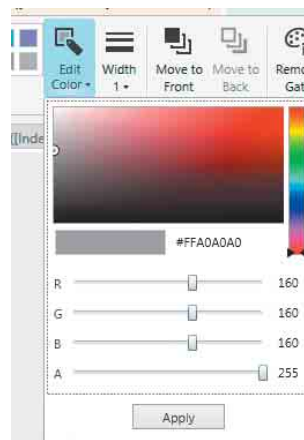

[Width]: Selects the line width of the selected gate.

[Color Order] group


Contains tool buttons for moving the gate with the selected color toward the front or back.

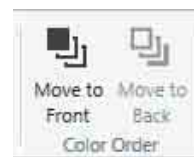

- [Move to Front]:** Moves the gate with the selected color one step toward the front.
- [Move to Back]:** Moves the gate with the selected color one step toward the back.

[Tools] group

Contains a tool button for copying and exporting gates.

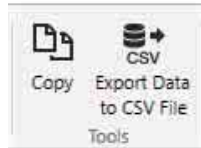

[Copy]: Copies the selected gate.

[Export to CSV File]: Exports the event data in the selected gate as a CSV-format file.

[Gates and Statistics] Table

The [Gates and Statistics] table displays the following information.

1 Name	2 Events	3 %Parent	4 %Total	5 Mean
All Events	0	0.00%	0.00%	
A	0	0.00%	0.00%	
B	0	0.00%	0.00%	
Brilliant V...				0
FTC-A				0
PE-A				0
APC-A				0
PerCP-Cy...				0
PE-Cy7-A				0

1 [Name]

Displays the names of gates and parameters.

2 [Events]

Displays the number of events.

3 [% Parent]

Displays the proportion of events of each gate relative to its parent.

4 [% Total]

Displays the proportion of events of each gate relative to the total of all events.

5 Other statistics values

Displays the values of statistics configured in the [Statistics Editor] dialog.

- [Mean]: Average value of events in population
- [Minimum]: Minimum value in a population
- [Maximum]: Maximum value in a population
- [SD]: Standard deviation of events from the mean value in a population
- [CV]: Coefficient of variation of events in a population
- [Median]: Median value (equal number of events above and below)
- [Mode]: Value with highest event count (corresponds to the highest peak on a histogram plot)
- [HPCV]: Half-peak coefficient of variation (coefficient of variation from width of distribution at half peak height)


For details, see “Displaying Statistics” (page 104).

6 Gate hierarchy

Displays the hierarchical relationship between gates. The color to the left of the gate name shows the color of the gates on plots.

[Acquisition] Tab (control pane)

The [Acquisition] tab controls data acquisition and recording.

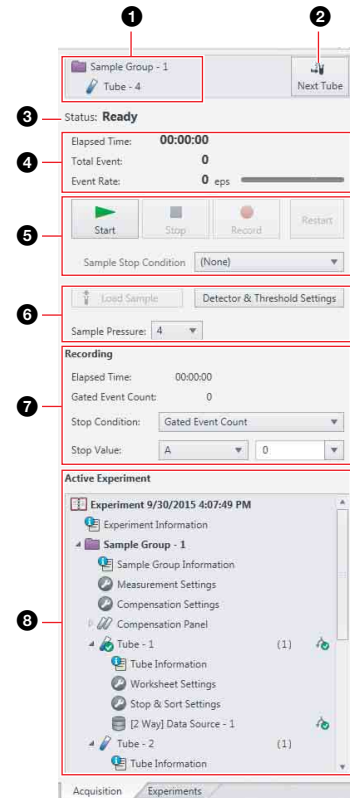

1 Currently assigned sample tube and sample group

2 [Next Tube]

Assigns the next tube.

If there are no further tubes, it duplicates the current tube and assigns the new tube.

3 Instrument status

Displays the status of the instrument.

The following states are displayed.

- Initializing
- Standby
- Ready
- Pause
- Measurement
- Probe Wash
- No Chip
- Chip De-bubble
- Ethanol Cleaning
- Error
- Waiting
- Bleach Cleaning
- DI Rinse
- Shutdown

4 Data acquisition status

Displays the status of data acquisition.

[Elapsed Time]: Displays the cumulative elapsed time since the start of data acquisition.

[Total Events]: Displays the total number of events detected during data acquisition.

[Event Rate]: Displays the rate events are detected during data acquisition in events per second (eps).

Range: 0 eps to 20,000 eps.

The event rate indicator on the right hand side shows the current event rate relative to the guaranteed sorting performance rate of 10,000 eps.



Tip

It is recommended that the event rate be reduced if the indicator exceeds 10k when sorting.

5 Data acquisition controls



(Start): Starts/resumes data acquisition.

It does not start data recording.



(Pause): Pauses data acquisition.

The elapsed time continues counting while acquisition is paused.



(Stop): Stops data acquisition manually.



(Record): Records the acquired data.



(REC stop): Stops data recording manually.



(Restart): Resets the total events and elapsed time counters to 0 and restarts data acquisition.

Note

Avoid keeping paused condition for a prolonged time, it may cause aggregates of sample and clog in the fluidics line.

Tip

If an automatic stop condition has been specified, data acquisition and/or recording stops automatically when the corresponding stop condition has been satisfied.

[Sample Stop Condition]: Selects the stop condition for stopping data acquisition automatically.

- None (data acquisition will stop automatically when the number of detected events exceeds 100,000,000 or the elapsed time exceeds 12:00:00, whichever occurs first)
- Recording (stops data acquisition automatically when the recording stop condition is satisfied)
- Sorting (stops data acquisition automatically when the sorting stop condition is satisfied)
- Recording and Sorting (stops data acquisition automatically when both the recording and sorting stop conditions are satisfied)

Tip

[Recording] cannot be selected for index sorting.

For details about index sorting, see “Index Sorting” (page 114).

6 Instrument controls

[Load Sample] / [Unload Sample]: Loads/unloads the sample tube.

These operations are enabled only when data acquisition is paused.

[Detector & Threshold Settings]: Displays the [Detector & Threshold Settings] dialog.

For details about the [Detector & Threshold Settings] dialog, see “Configuring Detector Settings” (page 86).

[Sample Pressure]: Sets the sample fluid pressure (range of 1 to 10).

7 Recording controls

[Elapsed Time]: Displays the cumulative elapsed time since the start of data recording.

[Event Count]: Displays the number of recorded events.

[Stop Condition]: Selects the stop condition for stopping data recording automatically.

- None
- Elapsed Time
Specifies elapsed time condition to stop recording automatically.
- Event Count
Specifies the event count condition to stop recording automatically.
- Gated Event Count
Specifies the event count condition of the selected gate to stop recording automatically. Specify the target gate of the stop condition, and set the number of events.

Tip

Recording control items cannot be configured for index sorting.

For details about index sorting, see “Index Sorting” (page 114).


8 Experiment Explorer (Active Experiment)

Displays the structure of the active experiment enabled for data acquisition, recording, and cell sorting.

For details about the Experiment Explorer, see “Experiment Explorer” (page 151) on the [Experiments] tab (control pane).

[Experiments] Tab (control pane)

The [Experiments] tab displays all current public (shared) and private experiments.

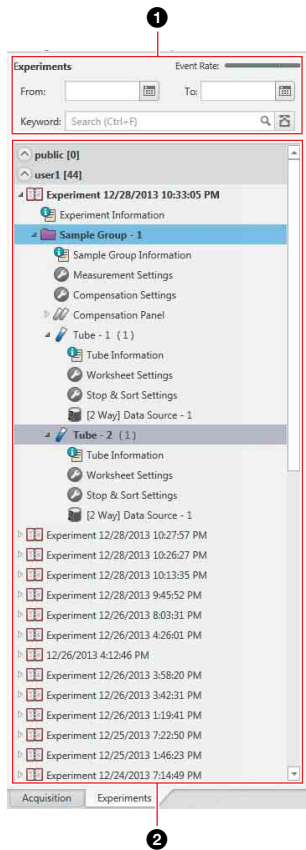

1 Search criteria

You can search experiments by date and keyword. The search results are displayed in the Experiment Explorer.

For details about searching, see “Searching Experiments” (page 84).

2 Experiment Explorer

Displays a list of all current experiments. The internal structure of each experiment is displayed in a hierarchical list view.

[public]: Lists shared experiments. Shared experiments can be viewed by all users.

User_name: Lists private experiments created by the logged-in user. Private experiments cannot be viewed by other users.

An experiment contains the following structural components.

Experiment

[Experiment Information]: Contains information about the experiment. Double-clicking this component displays the [Experiment Information] dialog (page 152).

Sample group: Displays the name of a sample group. A sample group contains the following components.

[Sample Group Information]: Contains information about the sample group. Double-

clicking this component displays the [Sample Group Information] dialog (page 152).

[Measurement Settings]: Contains instrument settings. Double-clicking this component displays the [Measurement Settings] dialog (page 80).

[Compensation Settings]: Contains the fluorescence compensation settings for the fluorochromes used. Double-clicking this component displays the [Compensation Settings] dialog (page 142).

[Compensation Panel]: Contains one negative (unstained) control tube and up to six positive (fluorochrome stained) control tubes. A compensation control tube is required for each fluorochrome used to mark samples in the sample group.

The icon indicates sample tubes for which data has been recorded or which have been used for sorting. These tubes cannot be reassigned as the active tube.

Double-clicking a tube displays the worksheet for the tube.

[Negative control]: Tube for recording data for the negative control sample used to compensate for cellular autofluorescence.

Positive control tubes: Tubes for recording single-stained positive control samples used to calculate the fluorescence compensation spillover matrix, one tube for each fluorochrome used to mark samples.

Sample tube: Tubes containing sample for data acquisition, recording, and cell sorting. Each tube also has worksheet settings and sort & stop settings associated with it.

Double-clicking a tube displays the worksheet for the tube.

To make a tube the active tube for data acquisition, select the tube and click [Assign Tube] in the [Target] group on the [Experiment] tab of the ribbon.

During acquisition the tube icon changes to .

The icon indicates sample tubes for which data has been recorded or which have been used for sorting. These tubes cannot be reassigned as the active tube.

The following icons are displayed for each tube on the right side in the Experiment Explorer.

: Indicates data is being recorded.



: Indicates data has been recorded.

: Indicates tube is being sorted.

: Indicates that the tube has been used for sorting into the collection tubes.

[Tube Information]: Contains information about the tube. Double-clicking this component displays the [Tube Information] dialog (page 152).

[Worksheet Settings]: Contains information about the plots, gates, and statistics on the worksheet. Double-clicking this component displays the [Worksheet Settings] dialog (page 153).

-  **[Stop & Sort Settings]:** Contains the sort settings, and the stop conditions for data acquisition, recording, and sorting for the tube. Double-clicking this component displays the [Stop & Sort Settings] dialog (*page 153*).
-  **Data source:** Indicates the presence of saved data and type of sort results.

Data source naming convention


The sorting method is prefixed to the name of the data source.

Sorting method	Data source name
2 Way Tubes	[2 Way] Data Source - #
8 Well Strips	[8 Well] Data Source - #
Slides	[Slide] Data Source - #
6 Well Plate	[6 Well] Data Source - #
12 Well Plate	[12 Well] Data Source - #
24 Well Plate	[24 Well] Data Source - #
48 Well Plate	[48 Well] Data Source - #
96 Well Plate	[96 Well] Data Source - #
384 Well Plate	[384 Well] Data Source - #
Index sorting Example: 96 Well Plate	[Index 96 Well] Data Source - # * The data source name has an "Index" prefix.
Imported data created using Cell Sorter Software version 1.3.1 or earlier Ex: 8 Well Strips	[Old 8 Well] Data Source - # * The data source name has an "Old" prefix.

[Experiment Information] dialog

The [Experiment Information] dialog is displayed by double-clicking [Experiment Information] for an experiment in the Experiment Explorer.

The [Experiment Information] dialog is used to view and edit information about the experiment.

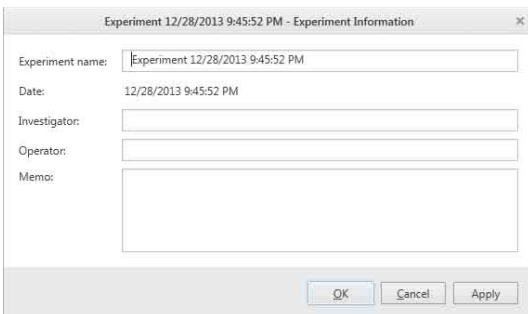

[Experiment Name]: Name of the experiment (required item).

[Date]: Displays the date the experiment was created. This parameter cannot be modified.

[Investigator]: Name of the lead researcher.

[Operator]: Name of the operator.

[Memo]: Descriptive comments.

[OK]: Applies the settings and closes the dialog.


[Cancel]: Cancels settings and closes the dialog.

[Apply]: Applies the settings without closing the dialog.

[Sample Group Information] dialog

The [Sample Group Information] dialog is displayed by double-clicking [Sample Group Information] in an experiment in the Experiment Explorer.

The [Sample Group Information] dialog is used to view and edit information about the sample group.

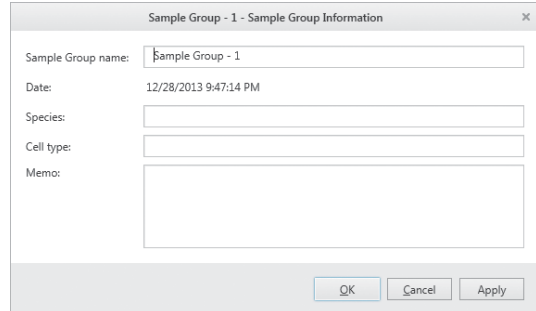

[Sample Group Name]: Unique name of the sample group (required item).

[Date]: Displays the date the sample group was created. This parameter cannot be modified.

[Species]: Name of the target biological species.

[Cell type]: Name of the cell type within the biological species.

[Memo]: Descriptive comments.

[OK]: Applies the settings and closes the dialog.


[Cancel]: Cancels settings and closes the dialog.

[Apply]: Applies the settings without closing the dialog.

[Tube Information] dialog

The [Tube Information] dialog is displayed by double-clicking [Tube Information] for a sample tube in the Experiment Explorer.

The [Tube Information] dialog is used to view and edit information about the sample tube.

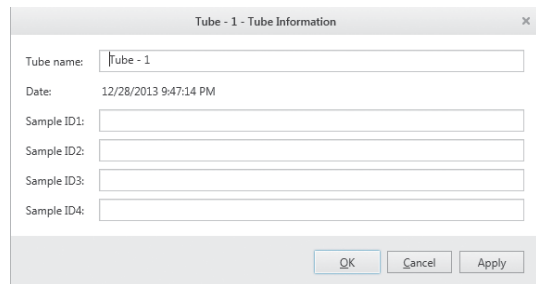

[Tube Name]: Unique name of the sample tube (required item).

[Date]: Displays the date the sample tube was created. This parameter cannot be modified.

[Sample ID1] to [Sample ID4]: ID labels for identifying the sample tube.

[OK]: Applies the settings and closes the dialog.

[Cancel]: Cancels settings and closes the dialog.

[Apply]: Applies the settings without closing the dialog.


[Worksheet Settings] dialog

The [Worksheet Settings] dialog is displayed by double-clicking [Worksheet Settings] for a sample tube in the Experiment Explorer.

The [Worksheet Settings] dialog is used to save the current worksheet settings (number of plots, type of plots, statistics settings, etc.) as the favorite settings. Saving current worksheet settings as the favorite settings allows you to restore the settings at a later time or to apply the settings to a different sample tube.

Tip

Data that is acquired and recorded for a tube is not included in the worksheet settings.

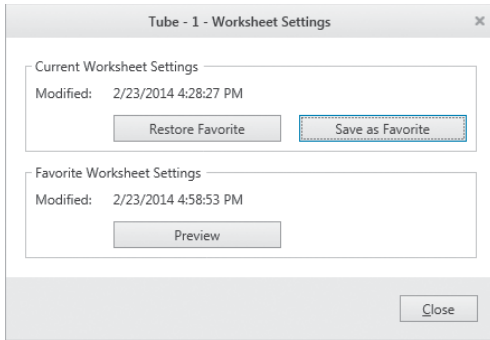

[Current Worksheet Settings]

[Modified]: Displays the date and time that the current worksheet settings were modified.

[Restore Favorite]: Restores the saved favorite worksheet settings as the current worksheet settings.

[Save as Favorite]: Saves the current worksheet settings as the favorite (default) worksheet settings.

[Favorite Worksheet Settings]

Modified: Displays the date and time that the favorite worksheet settings were saved.

[Preview]: Displays a preview of the worksheet with the saved favorite worksheet settings.


[Close]: Closes the dialog.

[Stop & Sort Settings] dialog

The [Stop & Sort Settings] dialog is displayed by double-clicking [Stop & Sort Settings] for a sample tube in the Experiment Explorer.

The stop and sort settings can be viewed in the [Stop & Sort Settings] dialog.

Example: 2-way tube sorting

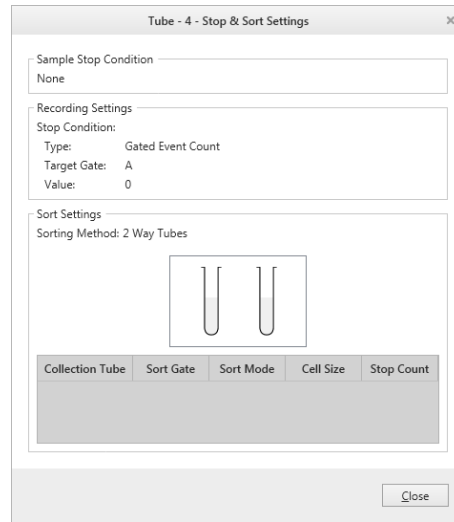

[Sample Stop Condition]: Displays the data acquisition stop condition.

[Recording Settings]

[Stop Condition]: Displays the recording stop condition.

[Sort Settings]

[Sorting Method]: Displays the sorting method for the sample tube.

You can change the order items are displayed by dragging and dropping the title cells.

[Close]: Closes the dialog.


[Sort Control] Pane

The [Sort Control] pane is used to configure and control cell sorting.

For details, see “Sorting” (page 58).

The items displayed in the [Sort Control] pane vary depending on the selected sorting method. This section describes the displayed items using two-way tube sorting and 96-well plate sorting as examples.

Tip

You can show/hide the [Sort Control] pane by clicking  in the top left corner of the pane. In analyzer mode, the [Sort Control] pane is automatically hidden.

For 2-way tube sorting

The following items are displayed when [2 Way Tubes] is selected in [Sorting Method].



1 [Sort Control] pane

Configures and controls droplet sorting.

[Sort Start] / [Sort Stop]: Starts/stops sorting.

[Load Collection] / [Unload Collection]: Loads/unloads the collection stage.

[Method]: Selects the sorting method. Set the sorting method to match the collection devices loaded on the collection stage.

The sorting parameters vary depending on the selected sorting method.

[Mode]: Selects the sorting mode.

Sorting mode

- Ultra Purity
- Purity
- Semi Purity
- Normal
- Semi Yield
- Yield
- Ultra Yield
- Single Cell
- Single Cell (3 drops) (100 μm /70 μm sorting chip only)
- Custom: Up to five user-defined modes

Tip

To adjust the position for sorting of target cells within a droplet, you can adjust the settings using [Edit setting].

For details, see “[Sort Mode Setting] dialog” (page 156).

Cell size


- Regular Cell
- Large Cell

For details about each sorting mode, see “Sorting Mode” (page 106).

Tips

- When sorting larger cells using a 100 μm /70 μm sorting chip, selecting [Single Cell (3 drops)] may improve the sorting performance. The cell size is fixed to [Large Cell] when [Single Cell (3 drops)] is selected.
- When a “Large cell” option is selected, the side stream deflection angle is adjusted for sorting of cells of larger size.
- The “Large cell” options are not displayed if sorting calibration is not performed. Select a “Large cell” option if droplet formation is irregular, or the droplet stream is hitting the side wall of the collection tubes or is missing the tubes when sorting cells larger than 15 μm in Normal mode.
- The side streams may become unstable when sorting cells of size 20 μm or larger.
- If droplet marks are visible on the collection tube holder after sorting, wipe clean with a lint-free, soft, clean cloth sprayed with ethanol and then another moistened with water.

Side stream monitor: Displays the image from the side stream monitor camera.

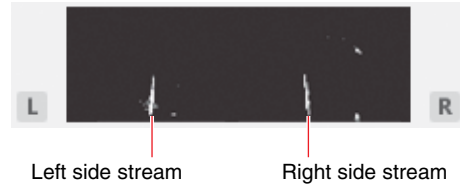

For details about adjusting the deflection angle of side streams, see “[Advanced Settings] dialog” (page 139).

[To Sort]: Selects the gates for the target events to sort into the left and right collection tubes.

[Stop Value]: Specifies the event count conditions to stop recording automatically for each of the left and right collection tubes.

You can select a value or enter a value from the keyboard.

Entering a value of “0” disables the stop condition.

2 [Sort Statistics] pane

Displays the statistics about events during sorting.

[Elapsed Time]: Displays the cumulative elapsed time since the start of sorting.

[Remaining Time]: Displays the estimated remaining sorting time.

[Sort Count]: Displays the number of events sorted.

[Sort Rate]: Displays the rate at which events are sorted in units of events per second (eps).

[Sort Efficiency]: Displays the number of sorts attempted as a percentage of targeted cells in the sample.

[Abort Count]: Displays the number of target events in the sort gate that do not match the sort mode criteria and which are aborted during sorting.

[Abort Rate]: Displays the rate at which events are aborted during sorting in units of events per second (eps).

3 [Droplet] pane

Displays the droplet stream immediately beneath the sorting chip nozzle.

The droplet stream is used to monitor droplet formation and to adjust the breakoff point.

Double-clicking the image displays the side streams in the [Droplet Viewer] window (page 156) for monitoring droplets.

4 [Hardware Status] pane

Displays the status of the interior lights in the sample loader and collection stage area, and the level in each of the tanks in the fluidics cart. It also displays the lasers enabled for the experiment.

[Light]

[Collection]: Turns the collection area light on/off.

[Sample]: Turns the sample injection chamber light on/off.


Tank: Indicates the fluid level in the sheath tank, ethanol tank, waste tank, and DI water tank.

A warning icon appears when the waste tank is close to full, or when the sheath, ethanol, or DI water tank is close to empty.

[Laser]: Indicates the lasers used for data acquisition.

For 8-well strips or slide glass sorting

The following items are displayed when [8 Well Strips] or [Slides] is selected in [Method].

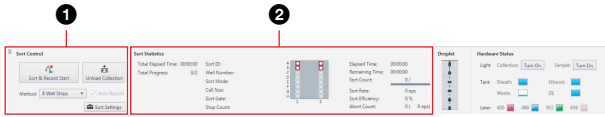

The [Droplet] and [Hardware Status] panes are the same as for two-way tube sorting.

1 [Sort Control] pane

Configures and controls droplet sorting.

[Sort Start] / [Sort Stop]: Starts/stops sorting.

[Load Collection] / [Unload Collection]: Loads/unloads the collection stage.

[Method]: Selects the sorting method. Set the sorting method to match the collection devices loaded on the collection stage.

The sorting parameters vary depending on the selected sorting method.

[Sort Settings]: Specifies the default sorting configuration.

Clicking this button displays the [Sort Settings] dialog.

For details about the [Sort Settings] dialog, see “Sorting” (page 58) and “Adjusting the Sort Position (SH800SP and SH800ZP)” (page 113).

2 [Sort Statistics] pane

Displays the statistics about events during sorting.

[Total Elapsed Time]: Displays the cumulative elapsed time since the start of sorting of the whole plate.

[Total Progress]: Displays the ratio of sorted wells to the total number of wells.

[Sort ID]: Displays the ID of the well currently being sorted.

[Well Number]: Displays the number of the well currently being sorted.

[Sort Mode]: Displays the sorting mode of the well currently being sorted.

[Cell Size]: Displays the current cell size setting.

[Sort Gate]: Displays the sort gate of the well currently being sorted.

[Stop Count]: Displays the number of events for the automatic stop condition of the well currently being sorted.

Plate (image): Displays the layout of the wells.

Double-clicking the plate image or right-clicking and selecting [Show Sorting Report] from the context menu displays the plate image in a separate window for monitoring progress.

[Elapsed Time]: Displays the cumulative elapsed time since the start of sorting.

[Remaining Time]: Displays the estimated remaining sorting time.

[Sort Count]: Displays the number of events sorted.

[Sort Rate]: Displays the rate at which events are sorted in units of events per second (eps).

[Sort Efficiency]: Displays the number of sorts attempted as a percentage of targeted cells in the sample.


[Abort Count]: Displays the number of target events in the sort gate that do not match the sort mode criteria and which are aborted during sorting.

[Abort Rate]: Displays the rate at which events are aborted during sorting in units of events per second (eps).

For multi-well plate sorting

The following items are displayed when a multi-well plate is selected in [Method].

Example: When [96 Well Plate] is selected

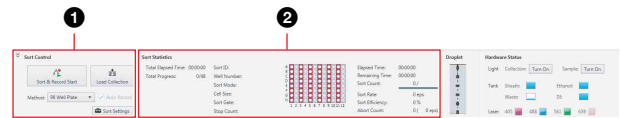

The [Droplet] and [Hardware Status] panes are the same as for two-way tube sorting.

1 [Sort Control] pane

Configures and controls droplet sorting.

[Sort Start] / [Sort Stop]: Starts/stops sorting.

[Load Collection] / [Unload Collection]: Loads/unloads the collection stage.

[Method]: Selects the sorting method. Set the sorting method to match the collection devices loaded on the collection stage.

The sorting parameters vary depending on the selected sorting method.

[Sort Settings]: Specifies the default sorting configuration.

Clicking this button displays the [Sort Settings] dialog.

For details about the [Sort Settings] dialog, see “Sorting” (page 58) and “Adjusting the Sort Position (SH800SP and SH800ZP)” (page 113).

2 [Sort Statistics] pane

Displays the statistics about events during sorting.

[Total Elapsed Time]: Displays the cumulative elapsed time since the start of sorting of the whole plate.

[Total Progress]: Displays the ratio of sorted wells to the total number of wells.

[Sort ID]: Displays the ID of the well currently being sorted.


[Well Number]: Displays the number of the well currently being sorted.

- [Sort Mode]:** Displays the sorting mode of the well currently being sorted.
- [Cell Size]:** Displays the current cell size setting.
- [Sort Gate]:** Displays the sort gate of the well currently being sorted.
- [Stop Count]:** Displays the number of events for the automatic stop condition of the well currently being sorted.
- Plate (image):** Displays the layout of the wells.
Double-clicking the plate image or right-clicking and selecting [Show Sorting Report] from the context menu displays the plate image in a separate window for monitoring progress.
- [Elapsed Time]:** Displays the cumulative elapsed time since the start of sorting.
- [Remaining Time]:** Displays the estimated remaining sorting time.
- [Sort Count]:** Displays the number of events sorted.
- [Sort Rate]:** Displays the rate at which events are sorted in units of events per second (eps).
- [Sort Efficiency]:** Displays the number of sorts attempted as a percentage of targeted cells in the sample.
- [Abort Count]:** Displays the number of target events in the sort gate that do not match the sort mode criteria and which are aborted during sorting.
- [Abort Rate]:** Displays the rate at which events are aborted during sorting in units of events per second (eps).

[Sort Mode Setting] dialog

The [Sort Mode Setting] dialog is displayed by selecting [Edit setting] from [Custom] under [Sort Mode]. The [Sort Mode Setting] dialog is used to adjust the position of target cells within a droplet for sorting. The adjustment values are valid for the currently selected tube only. To apply the value to other tubes, register the value in [User Preference].

To display [User Preference], see the tip (page 114) in step 3 in “Adjusting the Sort Position (SH800SP and SH800ZP)” (page 113).

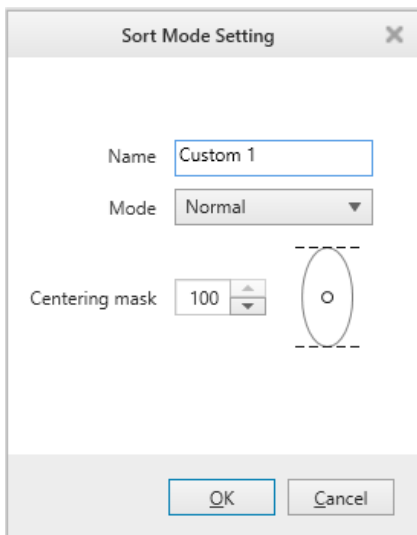

- [Name]:** Enter a name for the custom mode (required item).
- [Mode]:** Select the base sorting mode.
When a sorting mode is selected, the default value for that mode is entered in [Centering mask].
- [Centering mask]:** Specify the position of target cells within a droplet for sorting.


Tips

- The [Name] field cannot be left blank. Always specify a name.
- Specifying the same name as a preset sorting mode is not possible.

[Droplet Viewer] window

The [Droplet Viewer] window is displayed by double-clicking the image in the [Droplet] pane at the bottom of the main window.

The [Droplet Viewer] window is used to monitor droplet formation and as a guide when adjusting the breakoff point.

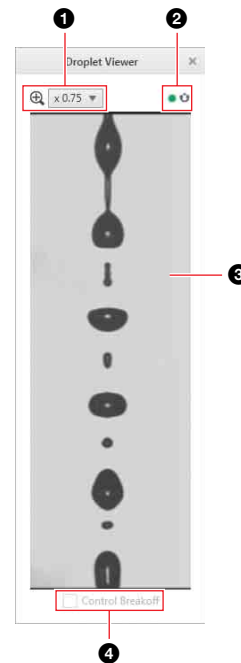

1 Zoom factor

Selects the zoom factor (×1.00, ×0.75, or ×0.50)

2 Status icon

●: Indicates the stability of droplet formation.

Indication	Status
Flashing slowly	Waiting for stream to stabilize
Flashing rapidly	Adjustment in progress
On	Stable

🔘: Indicates that the droplet control breakoff function is active.

No icon: Indicates unstable or absence of droplet formation.

③ Droplet monitor

Displays the droplet stream.

④ [Control Breakoff]

The control breakoff setting uses a feedback system to control the breakoff position automatically.

When enabled, the amplitude and frequency settings are adjusted automatically to maintain a stable breakoff point position with coherent side streams. When adjusting the breakoff point, first set the amplitude and frequency settings to obtain a stable breakoff point. After establishing a stable breakoff point, click the feedback control checkbox to enable the breakoff feedback control.

Tips

- The manual amplitude and frequency controls are disabled in this mode.
- When breakoff feedback control is enabled, if the breakoff point moves and cannot be recovered within a predetermined time, sorting stops and an error message is displayed on the screen.

Context Menu

The context menu displays common operations that can be accessed by right-clicking components in the Experiment Explorer or items on the worksheet.

Experiment Explorer Context Menu

This section describes the context menu items displayed for components in the Experiment Explorer.

Experiments

The following menu commands are displayed when right-clicking an experiment icon in the Experiment Explorer.

[New Sample Group]: Creates a blank sample group.

[Copy] / [Paste]: Copies the experiment and pastes a new experiment into the Experiment Explorer.

[Duplicate]: Duplicates the selected experiment.

[Save as Experiment Template]: Saves the selected experiment as a private template.

[Export FCS File]: Displays the [FCS File Export] dialog (*page 134*) for exporting the tubes in the selected experiment in FCS 3.0 or FCS 3.1 format.

[Send to Public]: Creates a duplicate of the selected experiment and places the duplicate under the [Public] node for sharing with other users.

[Delete]: Deletes the selected experiment.

[Rename]: Renames the selected experiment.

Compensation Panel

The following menu commands are displayed when right-clicking a compensation panel icon in the Experiment Explorer.

[Compensation Wizard]: Displays the [Compensation Wizard] (*page 67*) for recording control tubes in a compensation panel.

[Show Matrix]: Displays the [Compensation Settings] dialog (*page 142*) for viewing the spillover matrix, and loading/saving the spillover matrix.

[Calculate Matrix] Displays the [Calculate Compensation Settings] dialog (*page 142*) for calculating the coefficients of the fluorescence compensation spillover matrix based on the compensation panel.

[Export FCS File]: Displays the [FCS File Export] dialog (*page 134*) for exporting the tubes in the selected compensation panel in FCS 3.0 or FCS 3.1 format.

Control tubes in a compensation panel

The following menu commands are displayed when right-clicking a control tube in a compensation panel icon in the Experiment Explorer.

[Assign]: Assigns the selected control tube as the active tube for data acquisition and recording.

[Open Worksheet]: Opens the worksheet tab for the selected control tube.

[Export FCS File]: Displays the [FCS File Export] dialog (*page 134*) for exporting the selected control tube in FCS 3.0 or FCS 3.1 format.

[Reset Data]: Deletes the data recorded for the control tube.

Sample groups

The following menu commands are displayed when right-clicking a sample group icon in the Experiment Explorer.

[New Tube]: Creates a new sample tube.

Selecting this item creates a new sample tube in the sample group.

[Copy] / [Paste]: Allows you to copy and paste all settings (including sample tubes) in the selected sample group. However, the recorded data of the sample tubes will not be pasted.

[Paste to All Tubes in Sample Group]: Pastes copied worksheet settings to all sample tubes in the selected sample group.

This item is available when you copy the worksheet settings of a sample tube, and then right-click another sample group.

[Duplicate]: Duplicates the selected sample group.

[Save as Sample Group Template]: Displays the [Save as Sample Group Template] dialog (*page 131*) for saving the selected sample group as a private template.

[Apply Template]: Displays the [Apply Template] dialog (*page 132*) for selecting a template to apply to the selected sample group.

[Export FCS file]: Displays the [FCS File Export] (*page 134*) dialog for exporting the recorded data for the sample tubes included in the sample group as an FCS 3.0 or FCS 3.1 format file.

[Delete]: Deletes the selected sample group.

[Rename]: Renames the selected sample group.

Measurement settings

The following menu commands are displayed when right-clicking a [Measurement Settings] icon for a sample group in the Experiment Explorer.

[Copy] / [Paste]: Copies the measurement settings for the selected sample group and pastes them into the measurement settings for another sample group.

[Show Settings]: Displays the [Measurement Settings] dialog (*page 80*) for viewing measurement settings.

Compensation settings

The following menu commands are displayed when right-clicking a [Compensation Settings] icon for a sample group in the Experiment Explorer.

[Copy] / [Paste]: Copies the compensation settings for the selected sample group and pastes them into the compensation settings for another sample group.

[Show Settings]: Displays the [Compensation Settings] dialog (*page 142*) for viewing compensation settings.

Sample tubes

The following menu commands are displayed when right-clicking a sample tube icon in the Experiment Explorer.

[Assign]: Assigns the selected sample tube as the active tube for data acquisition, recording, and cell sorting.

[Open Worksheet]: Opens the worksheet tab for the selected sample tube.

[Apply Compensation]: Applies fluorescence compensation to the plots on the worksheet for the selected sample tube.

[Copy] / [Paste]: Copies the sample tube and pastes it into another sample tube.

[Duplicate]: Duplicates the selected sample tube.

[Save as Tube Template]: Saves the selected sample tube as a private template.

[Apply Template]: Displays the [Apply Template] dialog (*page 132*) for selecting a template to apply to the selected sample tube.

[Export FCS File]: Displays the [FCS File Export] dialog (*page 134*) for exporting the selected sample tube in FCS 3.0 or FCS 3.1 format.

[Delete]: Deletes the selected sample tube.

[Rename]: Renames the selected sample tube.

[Show Plate Sorting Monitor]: Displays the [Plate Sorting Monitor] dialog (*page 66*) for monitoring the status when sorting onto multi-well plates.

[Show Results]: Displays the [Tube Results] dialog (*page 132*) for displaying the recording results and sorting results for the selected sample tube.

Tube worksheet settings

The following menu commands are displayed when right-clicking a [Worksheet Settings] icon for a sample tube in the Experiment Explorer.

[Copy] / [Paste]: Copies the worksheet sheet settings for the selected sample tube and pastes them into the worksheet settings for another sample tube.

[Save as Favorite]: Saves the current worksheet settings as the favorite (default) worksheet settings.

[Restore Favorite]: Restores the saved favorite worksheet settings as the current worksheet settings.

[Apply to All Tubes In Same Sample Group]: Applies the worksheet settings for the selected sample tube to all sample tubes within the sample group.

[Show Settings]: Displays the [Worksheet Settings] dialog (*page 153*) for displaying, saving, and loading worksheet settings.

Data sources

The following menu commands are displayed when right-clicking a data source icon in the Experiment Explorer.

[Open Data Source]: Displays the data source on the worksheet.

[Delete]: Deletes the selected data source.

[Rename]: Renames the selected sample tube.

[Export FCS File]: Displays the [FCS File Export] dialog (*page 134*) for exporting the selected data source in FCS 3.0 or FCS 3.1 format.

[Show Results]: Displays the [Data Source Results] dialog (*page 133*) for displaying the recording results and sorting results for the selected data source.

Worksheet Context Menu

This section describes the context menu items displayed for items displayed on the worksheet and the worksheet itself.

Worksheet

The following menu commands are displayed when right-clicking on the worksheet.

[Refresh]: Updates the information displayed on the worksheet.

[New Density]: Adds a new density plot to the worksheet.

[New Dot Plot]: Adds a new dot plot to the worksheet.

[New Histogram]: Adds a new histogram plot to the worksheet.

[Show Table]: Shows/hides the [Gates and Statistics] table.

[Show Grid]: Displays the grid on the worksheet.

[Snap to Grid]: Aligns the plots and [Gates and Statistics] table on the worksheet with the grid when they are moved or added.

[Auto Arrange]: Automatically aligns the plots and [Gates and Statistics] table on the worksheet to the grid.

[Fit to Screen]: Expands the worksheet to fill the entire screen.

[Edit Gate]: Displays the [Gate Editor] dialog (*page 145*) to edit gates.

[Edit Statistics]: Displays the [Statistics Editor] dialog (*page 104*) to configure the statistics displayed in the [Gates and Statistics] table.

[Overlay Gate List]: Displays the [Overlay Gate List] dialog (*page 104*) for removing gates used for overlay display.

[Favorite Settings] >

[Save as Favorite]: Saves the current worksheet settings as the favorite (default) worksheet settings.

[Restore Favorite]: Restores the saved favorite worksheet settings as the current worksheet settings.

[Copy Worksheet Picture]: Copies the worksheet to the clipboard for pasting in other software, for example to produce reports.

[Print]: Displays the [Print] window (*page 124*) for printing the analysis results displayed on the worksheet.

[Custom Print]: Displays the [Custom Print] dialog (*page 146*) for selecting the items to print.

[Properties]: Displays the [Property Window] dialog (*page 147*).

Plots

The following menu commands are displayed when right-clicking a plot.

[Refresh]: Updates the information displayed on the worksheet.

[Auto Adjust]: Automatically adjusts the ranges of the X axis and Y axis of all plots to their minimum and maximum measured values.

[Copy Plot for Overlay] / [Paste as Overlay]: Copies and pastes an overlay source plate onto an overlay destination plot.

[Clear overlay plots]: Clears the overlay display of plots.

[Remove]: Removes the selected plot.

[Duplicate]: Duplicates the selected plot.

[Copy All Regions] / [Paste Region]: Copies all gates on a plot and pastes them onto another plot.

[Create Gate] >

[Rectangle]: Creates a rectangular gate on a dot plot or density plot.

[Ellipse]: Creates an elliptical gate on a dot plot or density plot.

[Polygon]: Creates a polygonal gate on a dot plot or density plot.

[Quadrant]: Creates a quadrant gate on a dot plot or density plot.

[Linear]: Creates a linear gate on a histogram plot.

[Plot Type] >

[Density]: Changes the type of the selected plot to a density plot.

[Dot Plot]: Changes the type of the selected plot to a dot plot.

[Histogram]: Changes the type of the selected plot to a histogram plot.

Click [Histogram] repeatedly to toggle between filled area view and outline view.

[Zoom]: Zooms in on the area selected by dragging the mouse on a plot.

[Open Gates Editor]: Displays the [Gate Editor] dialog (*page 145*) to edit gates.

[Open Statistics Editor]: Displays the [Statistics Editor] dialog (*page 104*) to configure the statistics displayed in the [Gates and Statistics] table.

[Copy Picture]: Copies the selected plot to the clipboard for pasting in other software, for example to produce reports.

[Properties]: Displays the [Property Window] dialog (*page 147*).

Gates

The following menu commands are displayed when right-clicking a gate.

[Create Density Plot]: Adds a density plot for the events within the selected gate.

[Create Dot Plot]: Adds a dot plot for the events within the selected gate.

[Create Histogram Plot]: Adds a histogram plot for the events within the selected gate.

[Move Color Order to Front]: Moves the gate with the selected color one step toward the front.

[Move Color Order to Back]: Moves the gate with the selected color one step toward the back.

[Convert to] >

[Rectangle]: Converts the selected gate to a rectangle gate.

[Ellipse]: Converts the selected gate to an ellipse gate.

[Polygon]: Converts the selected gate to a polygon gate.

[Copy Region] / [Paste Region]: Copies the selected gate and pastes it into another plot.

[Visible Gate]: Shows/hides the selected gate.

[Remove]: Deletes the selected plot.

[Send to Back]: Moves the selected gate to the back.

[Properties]: Displays the [Property Window] dialog (*page 147*).

[Gates and Statistics] table

The following menu commands are displayed when right-clicking within the [Gates and Statistics] table.

[Show Table]: Shows/hides the [Gates and Statistics] table.

[Save as CSV File]: Displays the [Exporting statistics table] dialog (*page 145*) for exporting the data in the [Gates and Statistics] table as a CSV-format file.

[Copy Picture]: Copies the [Gates and Statistics] table to the clipboard for pasting in other software, for example to produce reports.

[Open Gates Editor]: Displays the [Gate Editor] dialog (*page 145*) to edit gates.

[Open Statistics Editor]: Displays the [Statistics Editor] dialog (*page 104*) to configure the statistics displayed in the [Gates and Statistics] table.

[Gates and Statistics] table cells

The following menu commands are displayed when right-clicking within a [Gates and Statistics] table cell.

[Create Density Plot]: Adds a density plot for the events within the selected gate.

[Create Dot Plot]: Adds a dot plot for the events within the selected gate.

[Create Histogram Plot]: Adds a histogram plot for the events within the selected gate.

[Create Parent Plot]: Adds a parent plot to the worksheet.

[Copy Selected Gate]: Copies the selected gate.

[Cut Selected Gate]: Cuts the selected gate.

[Paste Selected Gate]: Pastes a previously copied or cut gate.

[Visible Gate]: Shows/hides the selected gate.

[Remove Selected Gate]: Deletes the selected gate.

[Show Table]: Shows/hides the [Gates and Statistics] table.

[Save as CSV File]: Displays the [Exporting statistics table] dialog (*page 145*) for exporting the data in the [Gates and Statistics] table as a CSV-format file.

[Open Gate Editor]: Displays the [Gate Editor] dialog (*page 145*) to edit gates.

[Open Statistics Editor]: Displays the [Statistics Editor] dialog (*page 104*) to configure the statistics displayed in the [Gates and Statistics] table.

[Refresh List]: Updates the content displayed in the [Gates and Statistics] table.

[Properties]: Displays the [Property Window] dialog (*page 147*).

This chapter describes the routine maintenance procedures required to maintain performance and longevity of the SH800.

DANGER

Fluids may contain biological, chemical, or other agents. Always wear gloves and other protective clothing, mask, and goggles, as required, when performing maintenance.

Maintenance Schedule

The maintenance procedures outlined in this manual should be performed according to the schedule below.

Preparations for the Day Before Measurement

Maintenance item	Schedule	See
Refilling the sheath tank	<ul style="list-style-type: none"> It is recommended to fill the tank the day before operation to allow air bubbles to escape. 	<i>"Refilling the Sheath Tank"</i> (page 164)

Maintenance for Each Startup Operation

Maintenance item	Schedule	See
Disposing of waste fluid Replacing the waste tank	<ul style="list-style-type: none"> Check that the tank is not full when starting up and shutting down Replace when nearly full or when prompted in Cell Sorter Software 	<i>"Emptying/ Changing the Waste Tank"</i> (page 166)
Checking that there is sufficient volume in the sheath tank	<ul style="list-style-type: none"> When starting up 	<i>"Refilling the Sheath Tank"</i> (page 164)
Refilling the sheath tank	<ul style="list-style-type: none"> When prompted in Cell Sorter Software. 	
Refilling the DI water tank	<ul style="list-style-type: none"> When starting up When prompted in Cell Sorter Software As required 	<i>"Refilling the DI Water Tank"</i> (page 167)

Maintenance for Each Shutdown Operation

Sample fluidics system cleaning

Maintenance item	Schedule	See
Sample fluidics system cleaning	<ul style="list-style-type: none"> When shutting down at day's end As required 	<i>"Cleaning the Sample Fluidics System using Bleach"</i> (page 170)

Parts cleaning

Maintenance item	Schedule	See
Cleaning the sample loader	• When shutting down	<i>"General Cleaning"</i> (page 172)
Cleaning the collection area and transparent protective covers	• When shutting down	
Cleaning the droplet camera windows	• When shutting down	
Cleaning the laser windows	• When shutting down	
Cleaning the collection tube holders, multi-well plate holder	• When shutting down • As required	
Cleaning the splash guard	• When shutting down • As required	

Monthly Maintenance

Maintenance item	Schedule	See
Changing the waste tank air filter	• Monthly	<i>"Changing Tank Air Filters"</i> (page 186)
Cleaning the DI water tank and DI water line	• Monthly • As required	<i>"Cleaning the DI Water Tank and DI Water Line"</i> (page 187)
Running ethanol cleaning	• Monthly • As required	<i>"Running ethanol cleaning"</i> (page 177)

Periodic Maintenance

Maintenance item	Schedule	See
Changing the PEEK sample line or the sample line	• Every 3 months • As required	<i>"Changing the PEEK Sample Line (PEEK sample line compatible models)"</i> (page 189)
Autoclaving the probe adapter and sample probe	• Every 3 months • As required	
Changing the sheath filter	• Every 3 to 12 months • Every 1 to 3 months (when ethanol cleaning of whole fluidics system) • As required	<i>"Changing the Sheath Filter"</i> (page 196) <i>"Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))"</i> (page 212)

Maintenance item	Schedule	See
Changing the DI water filter	• Every 3 to 12 months • Every 1 to 3 months (when ethanol cleaning of whole fluidics system) • As required	<i>"Changing the DI Water Filter"</i> (page 198) <i>"Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))"</i> (page 212)
Changing the sample loader O-ring	• Every 3 to 6 months	<i>"Changing the Sample Loader O-Ring"</i> (page 181)
Changing the sample probe	• Every 12 months	<i>"Changing the Sample Line and Probe (Non-PEEK sample line compatible models)"</i> (page 192)
Changing the sintered sheath line filter	• Every 12 months	<i>"Changing the Sintered Sheath Line Filter"</i> (page 200)
Changing the DI water tank and ethanol tank air filter	• Every 12 months	<i>"Changing Tank Air Filters"</i> (page 186)
Changing the waste catcher O-ring	• Every 12 months	<i>"Cleaning the Waste Catcher"</i> (page 176)

Miscellaneous

Maintenance item	Schedule	See
Cleaning the collection tube holder with embedded cooling agent (5 ml)	• As required	<i>"Cleaning the Collection Tube Holder with Embedded Cooling Agent (5 ml)"</i> (page 171)
Cleaning the deflection plates	• As required (when dirty)	<i>"Cleaning the Deflection Plates"</i> (page 174)
Cleaning the waste catcher	• As required (when dirty)	<i>"Cleaning the Waste Catcher"</i> (page 176)
Running Waste Line A maintenance	• After a week or more of non-use • As required	<i>"Running Waste Line A Maintenance"</i> (page 178)
Running partial ethanol cleaning or ethanol cleaning of all fluidics lines	• When a cleaning reminder is displayed on the login screen	<i>"When a Cleaning Reminder is Displayed"</i> (page 180)

Maintenance item	Schedule	See
Cleaning the sample loader O-ring	• As required	<i>“Changing the Sample Loader O-Ring” (page 181)</i>
Cleaning the sheath tank and sheath line	• As required	<i>“Cleaning the Sheath Tank and Sheath Line” (page 183)</i>
Releasing air in the sheath filter and sorting chip	• As required	<i>“Releasing Air in the Sheath Filter and Sorting Chip” (page 202)</i>
Releasing air in the DI water filter	• As required	<i>“Releasing Air in the DI Water Filter” (page 203)</i>
Autoclaving the sheath filter and DI water filter	• As required	<i>“Autoclaving the Sheath Filter and DI Water Filter” (page 204)</i>
Ejecting the sorting chip manually	• As required	<i>“Ejecting the Sorting Chip Manually” (page 206)</i>
Cleaning the optical filters	• As required	<i>“Cleaning and Handling of Optical Filters” (page 207)</i>
Disconnecting and reconnecting the fluidics cart	• As required	<i>“Disconnecting and Reconnecting the Fluidics Cart” (page 208)</i>
Cleaning the internal sheath and DI water line with ethanol	• As required	<i>“Cleaning the Internal Sheath Line and DI Water Line using Ethanol” (page 209)</i>

Fluidics Cart


Status of fluidics cart tank

The status of the three fluidics tanks in the fluidics cart is displayed on the LCD monitor, in addition to the Hardware Status pane in the bottom right corner of the main window in Cell Sorter Software.



Fluidics cart lock mechanism

The fluidics cart is fitted with a simple lock to prevent the tray from being withdrawn during operation. The cart should only be unlocked when performing fluidics cart maintenance. The cart can be locked/unlocked using a coin, or similar flat object.

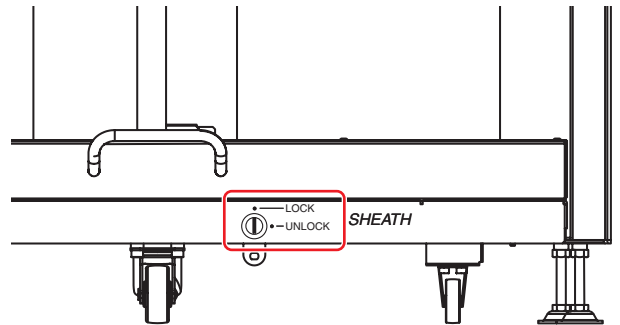

Refilling the Sheath Tank

The sheath tank should always be checked before starting operation. Refill the tank when the level is insufficient. Also, the tank must be refilled when prompted to do so in Cell Sorter Software.

Notes

- Sheath fluid is a sterile solution. Avoid contact with the contents of the tank to maintain sheath fluid sterility.
- Never use the supplied sheath tank for any other purpose.
- When opening the sheath tank cap, check that “Standby” is displayed on the LCD monitor or that the power to the main unit is turned off. If “Standby” is not displayed, place the unit in Standby mode using the following procedure.
 - ① Click [Settings] on the [Cytometer] tab of the ribbon.
The [Cytometer Settings] dialog appears.
 - ② Click [Advanced Settings].
The [Advanced Settings] dialog appears.
 - ③ Click [Standby] on the [Pressure Options] tab.
 - ④ Check that “Standby” is displayed on the LCD monitor.
- Do not force the lid down or place objects on the lid when the tank is pressurized.
- Store replacement sheath fluid in a room with the same temperature environment as the instrument, or allow the sheath fluid to acclimatize in the same room as the instrument before refilling the sheath tank. Droplet formation may become unstable if the sheath fluid added to the tank is a different temperature to the sheath fluid remaining in the tank.

Note

It is recommended that the sheath tank be filled one day before measurement to ensure all air bubbles are removed from the sheath fluid.

- 1 Pull out the fluidics cart tray.
- 2 Disconnect the sheath air line (clear) from the top of the sheath tank.



Note

It is important to disconnect the sheath air line before disconnecting the sheath fluid line.

- 3 Release the residual air pressure from the tank using the ring-pull air relief valve, then remove the tank from the tray.



- 4 Disconnect the sheath fluid line (blue) from the top of the tank.



- 5 Lift the lever securing the lid of the tank to open the lid.




- 6 Lower the lid slightly into the tank, turn it 90 degrees, and then remove the lid from the tank.



- 7 Refill the tank with sheath fluid.

Notes

- Do not touch the surfaces of the sheath tank lid that come into contact with sheath fluid. Take care not to spill any fluid.
- Do not use a pump or other object that has not been sterilized for filling the tank. It is recommended that the tank be placed immediately beneath the tap of the source sheath fluid container so that the tank is filled using gravity alone.
- Overfilling the sheath tank may damage the unit. Do not fill the tank beyond the full line inside the tank.

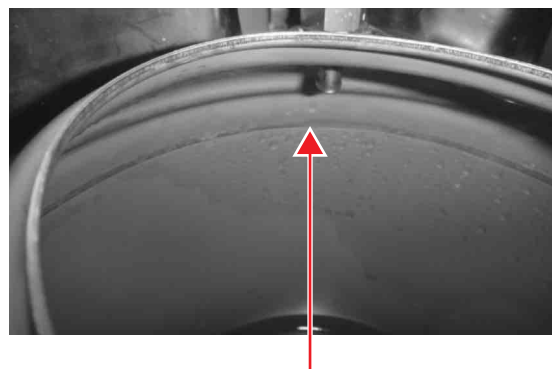

Fill with sheath fluid to below this line.

- 8 Attach the lid, and place the tank back on the fluidics cart tray.

Note

When inserting the sheath tank lid, make sure that the black rubber seal is attached to the lid.

- 9 Wipe off any residue liquid or dirt attached to the connector, and connect the sheath air line and sheath fluid line, in the reverse order of removal, until they lock into place.

- 10 Carefully clean any residue fluid from the outside surfaces of the tank and stow the tray.

Note

Always place the tank in the center of the tray area. The fluid level detector may not operate correctly if the tank is placed off-center.

- 11 If the sheath tank was refilled while the instrument is in use, click [Settings] on the [Cytometer] tab of the ribbon to display the [Advanced Settings] dialog, and then click [Ready] on the [Pressure Options] tab.

Note

To continue sorting, click [Sort Calibration] on the [Cytometer] tab of the ribbon to run auto calibration.

For details, see “Automatic Calibration” (page 50).

Emptying/Changing the Waste Tank

Check the fluid level in the waste tank daily before starting the unit. If it is full, dispose of the waste fluid or replace the tank with a new tank. Also, the waste fluid must be disposed or the waste tank must be replaced when prompted to do so in Cell Sorter Software.

Note

When handling the waste tank, always follow the laboratory rules for the handling of waste fluids.

DANGER

Waste fluid may contain virulent or hazardous substances that may be highly contagious or toxic, depending on the sample under test. Always wear gloves and other protective clothing, mask, and goggles, as required, when handling waste fluid.

- 1 Check that the SH800 main unit is turned off, or check that “Standby” is displayed on the LCD monitor.
- 2 Pull out the fluidics cart tray.
- 3 Unscrew the cap on the top of the waste tank and remove the waste tank nozzle.




- 4 Place the cap (with nozzle, connector, and line attached) in a beaker or other container to prevent waste fluid spills from the nozzle.



Notes

- Do not touch the surfaces of the waste tank cap and tank nozzle that come into contact with waste fluid. Take care not to spill any fluid.
 - If the beaker is moved or knocked over, the waste liquid will spill out and contaminate the area. Place the beaker in a safe location where it cannot be moved or knocked over.
- 5 Attach the inner cap on the tank, taking care not spill any waste fluid, then attach the outer cap and securely tighten.

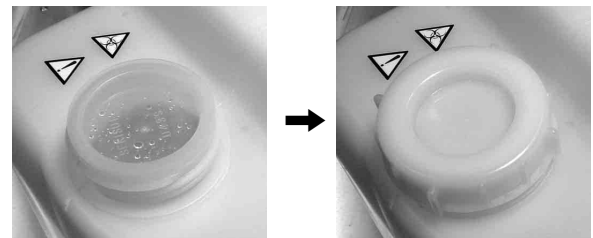

The inner and outer stopper caps are supplied with the tank. Always attach both the inner and outer caps to ensure the tank is sealed.

- 6 Remove the waste tank from the fluidics cart tray.
- 7 Empty the tank or prepare a new waste tank.

Note

Dispose of the waste tank and fluid in accordance with the laboratory rules and local ordinances and regulations.

- 8 Place an empty waste tank or a new tank on the fluidics cart tray.

- 9 Insert the waste tank nozzle from the beaker into the tank, securely tighten the cap, and stow the tray.

Notes

- Make sure that the waste fluid line is connected securely to the connector on the cap of the waste tank, and that the cap is securely fastened on the waste tank to prevent leakages.
- Always place the tank in the center of the tray area. The fluid level detector may not operate correctly if the tank is placed off-center.
- Do not autoclave fluid tanks. The shape of the fluid tanks will deform if placed in an autoclave.
- If the air filter attached to the waste tank becomes blocked, air cannot escape the waste tank as it fills with waste fluid.
- If the waste tank air filter gets wet, the filter may become clogged, causing the pressure inside the tank to rise because the air cannot escape, and may cause malfunction of the unit. If the waste tank is full, do not tilt or otherwise jostle the waste tank to prevent the filter from getting wet.

Refilling the DI Water Tank


The DI water tank should be refilled daily before starting the unit. The tank should also be refilled as required if the level becomes too low during operation and before cleaning with ethanol.

Tip

The DI water tank can be refilled at any time, except during auto calibration, measurement, and cleaning.

Precautions when handling DI water



- Refill the DI water tank with fresh DI water every time when starting up, replacing any unused DI water in the tank.
- Regularly clean the DI water tank with ethanol to suppress the growth of bacteria (*page 187*).
- Use ultra pure water or distilled water with extremely little or no bacteria or biological content to further suppress the growth of bacteria.

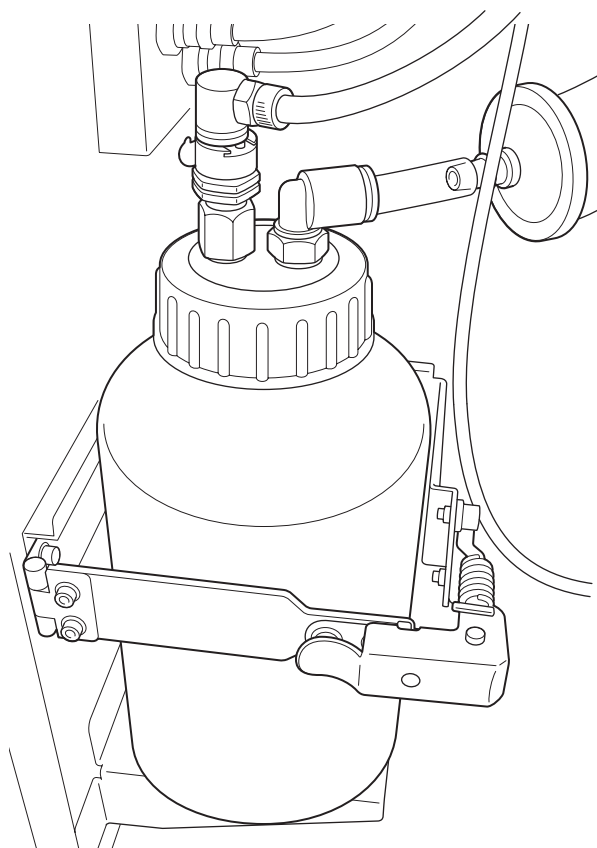

- 1 Check to make sure that auto calibration, data acquisition, or cleaning are not in progress.

- 2 Open the fluidics maintenance door, then disconnect the line from the top of the DI water tank.

Press the metal release catch to withdraw the connector.

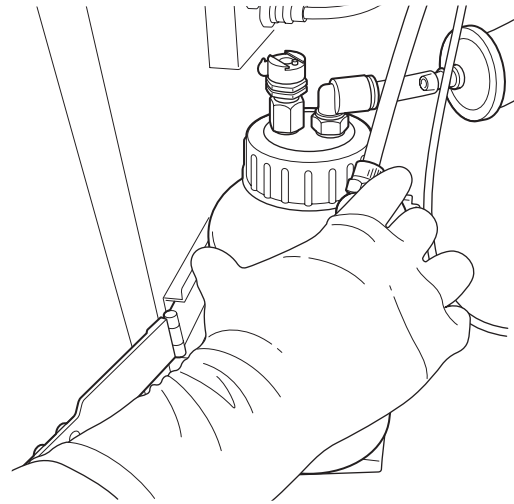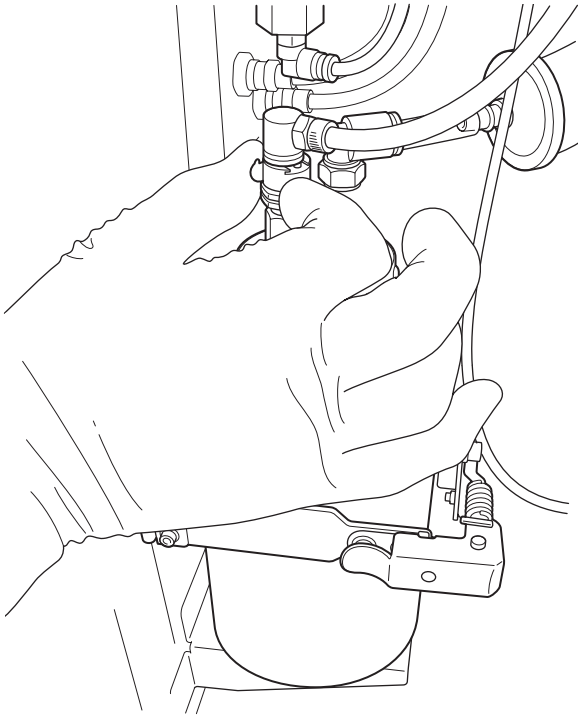

Note


Contact with skin cells and hair can cause bacterial contamination. When handling fluidics system components, always wear biology laboratory gloves and other protective clothing, mask, and goggles to prevent contamination.

- 4 Remove the cap, and withdraw the tank probe.
- 5 Discard any remaining DI water.


Note

Dispose of DI water in accordance with laboratory rules, and with local and federal ordinances and regulations.

- 3 Undo the metal restraining clamp and remove the DI water tank.

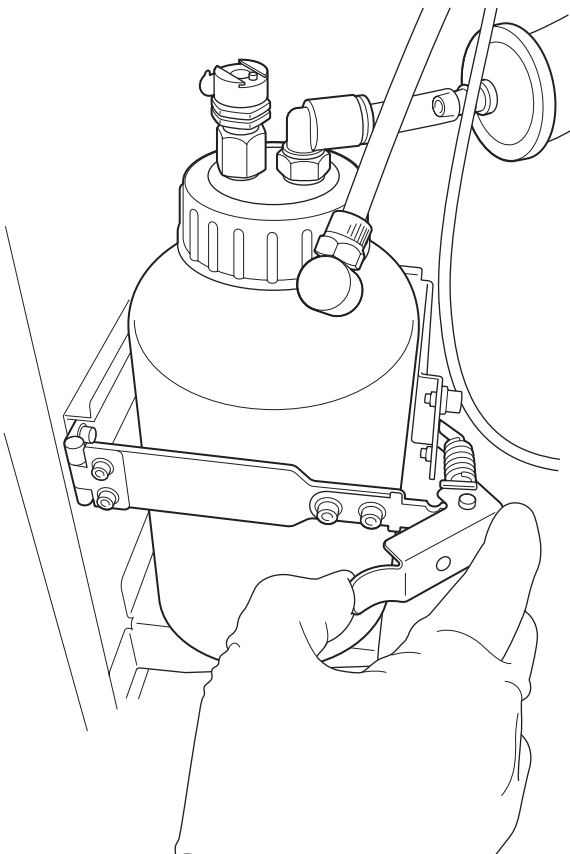

- 6 Refill the tank with fresh DI water up to the opening of the tank, and reattach the cap.
- 7 Re-install the DI water tank.
Install in the reverse order of removal.
- 8 Check the drip tray for any fluid spills or leakages, and clean as necessary.

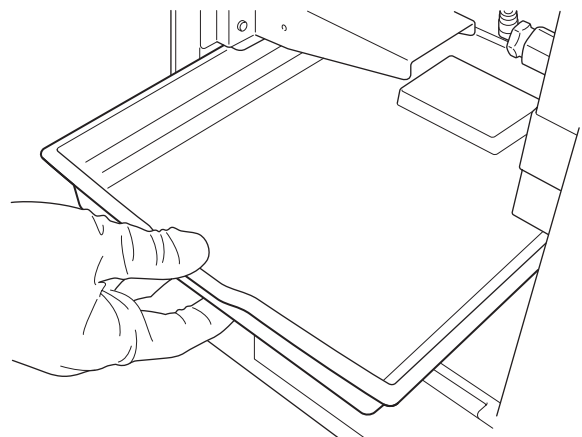

9 Close the door, and turn on the power supply.

Note

The DI water tank is fitted with a filter that should be replaced every 12 months. For details, see “*Changing Tank Air Filters*” (page 186).

Refilling the Ethanol Tank

The level of fluid in the ethanol tank can be monitored in the main window in Cell Sorter Software and on the LCD monitor.

Precautions when handling ethanol

- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Never pour ethanol into a tank other than the ethanol tank, and conversely, never pour other liquids into the ethanol tank.
- Make sure that the cap on the ethanol tank is correctly and securely fastened at all times.
- Make sure that the ethanol line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the ethanol tank itself.
- When the ethanol tank is not in use, remove all ethanol from the tank.

Note

The Cell Sorter is compatible with standard disinfectant ethanol (70%).

- 1** Pull out the fluidics cart tray.
- 2** Disconnect the ethanol line from the cap on the ethanol tank, and place it in a clean beaker or other container.



Press the metal release catch to withdraw the connector.

Note

Do not touch the surfaces of the ethanol connector that come into contact with ethanol. Take care not to spill any fluid.

- 3 Remove the ethanol tank from the fluidics cart for easy access to the cap of the tank.
- 4 Remove the cap and tank probe from the ethanol tank and place in a clean, safe location.



- 5 Refill the ethanol tank with the specified volume of ethanol, and reattach the cap.
- 6 Place the tank in the fluidics cart tray, connect the ethanol line, and stow the tray.

Insert the line connector and push until it locks into place. Carefully Wipe off any residue fluid from the connector and outside surfaces of the tank.

Cleaning the Sample Fluidics System using Bleach

The sample line and probe should be cleaned regularly using a protein-dissolving bleach to disinfect the lines that carry the sample fluid. The fluidics system should be disinfected on a regular basis, or as required, especially if running samples that contain, for example, infectious substances.

The sorting chip, PEEK sample line, sample line, and sample probe can be disinfected using a 1 to 3% Sodium Hypochlorite (bleach) solution.

Note

Cleaning the sample probe using a high-concentration bleach solution (5% or higher) may cause rusting or discoloration. Use normal concentration, aqueous solution.

- 1 In Cell Sorter Software, click [Bleach Cleaning] on the [Cytometer] tab of the ribbon.

Note

The bleach cleaning sequence must be followed to completion once it is started.

- 2 Load a 15 ml conical sample tube containing 10 ml of 1 to 3% Sodium Hypochlorite (bleach) solution in the sample loader.

Tip

When performing bleach cleaning during measurement, perform the cleaning sequence once using a 15 ml conical sample tube. When performing bleach cleaning before shutdown at the end of a day, perform the cleaning sequence three times using a 15 ml conical sample tube or once using a 30 ml cleaning tube.

- 3 Click [Start] to begin the cleaning process.

The solution passes through the sample probe, sample line, and sorting chip.

A progress bar is displayed during cleaning. Abort the process, if necessary, before the level of bleach in the sample tube falls below the tip of the sample probe.

- 4 When finished, remove the tube.
- 5 Click [DI Rinse] on the [Cytometer] tab of the ribbon.

- 6 Load a 15 ml conical sample tube containing 12 ml of DI water in the sample loader.

To prevent the accumulation of excess bleach, the level of DI water must exceed the level of bleach added in step 2.

- 7 Click [Start] to begin the rinse process.

Stop the DI rinse process, if necessary, before the level in the sample tube falls below the tip of the sample probe.

- 8 When finished, remove the tube.

- 9 Click [Probe Wash] on the [Cytometer] tab of the ribbon to flush the sample line and probe.

Notes

- Always use tubes with the specified fluid volume when using the Bleach Cleaning function. If the volume of fluid is insufficient, the cleaning fluids may be ejected from the sorting chip in a spray, contaminating the optical lenses and other components.
- If the main unit shuts down during waste line cleaning, always restart the main unit and run [DI water rinse]. If bleach solution remains in the instrument, it may damage the main unit.

Cleaning the Collection Tube Holder with Embedded Cooling Agent (5 ml)

Use the following procedure to clean the collection tube holder with embedded cooling agent (5 ml) if the tube holder is used and sample fluid spills onto the tube holder.


WARNING

Waste fluid may contain virulent or hazardous substances that may be highly contagious or toxic, depending on the sample under test. When handling waste fluid, always wear biology laboratory gloves and other protective clothing, mask, and goggles.


Note

Do not sort sample fluid above the permitted volume into the tube and do not overfill the tube with buffer solution to prevent sample fluid overflow from the tube.

- 1 Remove the screws (four locations) on both sides of the collection tube holder with embedded cooling agent (5 ml).

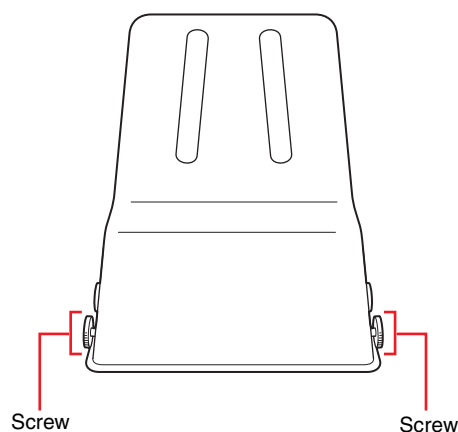

- 2 Remove the cover.

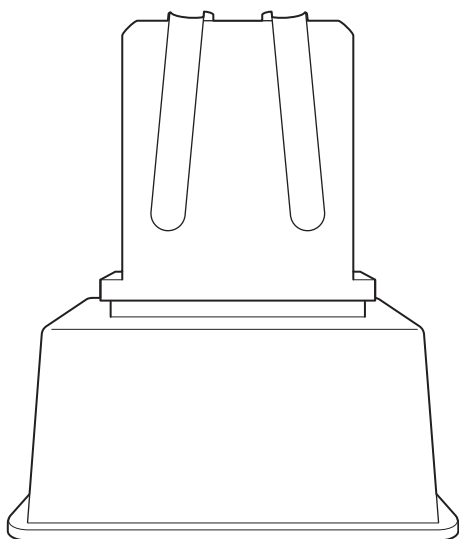

- 3 Wipe off any contamination.
- 4 Reattach the cover, and secure using the screws.

Note

Reattach the cover so that the grooves for the sample tubes align with the cutouts in the cover.

General Cleaning

The collection area and sample loader should be cleaned regularly to prevent the precipitation of salts from the sheath fluid and to prevent the spread of biological matter. Cleaning the instrument regularly also prevents trouble in the cultivation process after sorting.

Clean the side stream monitor window, laser windows, collection stage, collection tube holders, and sample loader using water and ethanol.

Notes

- Be aware of the door edges when doors are open.
- Always wear gloves and other protective clothing, mask, and goggles.

- 1 Turn on the power supply using the POWER/STANDBY button on the front panel.
- 2 Press the PUSH OPEN collection area door button on the front panel to access the collection area.
- 3 Moisten a clean, soft, lint-free laboratory-use cloth with clean water, and wipe with the cloth.


Side stream monitor window and laser windows

Carefully clean the side stream monitor window and laser windows to prevent reflections entering the side stream monitor. If the windows become dirty, an error may occur during automatic calibration.

Tip

The number of laser windows may vary, depending on the model.

Side stream monitor window

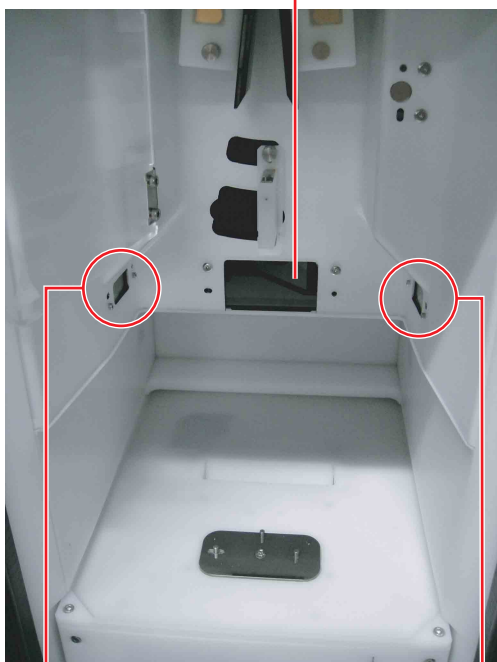

Laser windows

Laser windows

Collection stage and collection tube holders

Clean the collection stage and collection tube holders carefully, especially the upper portion of the collection stage, where the unit is most exposed to salts and biological matter.



Transparent protective covers in the collection area

Wipe the inner surfaces of the transparent protective covers.

Moisten a clean, soft, lint-free cloth with water, and wipe the covers.



Transparent protective covers

Note

Wiping with a dry cloth can generate static electricity, which may adversely affect the side streams during auto calibration. Never use a dry cloth.

Sample loader

Clean carefully to remove all salt products.



4 Moisten a clean, soft, lint-free laboratory-use cloth with clean water, and wipe the sample tube loader.

5 Spray a clean, soft, lint-free laboratory-use cloth with ethanol 70% from a spray bottle, and wipe with the cloth.

Notes


- Do not operate any hardware or software controls while cleaning the sample loader or the collection area.
- When cleaning the sample loader, remove the sample tube holder and clean separately.
- When cleaning the collection area, pull the collection stage out to its maximum travel and remove the collection tube holder and clean separately.
- When cleaning the collection area, also remove the deflection plates (*page 174*) and waste catcher (*page 176*), and clean separately.

Checking for leaks

- 1 Check that there are no fluid leaks in the drip tray inside the fluidics maintenance area.
- 2 Check that there are no leaks in the fluidics cart.

Cleaning the splash guard


- 1 Moisten a soft, lint-free laboratory use cloth with water and wipe the splash guard.
Take care to clean right into the edges and corners.

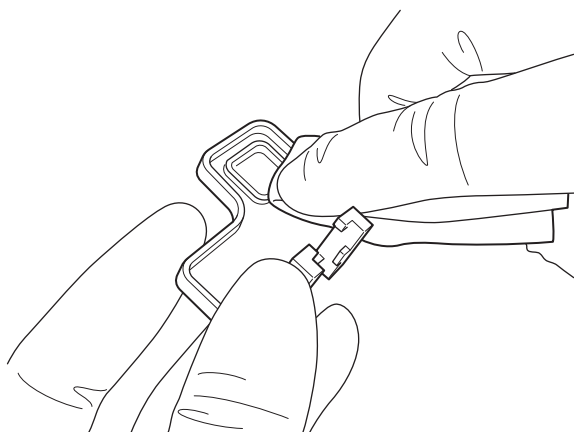

- 2 Moisten a clean, soft, lint-free cloth with water, and wipe the guard.
- 3 Spray a clean, soft, lint-free cloth with ethanol 70% from a spray bottle, and wipe the guard.

Cleaning the Deflection Plates


The deflection plates in the collection area should be removed and cleaned when cleaning the collection area to maintain sorting performance.

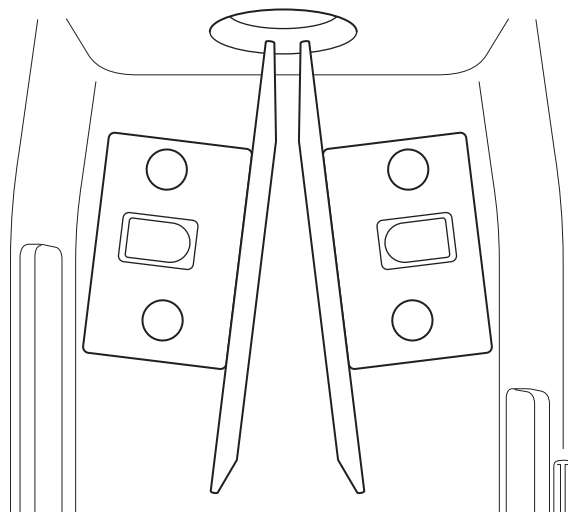


- 1 Check that the main unit is shut down.
- 2 Open the collection area door on the front panel.
- 3 Pull open the transparent safety cover to access the deflection plates.

The cover is held closed by a magnetic latch.


- 4** Remove the two plastic bolts securing each deflection plate by hand, and remove the deflection plates.

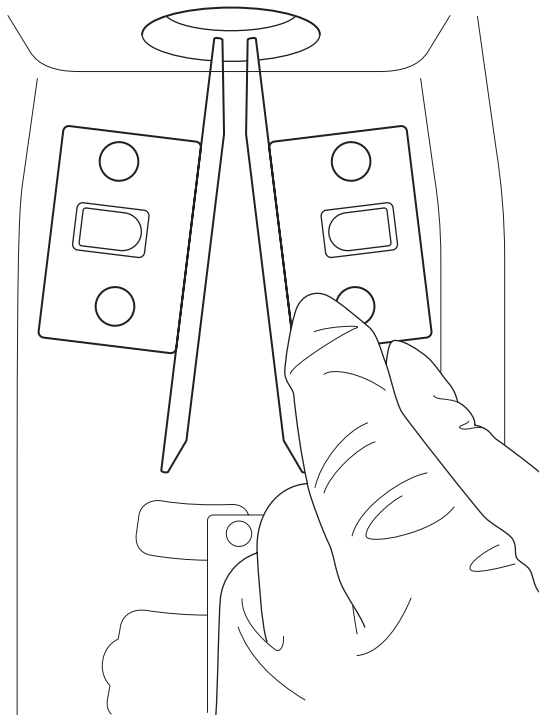

- 5** Clean the deflection plate mounting plate using a lint-free laboratory-use cloth moistened with water.

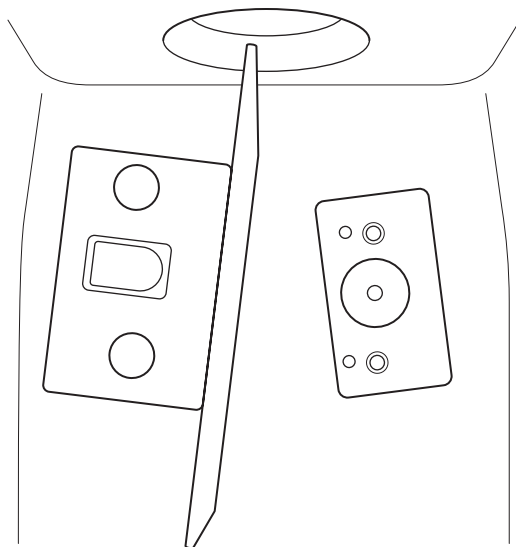


- 6** Clean the deflection plate.

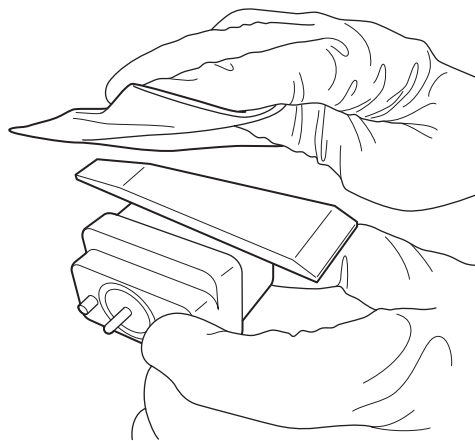

Note

The electrode surfaces of the deflection plates have been treated with a water-repellent coating. Wiping salt crystal deposits with a cloth may damage the coating. Wash the electrode surfaces with water to dissolve salt crystals, then wipe clean.

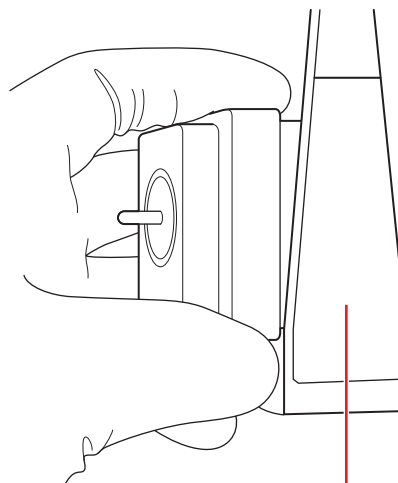

Electrode surface

- 7** Clean the deflection plate plastic bolts.
- 8** Spray a lint-free laboratory-use cloth with ethanol and clean the items in steps **5** to **7**.
- 9** Insert the deflection plate in the correct orientation.
- 10** Insert the two plastic bolts and tighten by hand.


Note

Do not overtighten using tools.


- 11** Close the transparent safety cover.

Cleaning the Waste Catcher


The waste catcher in the collection area should be removed and cleaned when cleaning the collection area to ensure the area is not contaminated between sorts.

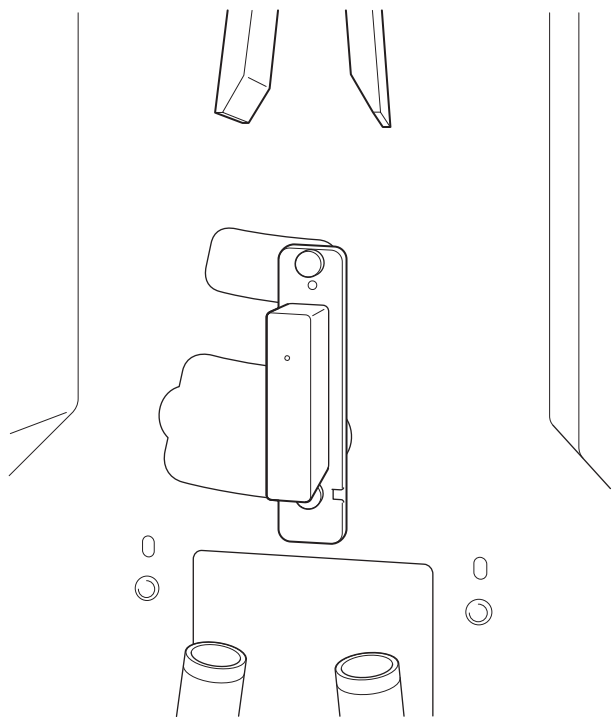


- 1 Check that the main unit is shut down.
- 2 Open the collection area door on the front panel.
- 3 Remove the two screws securing the waste catcher, and remove the catcher.

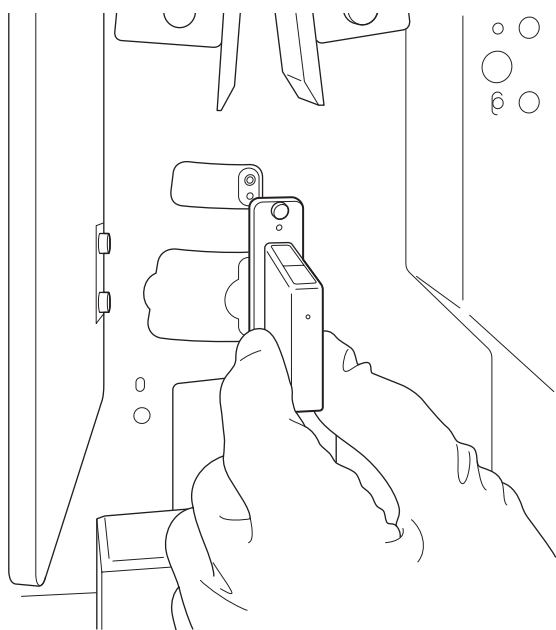

- 4 Clean the waste catcher mounting bracket, holes, and screws using a lint-free laboratory-use cloth moistened with water.

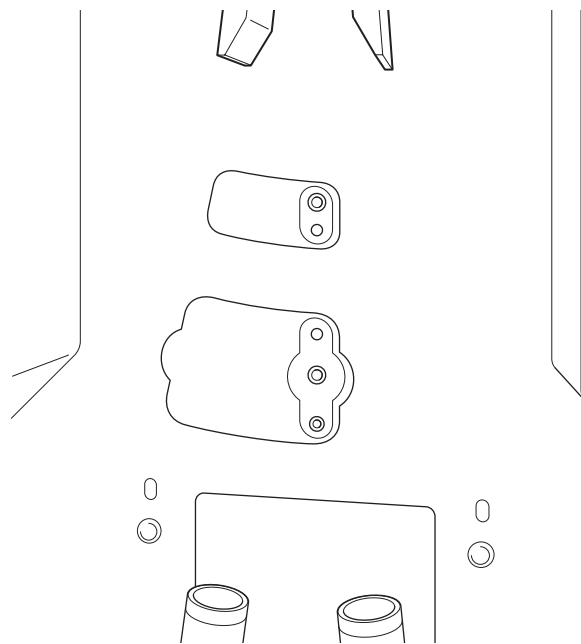

- 5 Clean the waste catcher.

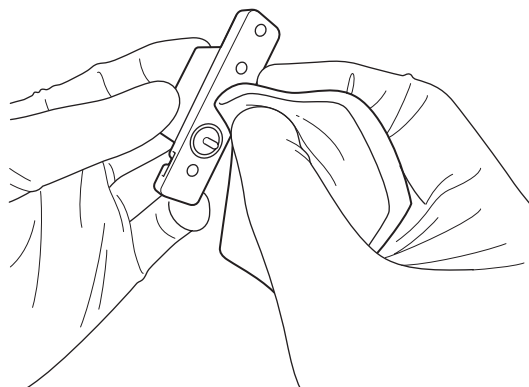

- 6 Spray a lint-free laboratory-use cloth with ethanol and clean the items in steps 4 and 5.

Wipe any moisture off the screws using a dry, lint-free laboratory-use cloth.



- 7 Reinstall the waste catcher.

Notes

- Electrostatic charge may accumulate in the collection area after cleaning the plastic covers with a cloth that is too dry. This charge can adversely affect the side stream position, causing an error in the automatic side stream calibration process. If this occurs, wipe the areas around the deflection plates with a cloth soaked in DI water to dissipate any accumulated electrostatic charge.
- Make sure the waste fluid outlet and O-ring of the waste catcher are correctly aligned with the port on the mounting plate, and the waste catcher is securely fastened to the mounting plate. There is a risk of waste

fluid leakage if not fitted correctly (especially, the O-ring), or risk of overflow if the port is obstructed.

- The O-ring on the rear of the waste catcher must be replaced every 12 months.
- Check that the waste catcher screw is not wet when tightening the screw. A malfunction may occur if the screw is wet when installed.

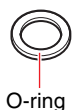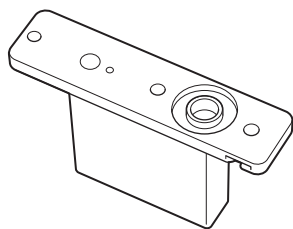

Running Maintenance Mode

The following maintenance can be performed in maintenance mode.

- Ethanol cleaning
- Ethanol cleaning of all fluidics lines
- Waste A maintenance (bleach cleaning)
- Changing the sample line

You can run maintenance mode from the Cell Sorter Software login screen.

Precautions when cleaning with ethanol


- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Do not pour any fluid other than the specified fluid into the fluidics system tanks.
- Make sure that the cap on the ethanol tank is correctly and securely fastened at all times.
- Make sure that the ethanol line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the ethanol tank itself.
- Make sure that the cap on the waste tank is correctly and securely fastened at all times.
- Make sure that the waste fluid line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the waste tank itself.

Note

The SH800 is compatible with standard disinfecting agents (ethanol 70%).


Running ethanol cleaning

- 1 Click [Maintenance] on the bottom right of the Cell Sorter Software login screen.

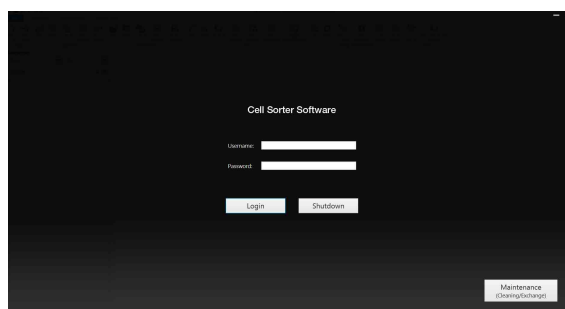

The [Maintenance] wizard appears.


- 2 Select [Ethanol cleaning], then click [Start].

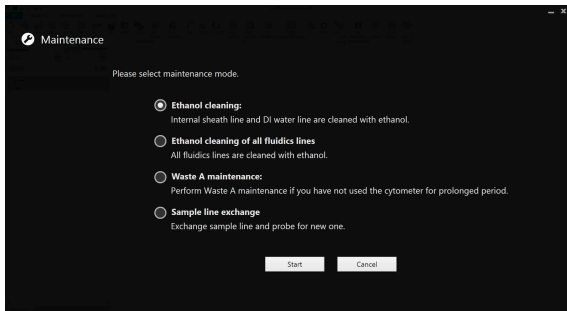

The [Ethanol Cleaning] wizard appears.

- 3 Start ethanol cleaning.

Follow the on-screen instructions.

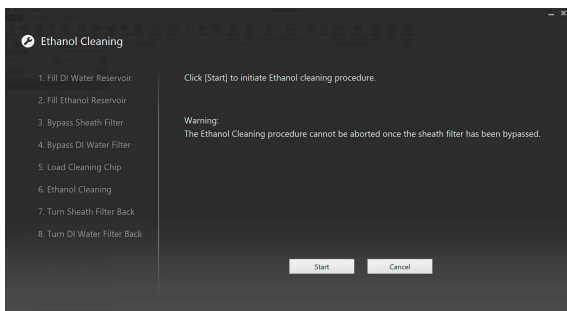


For details about ethanol cleaning, see “Cleaning the Internal Sheath Line and DI Water Line using Ethanol” (page 209).

When ethanol cleaning finishes, the login screen reappears.


Running ethanol cleaning of all fluidics lines

- 1 Click [Maintenance] on the bottom right of the Cell Sorter Software login screen.

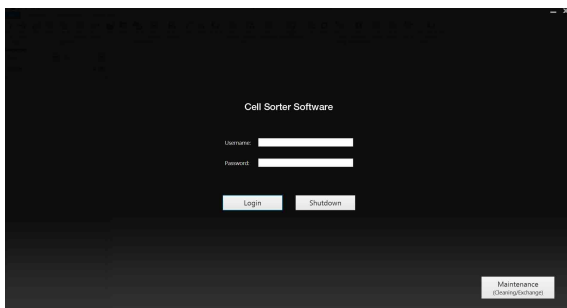

The [Maintenance] wizard appears.


- 2 Select [Ethanol cleaning for all fluidics line], then click [Start].

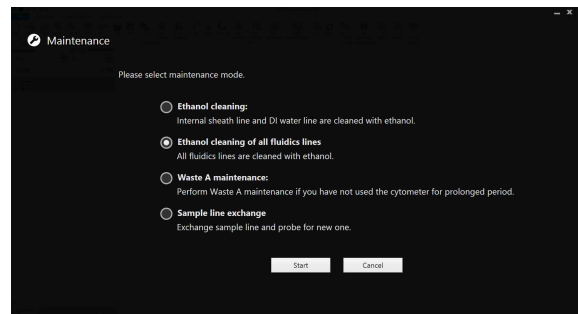

The [Ethanol cleaning for all fluidics line] wizard appears.

- 3 Start ethanol cleaning of all fluidics lines.

Follow the on-screen instructions.

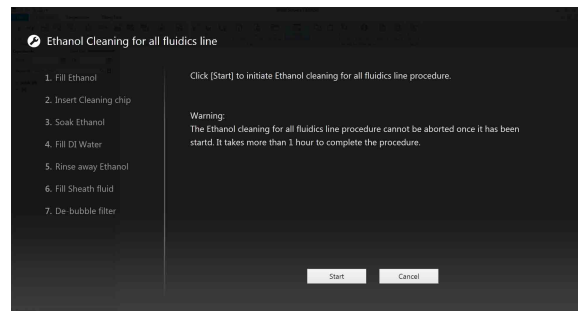

For details about ethanol cleaning of all fluidics lines, see “Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))” (page 212).

When ethanol cleaning of all fluidics lines finishes, the login screen reappears.


Tip

The sheath filter used for ethanol cleaning of all fluidics lines should be replaced once a month.

Running Waste Line A Maintenance


The instrument should be started in maintenance mode to prevent the buildup of salts and clean waste lines (Waste Line A).

- 1 Click [Maintenance] on the bottom right of the Cell Sorter Software login screen.

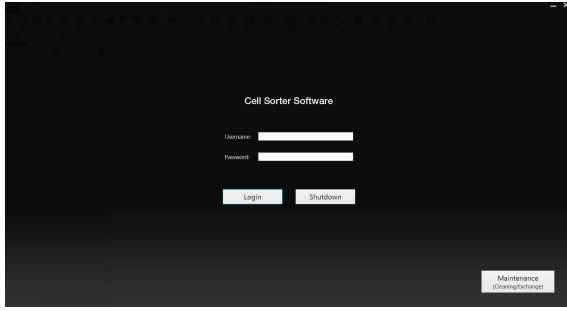

The [Maintenance] wizard appears.


- 2 Select [Waste A maintenance], then click [Start].

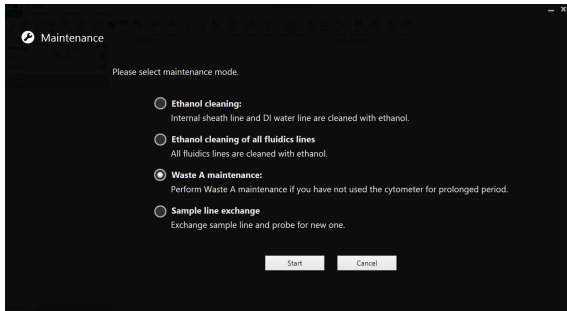

The [Waste A maintenance] wizard appears.

- 3 Start Waste A maintenance.

Follow the on-screen instructions. In Waste A maintenance, the fluidics system is cleaned using a sodium hypochlorite solution (bleach cleaning), the DI water lines are cleaned, and the internal sheath lines are cleaned, in that order.

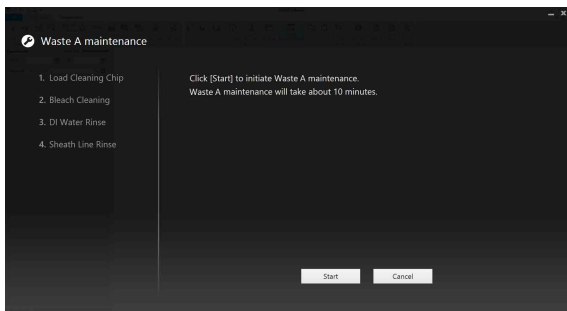


For details about cleaning using Sodium Hypochlorite solution, see “Cleaning the Sample Fluidics System using Bleach” (page 170).

When Waste A maintenance finishes, the login screen reappears.


Changing the Sample Line

- 1 Click [Maintenance] on the bottom right of the Cell Sorter Software login screen.

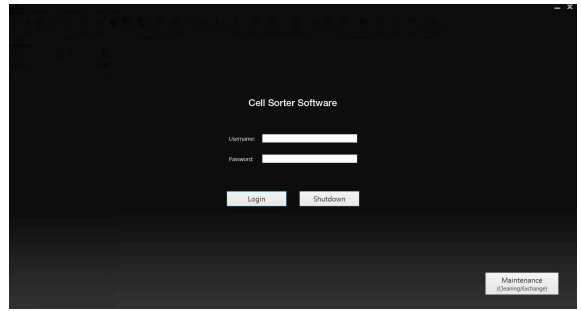

The [Maintenance] wizard appears.


- 2 Select [Sample line exchange], then click [Start].

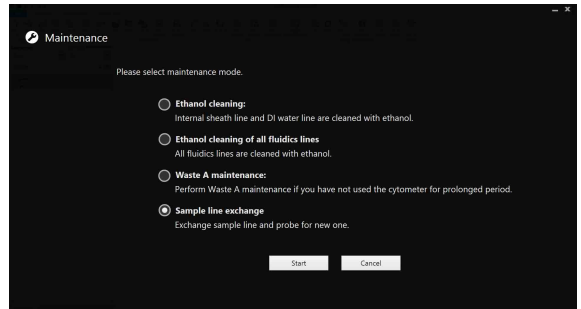

The [Sample Line Exchange] wizard appears.

- 3 Start changing the sample line.

Follow the on-screen instructions.

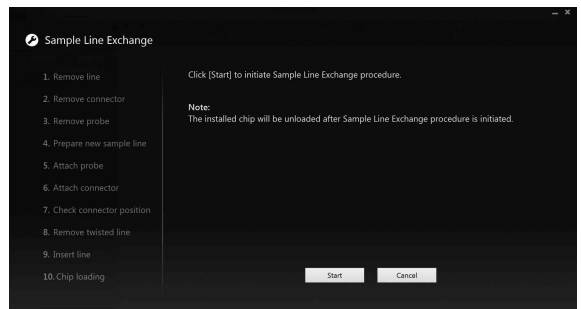


For details about changing the sample line, see “Changing the PEEK Sample Line (PEEK sample line compatible models)” (page 189) or “Changing the Sample Line and Probe (Non-PEEK sample line compatible models)” (page 192).

When the sample line has been changed, the login screen reappears.

When a Cleaning Reminder is Displayed

The instrument must be cleaned periodically using ethanol to reduce contamination and side stream instability resulting from poor system maintenance. You can set to display a message (cleaning reminder) on the login screen after a certain preset time has elapsed without performing ethanol cleaning.

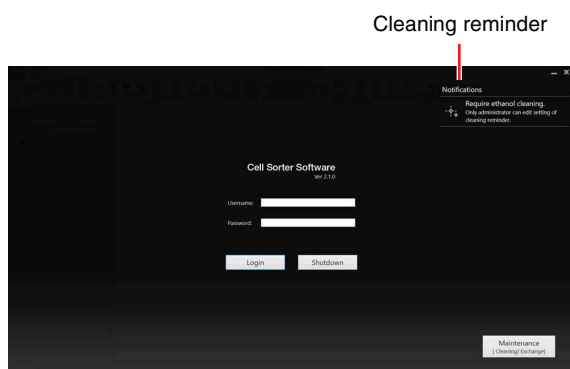

Tip

Only an administrator can set the ethanol cleaning intervals for sending cleaning reminder notifications.

For details about configuration, see “Setting a Cleaning Reminder” (page 42).


Precautions when cleaning with ethanol

- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Do not pour any fluid other than the specified fluid into the fluidics system tanks.
- Make sure that the cap on the ethanol tank is correctly and securely fastened at all times.
- Make sure that the ethanol line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the ethanol tank itself.
- Make sure that the cap on the waste tank is correctly and securely fastened at all times.
- Make sure that the waste fluid line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the waste tank itself.

Note


The SH800 is compatible with standard disinfecting agents (ethanol 70%).

- 1 Click the cleaning reminder.

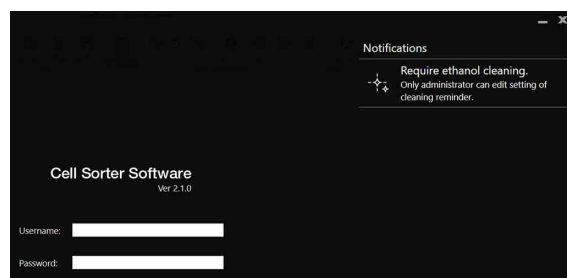

The [Cleaning Reminder] screen appears.

- 2 Select [Ethanol cleaning] or [Ethanol cleaning for all fluidics lines], then click [Start].

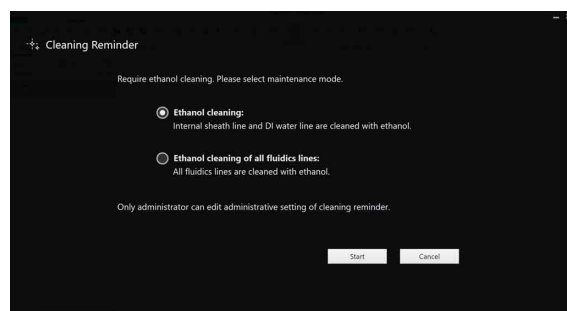

The [Ethanol cleaning for all fluidics lines] wizard appears.

Follow the on-screen instructions to clean the sheath line/DI water line or all fluidics lines with ethanol.

For details about ethanol cleaning, see “Cleaning the Internal Sheath Line and DI Water Line using Ethanol” (page 209).

For details about ethanol cleaning of all fluidics lines, see “Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))” (page 212).

When ethanol cleaning of the sheath line/DI water line or all fluidics lines finishes, the login screen reappears.


Tip

The sheath filter used for ethanol cleaning of all fluidics lines should be replaced once a month.

Changing the Sample Loader O-Ring


An O-ring is fitted into the base of the sample injection chamber that seals the chamber when the sample tube is lifted into the chamber. The O-ring should be cleaned regularly, and replaced every 3 to 6 months to ensure the integrity of the seal.

- 1 Check that the main unit is shut down.
- 2 Check that the compressed air supply is turned on, and open the flip-up door to access the sample probe.

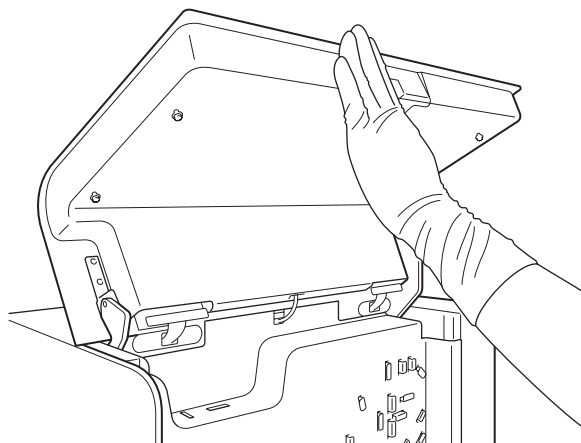

- 3 Turn off the compressed air supply.
Have an assistant help if you cannot reach the air supply control valve of the compressor supplying compressed air while holding the flip-up door open.

- 4 Lower the flip-up door slowly until it can support its own weight.

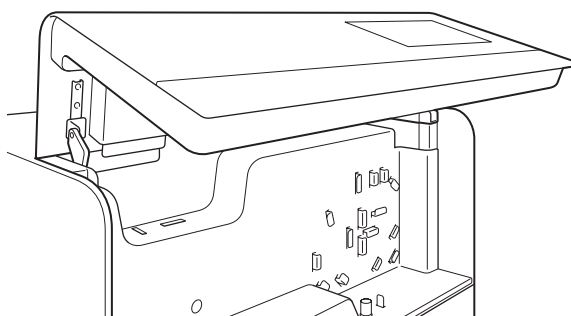

- 5 Remove the sample probe.

The sample probe is a precision component. It should be removed so that it cannot be damaged accidentally while removing the O-ring. For details, see “*Changing the PEEK Sample Line (PEEK sample line compatible models)*” (page 189) or “*Changing the Sample Line and Probe (Non-PEEK sample line compatible models)*” (page 192).

- 6 Push the bottom of the sample loader door in slightly, and then raise the door and hold it open.



If a sample tube is currently in the sample loader, remove it.

- 7 Push the sample loader waste tray (white) into its recess to provide access to the bottom of the injection chamber.



- 8 Pry out the O-ring using a small flat-blade screwdriver or a pair of tweezers.

Note


Take care not to damage the O-ring, if removing it for cleaning, or the O-ring groove.



- 9 Withdraw the O-ring.

To clean the O-ring, first wipe with a clean, soft, lint-free cloth that has been slightly moistened with water, then wipe with a similar cloth moistened with ethanol.

Also, clean the groove in the sample injection chamber and the top surface and circumference of the sample tube agitation unit that forms the seal with the O-ring.

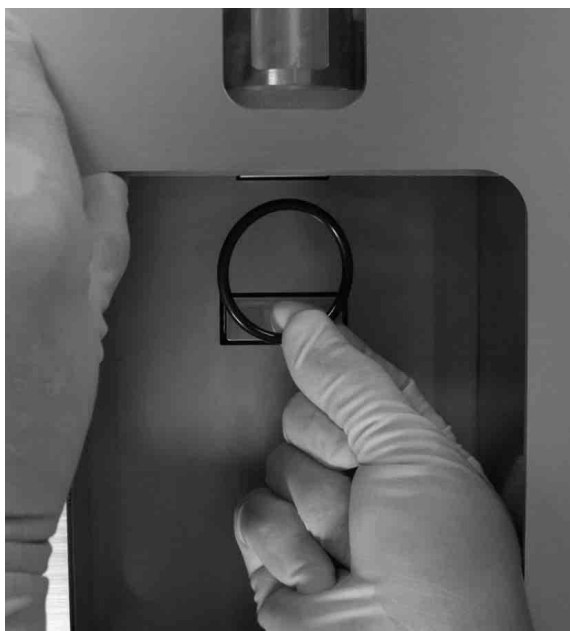

- 10 Insert a new or cleaned O-ring into the O-ring groove, and run your finger around the O-ring to ensure it is seated properly.



- 11 Lower the sample loader door and reinsert the sample probe.

- 12 Open the air supply valve on the compressor supplying compressed air.

- 13 Close the flip-up door.

- 14 Turn on the main power supply.

The sample loader waste tray will pop back out by itself.

Cleaning the Sheath Tank and Sheath Line

The sheath tank and sheath line should be cleaned regularly with ethanol to suppress the growth of bacteria.

Precautions when cleaning with ethanol

- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Never pour ethanol into a tank other than the ethanol tank, and conversely, never pour other liquids into the ethanol tank.
- Make sure that the cap on the ethanol tank is correctly and securely fastened at all times.
- Make sure that the ethanol line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the ethanol tank itself.
- Make sure that the lid on the waste tank is correctly and securely fastened at all times.
- Make sure that the waste fluid line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the waste tank itself.

Note

The SH800 is compatible with standard disinfecting agents (ethanol 70%).

- 1 Replace or empty the waste tank.

For details, see “Emptying/Changing the Waste Tank” (page 166).


- 2 Launch Cell Sorter Software.

On the Auto Calibration screen, click [Skip Auto Calibration] to bypass automatic calibration.

- 3 Click [Settings] on the [Cytometer] tab of the ribbon.


The [Cytometer Settings] dialog appears.

- 4 On the [Control] tab, click [Advanced Settings].

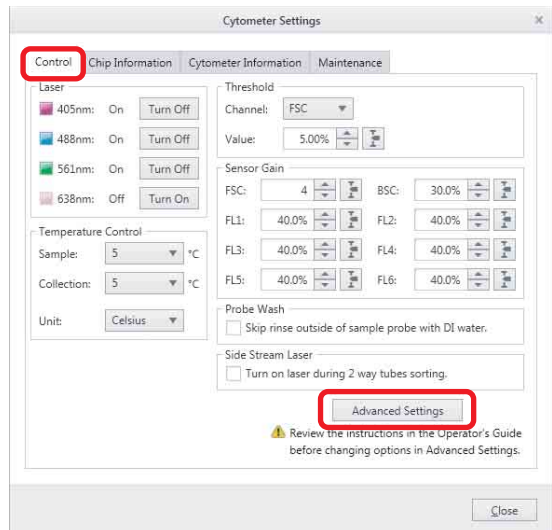

The [Advanced Settings] dialog appears.

- 5 On the [Pressure Options] tab, click [Standby] to stop the sheath fluid flow.

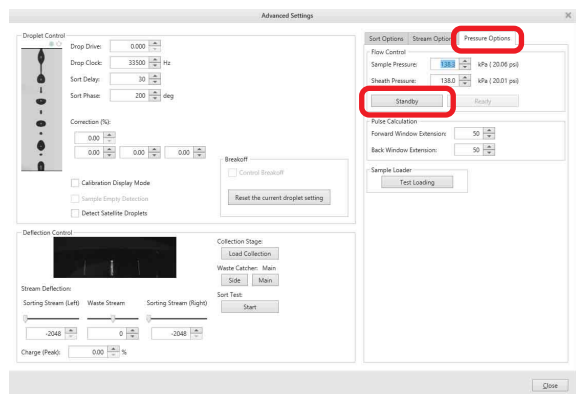

- 6 Perform steps 3 to 6 in “Refilling the Sheath Tank” (page 164) to remove the lid of the sheath tank.



7 Dispose of the remaining sheath fluid in the tank.

Note


Dispose of the sheath fluid in accordance with the laboratory rules and local ordinances and regulations.

8 Clean the sheath tank using DI water, and then pour 3 liters (0.8 gallons US) of ethanol into the sheath tank and reattach the lid.

Notes

- Do not touch the inner surface of the lid of the sheath tank. Also, take care not to spill any ethanol.
- Do not use a pump or other object that has not been sterilized for pouring ethanol into the tank.
- When inserting the sheath tank lid, make sure that the black rubber seal is attached to the lid.

9 Grasp the handles and shake the sheath tank to clean the inside of the tank (if necessary).

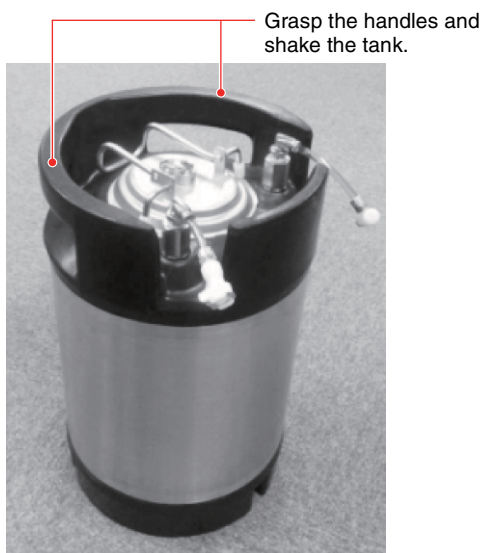

10 Place the tank in the cart tray, and connect the sheath fluid line (blue) and sheath air line (clear) in the reverse order of removal, until they lock into place.


Note

Always connect the sheath fluid line (blue) and sheath air line (clear) in that order. Connecting the sheath air line first will pressurize the sheath tank, causing ethanol to seep out of the sheath fluid line.

11 Clean any fluid from the connectors and outside surfaces of the tank and stow the tray.

Insert the line connector and push until it locks into place. Clean any fluid from the connector and outside surfaces of the tank.

12 Disconnect the sheath filter lines in the main unit, and insert the filter bypass line.

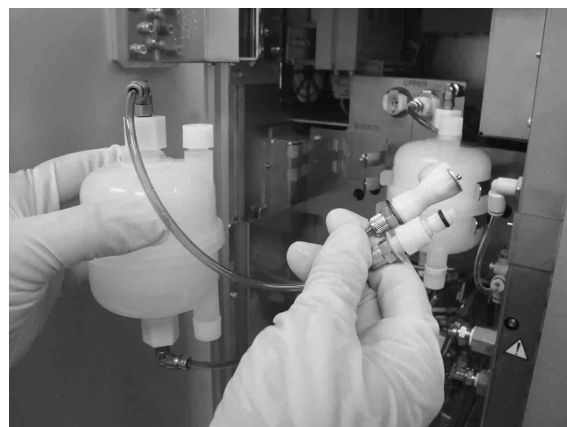

13 In Cell Sorter Software, click [Sheath Filter] in the [De-bubble] group on the [Cytometer] tab of the ribbon.

A popup window appears.

- 14 Click [Start] to begin cleaning, then click [Close] when finished.
- 15 Click [Ethanol Cleaning] in the [Shutdown] group on the [Cytometer] tab of the ribbon.

Follow the on-screen instructions to clean the instrument with ethanol.
When cleaning is completed, both the main unit and Cell Sorter Software automatically shut down.

Note

A prompt to reconnect the sheath filter appears on the Ethanol Cleaning screen. However, when cleaning the sheath tank and sheath line, leave the bypass line in place and proceed to the next step.

- 16 Dispose of the ethanol in the tank, using the same procedure in steps 5 and 6.

Note

Dispose of the ethanol in accordance with the laboratory rules and local ordinances and regulations.

- 17 Pour 4 liters (1.1 gallons US) of DI water into the sheath tank and reattach the lid.

Notes

- Do not touch the inner surface of the lid of the sheath tank. Also, take care not to spill any DI water.
- Do not use a pump or other object that has not been sterilized for pouring DI water into the tank.
- When inserting the sheath tank lid, make sure that the black rubber seal is attached to the lid.

- 18 Stow the sheath tank in the tray using the same procedure as in steps 9 and 10.
- 19 Turn on the POWER/STANDBY button on the front panel of the main unit.
- 20 Launch Cell Sorter Software.

On the Auto Calibration screen, click [Skip Auto Calibration] to bypass automatic calibration.
- 21 Click [Sheath Filter] in the [De-bubble] group on the [Cytometer] tab of the ribbon.

A popup window appears.
- 22 Click [Start] to begin cleaning, then click [Close] when finished.

Repeat this procedure at least five times to remove all traces of ethanol in the sheath fluid line.

- 23 Disconnect the filter bypass line, and reconnect the sheath filter lines.

- 24 Perform steps 2 to 6 in “Refilling the Sheath Tank” (page 164) to remove the sheath tank from the tray.

- 25 Grasp the handles and shake the sheath tank to clean the inside of the tank, as described in step 8.

- 26 Remove the lid of the sheath tank, and dispose of the DI water.

Note

Dispose of the DI water in accordance with the laboratory rules and local ordinances and regulations.

- 27 Refill the tank with sheath fluid.


For details, see step 7 and later in “Refilling the Sheath Tank” (page 165).

Changing Tank Air Filters

The DI water tank in the main unit, and the ethanol tank and waste tank in the fluidics cart are each fitted with an air filter. The filter allows air to flow into or out of the tank as the fluid level in the tank changes.

The air filters on the three tanks are connected using the same type of connector. This section describes the replacement procedure using the DI water tank as an example.

- 1** Grasp the short length of tubing with one hand, and push the quick release collet into the connector using your other hand.

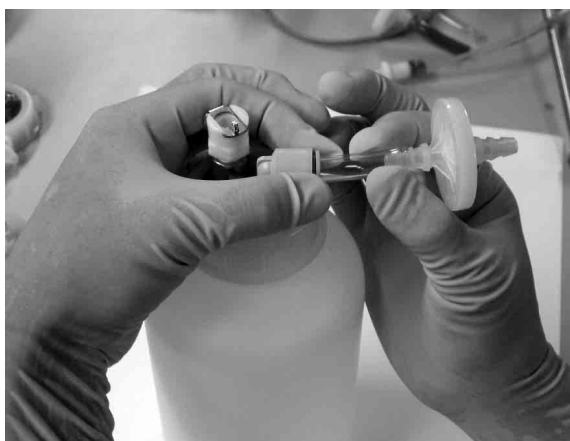

- 2** Pull the tubing out of the connector.



- 3** Insert the replacement air filter tubing into the connector and push firmly into place.

Note

The filter may become clogged if fluid gets inside the filter, adversely affecting unit operation. If any fluid enters the filter while emptying, replacing, or refilling tanks, replace the filter as soon as possible.

Cleaning the DI Water Tank and DI Water Line

The DI water tank and line should be cleaned regularly with ethanol to suppress the growth of bacteria.


To suppress the growth of bacteria, use sterile distilled water or sterile purified water as the DI water supply for normal use and cleaning.

Precautions when cleaning with ethanol

- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Do not place any fluid other than the specified fluids in the DI water tank.
- Make sure that the cap on the waste tank is correctly and securely fastened at all times.


Note

The SH800 is compatible with standard disinfecting agents (ethanol 70%).

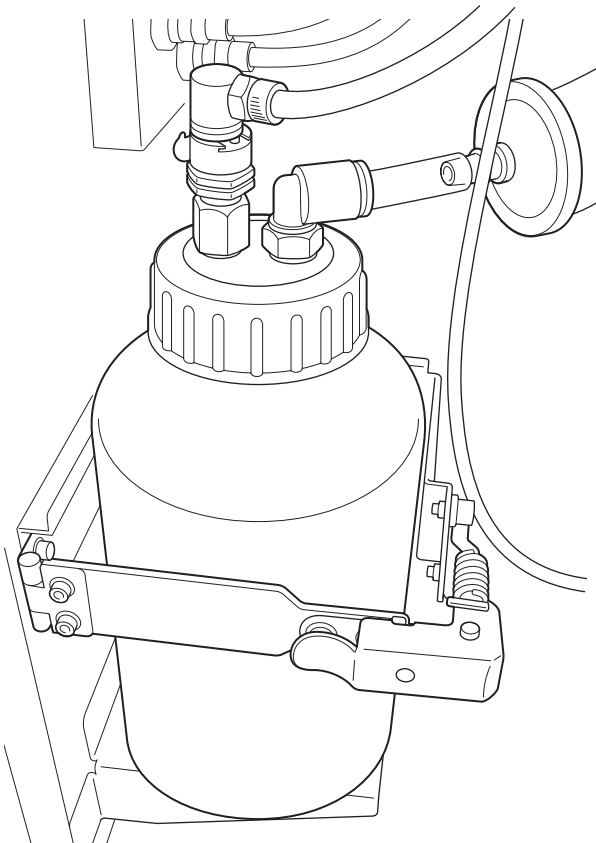



- 1 Shut down the main unit.
- 2 Open the fluidics maintenance door, then disconnect the line from the top of the DI water tank.

Press the metal release catch to withdraw the connector.

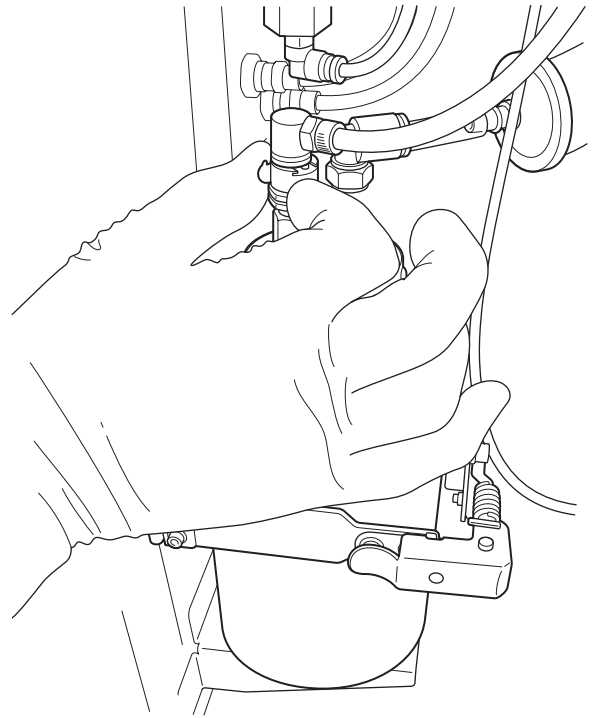

- 3 Undo the metal restraining clamp and remove the DI water tank.

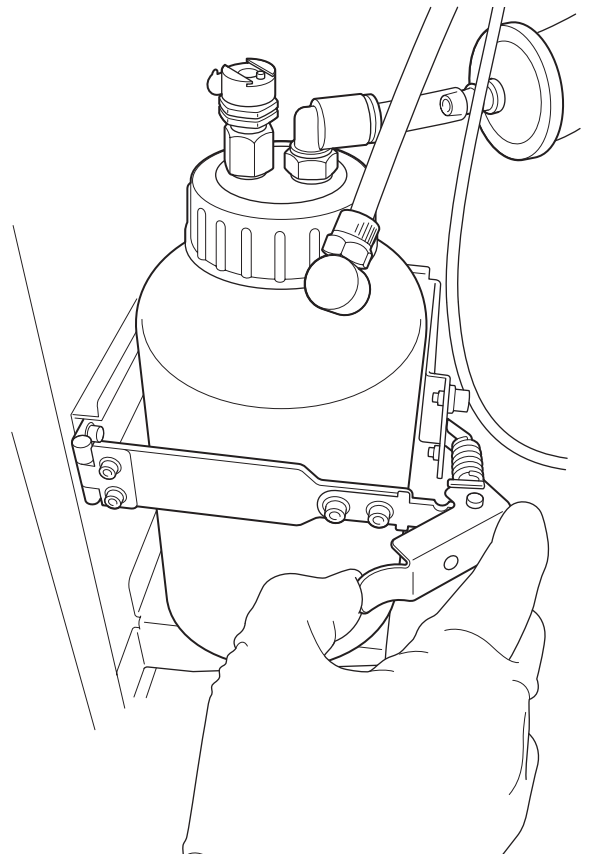

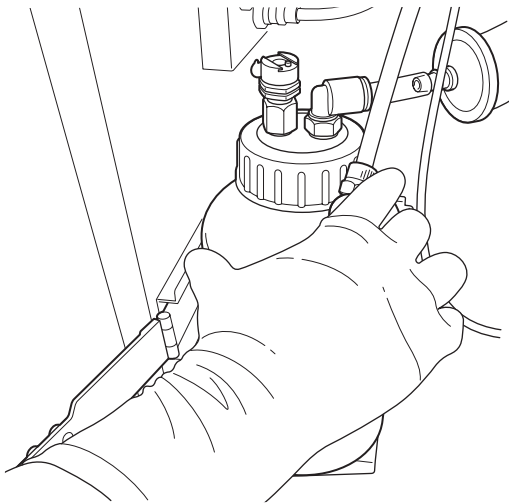

Note

Always wear gloves and other protective clothing, mask, and goggles, when handling fluidics system components. Skin cells and hair are the most common cause of bacterial contaminations.


- 4 Remove the cap, and withdraw the tank probe.
- 5 Disconnect the quick-release connector from the filter.
- 6 Discard any remaining DI water.

Note

Dispose of DI water in accordance with laboratory rules, and with local and federal ordinances and regulations.

- 7 Pour 0.5 liters (0.15 gallons US) of ethanol into the DI water tank, and reattach the cap.
- 8 Shake the tank while holding your finger over the filter connection port to clean the tank.

- 9 Remove the DI water filter and connect the DI water filter bypass line (green).

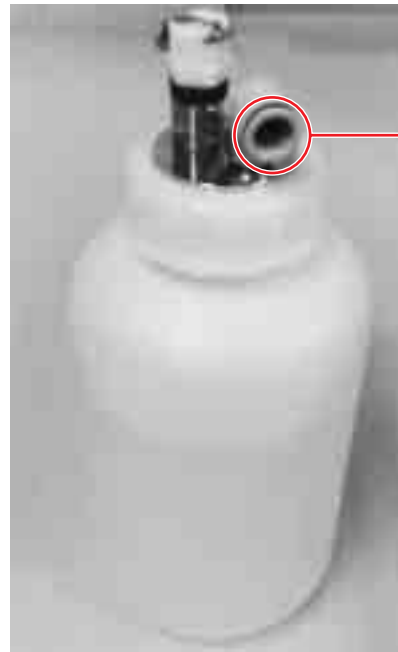

Shake gently while holding here

- 10 Turn on the power supply to the main unit, and start Cell Sorter Software.
- 11 Click [Probe Wash] on the [Cytometer] tab of the ribbon to wash the sample probe.
Repeat ten times to fill the DI water line with ethanol.
- 12 Let stand for 20 minutes, then remove the cap of the DI water tank and dispose of the ethanol.

Note

Dispose of ethanol in accordance with laboratory rules, and with local and federal ordinances and regulations.


- 13 Fill the DI water tank with DI water, then shake the tank with your finger over the air filter connection port to rinse the tank.
- 14 In Cell Sorter Software, click [Probe Wash] on the [Cytometer] tab of the ribbon to wash the sample probe.
Repeat ten times to remove all traces of ethanol from the DI water line.
- 15 Remove the cap from the DI water tank and dispose of the DI water.

Note

Dispose of DI water in accordance with laboratory rules, and with local and federal ordinances and regulations.

- 16 Connect the quick-release connector to the filter.

- 17 Refill the tank with fresh DI water up to the opening of the tank, and reattach the cap.
- 18 Re-install the DI water tank.
Install in the reverse order of removal.
- 19 Disconnect the DI water filter bypass line (green) and reattach the DI water filter.
- 20 Check the drip tray for any fluid spills or leakages, and clean as necessary.

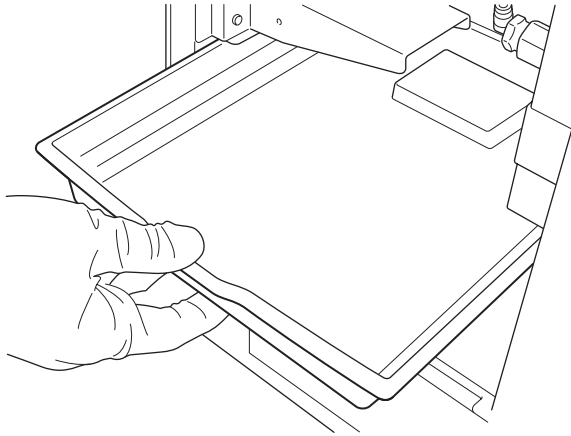

- 21 Close the fluidics maintenance door.
- 22 In Cell Sorter Software, click [Probe Wash] on the [Cytometer] tab of the ribbon, as required, to fill the DI water line with DI water.

Changing the PEEK Sample Line (PEEK sample line compatible models)

The PEEK sample line must be replaced regularly to prevent the buildup of salts restricting the sample flow.


Notes

- The PEEK sample line should be checked for any signs of damage or wear. A damaged PEEK sample line may split and cause sample fluid leakage.
- The PEEK sample line does not support autoclaving (however, the probe adapter does support autoclaving).
- Always leave the PEEK sample line connected when the unit is not in use for protection against dust.
- Use the following procedure to attach the PEEK sample line. There is a risk of fluid leakage if the sample line is not attached correctly.

For details, see “Changing the Probe Adapter (PEEK sample line compatible models)” (page 195).

Tip

Sample fluid may contain biological, chemical, or other agents. Always wear biology laboratory gloves and other protective clothing, mask, and goggles when changing the PEEK sample line.

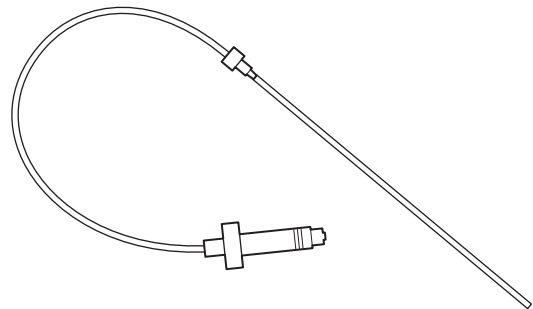


- 1 In Cell Sorter Software, click [Sample Line] in the [Exchange] group on the [Cytometer] tab of the ribbon.
The [Sample Line Exchange] wizard appears.
- 2 Select [PEEK sample line] and click [Start] to begin changing the PEEK sample line.

Notes

- Always follow the instructions displayed in the wizard when changing the PEEK sample line.
- The sorting chip must be replaced or reinserted after changing the PEEK sample line.


The PEEK sample line can also be changed using [Sample line exchange] on the Initial Instrument Setup screen or using [Sample line exchange] in maintenance mode. For details, see “Logging In” (page 45) and “Running Maintenance Mode” (page 177).

3 Open the flip-up door.

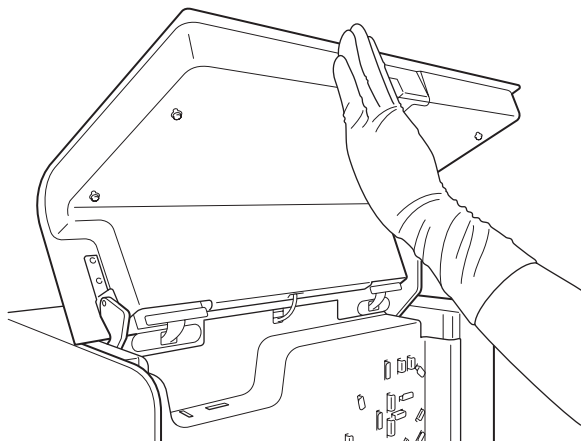


4 Remove the sample line from the pinch valve, then click [Next].

To remove, grasp the sample line on both sides of the pinch valve and pull the sample line out towards you.

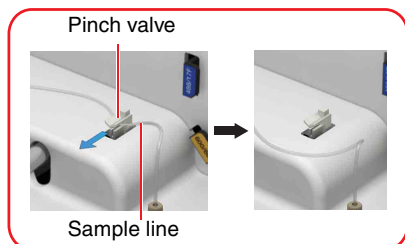

5 Grasp the T-shaped portion of the sample line connector, and gently unscrew it counterclockwise to remove it from the inlet port on the chip loader, then click [Next].

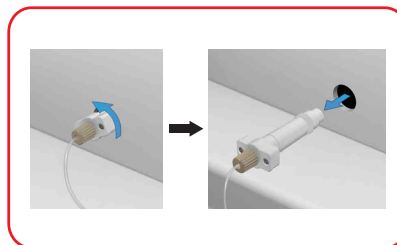


Note

Take care when removing the sample line. Sample fluid can damage electronic circuits if any droplets fall down between the gaps in the sample loader cover and come into contact with internal components.

6 Grasp the black fitting on the PEEK sample line and gently unscrew it counterclockwise to remove it, then click [Next].

Notes

- Unscrew the black fitting while holding the probe adapter down.
- If the sample probe cannot be pulled out, the black fitting may be turning in the reverse direction due to twists in the sample line. Unscrew slowly without releasing the black fitting.

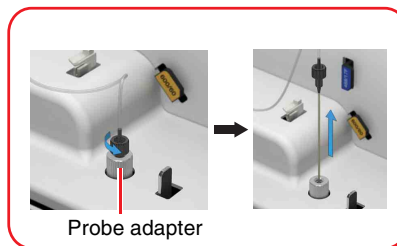


Notes


- Take care when handling the sample probe. It is a high-precision component, which can be damaged easily if handled carelessly.

- Discard the sample line in accordance with the laboratory rules and local ordinances and regulations.

7 Prepare a new PEEK sample line, then click [Next].

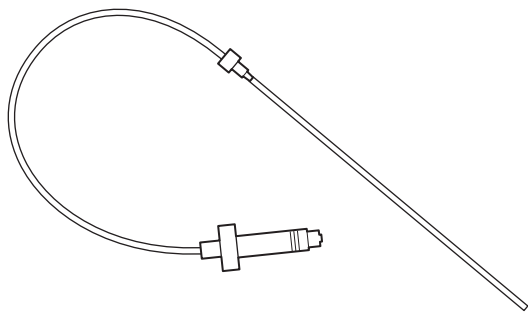

8 Insert the PEEK sample line into the probe adapter, grasp the black fitting, and gently screw it clockwise to secure it, then click [Next].

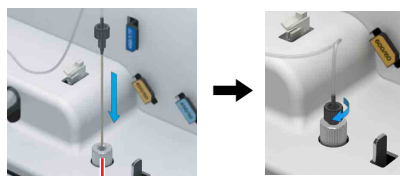


Probe adapter

Note

Fully insert and tighten the threaded portion of the fitting. If the fitting is not fully secured, the correct sample pressure may not be applied.

9 Attach the new sample line to the inlet port on the front of the chip loader, then click [Next].

Screw part A until it clicks into place, making sure it is securely attached.

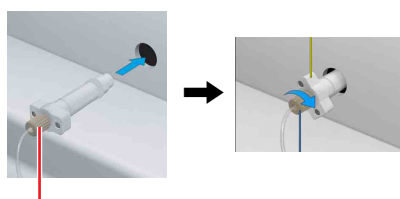

Part A

Note


When attaching the sample line, always hold part A of the connector, not the T-shaped portion. Holding the connector by the T-shaped portion may cause tightening that exceeds the appropriate tightening torque, and adversely affect instrument performance.

Tips

- When attaching the sample line connector, always screw in the connector until you hear a clicking sound.
- Check the sample line connector regularly for any signs of looseness.

10 Check that the sample line connector is attached correctly, then click [Next].

When installed correctly, the stepped portion of the sample line connector fits under the cover. View the connector from the side and check the stepped portion is not visible.

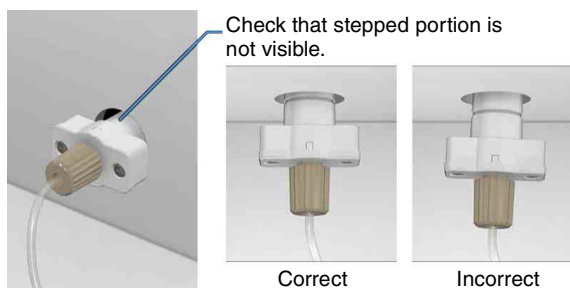


Check that stepped portion is not visible.

Correct

Incorrect


11 Turn the probe adapter, removing any twists in the sample line, then click [Next].

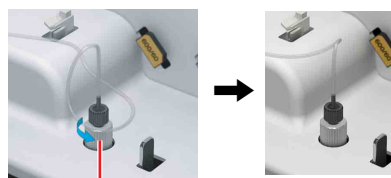

Probe adapter

12 Insert the sample line into the pinch valve, then click [Next].

Make sure the sample line is inserted all the way to the back of the pinch valve.

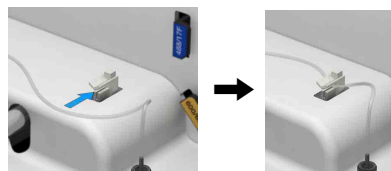

Notes

- Insert the sample line securely into the pinch valve. If the sample line is not correctly inserted in the pinch valve, sample fluid may leak.
- Make sure the sample line connector is securely fastened and that the sample line is correctly inserted in the pinch valve. If the sample line is not inserted correctly, sheath fluid may flow in the reverse direction, potentially causing serious damage to internal components and increasing the risk of biohazard contamination.

13 Click [OK] to exit the [Sample Line Exchange] wizard.

Note

When changing the sample line in maintenance mode, the chip loading screen appears before exiting the [Sample Line Exchange] wizard. Reinsert the unloaded sorting chip, then click [Next].

When the sample line has been changed, the chip exchange screen appears.

Changing the Sample Line and Probe (Non-PEEK sample line compatible models)


The sample line must be replaced regularly. The sample line and sample probe can be removed independently. The sample probe can be removed and sterilized using an autoclave.

Notes

- The sample line should be checked regularly for any signs of damage or wear. A damaged sample line may split and cause sample fluid leakage.
- Always leave the sample line connected when the unit is not in use for protection against dust.
- Use the following procedure to attach the sample line. There is a risk of fluid leakage if the sample line is not attached correctly.

WARNING

Sample fluid may contain biological, chemical, or other agents. Always wear gloves and other protective clothing, mask, and goggles, when changing the sample line or sample probe.

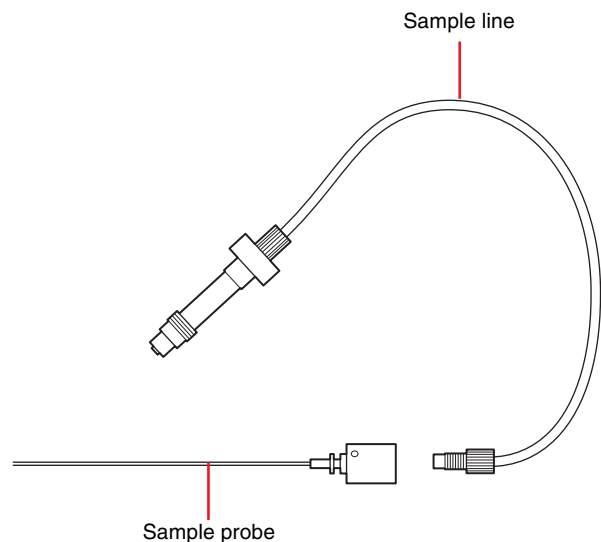


- 1** In Cell Sorter Software, click [Sample Line] in the [Exchange] group on the [Cytometer] tab of the ribbon.
The [Sample Line Exchange] wizard appears.
- 2** Click [Start] to begin changing the sample line and sample probe.

Notes

- Always follow the instructions displayed in the wizard when changing the sample line and sample probe.
- The sorting chip must be replaced or reinserted after changing the sample line and sample probe.


The sample line and sample probe can also be changed using [Sample line exchange] on the Initial Instrument Setup screen or using [Sample line exchange] in maintenance mode. For details, see “Logging In” (page 45) and “Running Maintenance Mode” (page 177).

3 Open the flip-up door.

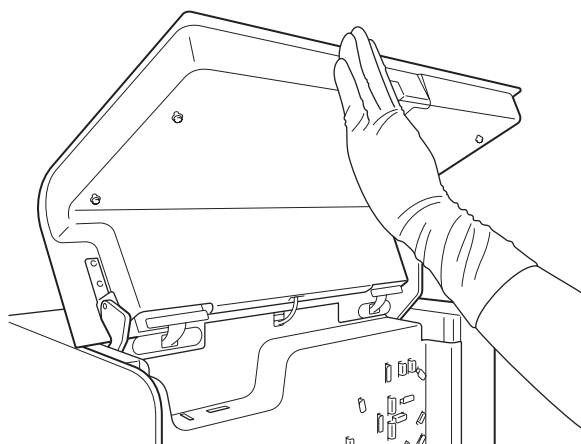


4 Remove the sample line from the pinch valve, then click [Next].

To remove, grasp the sample line on both sides of the pinch valve and pull the sample line out towards you.

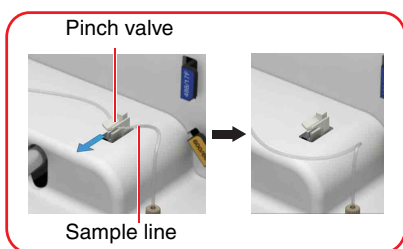


5 Grasp the T-shaped portion of the sample line connector, and gently unscrew it counterclockwise to remove it from the inlet port on the chip loader, then click [Next].

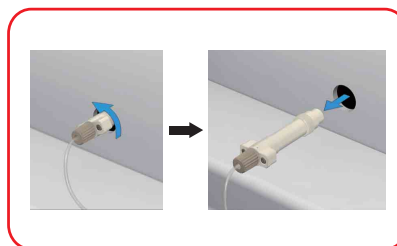

Note

Take care when removing the sample line. Sample fluid can damage electronic circuits if any droplets fall down between the gaps in the sample loader cover and come into contact with internal components.

6 Push and hold the sample probe release catch to the right, and pull the probe straight up and out, then click [Next].

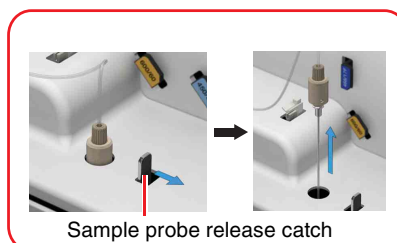


Notes

- Take care when handling the sample probe. It is a high-precision component, which can be damaged easily if handled carelessly.
- Discard the sample line as biohazard waste in accordance with the laboratory rules and local ordinances and regulations.

- 7** Prepare a new sample line and sample probe, attach the sample line to the sample probe, then click [Next].

Hold part B of the sample probe, and screw part C on the sample line until it clicks into place, making sure it is securely attached.

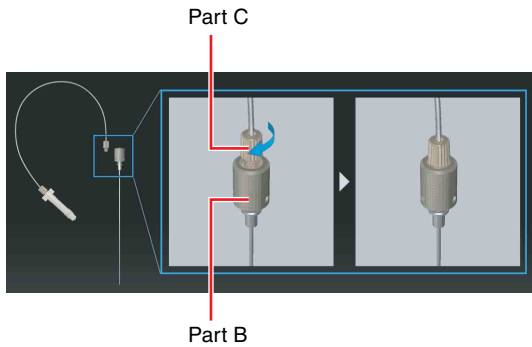


Note

If reattaching the removed sample line and sample probe, disinfect the components using the following procedure, as required.

- Place the sample line in a beaker (or other container) of ethanol and let sit for about 30 minutes.
- Autoclave the sample probe under general autoclave conditions (121 °C (250 °F), 2 atmospheres, approx. 20 minutes).

- 8** Insert the sample probe while pressing its metal release catch to the right.

Press down on the sample probe to insert it securely, and release the metal release catch. Verify that the metal release catch has returned to the left side.

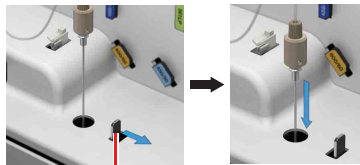


Sample probe release catch



- 9** Open the flip-up door, and attach the new sample line to the inlet port on the front of the chip loader, then click [Next].

Screw part A until it clicks into place, making sure it is securely attached.

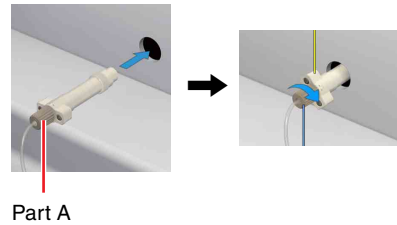

Part A

Note


When attaching the sample line, always hold part A of the connector, not the T-shaped portion. Holding the connector by the T-shaped portion may cause tightening that exceeds the appropriate tightening torque, and adversely affect instrument performance.

Tips

- When attaching the sample line connector, always screw in the connector until you hear a clicking sound.
- Check the sample line connector regularly for any signs of looseness.

- 10** Check that the sample line connector is attached correctly, then click [Next].

When installed correctly, the stepped portion of the sample line connector fits under the cover. View the connector from the side and check the stepped portion is not visible.

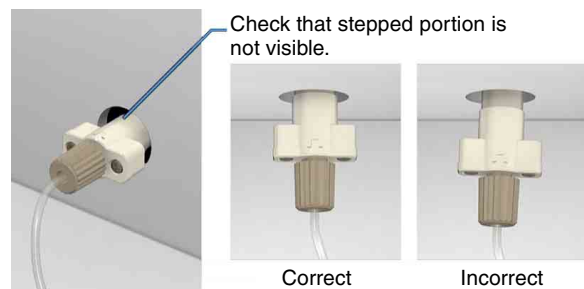


Check that stepped portion is not visible.

Correct

Incorrect


- 11** Turn the probe adapter, removing any twists in the sample line, then click [Next].

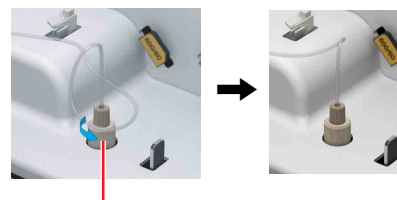

Probe adapter

- 12** Insert the sample line into the pinch valve, then click [Next].

Make sure the sample line is inserted all the way to the back of the pinch valve.

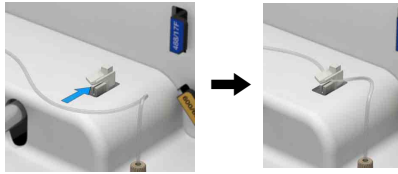

Notes

- Insert the sample line securely into the pinch valve. If the sample line is not correctly inserted in the pinch valve, sample fluid may leak.
- Make sure the sample line connector is securely fastened and that the sample line is correctly inserted in the pinch valve. If the sample line is not inserted correctly, sheath fluid may flow in the reverse direction, potentially causing serious damage to internal components and increasing the risk of biohazard contamination.

- 13** Click [OK] to exit the [Sample Line Exchange] wizard.

Note

When changing the sample line in maintenance mode, the chip loading screen appears before exiting the [Sample Line Exchange] wizard. Reinsert the unloaded sorting chip, then click [Next].

When the sample line has been changed, the chip exchange screen appears.

Changing the Probe Adapter (PEEK sample line compatible models)

The probe adapter can be removed, disinfected using autoclaving, and then reinserted back into the unit.

WARNING

Sample fluid may contain biological, chemical, or other agents. Always wear biology laboratory gloves and other protective clothing, mask, and goggles when changing the probe adapter.

Note


Change the probe adapter using the following procedure after removing the PEEK sample line. In particular, when disinfecting the probe adapter for reuse, always follow the procedure correctly to change the adapter.

- 1** Remove the PEEK sample line.

For details about removal, see steps 1 to 6 in “Changing the PEEK Sample Line (PEEK sample line compatible models)” (page 189).

- 2** Remove the probe adapter.

Push and hold the sample probe release catch to the right, then pull the probe straight up and out.

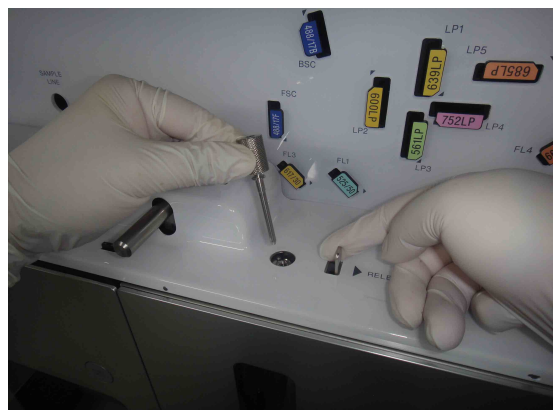

3 Disinfect the probe adapter, as required.

Autoclave the probe adapter under general autoclave conditions (121 °C (250 °F), 2 atmospheres, approx. 20 minutes).

Note

Make sure the probe adapter is autoclaved if reusing the probe adapter.

To reinstall

1 Install the probe adapter.


Push and hold the sample probe release catch to the right, insert the probe straight down, and then let go of the release catch. Check that the release catch returns all the way to the left.

2 Install the PEEK sample line.

For details about installation, see steps **7 to 13** in “Changing the PEEK Sample Line (PEEK sample line compatible models)” (page 189).

Changing the Sheath Filter


The sheath filter in the fluidics module should be replaced regularly to prevent small particles of dust that can cause fluctuations in the fluid flow and bacteria that can cause contamination of the sample.

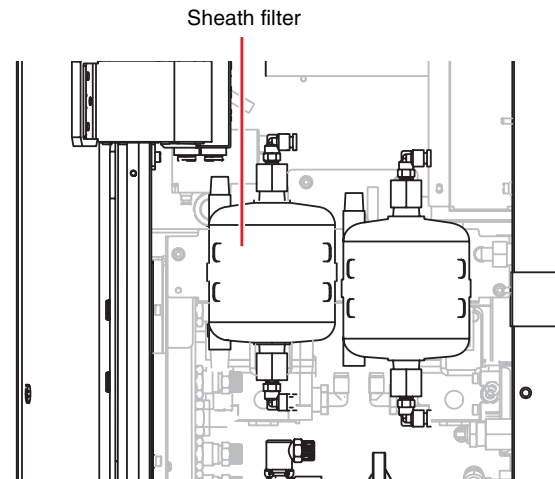

Perform the following procedure by clicking [Skip Auto Calibration] in “Initial Instrument Setup” (page 46) or “Automatic Calibration” (page 50) with the main window displayed.

- 1 In Cell Sorter Software, click [Settings] on the [Cytometer] tab of the ribbon.
The [Cytometer Settings] dialog appears.
- 2 Click [Advanced Settings].
The [Advanced Settings] dialog appears.
- 3 Click [Standby] in the [Pressure Options] tab of the [Advanced Settings] dialog.
The sheath fluid flow stops.

- 4** Open the fluidics maintenance door and remove the sheath filter from the retaining clamp without disconnecting the connectors.

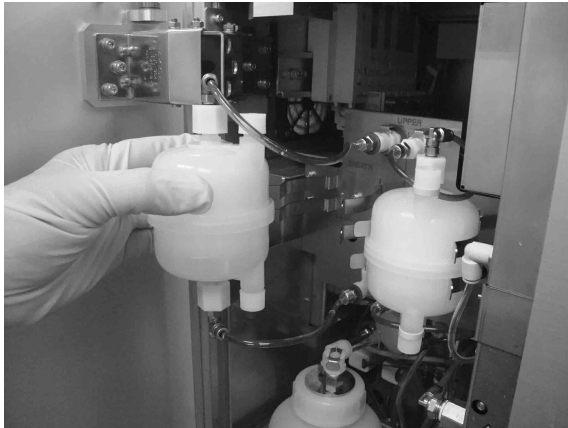


- 5** Disconnect the sheath filter lines from the “UPPER” and “LOWER” connectors.

Press the metal release catch and withdraw each connector.

Tip

The sheath filter connectors have stop valves that prevent leakage of fluid when disconnected.

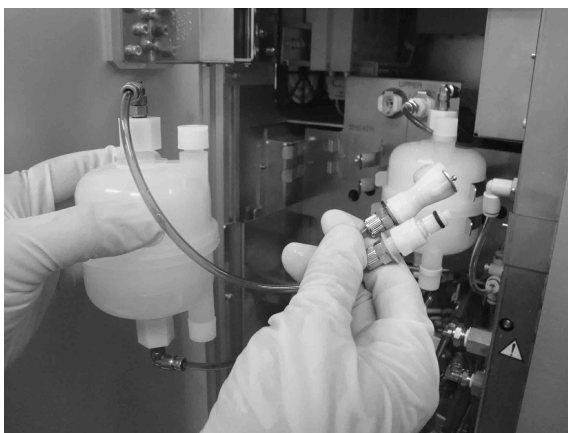


- 6** Discard the used sheath filter.

Note

Discard the sheath filter as biohazard waste in accordance with the laboratory rules and local ordinances and regulations.

- 7** Hold a new sheath filter with the arrow on the side of the filter facing up.

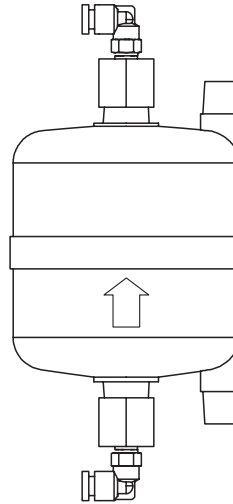

- 8** Connect the two fluidics lines from the sheath filter to the corresponding connectors.

Connect the line from the top of the filter to the “UPPER” connector, and the line from the bottom to the “LOWER” connector. Insert each connector and push it until it locks into place. A click sound occurs when the connectors lock into place.

- 9** Insert the sheath filter into the retaining clamp.
- 10** In Cell Sorter Software, click [Sheath Filter] in the [De-bubble] group on the [Cytometer] tab of the ribbon.

The [Sheath Filter De-bubble] dialog appears.

- 11** Click [Start]. When the de-bubble function is completed, click [Close].

Repeat two more times to remove residual air from the sheath filter.

Note

The sheath filter de-bubble process takes a short time to finish. Run the sheath filter de-bubble function at least three times to make sure that all air and small particles have been removed from the filter and sheath line. Failing to do so may cause any solid matter not removed from accumulating and clogging the sorting chip.

- 12** Perform steps **4** to **8** as described in “To release air trapped in the sheath filter manually” (page 202).

- 13** Click [Ready] in the [Pressure Options] tab of the [Advanced Settings] dialog.


The sheath fluid flow restarts.

- 14** Load and run a sample tube containing sheath fluid, and check for leaks around the connectors.

Check the drip tray for any fluid spills or leakages, and clean as necessary.

Changing the DI Water Filter

The DI water filter in the fluidics module should be replaced regularly to prevent small particles of dust that can cause fluctuations in the fluid flow and bacteria that can cause contamination of the sample.

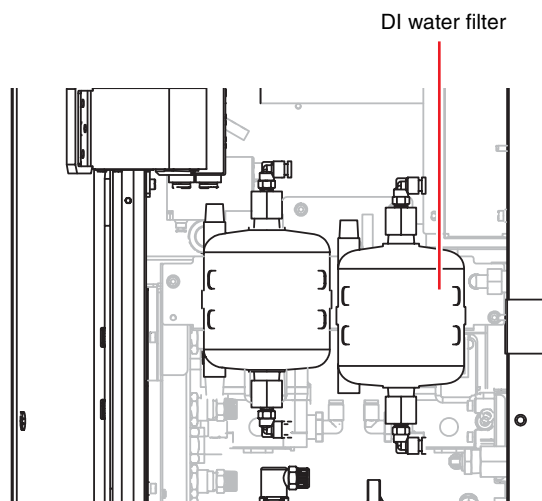

- 1** Shut down the main unit.
- 2** Open the fluidics maintenance door and remove the DI water filter from the retaining clamp without disconnecting the connectors.




- 3 Disconnect the DI water filter lines from the “UPPER” and “LOWER” connectors.

Press the metal release catch and withdraw each connector.

Tip

The DI water filter connectors have stop valves that prevent leakage of fluid when disconnected.

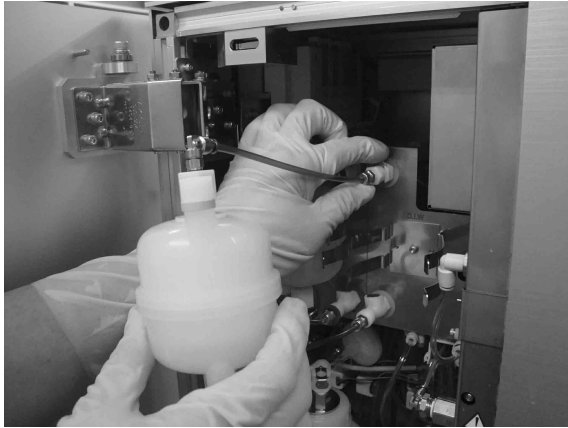


- 4 Discard the used DI water filter.

Note

Discard the DI water filter as biohazard waste in accordance with the laboratory rules and local ordinances and regulations.

- 5 Hold a new DI water filter with the arrow on the side of the filter facing up.

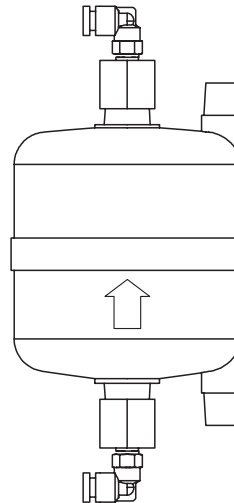

- 6 Connect the two fluidics lines from the DI water filter to the corresponding connectors.

Connect the line from the top of the filter to the “UPPER” connector, and the line from the bottom to the “LOWER” connector. Insert each connector and push it until it locks into place. A click sound occurs when the connectors lock into place.

- 7 Insert the DI water filter into the retaining clamp.
- 8 Close the fluidics maintenance door and turn on the power supply and log in.
- 9 Open the fluidics maintenance door, remove the DI water filter from the retaining clamp (without disconnecting the connectors), and check the fluid level.



- 10 Turn on the power supply to the main unit, and start Cell Sorter Software.

- 11 Click [Settings] on the [Cytometer] tab of the ribbon to display the [Cytometer Settings] dialog. Click [DI Filter de-bubble] on the [Maintenance] tab.

The [DI Filter de-bubble] dialog appears.

- 12 Click [Start].


13 Release the air trapped in the DI water filter.

Slowly open the air release port on the DI water filter, within five minutes of startup, to release the trapped air.

If air is not released within five minutes after startup, turn the power supply to the main unit off and then on again, then repeat the procedure.

Notes

- It is recommended that the drip tray be pulled out partway before proceeding. Slowly open the air release port while holding a soft, clean cloth ready to wipe away any DI water overflow.
- Keep the DI water filter clear of all other objects when the air release port is open.

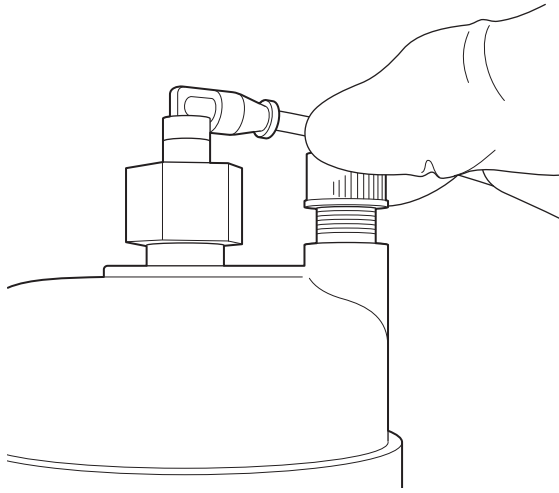

The trapped air escapes and the fluid level in the DI water filter begins to rise.

14 Close the air release port when the fluid level nears the top of the filter.

Make sure that the air release port is tightened and that there is no leakage of fluid.

Note

If the air release port is overtightened, the filter may become damaged. Do not apply excess force when tightening.

15 Click [Stop].

16 Place the DI water filter in the retaining clamp, and close the fluidics maintenance door.

17 When finished, click [Close].

Changing the Sintered Sheath Line Filter

A sintered sheath line filter is placed in the sheath line to filter residue or small particles from the sheath fluid supply. The filter should be replaced every 12 months.



Sintered sheath filter

1 Shut down the main unit.

Make sure the main unit is completely shut down before proceeding.

2 Turn off the compressed air supply.

3 Release the residual air pressure from the sheath tank using the ring-pull air relief valve.

4 Hold the sintered sheath line filter with one hand, push and hold in the quick release collet with the other hand, and pull out the sheath fluid line (blue) from the side furthest from the sheath tank.

Exert gentle force if the tube feels difficult to pull out from the connector.



- 5** In the same way, remove the white connector on the sheath tank side, and discard the filter.



Note

Discard the sintered sheath line filter as biohazard waste in accordance with the laboratory rules and local ordinances and regulations.

- 6** Insert a new sintered sheath line filter in the sheath fluid line, pushing the line all the way into the filter, then connect the white connector to the sheath tank.

Releasing Air in the Sheath Filter and Sorting Chip

During operation, air may become trapped in the sheath filter, interrupting the smooth flow of sheath fluid to the sorting chip. Air may also become trapped in the sorting chip itself, preventing the formation of regular-shaped droplets.

To release air trapped in the sheath filter using the de-bubble function

- 1 Click [Sheath Filter] in the [De-bubble] group on the [Cytometer] tab of the ribbon.

The [Sheath Filter De-bubble] dialog appears.

- 2 Click [Start], and then click [Close] when finished.

The sheath filter de-bubble process takes a short time to finish.

If changing the sheath filter, run the sheath filter de-bubble process at least three times to make sure that all air and small particles have been removed from the filter and the sheath line.

To release air trapped in the sorting chip using the de-bubble function

If the formation of droplets is irregular even though there is no air trapped in the sheath filter, de-bubble the sorting chip. Click [Chip] in the [De-bubble] group on the [Cytometer] tab of the ribbon.

To release air trapped in the sheath filter manually

If the volume of air trapped in the sheath filter is large, it may be more efficient to release the trapped air manually before using the de-bubble function.

This procedure is performed with the power supply turned on and Cell Sorter Software up and running.

- 1 In Cell Sorter Software, click [Settings] on the [Cytometer] tab of the ribbon.

The [Cytometer Settings] dialog appears.

- 2 Click [Advanced Settings].

The [Advanced Settings] dialog appears.

- 3 Click [Standby] in the [Pressure Options] tab of the [Advanced Settings] dialog.

The sheath fluid flow stops.

- 4 Open the fluidics maintenance door, remove the sheath filter from the retaining clamp (without disconnecting the connectors), and check for the presence of air bubbles.

- 5 If there are air bubbles, slowly open the air release port on the top of the sheath filter.

Notes

- It is recommended that the drip tray be pulled out partway before proceeding. Slowly open the air release port while holding a soft, clean cloth ready to wipe away any sheath fluid overflow.
- Keep the sheath filter clear of all other objects when the air release port is open.



The trapped air escapes and the fluid level in the sheath filter begins to rise.

- 6 Close the air release port when the fluid level nears the top of the filter.


Make sure that the air release port is tightened and that there is no leakage of fluid.

Note

If the air release port is overtightened, the filter may become damaged. Do not apply excess force when tightening.

- 7 Tap the side of the sheath filter with your hand 2 or 3 times.

This dislodges any trapped air in the lines, which then collects in the sheath filter.

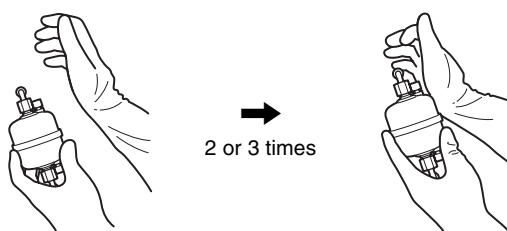

Use the following procedure if there are air bubbles in the sheath filter.

- ① Open the air release port to expel the trapped air.
- ② Check that the air has been released and then close the air release port.
- ③ Tap the filter again and check that there are no air bubbles being formed.

Repeat steps ① to ③ until all air has been released.

- 8 Place the sheath filter in the retaining clamp, then close the fluidics maintenance door.
- 9 Click [Ready] in the [Pressure Options] tab of the [Advanced Settings] dialog.

The sheath fluid flow restarts.

Note

Check that there are no air bubbles in the sheath filter before running samples for measurement. It is recommended that you use this procedure if there are any air bubbles in the sheath filter.

Releasing Air in the DI Water Filter

During operation, air may become trapped in the DI water filter, interrupting the smooth flow of DI water.

To release air trapped in the DI water filter using the de-bubble function

- 1 Click [Settings] on the [Cytometer] tab of the ribbon.
The [Cytometer Settings] dialog appears.
- 2 Click [DI Water de-bubble] on the [Maintenance] tab
The [DI Water de-bubble] dialog appears.
- 3 Click [Start].
- 4 Open the fluidics maintenance door, remove the DI water filter from the retaining clamp (without disconnecting the connectors).
- 5 Slowly open the air release port on the top of the DI water filter.

Notes

- It is recommended that the drip tray be pulled out partway before proceeding. Slowly open the air release port while holding a soft, clean cloth ready to wipe away any DI water overflow.
- Keep the DI water filter clear of all other objects when the air release port is open.



The trapped air escapes and the fluid level in the DI water filter begins to rise.

- 6 Close the air release port when the fluid level nears the top of the filter.

Make sure that the air release port is tightened and that there is no leakage of fluid.

Note

If the air release port is overtightened, the filter may become damaged. Do not apply excess force when tightening.

- 7 Click [Stop].
- 8 Place the DI water filter in the retaining clamp, and close the fluidics maintenance door.
- 9 When finished, click [Close].

Autoclaving the Sheath Filter and DI Water Filter


You can autoclave the sheath filter and DI water filter if using autoclavable parts to sterilize the filters.

Notes

- Only the filters are autoclavable.
- Each filter can be autoclaved up to ten times.
- After removing the filter from the main unit, perform work over a sink to avoid losing the parts.

Autoclaving the Sheath Filter and DI Water Filter

If necessary, you can autoclave the sheath filter and DI water filter.

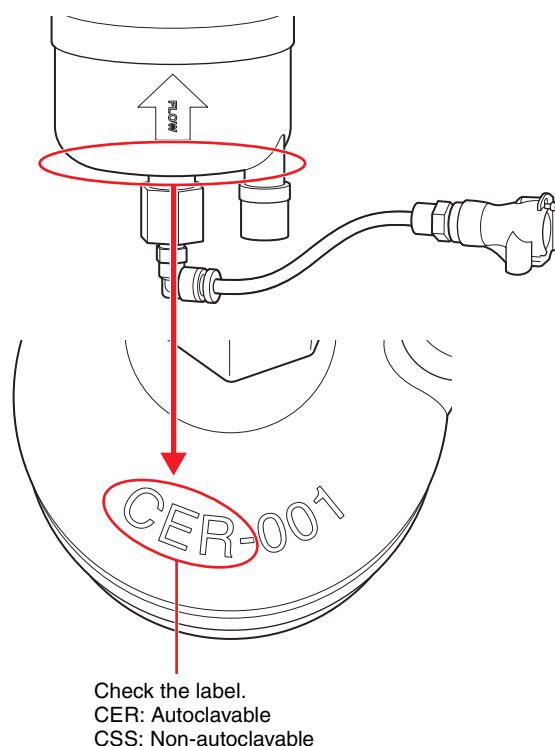


- 1 Remove the filter from the main unit.
For details about removing filters, see “Changing the Sheath Filter” (page 196) and “Changing the DI Water Filter” (page 198).
- 2 Remove the top and bottom caps, and dispose of any fluid inside the filter.
Drain the fluid from the filter until there is no remaining fluid.


- 3** Remove the IN and OUT connectors from the filter, and remove the spacer and sealing from inside each connector.

Notes

- Do not lose the spacers and sealings after removing them.
- You can identify the orientation of the filter by the FLOW arrow on the filter's body.

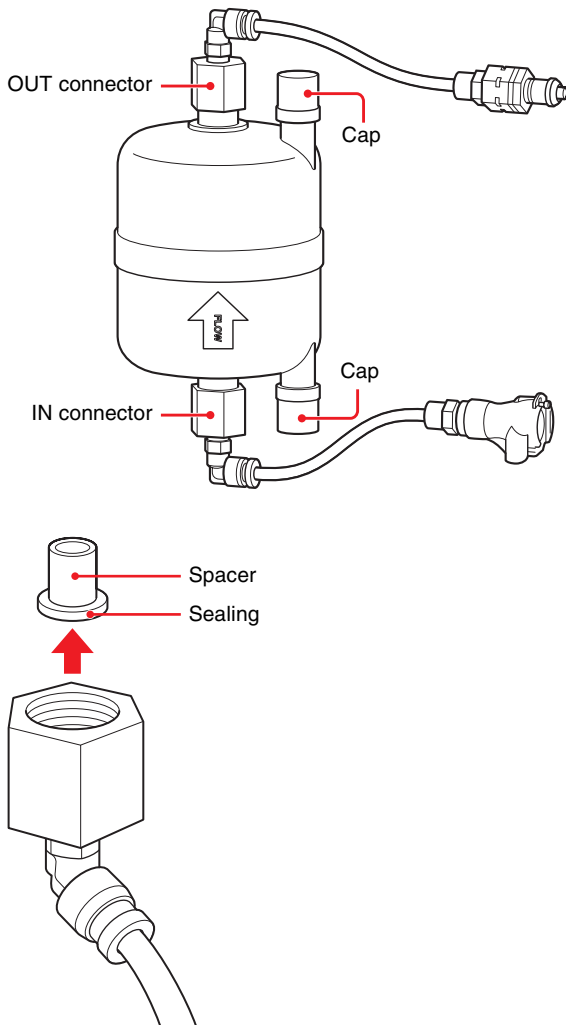

The filter should have external parts removed.

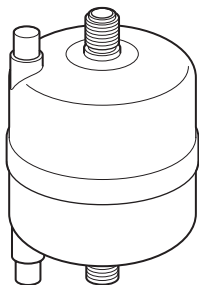

- 4** Autoclave the filter under the following conditions.

Temperature: 121 °C (250 °F)
 Pressure: 2 atmospheres
 Duration: 30 minutes

- 5** After autoclaving, attach the couplers and reinsert into the main unit.


Notes

- The IN and OUT connectors differ at their tube ends. Be sure to verify the orientation of the filter and identify the IN and OUT connectors beforehand to ensure proper attachment.
- Make sure there is no dust or particles on the spacers and sealings to ensure leak prevention.

- ① Place the sealing on the spacer. Make sure the head of the spacer protrudes slightly from the sealing.



- ② Insert the spacer into the connecting point on the filter.

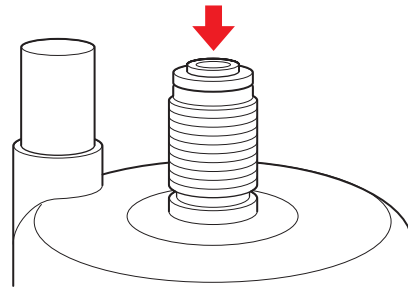


Note


Do not press on the spacer portion. Insert the spacer gently without altering the state prepared in step ①.

- ③ Attach the IN connector to the filter.

- ④ Repeat the procedure to attach the OUT connector to the filter.

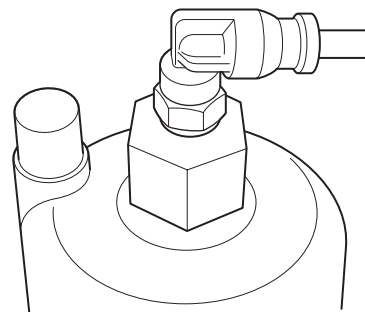

Tighten the connectors by hand, and then tighten them further by 1/6 turn using a wrench.

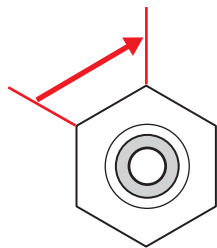


Notes

- Excessive tightening may damage the filter (recommended tightening torque: 0.5 ± 0.1 N·m).
- After attaching the filter, release the air from inside the filter by performing the procedures in the “Releasing Air in the Sheath Filter and Sorting Chip” - “To release air trapped in the sheath filter manually” (page 202) and “Releasing Air in the DI Water Filter” (page 203).


Ejecting the Sorting Chip Manually

If for any reason the chip will not eject normally, the chip can be removed manually.


- 1 Check that the compressed air supply is turned on, and open the flip-up door to access the sorting chip.

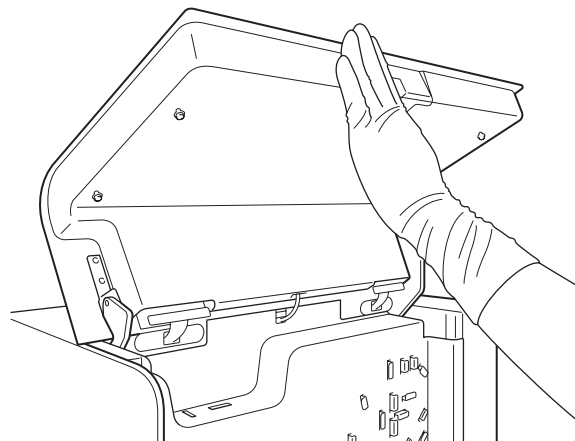

If the sample line connector is out at the front, go to step 6.

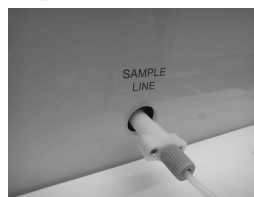


If the sample line connector is not out at the front, go to step 2.

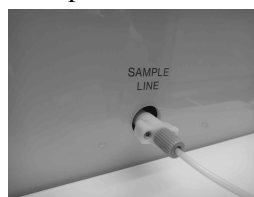


- 2 Turn off the compressed air supply.

Have an assistant help if you cannot reach the air supply control valve of the compressor supplying compressed air while holding the flip-up door open.

- Lower the flip-up door slowly until it can support its own weight.

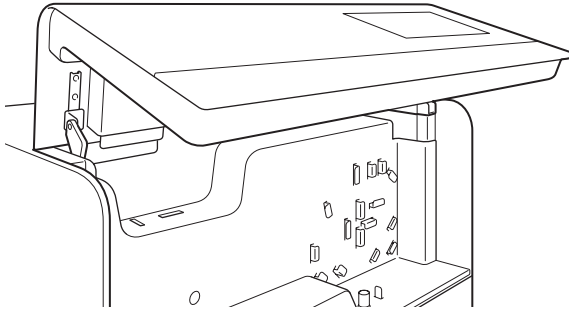


- Shut down the main unit.
- Grasp the sample line connector firmly and pull it out towards you to the position, if not in position already.

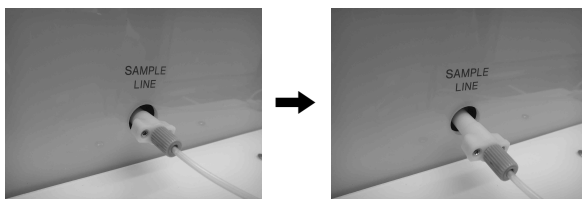

Pulling the sample line connector withdraws the sorting chip connector interface, allowing the chip to be removed.

- Eject the chip by pushing the top of the thumbwheel from front to back.

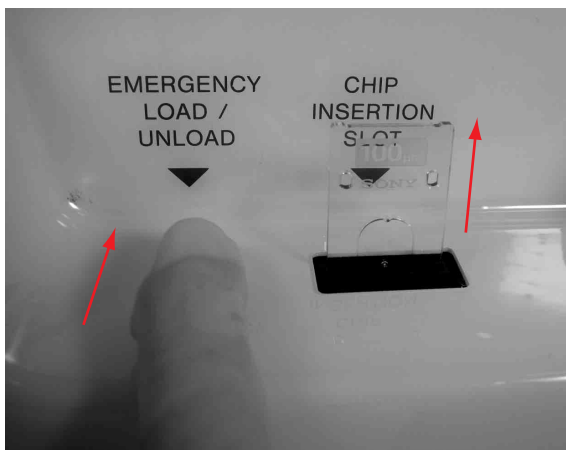

Note

Touching the thumbwheel or the sample line will directly affect the chip alignment. If you touch either and then opt to leave the chip inserted, the chip must be realigned. To align the chip, click [Chip Alignment] on the [Cytometer] tab of the ribbon.

- When finished, turn on the compressed air supply.
- Close the flip-up door.
- Turn on the main power supply.

Cleaning and Handling of Optical Filters

The bandpass optical filters and longpass optical filters in the detection module require cleaning to remove any dust on the surface of the filters. The filters should be checked and cleaned every 6 to 12 months, or more often if the optical filter pattern is frequently changed.

Note

The optical filters are high-precision components that require the utmost care. If an optical filter is scratched or damaged in any way, it must be replaced.


- Remove an optical filter by placing your finger under the lip of the retaining holder and pulling directly out.



- Grip the sides of the optical filter and withdraw it from the filter slot.

Note

Do not touch the surfaces of the optical filters.

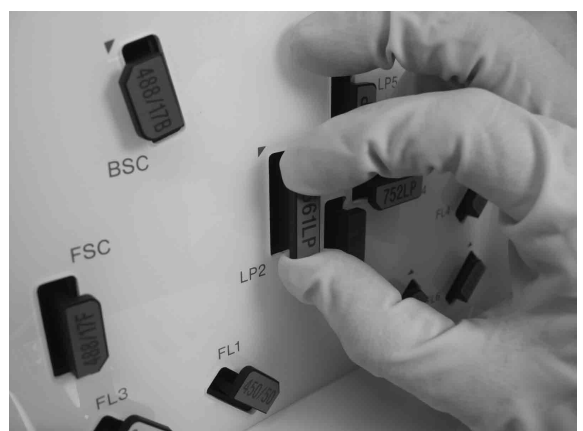

It is recommended that only one optical filter be withdrawn at any one time to ensure that it is returned to the original slot.

- 3 Remove loose dust or particles using a canned air blower or a bulb puffer hand duster, similar to those used in photography.

Direct the airflow across the surface at an oblique angle. If necessary, gently wipe the surface using lens paper. Use a fresh surface of the cloth for each wipe. Maintain a continuous wiping motion at a constant speed until the cloth passes the edge of the optical filter.

- 4 After cleaning, insert the optical filter into the slot and push the lip of the optical filter until it clicks into place.



Notes

- Avoid touching or wiping coated or metal mirror surfaces at all times, except when cleaning.
- Avoid handling exposed coatings with bare fingers.
- Always wear gloves, and also wear a face mask, goggles, and other protective clothing as required to prevent oils from your hands contaminating the optical glass, and to prevent solvents and chemicals coming into contact with your skin or eyes.

Disconnecting and Reconnecting the Fluidics Cart

After installation, it may be necessary to disconnect and then reconnect the fluidics cart if you wish to move the fluidics cart and need to re-route the path of the fluidics lines, sheath air line, and cart connection cable between the fluidics cart and the main unit.

Note

If you wish to move the SH800 main unit, contact your Sony distributor. Do not attempt to relocate it yourself. The main unit contains high-precision optical and laser instruments that require alignment after installation that only your Sony distributor can provide.

- 1 With Cell Sorter Software shut down, turn off the power supply using the MAIN POWER switch on the rear panel of the main unit.
- 2 Disconnect the three fluidics lines (sheath fluid, waste fluid, and ethanol) and sheath air line from the rear of the fluidics cart.

Press the metal release catch to withdraw each connector.



Notes

- Do not touch the surfaces of the connectors that come into contact with fluids. Take care not to spill any fluid.
- Always wear gloves and other protective clothing, mask, and goggles, as required, when handling waste fluid.

- 3 Disconnect the connection cable from the rear of the fluidics cart.

Unscrew the retaining screws and pull the connector out.

Cart connection cable



- 4 Move the fluidics cart and re-route fluidics lines, sheath air line, and cable, as required, then reconnect the lines, air line, and cable.

Insert each connector and push until it locks into place.

- 5 Turn on the power supply.

Check for any fluid spills, and clean as necessary.

- 6 Check the operation of the fluid level sensors on the LCD monitor.

Check the electrical cable connection if the status information does not appear.

Cleaning the Internal Sheath Line and DI Water Line using Ethanol

The SH800 internal sheath line and DI water line should be cleaned with ethanol regularly to prevent the growth of bacteria and other agents. In addition, the fluidics system should also be cleaned as required after running samples that contain infectious substances.

Precautions when cleaning with ethanol

- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Never pour ethanol into a tank other than the ethanol tank, and conversely, never pour other liquids into the ethanol tank.
- Make sure that the cap on the ethanol tank is correctly and securely fastened at all times.
- Make sure that the ethanol line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the ethanol tank itself.
- Make sure that the lid on the waste tank is correctly and securely fastened at all times.
- Make sure that the waste fluid line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the waste tank itself.

Note

The SH800 is compatible with standard disinfecting agents (ethanol 70%).

- 1 In Cell Sorter Software, click [Ethanol Cleaning] on the [Cytometer] tab of the ribbon.

Notes

- The ethanol cleaning sequence cannot be aborted once it is started.
- The SH800 hardware and software are automatically shut down after cleaning with ethanol.

You can also run ethanol cleaning in maintenance mode. For details, see “Running Maintenance Mode” (page 177).

- 2 Click [Start] to begin the cleaning process and follow the on-screen instructions.

- 3** Refill the DI water tank in the main unit with distilled, deionized water, then click [Next].

Make sure that there is sufficient DI water in the tank. An alarm will be raised if there is insufficient volume of DI water in the DI water tank before cleaning.

See “Refilling the DI Water Tank” (page 167).

- 4** Refill the ethanol tank in the fluidics cart with ethanol, then click [Next].

Make sure that there is at least 0.5 liters (18 fl oz.) in the ethanol tank. An alarm will be raised if there is insufficient volume of ethanol in the ethanol tank before cleaning.

See “Refilling the Ethanol Tank” (page 169).

- 5** Remove the sheath filter.


- ① Remove the sheath filter from the retaining clamp without disconnecting the connectors.




- ② Disconnect the sheath filter lines from the “UPPER” and “LOWER” connectors. Press the metal release catch and withdraw each connector.

Tip


The sheath filter connectors have stop valves that prevent leakage of fluid when disconnected.

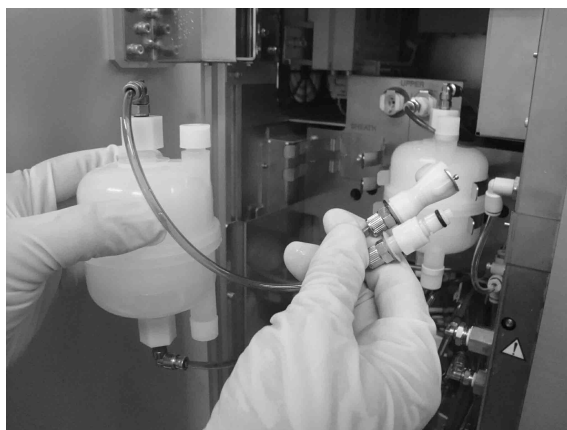

- 6** Remove the DI water filter in the same way.

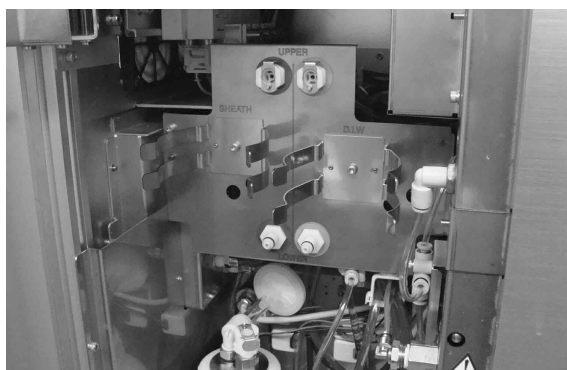

- 7** Connect the sheath filter bypass line (yellow) and DI water filter bypass line (green) to the corresponding “UPPER” and “LOWER” connectors.

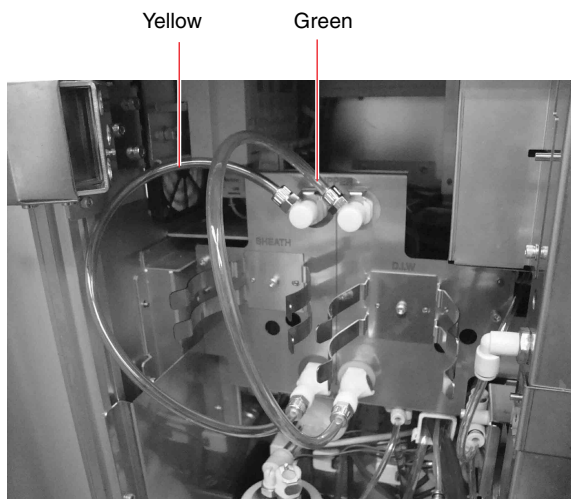

- 8** Hold the QR code on the packaging for the cleaning chip in front of the built-in camera on the PC.

The currently loaded sorting chip, if one is present, is automatically ejected.

- 9** Insert the cleaning chip in the insertion slot on the top of the chip loader, then click [Next].

- 10** Select the ethanol cleaning time, then click [Start].

The recommended ethanol cleaning time is 12 minutes.


A progress bar is displayed during cleaning.

The sheath fluid line is cleaned with ethanol for the specified duration. When the cleaning time has elapsed, the sheath fluid line is flushed with DI water to remove all traces of ethanol.

- 11** When cleaning is completed, click [Next].

- 12** Remove the sheath filter bypass line (yellow) and DI water filter bypass line (green), and reattach the sheath filter and DI water filter.

The DI water filter goes on the right, and the sheath filter on the left. Check that the filters are not connected incorrectly.

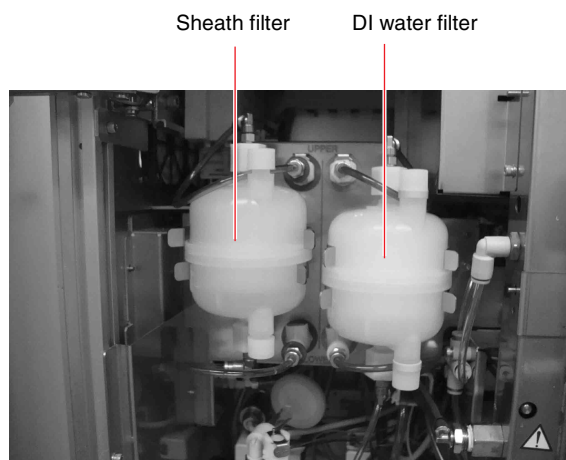

- 13** Click [Next].

The SH800 hardware and software shut down automatically.

The cleaning chip remains loaded in the chip loader after shutdown.

Cleaning All Fluidics Lines using Ethanol (SH800S, SH800SP, SH800Z, SH800ZP, SH800 (96-well plate models))

The entire fluidics system, including the sheath tank, sheath line, DI water tank, DI water line, and sample line, should be cleaned with ethanol regularly to prevent the growth of bacteria and other agents. The fluidics system should be cleaned on a regular basis, or as required, when running samples that contain, for example, infectious substances.

Tip

Prepare a cleaning chip before starting the cleaning process.

Precautions when cleaning with ethanol

- Ensure the laboratory is well ventilated during cleaning with ethanol. Inhaling ethanol vapor can cause irritation of the eyes, skin and respiratory tract, loss of coordination, drowsiness, and in sufficient concentration, unconsciousness.
- Never place a naked flame, embers, or other material that can emit sparks near the main unit or the fluidics cart. Ethanol vapor is highly flammable at normal room temperature.
- Never pour ethanol into a tank other than the ethanol tank, and conversely, never pour other liquids into the ethanol tank.
- Make sure that the cap on the ethanol tank is correctly and securely fastened at all times.
- Make sure that the ethanol line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the ethanol tank itself.
- Make sure that the lid on the waste tank is correctly and securely fastened at all times.
- Make sure that the waste fluid line is securely connected to the connectors on the rear panel of the main unit, rear panel of the fluidics cart, and the waste tank itself.

Note


The SH800 is compatible with standard disinfecting agents (ethanol 70%).

- 1 Replace or empty the waste tank.

For details, see “Emptying/Changing the Waste Tank” (page 166).

- 2 Launch Cell Sorter Software and enter maintenance mode using [Maintenance Mode] on the bottom right of the login screen, select [Ethanol cleaning of all fluidics lines], then click [Start].


For details, see “Running Maintenance Mode” (page 177).

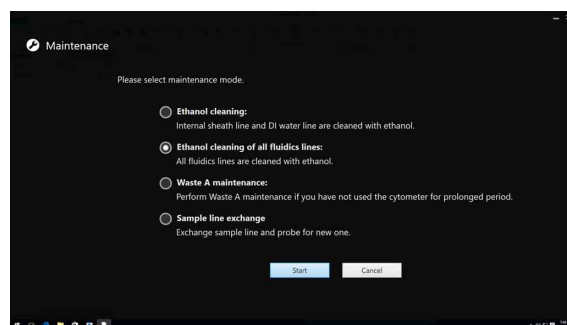

- 3 Click [Start] to start the [Ethanol Cleaning of all fluidics lines] wizard.

- 4 Change the sheath filter, as required, then click [Next].

For details, see “Changing the Sheath Filter” (page 196).

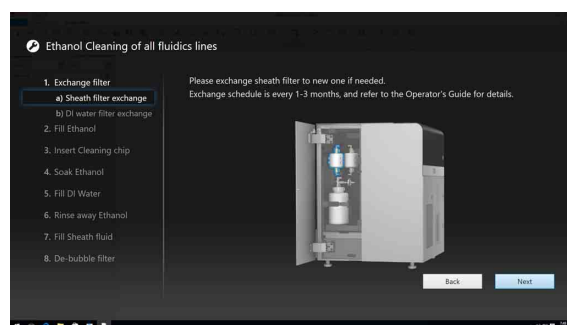


Note

The sheath filter used for ethanol cleaning of all fluidics lines should be replaced once a month or every 3 months according to the following conditions.

- If the cleaning time using ethanol is less than 24 hours: Every 3 months
- If the cleaning time using ethanol is more than 24 hours: Every month


- 5 Open the air release port on the sheath filter, to release the air in the sheath filter, and close the air release port, then click [Next].

For details, see steps 4 to 6 of “To release air in the sheath filter manually” in “Releasing Air in the Sheath Filter and Sorting Chip” (page 202).

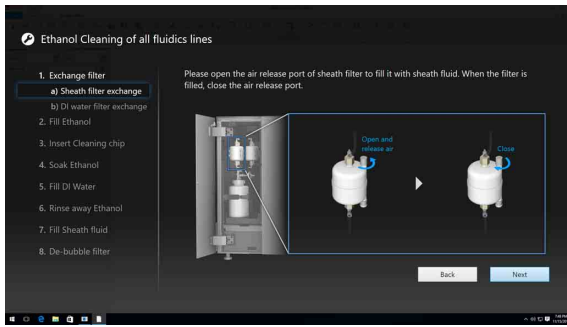

6 Change the DI water filter, as required, then click [Next].

For details, see steps 2 to 7 of “Changing the DI Water Filter” (page 198).

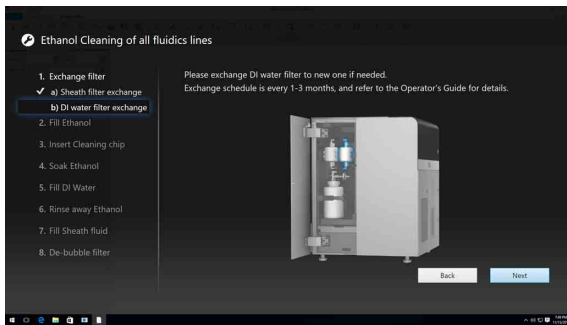


Note

The DI water filter used for ethanol cleaning of all fluidics lines should be replaced once a month or every 3 months according to the following conditions.

- If the cleaning time using ethanol is less than 24 hours: Every 3 months
- If the cleaning time using ethanol is more than 24 hours: Every month


7 Open the air release port on the DI water filter, releasing the air in the DI water filter, and close the air release port, then click [Next].

For details, see steps 4 to 8 (skipping step 7) in “Releasing Air in the DI Water Filter” (page 203).

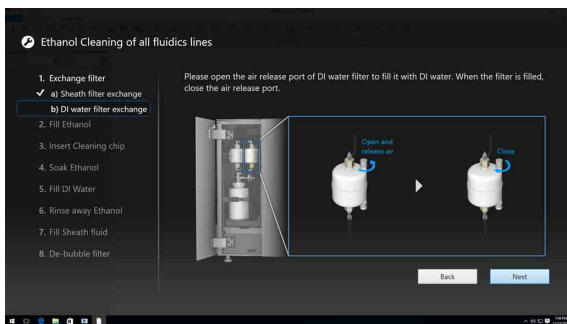

8 Remove the sheath tank and discard any remaining sheath fluid, then click [Next].

For details, see steps 6 to 8 in “Cleaning the Sheath Tank and Sheath Line” (page 183).

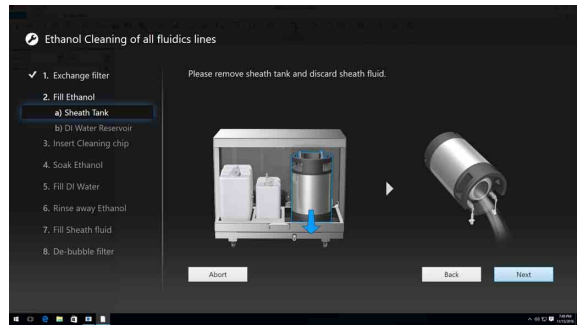


9 Pour 1 liter (0.25 gallons US) of ethanol into the sheath tank and reattach the lid.

Notes

- Do not touch the inner surface of the lid of the sheath tank. Also, take care not to spill any ethanol.
- Do not use a pump or other object that has not been sterilized for pouring ethanol into the tank.
- When inserting the sheath tank lid, make sure that the black rubber seal is attached to the lid.

10 Grasp the handles and shake the sheath tank to clean the inside of the tank, and place the sheath tank back in the fluidics cart, then click [Next].

For details, see steps 9 to 11 in “Cleaning the Sheath Tank and Sheath Line” (page 183).

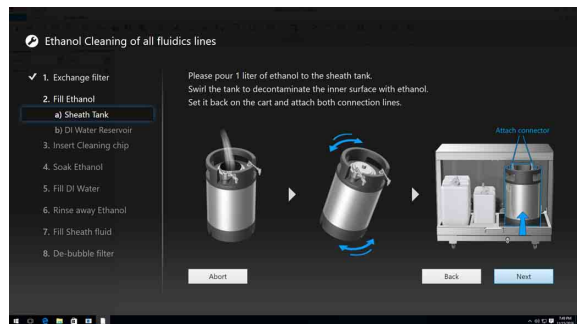

11 Remove the DI water tank and discard any remaining DI water, then click [Next].

For details, see steps 2 to 6 in “Cleaning the DI Water Tank and DI Water Line” (page 187).



12 Pour 600 ml (0.2 gallons US) of ethanol into the DI water tank and reattach the lid.


Notes

- Take care not to spill any ethanol.
- Do not use a pump or other object that has not been sterilized for pouring ethanol into the tank.

13 Remove the DI water tank filter from the DI water tank, and shake the tank while holding your finger over the filter connection port to clean the tank.


For details, see steps 5 and 8 in “Cleaning the DI Water Tank and DI Water Line” (page 187).

14 Attach the DI water tank filter to the DI water tank and place the DI water tank back in the main unit, then click [Next].

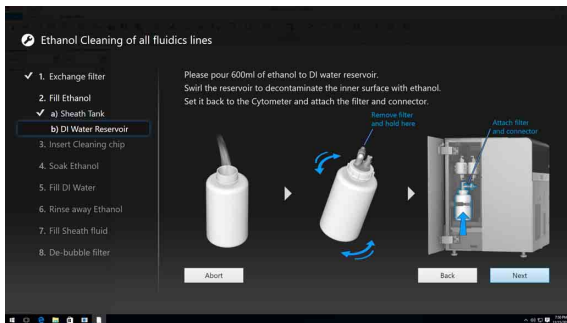


15 Hold the QR code on the cleaning chip packaging in front of the camera on the host computer.


The currently loaded sorting chip, if one is present, is ejected from the chip loader.

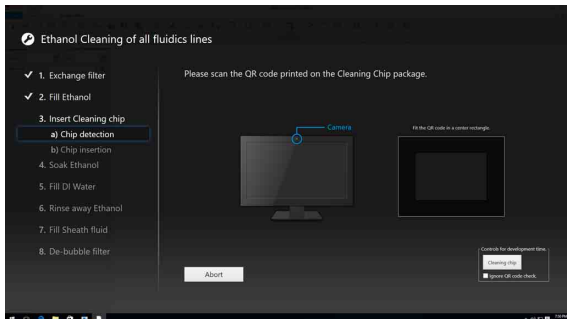

16 Insert the cleaning chip into the chip loading slot on the top of the chip loader, then click [Next].

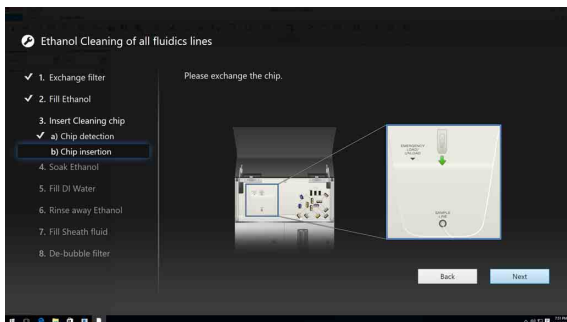


17 Load an empty 15 ml conical tube in the sample loader and place a check mark in the checkbox, then click [OK].

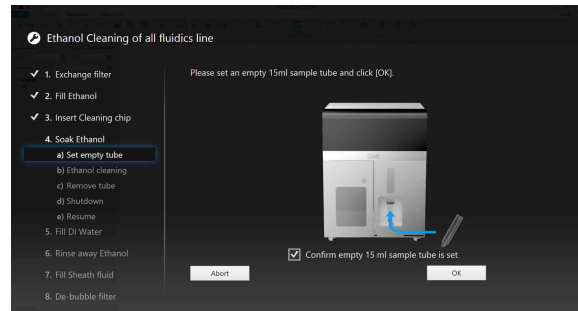

Note


Be sure to load an empty 15 ml conical tube in the sample loader before placing a check mark in the checkbox.

18 Select the ethanol cleaning time, then click [Start].

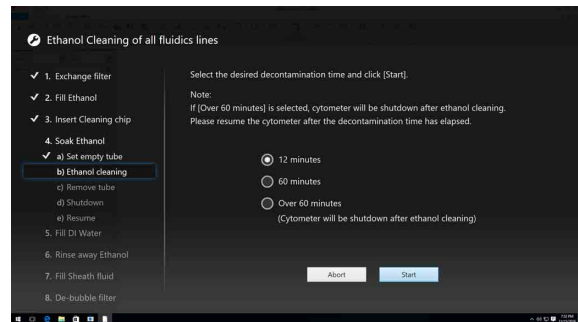

Ethanol cleaning starts.


During cleaning, a progress bar is displayed. Ethanol cleaning ends when the selected ethanol soak time elapses.

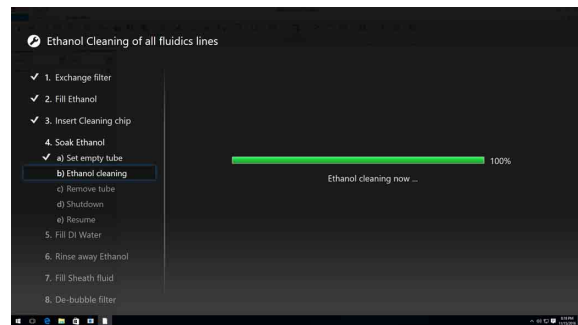

Note

If [Over 60 minutes] is selected, the main unit must be shut down after ethanol cleaning has finished for soak rinsing. Do not shut down the host computer. Leave the software running.

- 19** Remove the 15 ml conical tube and discard of the fluid in the tube, then click [Next].

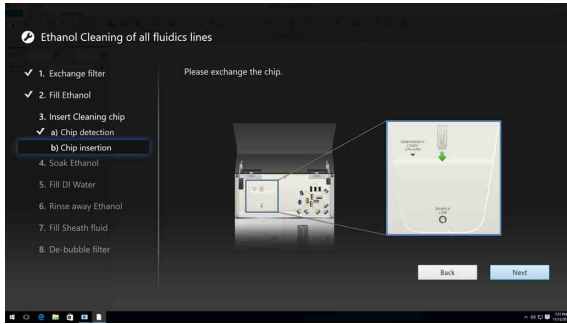


Tip

If [12 minutes] or [60 minutes] is selected in step **18**, proceed to step **22**. If [Over 60 minutes] is selected, proceed to step **20**.

Note

Dispose of the ethanol in accordance with the laboratory rules and local ordinances and regulations.

- 20** Click [Shutdown] to shut down the main unit and start the soak rinsing.

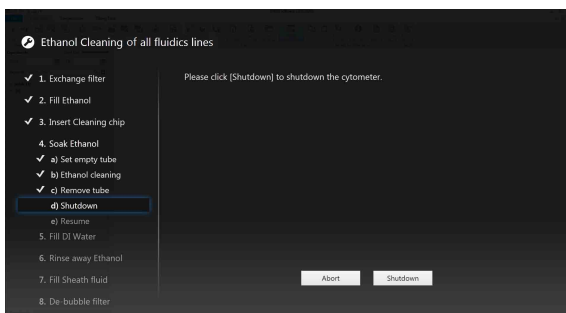


Tip

If [12 minutes] or [60 minutes] is selected in step **18**, this step is skipped.

- 21** After the soak rinsing time has elapsed, check that the compressed air supply is turned on, press the POWER button on the main unit to turn the unit on, then click [Resume].


Cell Sorter Software restarts when [Resume] is clicked.

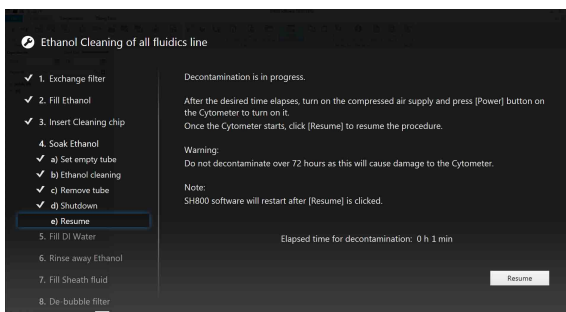

Tip

If [12 minutes] or [60 minutes] is selected in step **18**, this step is skipped.

- 22** Remove the sheath tank as in step **8** and discard any remaining ethanol, then click [Next].

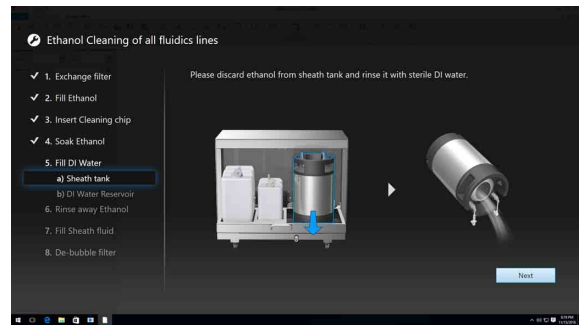

Note

Dispose of the ethanol in accordance with the laboratory rules and local ordinances and regulations.

- 23** Pour 1.5 liters (0.4 gallons US) of sterilized DI water into the sheath tank and reattach the lid.

Notes

- Do not touch the inner surface of the lid of the sheath tank. Also, take care not to spill any DI water.
- Do not use a pump or other object that has not been sterilized for pouring DI water into the tank.
- When inserting the sheath tank lid, make sure that the black rubber seal is attached to the lid.

- 24** Grasp the handles as in step **10** and shake the sheath tank to clean the inside of the tank, and place the sheath tank back in the fluidics cart, then click [Next].



25 Remove the DI water tank as in step **11** and discard any remaining ethanol, then click [Next].




Note


Dispose of the ethanol in accordance with the laboratory rules and local ordinances and regulations.

26 Pour 1 liter (0.25 gallons US) of sterilized DI water into the DI water tank and reattach the lid.

27 Clean the inside of the DI water tank as in steps **13** to **14**, and place the DI water tank back in the main unit, then click [Next].

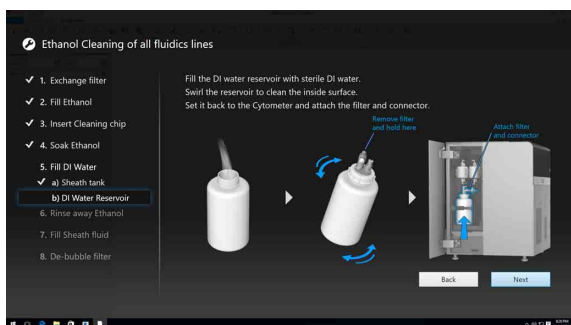


28 Load an empty 15 ml conical tube in the sample loader and place a check mark in the checkbox, then click [OK].

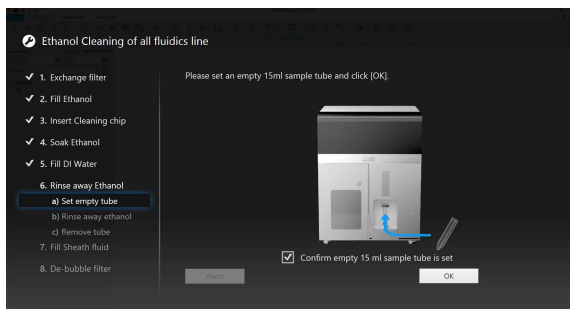


Note

Be sure to load an empty 15 ml conical tube in the sample loader before placing a check mark in the checkbox.


29 Click [Start].

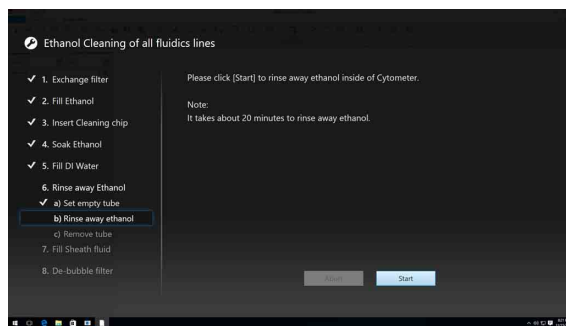

Ethanol removal rinsing starts. During rinsing, a progress bar is displayed.

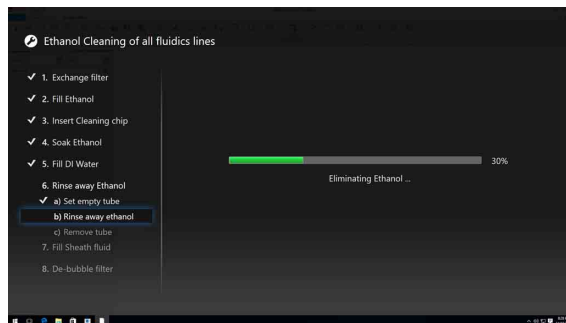

30 Remove the 15 ml conical tube and discard of the DI water in the tube, then click [Next].

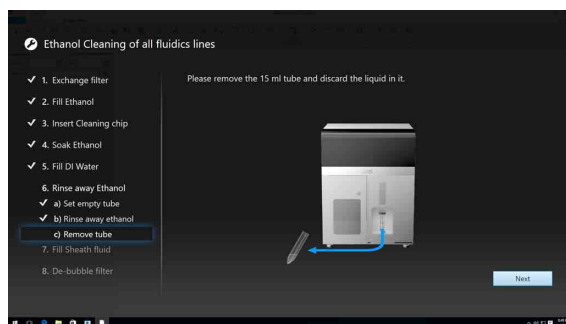

Note

Dispose of the DI water in accordance with the laboratory rules and local ordinances and regulations.

31 Remove the sheath tank as in step **8** and discard any remaining DI water, then click [Next].




Note

Dispose of the DI water in accordance with the laboratory rules and local ordinances and regulations.

- 32** Fill the sheath tank with sheath fluid and place the sheath tank back in the fluidics cart, then click [Next].


For details, see steps **10 to 11** in “Cleaning the Sheath Tank and Sheath Line” (page 183).

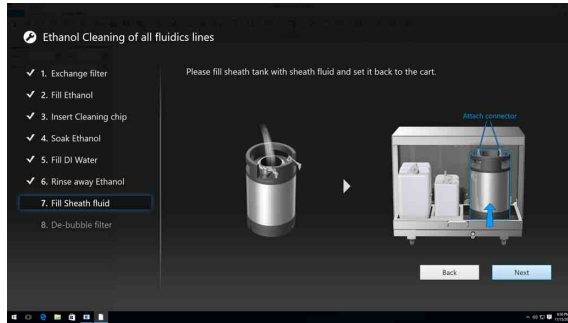


- 33** Remove the DI water tank and discard any remaining DI water.

- 34** Refill the DI water tank with DI water, place the tank back in the main unit, and then click [Next].

For details, see “Refilling the DI Water Tank” (page 167).

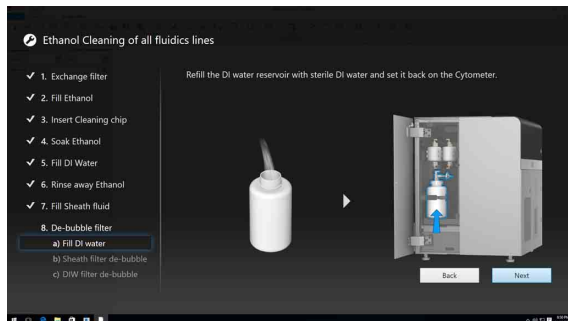


- 35** Click [Start] to begin de-bubbling the sheath filter. When de-bubble is completed, click [Next].

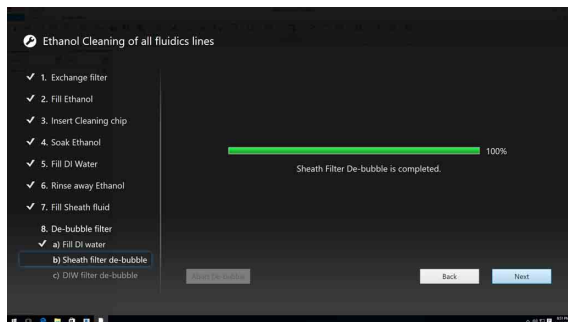


- 36** Follow the on-screen instructions to release air from the sheath filter, then click [Next].

For details, see steps **4 to 8** of “To release air in the sheath filter manually” in “Releasing Air in the Sheath Filter and Sorting Chip” (page 202).

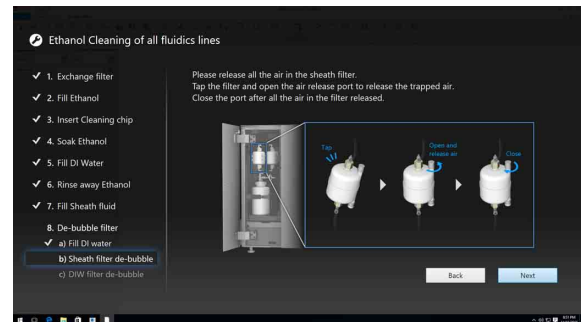


- 37** Check that all air has been released from the sheath filter, then click [Next].

If there is still air in the filter, click [De-bubble again] and repeat steps **35 to 36**.

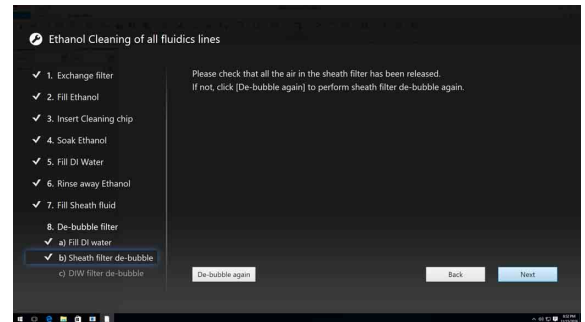


- 38** Follow the on-screen instructions to open the air release port on the DI water filter, releasing the air in the DI water filter, and close the air release port, then click [Next].

For details, see steps **4 to 8** (skipping step **7**) in “Releasing Air in the DI Water Filter” (page 203).

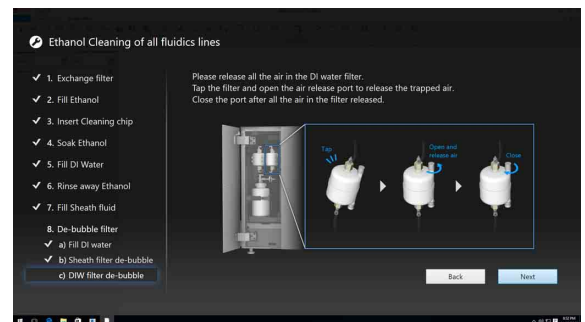


39 Click [OK] to exit the [Ethanol Cleaning for all fluidics line] wizard.

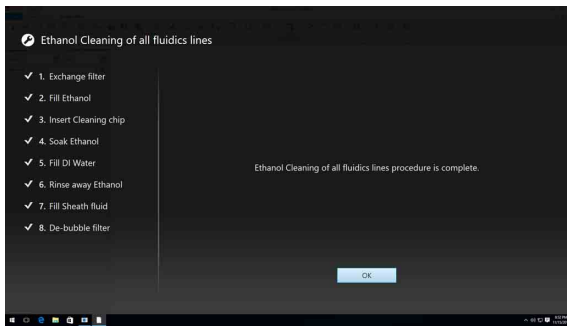

This chapter describes some of the basic concepts of flow cytometry and cell sorter operating principles of the SH800.

Fluid Flow

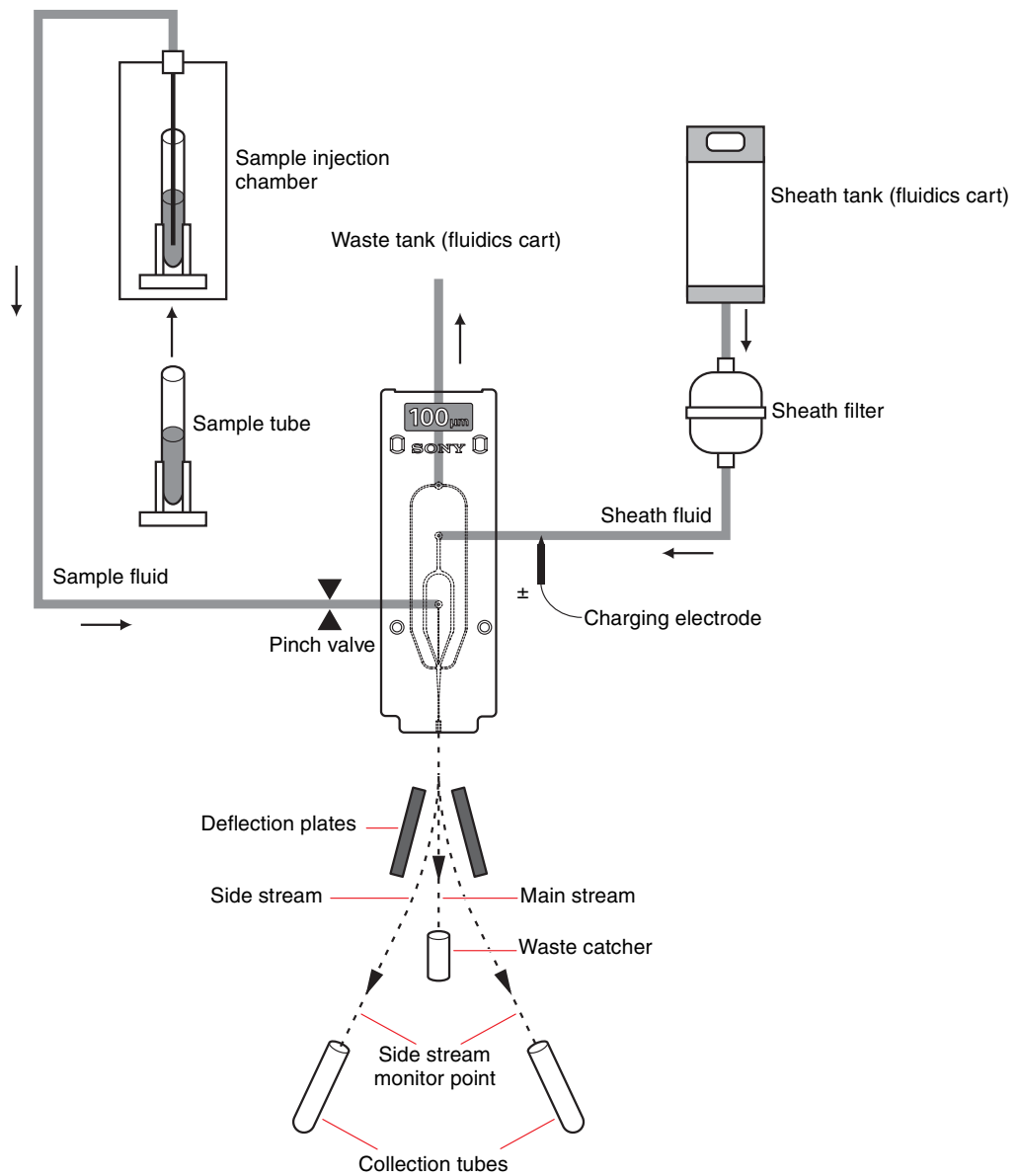

Appendix A Operating Principles

Sample Fluid

Sample preparation is an important part of the testing procedure in flow cytometry. Sample fluid must contain cells in a monodispersed suspension that has been filtered using an appropriate filter mesh to remove aggregates. This is required to maintain a consistent sample concentration. If the sample is not filtered, the sample may clog or adhere to the surfaces of fluidics components, and stable performance cannot be maintained. For more information about sample preparation, refer to reference sources.

Before running samples, check that the sorting chip nozzle size is appropriate for the size of cells in the sample fluid. As a generally accepted rule of thumb, the nozzle size should be approximately five times (min) the diameter of the largest cell in the sample fluid to maintain stable sample fluid flow without clogging the chip.

The standard sorting chip has a 100 μm nozzle, which is suitable for most cell sorting applications. Cells ranging from approximately 1 μm to 20 μm in diameter can be analyzed and sorted.

Note

The size of the sorting chip nozzle affects the droplet size, the gap between droplets, and the consumption of sample fluid and sheath fluid.

Sample fluid can be loaded in 0.5 ml micro tubes, 1.5 ml micro tubes, 5 ml round tubes or 15 ml conical tubes for use with the SH800. Each type of tube fits into a corresponding tube holder for the sample tube. The tube holder all fit into a D-cut slot in the base of the sample loader.

After the sample is loaded and [Start] is clicked in Cell Sorter Software, the sample door closes and the sample tube containing the sample is raised into the injection chamber. The chamber is hermetically sealed and then pressurized using regulated compressed air. The pressure inside the injection chamber forces the sample fluid up through the sample probe and into the sample line connected to the probe at the top of the chamber. Sample fluid flows through the sample line and into the sample fluid inlet port on the front of the sorting chip, via a pinch valve which opens/closes the sample line to control the sample flow.

The sample fluid flow rate is controlled by the [Sample Pressure] setting on the main window in Cell Sorter Software. The sample pressure determines the speed of the sample flow, which indirectly controls the distance between cells in the sample fluid flow as they are injected into the sheath fluid in the sorting chip.

An agitation unit is provided to maintain consistency of the sample fluid during testing. The unit is switched on/off from Cell Sorter Software.

Sheath Fluid


Sheath fluid is supplied from the sheath tank in the fluidics cart. A level sensor in the fluidics cart monitors the level in the tank. The sensor output is displayed on the LCD monitor and on the bottom right of the main window in Cell Sorter Software. A warning is displayed if the level gets close to empty. The tank is pressurized using regulated compressed air, which is used to force the sheath fluid through the fluidics system components at constant pressure, avoiding the pressure fluctuation problems common to pumps.

When [Start] is clicked in Cell Sorter Software to begin sorting, sheath fluid is forced under pressure through the sheath filter and into the sheath fluid inlet port on the front of the sorting chip. A charging electrode mounted near the inlet port is used to electrostatically charge the sheath fluid for sorting.


The sheath fluid flow rate is controlled by the [Sheath Pressure] setting in Cell Sorter Software. The sheath pressure determines the speed of the sheath fluid flow through the sorting chip.

Hydrodynamic Focusing

Sample fluid and sheath fluid are forced into the sorting chip using air pressure. The sample fluid is injected into the center of the sheath fluid stream, hydrodynamically focusing the sample fluid within the sheath fluid, according to the principles of fluid dynamics. The internal micro-channel structure of the sorting chip encases the sample fluid within the sheath fluid to ensure high-sensitivity, stable optical detection characteristics.

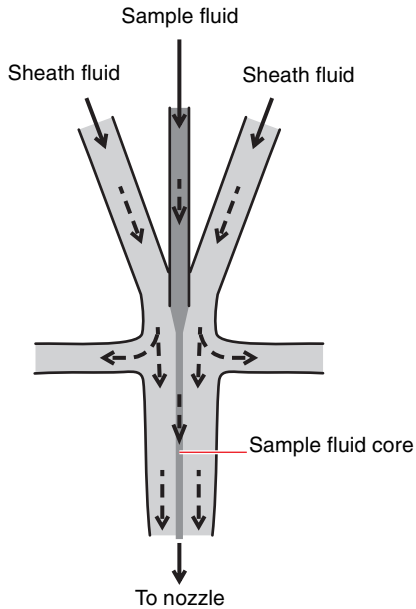

The relative difference in pressure between the sheath fluid flow and the sample fluid flow forces the diameter of the sample fluid core to vary. When the sample pressure is lowered, the diameter of the sample fluid core decreases, “focusing” the sample fluid into a narrow core stream. When the sample pressure is raised, the diameter increases with a corresponding increase in the sample event detection rate.

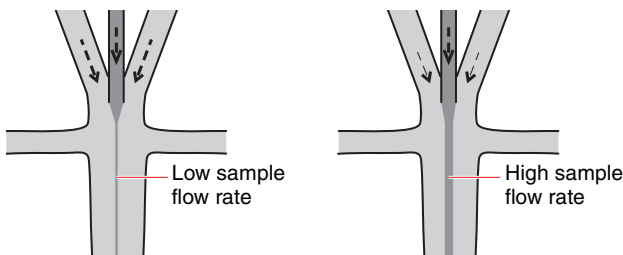


The flow of sample cells encased in sheath fluid passes the optical detection point where the fluid core is interrogated by up to four lasers of different wavelengths.

The highest measurement resolution is obtained when the sample fluid flow is focused into a single-file stream of cells moving past the optical detection point at low sample flow pressure, maximizing the exposure to excitation laser energy that each cell receives.

The diameter of the channel in the sorting chip narrows as it approaches the nozzle at the bottom of the chip, increasing the flow speed while maintaining the relative proportion between the sheath and sample fluid. Steady fluid flow ejected from the nozzle is critical for accurate sorting, hence vibrations of the main unit or the sheath line on the rear panel must be avoided during sorting.

Droplet Sorting

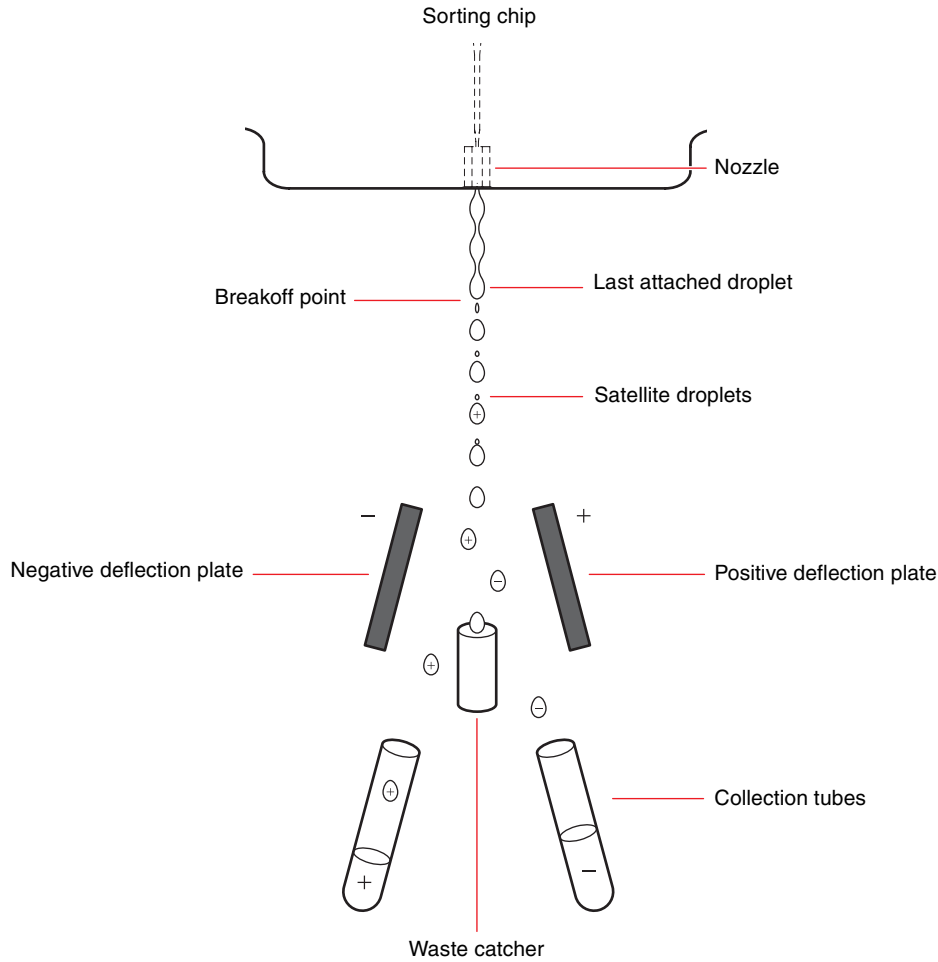

When the sorting chip is inserted in the chip loader, a transducer causes the chip to vibrate at an ultrasonic frequency as fluid flows through the chip. The vibrations cause the fluid flow to be ejected in a jet of uniform droplets from the nozzle at the bottom of the chip.

At a point a short distance beneath the sorting chip, called the breakoff point, the fluid flow breaks into a very regular stream of uniform droplets. The size of the droplets is determined by the nozzle size, frequency of vibration, and the speed of the fluid flow. The amplitude of vibration (drop drive voltage) controls the position of the breakoff point, and the frequency of vibration determines the distance between droplets.

If the last attached droplet satisfies the sorting criteria, the droplet containing the cell receives an electrostatic charge just before the fluid flow reaches the breakoff point from a charging electrode located near the sheath fluid inlet on the sorting chip. The last attached droplet retains the electrical charge after it breaks off from the fluid flow. The electrostatic charge is distributed on the periphery of the droplet and does not affect any sample cells contained within the droplet.



The polarity (positive or negative) of the charge applied to a droplet depends on the cell sorting criteria defined by the characteristics of the cell populations that you want to isolate for sorting using gates specified in Cell Sorter Software. This information is then used to drive the charging pulse for sorting. Unwanted droplets are left uncharged and are collected by the waste catcher.

In an ideal sort, all cells travel pass the optical detection point in single file, and are sorted one cell at a time in separate droplets, with an interval between droplets. In practice, however, cells can appear very close together at the optical detection point, which means they may be found within the same droplet or in adjacent droplets. The sorting mode selection in Cell Sorter Software determines whether to sort or abort droplets whenever these “coincidences” occur.

See “*Sorting Mode*” (page 106).

The time taken for droplets to traverse the distance between laser interrogation at the optical detection point and the breakoff point is called the sort delay. The sort delay controls the timing of the electrostatic charging of the target droplet for stable sorting.

The droplet camera monitors the droplet stream before and after the breakoff point. The stroboscopic output image of the droplet camera is displayed in the bottom of the main window in Cell Sorter Software. The output image is also used by control circuits to automatically adjust the amplitude and frequency of vibration of the sorting chip to control the position of the breakoff point and the timing of the electrostatic charging pulse. Manual droplet control is also supported.

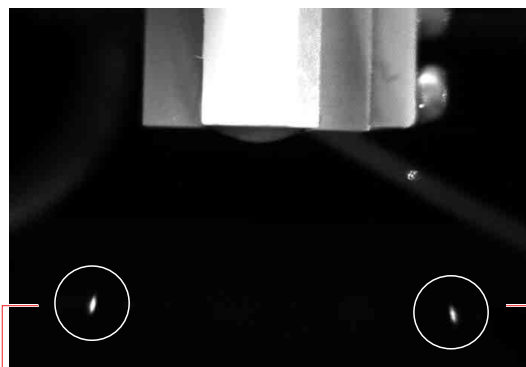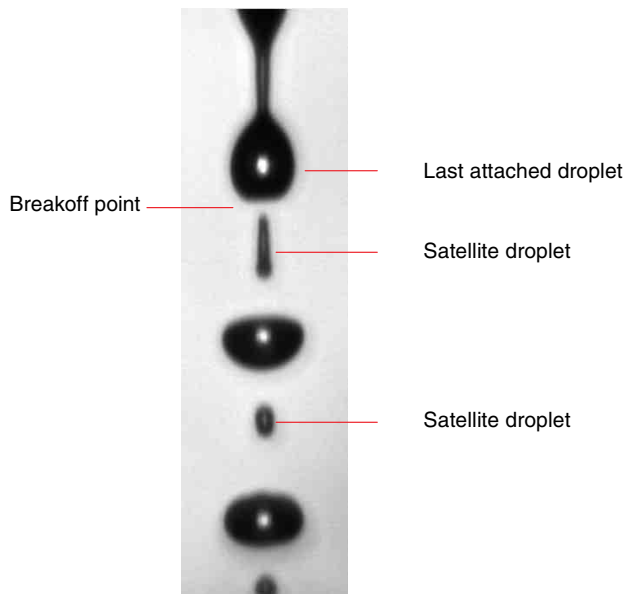

Left side stream

Right side stream

WARNING

The deflection plates are charged at extremely high voltages. Contact with charged deflection plates can result in serious injury or electrical shock. Do not touch the deflection plates while power is connected to the main unit.

The collection area is sealed during operation to prevent contamination of the air by aerosols generated by falling droplets and to prevent air circulation in the room affecting the accuracy of the droplet sorting. Negative pressure (vacuum suction) helps reduce the risks associated with aerosol generation.

WARNING

Waste fluid may contain biological, chemical, or other agents. Always wear gloves and other protective clothing, mask, and goggles, as required, when handling samples. The handling of waste fluid should be performed in accordance with biological hazard handling safety procedures.

The formation of droplets at the breakoff point and the angular deflection of side streams are calibrated at startup using automatic setup beads.

For details about calibration, see “Automatic Calibration” (page 50).

High-voltage deflection plates on either side of the droplet stream deflect the charged droplets, using electrostatic attraction and repulsion, to the left and right as the droplets fall between the deflection plates, sorting the droplets into two side streams (two-way sorting). The polarity of the charge on each droplet determines the direction of the stream (left or right). Uncharged droplets (main stream) are not deflected and fall into the waste catcher. The side streams are directed toward the collection tubes on the collection stage.

The side stream monitor detects the position of the side streams. The side stream laser irradiates the side streams, and an image captured by the side stream camera is displayed at the bottom of the main window in Cell Sorter Software. The output image is also used by control circuits to automatically adjust the angle of deflection of the side streams so that they are directed into the collection tubes. The collection stage supports 15 ml conical tubes and 5 ml round tubes inserted in collection tube holders.


Optical System

This section describes the signals that can be generated using the SH800.

Light and Signal Generation

Light Scatter


When laser light hits a cell passing the optical detection point in the sorting chip, light is scattered in all directions. Light scattered in the forward direction is called forward scatter (FSC). Light scattered to the sides and back toward the light source is called back scatter (BSC).

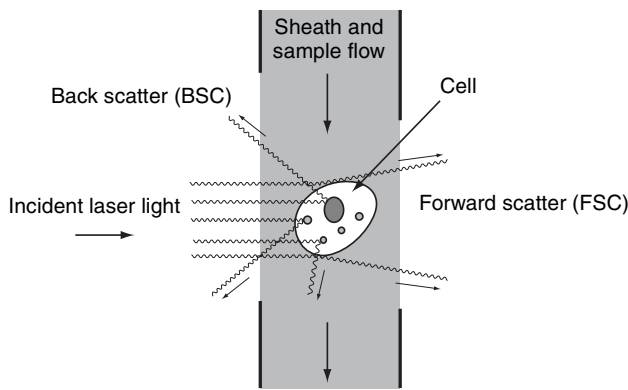


In general terms, forward scatter can provide information about the size of the cell or can indicate the state of living cells. Back scatter can provide information about the internal detail of the cell, indicating complexity, granularity, and irregularities in a cell.

Fluorescence

When laser light hits a cell passing the optical detection point in the sorting chip, the light excitation can cause fluorochromes in molecules or antibodies that have been labeled with fluorescent markers to fluoresce in the sample, producing fluorescent light (FL). Fluorochromes for markers emit high-level fluorescent light when excited by lasers at certain wavelengths and low-level fluorescent light at other wavelengths. The wavelength of the excitation lasers should be selected to match the fluorochromes for the markers.

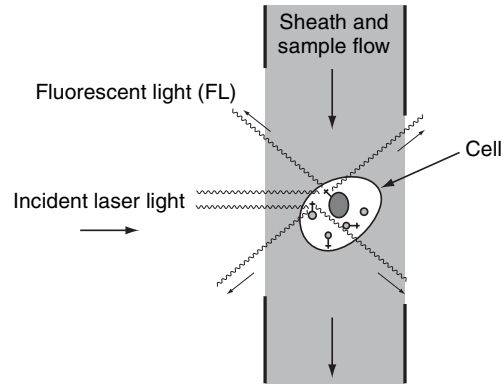


Fluorescence occurs when an electron in a molecule is raised to an excited state as the electron absorbs a photon of the incident excitation light. The electron loses some of this energy before the molecule emits a different photon and the electron returns to its ground state. The emitted photon has less energy than the absorbed photon, so the emitted light has a longer wavelength than the incident light and hence a different color. The difference between the peak excitation wavelength and peak emitted wavelength is called the Stokes shift. Some fluorochromes are available that exhibit a narrow Stokes shift while others exhibit a wide Stokes shift. Multicolor fluorescent light analysis can be used to provide a wealth of structural and other functional information about cells.

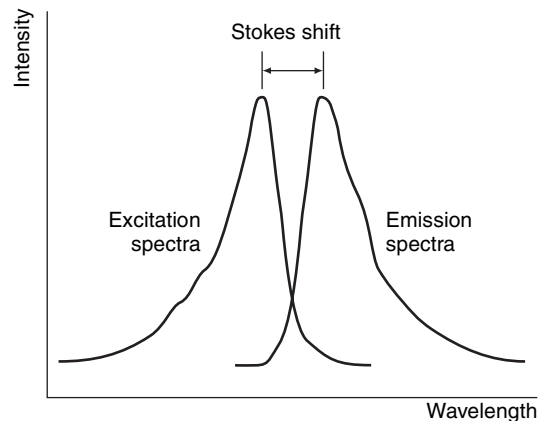

Laser Interrogation


Light source

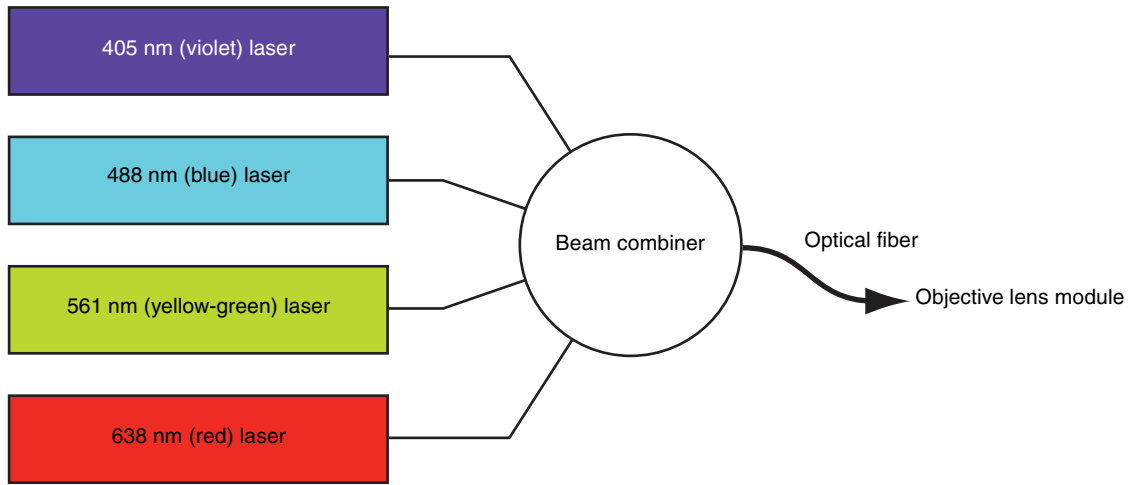

* 4-laser models

The excitation laser module supplies the laser light used to interrogate individual cells as they flow past the optical detection point. The excitation laser beams are combined into a single collinear beam and coupled to a single-mode fiber to the objective lens module. The excitation laser module uses Automatic Power Control (APC) to provide optical measurement accuracy and reproducibility.

Light collection

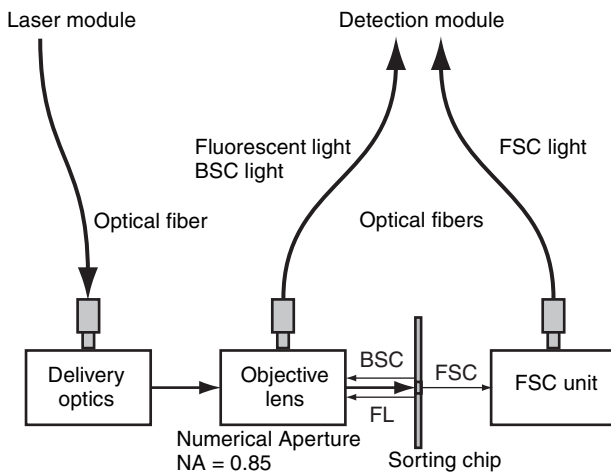


The objective lens module focuses the beam from the laser module onto the optical detection point in the sorting chip.



The laser light source from the excitation laser module is focused on the optical detection point in the sorting chip.

Detection

The detection module takes the light collected from the objective lens module and separates it into eight channels of light, comprising forward scatter light (FSC), back scatter light (BSC), and six channels of fluorescent color light (FL1 to FL6). The output from each channel is received by a detector (photodiode (PD) or photomultiplier tube (PMT)) that converts the light energy into an electrical signal for processing by the acquisition module.

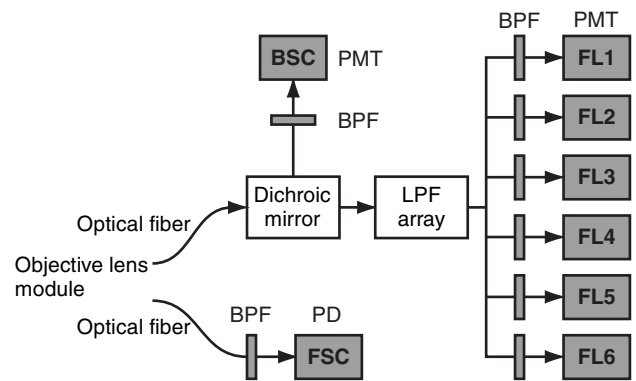


The detection module is comprised by an array of dichroic longpass optical filters (LPF) and bandpass optical filters (BPF). Dichroic longpass optical filters reflect light below a certain wavelength and transmit light above a certain wavelength. Bandpass optical filters transmit light in a narrow band of wavelengths and block light of all other wavelengths. For bandpass optical filters, the first number in the name indicates the center wavelength and the second indicates the bandpass width (e.g. 525/50 is a 525 nm bandpass optical filter with 50 nm bandwidth).

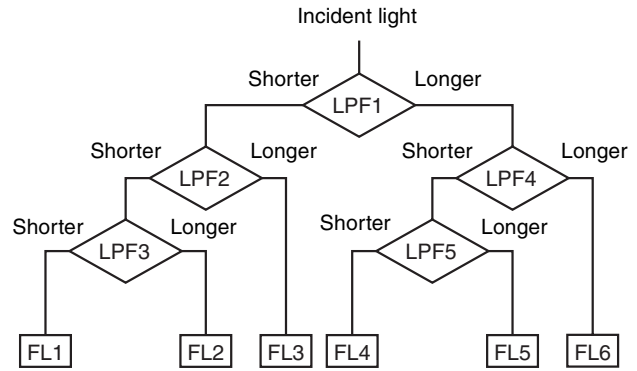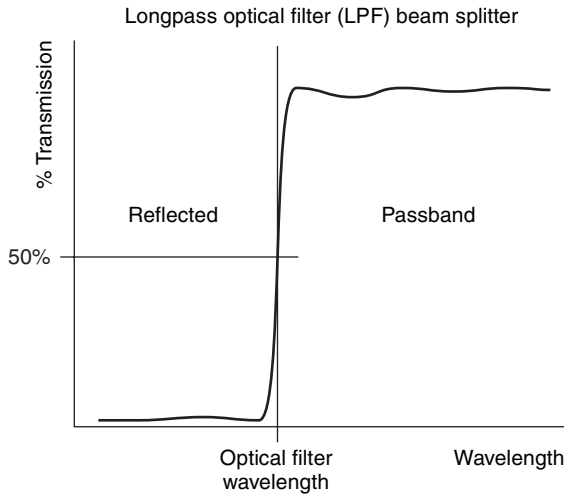


The SH800 can detect six channels of fluorescent light simultaneously. The laser configuration and the optical filter pattern determine the fluorescent markers that can be detected.

For a list of fluorochromes that can be detected for each model, see "Fluorochrome Detection Matrix" (page 232).

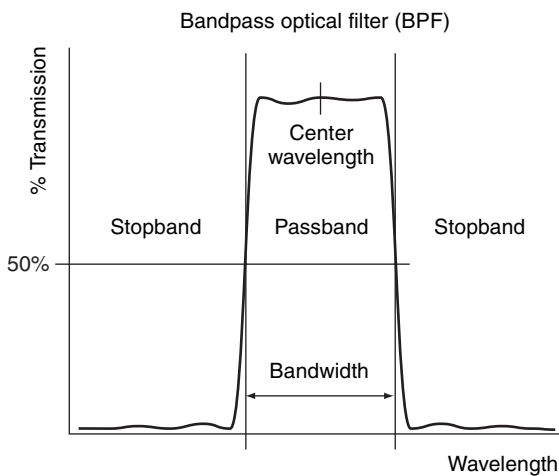

The output from the photodiode and PMT detectors is passed to the acquisition module. The pulse shape that is detected by each detector varies depending on the size and type of cell, and the type of light being collected. The SH800 measures three parameters per pulse (area, height, and width) to provide a wealth of information about cells. The width of the pulse provides an indication of the size of a cell, the height of the pulse provides an indication of the peak fluorescence intensity (brightness) of the cell, and the area provides an indication of the total fluorescent material contained in the cell. The pulse parameters can be selected independently for each channel.

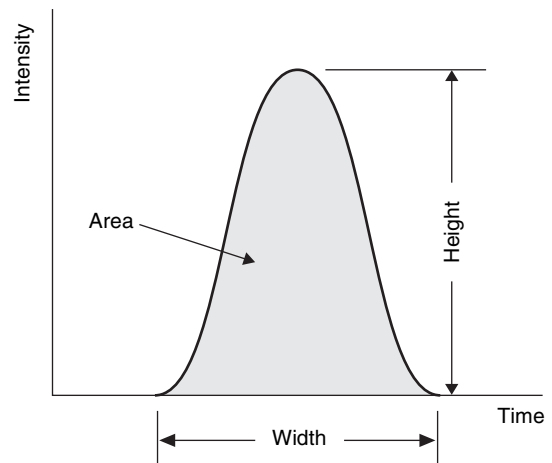

Forward scatter light passes through a collimator and a bandpass optical filter (BPF), and is detected by a photodiode (PD). The bandpass optical filter determines the wavelength of the laser source for which forward scatter is measured.

Note

The SH800 is configured by default to measure forward scatter light at a wavelength of 488 nm.

Back scatter light passes through a bandpass optical filter and is detected and amplified by a photomultiplier tube (PMT) detector. The bandpass optical filter determines the wavelength of the laser source for which back scatter is measured.

Note


The SH800 is configured by default to measure back scatter light at a wavelength of 488 nm.

Fluorescent lights passes through an array of longpass optical filters (LPFs), separating the fluorescent light into individual fluorescent channels. Each color then passes through a bandpass optical filter and is detected and amplified by a PMT detector.

When sorting or analyzing cells, you can set a threshold level for the pulse height for one of the channels (BSC, FSC, FL1 to FL6) to discriminate between real events and background noise. The threshold level for a channel sets the minimum light detector output that is required to record an event. When the trigger is set, data is acquired only for events for which the detector output exceeds the threshold level. Values below the threshold level are ignored. The FSC channel threshold level is the most

commonly used trigger, as this channel is less susceptible to the effects of background noise.

The pulse at the threshold level can also be effectively widened to increase the region of the pulse included in the detected signal. This is achieved by adding pulse data that was digitized just before the pulse rises above the threshold value and just after the pulse falls below the threshold value. These two additional components are called forward window extension (FWE) and back window extension (BWE), respectively. Window extensions are used to provide more accurate measurements of the area pulse parameter. When the window extensions are set to zero, only the area above the threshold level is included in the area calculation. By setting window extensions before and after the threshold crossing points, a more accurate determination of the total area under the pulse can be achieved.

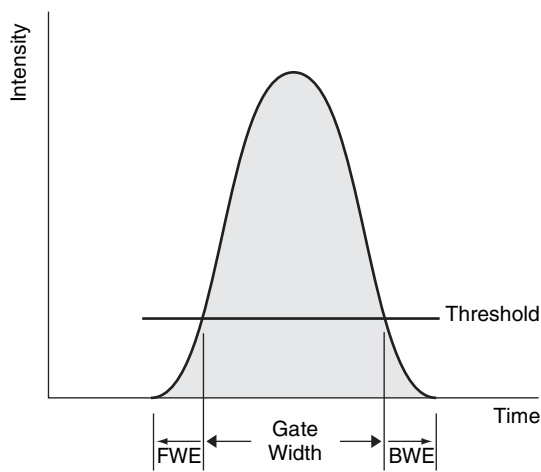


The optimum length of the window extensions depends on the rise height and width of the pulse. Hence, the optimum window extension will vary from pulse to pulse within the same sample, and from experiment to experiment. If the window extensions are too narrow, the measurement for the area pulse parameter will be smaller than its real value. If the window extensions are too wide, then the window extension may begin to overlap with other pulses, degrading the calculation.

When a cell is detected during sorting, it is generally not possible to know with great accuracy if the signal represents the passage of a large single cell, small multiple cells adhering to one another, or other non-cellular material. Accordingly, it is more common to refer to each signal detection not as a cell, but as an event.

Fluorescence Compensation


When detecting the emitted fluorescent light for a fluorochrome, components of the emitted light can appear in one or more of the fluorescent light detectors (FL1 to FL6) due to the wide emission spectrum of each fluorochrome. When detecting the emitted light for more than one fluorochrome, components of the spectrum of all fluorochromes can appear in multiple detectors, a

phenomenon called spillover. The problem then becomes one of working out the proportion of the light detected in each detector due to the individual fluorochromes. For example, in the following diagram, the output of detector FL1 is almost totally due to the spectra of fluorochrome A. The output of detector FL3 is predominantly due to the spectra of fluorochrome B, but it does contain a small component (approximately 10%) due to fluorochrome A. However, the output of FL2 contains significant components from the spectra of both fluorochromes A and B. This problem increases considerably as the number of fluorochromes being simultaneously detected increases.

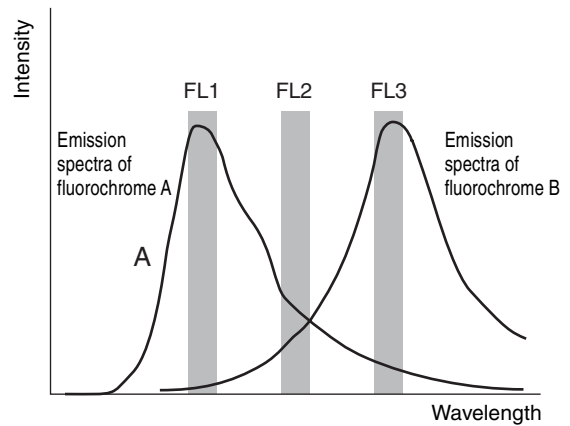

Fluorescence compensation is the process of calculating the components of the detected signals due to each fluorochrome in the sample and subtracting the “unwanted” components from each channel. This converts the channel detector output from signals of fluorescence intensity at a given wavelength to signals of fluorescence intensity due to each fluorochrome.

Compensation adjusts the mean value of fluorescence intensity of populations that are not stained with a particular fluorochrome to the mean value of a negative control sample, which by definition is not stained with any fluorochromes. When compensated, the populations effectively appear aligned into quadrants on 2D plots.

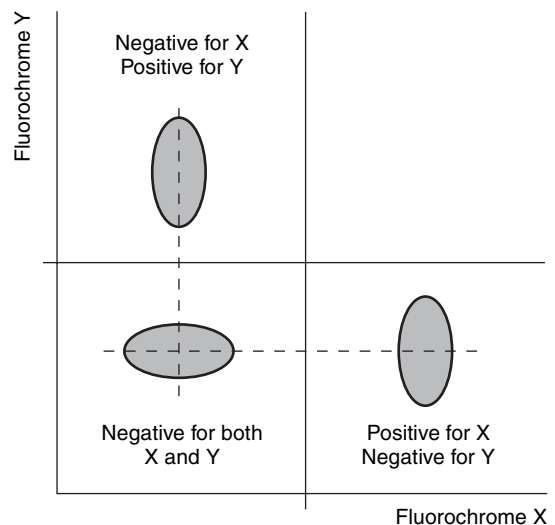

Notes

- The visual representation of the compensated data on a 2D plot will vary depending on the axis type (log, linear, or biexponential).
- The negative and positive populations may or may not appear as diverse populations on 2D plots, depending on the type of cells in the sample and the selected fluorochromes. For example, if samples contain cells that are actively acquiring or losing the antigen with the fluorochrome marker, the two populations may tend to merge into one another.
- Cell Sorter Software also supports manual specification of the fluorescent compensation parameters. See “*Adjusting Fluorescence Compensation Manually*” (page 89).



Series Lineup

The SH800 cell sorter is available in six different models equipped with a combination of one, two, three, or four excitation lasers of different wavelengths. The models are indicated by a two-letter suffix (AC to FC).

SH800S

Model	No. of lasers	Laser wavelengths (nm)
LE-SH800SA	1	488
LE-SH800SB	2	488, 638
LE-SH800SC	2	405, 488
LE-SH800SD	2	488, 561
LE-SH800SE	3	405, 488, 638
LE-SH800SF	4	405, 488, 638, 561
LE-SH800SG	3	405, 488, 561
LE-SH800SH	3	488, 638, 561

SH800SP

Model	No. of lasers	Laser wavelengths (nm)
LE-SH800SAP	1	488
LE-SH800SBP	2	488, 638
LE-SH800SCP	2	405, 488
LE-SH800SDP	2	488, 561
LE-SH800SEP	3	405, 488, 638
LE-SH800SFP	4	405, 488, 638, 561
LE-SH800SGP	3	405, 488, 561
LE-SH800SHP	3	488, 638, 561

SH800Z

Model	No. of lasers	Laser wavelengths (nm)
LE-SH800ZA	1	488
LE-SH800ZB	2	488, 638
LE-SH800ZC	2	405, 488
LE-SH800ZD	2	488, 561
LE-SH800ZE	3	405, 488, 638
LE-SH800ZF	4	405, 488, 638, 561
LE-SH800ZG	3	405, 488, 561
LE-SH800ZH	3	488, 638, 561

SH800ZP

Model	No. of lasers	Laser wavelengths (nm)
LE-SH800ZAP	1	488
LE-SH800ZBP	2	488, 638
LE-SH800ZCP	2	405, 488
LE-SH800ZDP	2	488, 561
LE-SH800ZEP	3	405, 488, 638
LE-SH800ZFP	4	405, 488, 638, 561
LE-SH800ZGP	3	405, 488, 561
LE-SH800ZHP	3	488, 638, 561

SH800

Model	No. of lasers	Laser wavelengths (nm)
LE-SH800AC	1	488
LE-SH800BC	2	488, 638
LE-SH800CC	2	488, 405
LE-SH800DC	2	488, 561
LE-SH800EC	3	488, 405, 638
LE-SH800FC	4	488, 405, 638, 561

- 488 nm blue laser (standard)
- 405 nm violet laser
- 561 nm yellow-green laser
- 638 nm red laser

The fluorochromes that can be excited and the resulting emission spectra that can be detected by each laser configuration depend on the wavelength of the lasers and the optical filter pattern. The fluorochrome detection matrix for each laser configuration is given in the “*Fluorochrome Detection Matrix*” (page 232).

Tip

On multi-laser models, you specify the lasers used for experiments when Cell Sorter Software launches.

Notes

- Do not use the 405 nm laser for extended periods of time. Extended use can damage the sorting chip.
- The use of optical adjustment procedures not specified in this document or installation/removal of optical components may result in hazardous radiation exposure. Do not remove or otherwise modify the excitation laser or detection module.

Optical Filter Patterns

The detection module takes the scattered and fluorescent light collected from cells and separates it into eight channels of light, comprising the forward scatter channel, back scatter channel, and six channels of fluorescent light of various wavelengths.

The SH800 uses a combination of longpass optical filters (in slots labeled LP1 to LP5) and bandpass optical filters (in slots labeled BSC, FSC, and FL1 to FL6) arranged in an optically optimized layout, according to the laser configuration for an experiment.

The SH800 uses two standard optical filter patterns.

Optical filters	Optical filter pattern 1*1	Optical filter pattern 2*2
LP1	639LP	639LP
LP2	600LP	561LP
LP3	561LP	487.5LP
LP4	752LP	752LP
LP5	685LP	685LP
FL1	525/50	450/50
FL2	585/30	525/50
FL3	617/30	600/60
FL4	665/30	665/30
FL5	720/60	720/60
FL6	785/60	785/60
FSC	488/17F	488/17F
BSC	488/17B	488/17B

*1. On the following models, dummy filters are installed in the RESERVE slots.

SH800AC, SH800BC, SH800DC, SH800ZA, SH800ZB, SH800ZD, SH800ZH, SH800ZAP, SH800ZBP, SH800ZDP, SH800ZHP, SH800SA, SH800SB, SH800SD, SH800SH, SH800SAP, SH800SBP, SH800SDP, SH800SHP

*2. On the following models, filters not used in the current optical filter pattern should be stored in the RESERVE slots. Filter pattern 1 can also be used when the 405 nm laser is turned off to make efficient use of the six fluorescence detection channels.

SH800CC, SH800EC, SH800FC, SH800ZC, SH800ZE, SH800ZF, SH800ZG, SH800ZCP, SH800ZEP, SH800ZFP, SH800ZGP, SH800SC, SH800SE, SH800SF, SH800SG, SH800SCP, SH800SEP, SH800SFP, SH800SGP

- Filter pattern 1 can also be used when the 405 nm laser is turned off to make efficient use of the six fluorescence detection channels.
- The shaded portion of the table indicates the filters that are different between optical filter patterns 1 and 2.


Tips

- Optic filter pattern 2 is used only on the following models equipped with a 405 nm laser.
SH800CC, SH800EC, SH800FC, SH800ZC, SH800ZE, SH800ZF, SH800ZG, SH800ZCP, SH800ZEP, SH800ZFP, SH800ZGP, SH800SC, SH800SE, SH800SF, SH800SG, SH800SCP, SH800SEP, SH800SFP, SH800SGP
Models other than the above use optic filter pattern 1 only.
- The longpass optical filters are bigger than the bandpass optical filters to prevent the accidental

insertion of optical filters in the wrong slots. The LP1 to LP5 longpass optical filters are labeled with the wavelength rating of the optical filter and an “LP” suffix (e.g. “639LP” for 639 nm longpass optical filter). The FL1 to FL6 bandpass optical filters are labeled with the center wavelength of the optical filter and the bandwidth of the optical filter (e.g. “525/50” for 525 nm bandpass optical filter 50 nm bandwidth). The FSC and BSC bandpass optical filter labels have an “F” and “B” suffix, respectively (e.g. “488/17F” for 488 nm bandpass optical filter with 17 nm bandwidth for forward scatter detection).


- Optical filters not used in the current filter pattern can be stored in the RESERVE 1 to 3 slots.

Optical Filter Pattern 1

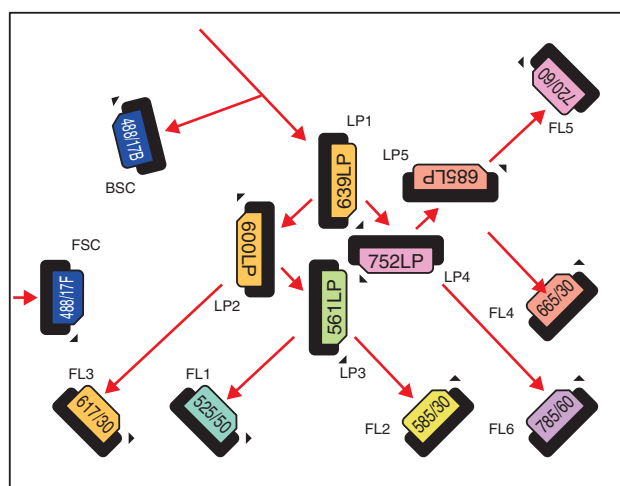

Optical filter label	Installed optical filter
LP1	639LP
LP2	600LP
LP3	561LP
LP4	752LP
LP5	685LP
FL1	525/50
FL2	585/30
FL3	617/30
FL4	665/30
FL5	720/60
FL6	785/60
FSC	488/17F
BSC	488/17B


Optical Filter Pattern 2

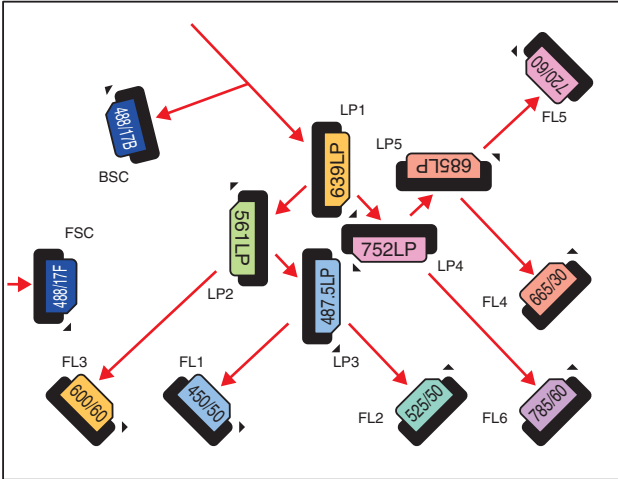

Optical filter label	Installed optical filter
LP1	550LP
LP2	525LP
LP3	495LP
LP4	725LP
LP5	605LP
FL1	450/50
FL2	510/20
FL3	535/30
FL4	585/40
FL5	640/30
FL6	785/60
FSC	488/17F
BSC	488/17B


Optical filter label	Installed optical filter
LP1	639LP
LP2	561LP
LP3	487.5LP
LP4	752LP
LP5	685LP
FL1	450/50
FL2	525/50
FL3	600/60
FL4	665/30
FL5	720/60
FL6	785/60
FSC	488/17F
BSC	488/17B

Filter Pattern for BV/PE/PI

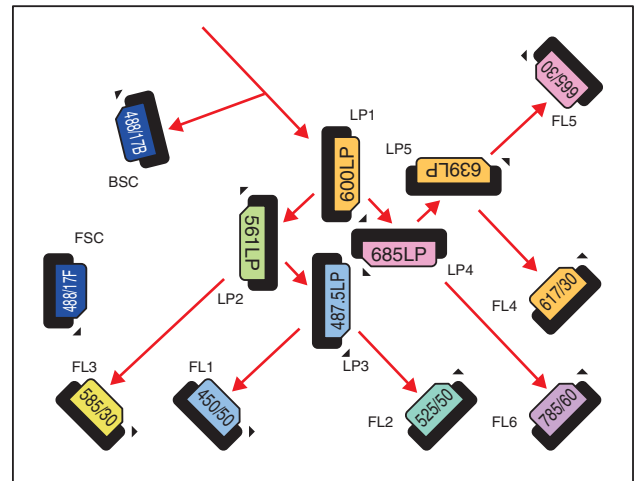

Optical filter label	Installed optical filter
LP1	600LP
LP2	561LP
LP3	487.5LP
LP4	685LP
LP5	639LP
FL1	450/50
FL2	525/50
FL3	585/30
FL4	617/30
FL5	665/30
FL6	785/60
FSC	488/17F
BSC	488/17B

Fluorescent Protein Optical Filter Pattern (Option)

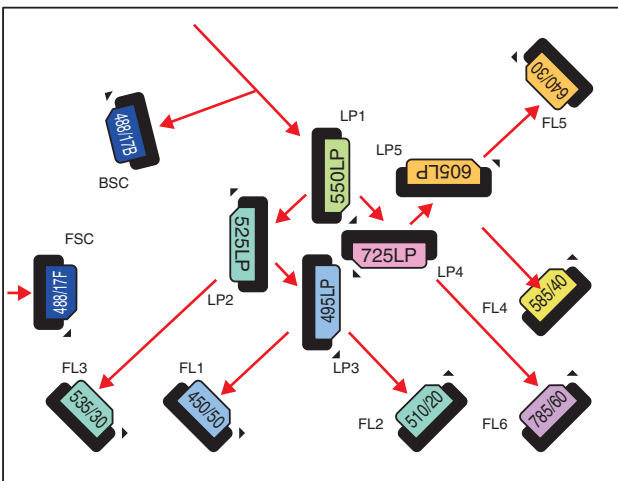

Optical filter label	Installed optical filter
LP1	600LP
LP2	561LP
LP3	487.5LP
LP4	685LP
LP5	639LP
FL1	450/50
FL2	525/50
FL3	585/30
FL4	617/30
FL5	665/30
FL6	785/60
FSC	488/17F
BSC	488/17B

Fluorochrome Detection Matrix

1 laser (488 nm), optical filter pattern 1

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
488 nm laser	EGFP (Enhanced GFP)	508	FL1 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL1 (525/50)
	FITC	518	FL1 (525/50)
	Alexa Fluor 488	519	FL1 (525/50)
	EYFP (Enhanced YFP)	527	FL1 (525/50)
	mCitrine	529	FL1 (525/50)
	PE (R-Phycoerythrin)	576	FL2 (585/30)
	PE-Texas Red	615	FL3 (617/30)
	PI (Propidium Iodide)	617	FL3 (617/30)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
	PE-Cy7	779	FL6 (785/60)

2 lasers (488 and 638 nm lasers), optical filter pattern 1

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
488 nm laser	EGFP (Enhanced GFP)	508	FL1 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL1 (525/50)
	FITC	518	FL1 (525/50)
	Alexa Fluor 488	519	FL1 (525/50)
	EYFP (Enhanced YFP)	527	FL1 (525/50)
	mCitrine	529	FL1 (525/50)
	PE (R-Phycoerythrin)	576	FL2 (585/30)
	PE-Texas Red	615	FL3 (617/30)
	PI (Propidium Iodide)	617	FL3 (617/30)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
	PE-Cy7	779	FL6 (785/60)
638 nm laser	APC (Allophycocyanin)	660	FL4 (665/30)
	Cy5	670	FL4 (665/30)
	Alexa Fluor 647	668	FL4 (665/30)
	APC-Cy5.5	695	FL5 (720/60)
	Alexa Fluor 700	719	FL5 (720/60)
	APC-Cy7	779	FL6 (785/60)
	APC-Alexa Fluor 750	775	FL6 (785/60)

2 lasers (405 and 488 nm lasers), optical filter pattern 2

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
405 nm laser	Brilliant Violet 421	421	FL1 (450/50)
	Alexa Fluor 405	421	FL1 (450/50)
	DAPI	455	FL1 (450/50)
	Pacific Blue	455	FL1 (450/50)
	mCFP	475	FL1 (450/50)
	Hoechst 33342	483	FL1 (450/50)
	AmCyan1	491	FL2 (525/50)
	T-Sapphire	511	FL2 (525/50)
	Qdot 525	525	FL2 (525/50)
	Qdot 545	545	FL2 (525/50)
	Pacific Orange	551	FL2 (525/50)
	Brilliant Violet 570	570	FL3 (600/60)
	Qdot 585	588	FL3 (600/60)
	Qdot 605	603	FL3 (600/60)
	Qdot 655	654	FL4 (665/30)
	Qdot 705	705	FL5 (720/60)
	Qdot 800	800	FL6 (785/60)
488 nm laser	EGFP (Enhanced GFP)	508	FL2 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL2 (525/50)
	FITC	518	FL2 (525/50)
	Alexa Fluor 488	519	FL2 (525/50)
	EYFP (Enhanced YFP)	527	FL2 (525/50)
	mCitrine	529	FL2 (525/50)
	PE (R-Phycoerythrin)	576	FL3 (600/60)
	PE-Texas Red	615	FL3 (600/60)
	PI (Propidium Iodide)	617	FL3 (600/60)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
PE-Cy7	779	FL6 (785/60)	

2 lasers (488 and 561 nm lasers), optical filter pattern 1

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
488 nm laser	EGFP (Enhanced GFP)	508	FL1 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL1 (525/50)
	FITC	518	FL1 (525/50)
	Alexa Fluor 488	519	FL1 (525/50)
	EYFP (Enhanced YFP)	527	FL1 (525/50)
	mCitrine	529	FL1 (525/50)
	PE (R-Phycoerythrin)	576	FL2 (585/30)
	PE-Texas Red	615	FL3 (617/30)
	PI (Propidium Iodide)	617	FL3 (617/30)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
	PE-Cy7	779	FL6 (785/60)
561 nm laser	mOrange	562	FL2 (583/30)
	DsRed-Monomer	586	FL2 (583/30)
	tdTomato	581	FL2 (583/30)
	mCherry	610	FL3 (617/30)
	mPlum	649	FL4 (665/30)

3 lasers (405, 488, and 638 nm lasers), optical filter pattern 2

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
405 nm laser	Brilliant Violet 421	421	FL1 (450/50)
	Alexa Fluor 405	421	FL1 (450/50)
	DAPI	455	FL1 (450/50)
	Pacific Blue	455	FL1 (450/50)
	mCFP	475	FL1 (450/50)
	Hoechst 33342	483	FL1 (450/50)
	AmCyan1	491	FL2 (525/50)
	T-Sapphire	511	FL2 (525/50)
	Qdot 525	525	FL2 (525/50)
	Qdot 545	545	FL2 (525/50)
	Pacific Orange	551	FL2 (525/50)
	Brilliant Violet 570	570	FL3 (600/60)
	Qdot 585	588	FL3 (600/60)
	Qdot 605	603	FL3 (600/60)
	Qdot 655	654	FL4 (665/30)
	Qdot 705	705	FL5 (720/60)
	Qdot 800	800	FL6 (785/60)
488 nm laser	EGFP (Enhanced GFP)	508	FL2 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL2 (525/50)
	FITC	518	FL2 (525/50)
	Alexa Fluor 488	519	FL2 (525/50)
	EYFP (Enhanced YFP)	527	FL2 (525/50)
	mCitrine	529	FL2 (525/50)
	PE (R-Phycoerythrin)	576	FL3 (600/60)
	PE-Texas Red	615	FL3 (600/60)
	PI (Propidium Iodide)	617	FL3 (600/60)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
PE-Cy7	779	FL6 (785/60)	
638 nm laser	APC (Allophycocyanin)	660	FL4 (665/30)
	Cy5	670	FL4 (665/30)
	Alexa Fluor 647	668	FL4 (665/30)
	APC-Cy5.5	695	FL5 (720/60)
	Alexa Fluor 700	719	FL5 (720/60)
	APC-Cy7	779	FL6 (785/60)
	APC-Alexa Fluor 750	775	FL6 (785/60)

3 lasers (405, 488, and 561 nm lasers), optical filter pattern 2

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
405 nm laser	Brilliant Violet 421	421	FL1 (450/50)
	Alexa Fluor 405	421	FL1 (450/50)
	DAPI	455	FL1 (450/50)
	Pacific Blue	455	FL1 (450/50)
	mCFP	475	FL1 (450/50)
	Hoechst 33342	483	FL1 (450/50)
	AmCyan1	491	FL2 (525/50)
	Brilliant Violet 510	510	FL2 (525/50)
	T-Sapphire	511	FL2 (525/50)
	Qdot 525	525	FL2 (525/50)
	Qdot 545	545	FL2 (525/50)
	Pacific Orange	551	FL2 (525/50)
	Brilliant Violet 570	570	FL3 (600/60)
	Qdot 585	588	FL3 (600/60)
	Qdot 605	603	FL3 (600/60)
	Brilliant Violet 605	605	FL3 (600/60)
	Brilliant Violet 650	650	FL4 (665/30)
	Qdot 655	654	FL4 (665/30)
	Qdot 705	705	FL5 (720/60)
	Brilliant Violet 711	711	FL5 (720/60)
Brilliant Violet 785	785	FL6 (785/60)	
Qdot 800	800	FL6 (785/60)	
488 nm laser	EGFP (Enhanced GFP)	508	FL2 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL2 (525/50)
	FITC	518	FL2 (525/50)
	Alexa Fluor 488	519	FL2 (525/50)
	SYBR Green I	522	FL2 (525/50)
	EYFP (Enhanced YFP)	527	FL2 (525/50)
	Vinus	528	FL2 (525/50)
	mCitrine	529	FL2 (525/50)
	PE (R-Phycoerythrin)	576	FL3 (600/60)
	PE-Texas Red	615	FL3 (600/60)
	PI (Propidium Iodide)	617	FL3 (600/60)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
	PE-Cy7	779	FL6 (785/60)
561 nm laser	mOrange	562	FL3 (600/60)
	DsRed-Monomer	586	FL3 (600/60)
	tdTomato	581	FL3 (600/60)
	mCherry	610	FL3 (600/60)
	mPlum	649	FL4 (655/30)

3 lasers (488, 561, and 638 nm lasers), optical filter pattern 1

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
488 nm laser	EGFP (Enhanced GFP)	508	FL1 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL1 (525/50)
	FITC	518	FL1 (525/50)
	Alexa Fluor 488	519	FL1 (525/50)
	SYBR Green I	522	FL1 (525/50)
	EYFP (Enhanced YFP)	527	FL1 (525/50)
	Venus	528	FL1 (525/50)
	mCitrine	529	FL1 (525/50)
	PE (R-Phycoerythrin)	576	FL2 (585/30)
	PE-Texas Red	615	FL3 (617/30)
	PI (Propidium Iodide)	617	FL3 (617/30)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (655/30)
	PE-Cy5	670	FL4 (655/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (655/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
	PE-Cy7	779	FL6 (785/60)
561 nm laser	mOrange	562	FL2 (585/30)
	DsRed-Monomer	586	FL2 (585/30)
	tdTomato	581	FL2 (585/30)
	mCherry	610	FL3 (617/30)
	mPlum	649	FL4 (655/30)
638 nm laser	APC (Allophycocyanin)	660	FL4 (655/30)
	Cy5	670	FL4 (655/30)
	Alexa Fluor 647	668	FL4 (655/30)
	APC-Cy5.5	695	FL5 (720/60)
	Alexa Fluor 700	719	FL5 (720/60)
	APC-Cy7	779	FL6 (785/60)
	APC-Alexa Fluor 750	775	FL6 (785/60)

4 lasers (405, 488, 561, and 638 nm lasers), optical filter pattern 2

Excitation laser	Fluorochrome		Channel (optical filter)
	Name	Peak emission wavelength (nm)	
405 nm laser	Brilliant Violet 421	421	FL1 (450/50)
	Alexa Fluor 405	421	FL1 (450/50)
	DAPI	455	FL1 (450/50)
	Pacific Blue	455	FL1 (450/50)
	mCFP	475	FL1 (450/50)
	Hoechst 33342	483	FL1 (450/50)
	AmCyan1	491	FL2 (525/50)
	T-Sapphire	511	FL2 (525/50)
	Qdot 525	525	FL2 (525/50)
	Qdot 545	545	FL2 (525/50)
	Pacific Orange	551	FL2 (525/50)
	Brilliant Violet 570	570	FL3 (600/60)
	Qdot 585	588	FL3 (600/60)
	Qdot 605	603	FL3 (600/60)
	Qdot 655	654	FL4 (665/30)
	Qdot 705	705	FL5 (720/60)
	Qdot 800	800	FL6 (785/60)
488 nm laser	EGFP (Enhanced GFP)	508	FL2 (525/50)
	CFSE (Carboxyfluorescein Succinimidyl ester)	517	FL2 (525/50)
	FITC	518	FL2 (525/50)
	Alexa Fluor 488	519	FL2 (525/50)
	EYFP (Enhanced YFP)	527	FL2 (525/50)
	mCitrine	529	FL2 (525/50)
	PE (R-Phycoerythrin)	576	FL3 (600/60)
	PE-Texas Red	615	FL3 (600/60)
	PI (Propidium Iodide)	617	FL3 (600/60)
	7-AAD (7-Aminoactinomycin D)	647	FL4 (665/30)
	PE-Cy5	670	FL4 (665/30)
	PerCP (Peridinin chlorophyll protein)	675	FL4 (665/30)
	PE-Cy5.5	695	FL5 (720/60)
	PerCP-Cy5.5	695	FL5 (720/60)
	PerCP-eFluor 710	710	FL5 (720/60)
	PE-Cy7	779	FL6 (785/60)
561 nm laser	mOrange	562	FL3 (600/60)
	DsRed-Monomer	586	FL3 (600/60)
	tdTomato	581	FL3 (600/60)
	mCherry	610	FL3 (600/60)
	mPlum	649	FL4 (665/30)
638 nm laser	APC (Allophycocyanin)	660	FL4 (665/30)
	Cy5	666	FL4 (665/30)
	Alexa Fluor 647	668	FL4 (665/30)
	APC-Cy5.5	695	FL5 (720/60)
	Alexa Fluor 700	719	FL5 (720/60)
	APC-Cy7	779	FL6 (785/60)
	APC-Alexa Fluor 750	775	FL6 (785/60)

Relationship between Sample Pressure and Sample Flow/Event Rate

100 µm sorting chip

[Sample Pressure]		1	2	3	4	5	6	7	8	9	10
Flow rate (µl/min)		6	11	16	21	27	37	47	58	68	89
Event rate (eps)	Sample concentration (1e6/ml)	60	160	260	350	450	650	840	1,000	1,200	1,600
	Sample concentration (1e7/ml)	600	1,300	2,100	2,800	3,600	5,100	6,600	8,100	9,600	12,600

130 µm sorting chip

[Sample Pressure]		1	2	3	4	5	6	7	8	9	10
Flow rate (µl/min)		1	3	7	10	16	22	28	34	40	50
Event rate (eps)	Sample concentration (1e6/ml)	10	30	100	150	250	400	500	600	700	900
	Sample concentration (1e7/ml)	100	600	1,200	2,000	2,800	3,600	4,500	5,400	6,500	8,000

70 µm sorting chip

[Sample Pressure]		1	2	3	4	5	6	7	8	9	10
Flow rate (µl/min)		8	18	28	42	62	80	100	120	140	168
Event rate (eps)	Sample concentration (1e7/ml)	1000	2200	3500	5300	7700	10000	12500	15000	17500	21000

The values above are reference values. Values may vary depending on the environment.

Troubleshooting

This section lists common problems that may occur and the actions to take to resolve each problem. Check these items before calling for service. If the problem still persists after carrying out the recommended actions, contact your Sony distributor.

Problem	Possible Cause/Recommended Solution
The event rate does not rise during data acquisition, sorting, or adjustment.	The sample line may be clogged. Click [Probe Wash] in the [Cleaning] group on the [Cytometer] tab of the ribbon to clean the sample line. If the problem persists, click [DI Rinse] in the [Cleaning] group on the [Cytometer] tab of the ribbon to flush the clog. If the problem persists, change the sample line (<i>page 192</i>). After DI Rinse or changing the sample line run automatic calibration again.
Measurement values cannot be acquired, or sorting performance cannot be maintained.	The sorting chip may be clogged or contain trapped air. Click [Chip] in the [De-bubble] group on the [Cytometer] tab of the ribbon several times as required to clear the chip. If the problem persists, click [DI Rinse] in the [Cleaning] group on the [Cytometer] tab of the ribbon to flush the clog or trapped air. If the problem persists, replace the sorting chip. After DI Rinse or replace sorting chip run automatic calibration again. If the problem persists, the optical module lens may require cleaning. Contact your Sony distributor for service.
Droplet image in droplet viewer shakes up and down, or flows in a curved path.	The sorting chip may be clogged or contain trapped air. Click [Chip] in the [De-bubble] group on the [Cytometer] tab of the ribbon several times as required to clear the chip. If the problem persists, click [DI Rinse] in the [Cleaning] group on the [Cytometer] tab of the ribbon to flush the clog or trapped air. The fluid consistency of the sample may be causing fluctuations in the sample flow. Click [Agitate] on the [Cytometer] tab of the ribbon to turn on sample agitation. Check that there is no air leaking from the sheath tank. If air is leaking, reattach the lid until it is correctly sealed. Check that there are no air bubbles in the sheath filter. If the problem persists, there may be air bubbles trapped in the sheath filter. Remove the air bubbles from the sheath filter. See <i>"Releasing Air in the Sheath Filter and Sorting Chip"</i> (<i>page 202</i>). If the problem persists, there may be salt deposits in the flow channels. Log off and restart in maintenance mode (<i>"Running Maintenance Mode"</i> (<i>page 177</i>)). The sorting chip is ejected automatically. Check that the surface of the chip is not dirty or smudged, then perform chip realignment. If the sorting chip is dirty or smudged, replace the chip. If the problem persists, contact your Sony distributor for service.
Shape of droplets is irregular.	The sorting chip may contain trapped air. Check that there is no air trapped in the sheath filter. Click [Chip] in the [De-bubble] group on the [Cytometer] tab of the ribbon several times as required to clear the chip. Check that there are no air bubbles in the sheath filter. If the problem persists, there may be air bubbles trapped in the sheath filter. Remove the air bubbles from the sheath filter. See <i>"Releasing Air in the Sheath Filter and Sorting Chip"</i> (<i>page 202</i>). If the problem persists, there may be salt deposits in the flow channels. Log off and restart in maintenance mode (<i>"Running Maintenance Mode"</i> (<i>page 177</i>)). The sorting chip is ejected automatically. Check that the surface of the chip is not dirty or smudged, then perform chip realignment. If the sorting chip is dirty or smudged, replace the chip. The sorting parameters may be set incorrectly. The sorting parameters affect the droplet shape, breakoff point, and electric charge. Increase the amplitude or decrease the frequency to make droplets larger. Decrease the amplitude or increase the frequency to make droplets smaller.
Sheath fluid is not flowing.	Check that the connectors for all fluidics lines in the fluidics cart and main unit are connected securely and that the lid of the sheath tank is closed tightly. Check that the sheath tank lid has a black rubber seal attached. Check that there are no air bubbles in the sheath filter. Check the pressure of the compressed air supply to the main unit.
Electrical discharge between deflection plates causes sorting errors.	Discharge may occur due to fluid droplets adhering to the surface of the plates. Turn off the power, open the collection area door, and wipe the deflection plates dry. If residue remains on the plates, remove and clean the plates.
The plot position does not change after increasing the gain.	Increasing the gain too much can cause output saturation. In this case, reduce the gain and readjust settings to find the optimum operating point.

Problem	Possible Cause/Recommended Solution
Sorting chip does not load correctly.	Log off and restart in maintenance mode (<i>“Running Maintenance Mode” (page 177)</i>). If the problem persists, replace the sorting chip. If the problem persists, contact your Sony distributor for service.
Chip does not eject.	Use the manual load/unload thumbwheel on the chip loader to remove the chip.
Sorting chip clogs after changing the sheath filter.	There may be some small particles in the replacement filter that have not been removed. Run the sheath filter de-bubble function at least three times.
Sample tube hits the sample probe.	Use only standard tubes and fully insert the tube in the sample tube holder when loading a sample. An ill-fitting or non-standard tube can damage the sample probe.
Fluid is leaking from the sample probe.	Check that the sample line is fully inserted in the pinch valve.
Fluid is leaking from the top in the collection area.	The sample fluid connector may not be securely fastened. Securely fasten the sample line connector.
Fluidics component (pump, fluid connectors, etc.) is leaking fluid.	The seals may have deteriorated due to age, causing ethanol or waste fluid to leak. Clean the area, and contact your Sony distributor for service. Note Always wear gloves and other protective clothing, mask, and goggles, as required, when cleaning spills and leakages.
Droplet camera image is erratic.	The droplet camera may require cleaning. Contact your Sony distributor for service.
Droplets are not being formed or the droplet camera image is not displayed.	The droplet camera may be defective. Contact your Sony distributor for service.
Data is not acquired in a given fluorescence detection channel or condition less than favorable.	The optical filters may not be inserted in the correct filter pattern or may require cleaning. Check that the filters are clean and inserted in the correct slots.
Gain settings appear to have deviated during measurement.	The optical filters may have been subjected to static electricity discharge. Pause or stop data acquisition, and then resume data acquisition to restore normal gain settings.
No output from the lasers (laser temperature error).	If the ambient operating temperature exceeds the maximum operating temperature of 35 °C (95 °F), the lasers shutdown automatically and an error message is displayed. To clear the error, shut down the main unit and wait for the ambient temperature to cool sufficiently before restarting the main unit. If the error persists even after allowing the ambient temperature to cool below 35 °C (95 °F), contact your Sony distributor for service. If the ambient temperature drops below the minimum operating temperature of 10 °C (50 °F), the lasers shut down automatically and an error message is displayed. To clear the error, shut down the main unit and wait for the ambient temperature to rise above 10 °C (50 °F) before restarting the main unit. If the error persists even after the temperature has risen above 10 °C (50 °F), contact your Sony distributor for service.
Waste fluid is not being collected in the waste tank.	Waste fluid may be leaking in the fluidics cart due to an unsecured waste line connector on the waste tank. Make sure the waste line connector is securely connected on the waste tank. Replace the waste tank air filter if more than one month has elapsed since the air filter was last replaced. If the problem persists, contact your Sony distributor for service.
Communication error between the SH800 and the host computer.	Shutdown and restart both the SH800 main unit and the host computer.
POWER/STANDBY button indicator is flashing.	If the doors can be opened and closed, remove the sample tube and collection tubes before performing the following procedure. 1 Press and hold the POWER/STANDBY button until the indicator stops flashing. 2 Press the POWER/STANDBY button again to start the main unit. If the POWER/STANDBY button indicator does not flash, continue normal operation. 3 If the POWER/STANDBY button indicator starts flashing again, press and hold the POWER/STANDBY button until the indicator stops flashing, then turn off the main power switch on the rear panel. 4 Contact your Sony distributor for service.

Problem	Possible Cause/Recommended Solution
An error occurs in automatic calibration or automatic calibration stops and does not proceed.	<ul style="list-style-type: none"> • There may be air bubbles trapped in the sheath filter. Remove the air bubbles from the sheath filter. See <i>“Releasing Air in the Sheath Filter and Sorting Chip” (page 202)</i>. • If an error occurs while adjusting the side streams, the sensor windows may be dirty, preventing normal side stream detection. Clean the windows using a soft, lint-free cloth lightly moistened with water. Clean the black sensor windows on the left, right, and back. • The temperature within the unit may have changed. With the main unit up and running, leave standing until the room temperature stabilizes, then run automatic calibration again. • If the problem persists, there may be salt deposits in the flow channels. Log off and run the unit in maintenance mode (<i>“Running Maintenance Mode” (page 177)</i>). • If the problem persists, change the sorting chip. • If the problem persists, contact your Sony distributor for service.
Side streams are not visible during automatic calibration.	The transparent protective covers in the collection area may be charged with static electricity. Clean the covers using a soft, lint-free cloth moistened with water. For details, see <i>“General Cleaning” (page 172)</i> .
The output pressure of the air compressor drops.	Connect the power cord for the air compressor directly to a power outlet. Do not connect using an extension cord.
The waste tank is bulging.	The waste tank air filter may be blocked. Replace the waste tank air filter. See <i>“Changing Tank Air Filters” (page 186)</i> .
Water splashes occur when cleaning the sample probe.	De-bubble the DI water filter. For details, see <i>“Changing the DI Water Filter” (page 198)</i> .
The instrument will not operate after adding a printer in [Control Panel] on the host computer.	Adding a printer in [Control Panel] while the SH800 main unit and host computer are connected may cause the SH800 main unit to stop functioning. Press and hold the POWER/STANDBY button on the front panel to turn off the power, then turn the power on again and restart Cell Sorter Software. To add a printer, first turn off the SH800 power supply or disconnect the network cable before adding the printer.
The main window in Cell Sorter Software does not update (e.g. the event rate does not update).	If the main window display does not update, click the SH800 icon in the task bar to minimize the main window. Then, click the SH800 icon again to restore the window. If pull-down menus still do not open in the main window, open the [Detector & Threshold Settings] dialog (<i>page 86</i>) and operate a [Threshold] pull-down menu, for example.
Button labels displayed for dialogs on the main window are transparent.	If the dialog display is corrupted, press the ESC key to close the dialog (clicking the [X] icon in the top right corner of the dialog has no affect). Then, reopen the dialog from the main window. If the problem persists, click the SH800 icon in the task bar to minimize the main window and then click it again to restore the window.

Error Messages

This section lists the error messages that may be displayed in Cell Sorter Software. When an error occurs, check these items for the actions to take to resolve each error. If the problem still persists after carrying out the recommended actions, make a note of the error message and then contact your Sony distributor.

Error message	Solution
Unload Chip Error. Discard and load a new chip. If the error reoccurs, please contact your service representative. (Error ID : 00 - 100)	Shut down the main unit and turn off the compressed air supply. Pull the sample line out from the chip connector, and remove the sorting chip manually.
Load Chip Error. Error on loading mechanism. Please reinsert the chip. If the error reoccurs, please contact your service representative. (Error ID : 00 - 101)	Insert a new sorting chip and try to load the chip again. If the problem persists, contact your Sony distributor for service.
Host Communication Error. 1 Shutdown the cytometer. 2 Make sure the cable is connected correctly and firmly and restart the cytometer and software. If the error reoccurs, please contact your service representative. (Error ID : 00 - 102)	Check the connection between the main unit and the host computer and restart Cell Sorter Software. If the problem persists, contact your Sony distributor for service.

Error message	Solution
<p>Unload Sample tube Error. If the error reoccurs, please contact your service representative. (Error ID : 00 - 104)</p>	<p>Use the following procedure.</p> <p>1 Check that the compressed air supply is supplying rated pressure.</p>
<p>Unload Sample tube Error. Failed to reduce the sample pressure. If the error reoccurs, please contact your service representative. (Error ID : 00 - 509)</p>	<p>2 Shut down the main unit, and turn off the compressed air supply.</p> <p>3 Clean or replace the sample loader O-ring.</p> <p>If the problem persists, contact your Sony distributor for service.</p>
<p>Unload Sample tube Error. Failed to down the sample loader. If the error reoccurs, please contact your service representative. (Error ID : 00 - 506)</p>	
<p>Unload Sample tube Error. Failed to move the support guide. If the error reoccurs, please contact your service representative. (Error ID : 00 - 507)</p>	
<p>Unload Sample tube Error. Failed to move the sample stage. If the error reoccurs, please contact your service representative. (Error ID : 00 - 508)</p>	
<p>Load Sample tube Error. Please reload the tube. If the error reoccurs, please contact your service representative. (Error ID : 00 - 105)</p>	
<p>Load Sample tube Error. Failed to move the waste tray. Please reload the tube. If the error reoccurs, please contact your service representative. (Error ID : 00 - 501)</p>	
<p>Load Sample tube Error. Failed to lift up the sample loader. Please reload the tube. If the error reoccurs, please contact your service representative. (Error ID : 00 - 502)</p>	
<p>Load Sample tube Error. Failed to move the support guide. Please reload the tube. If the error reoccurs, please contact your service representative. (Error ID : 00 - 503)</p>	<p>Use the following procedure.</p> <p>1 Check that the compressed air supply is supplying rated pressure.</p> <p>2 Shut down the main unit, and turn off the compressed air supply.</p> <p>3 Clean or replace the sample loader O-ring.</p> <p>If the problem persists, contact your Sony distributor for service.</p>
<p>Load Sample tube Error. Failed to move the sample stage. Please reload the tube. If the error reoccurs, please contact your service representative. (Error ID : 00 - 504)</p>	
<p>Sheath Fluid Tank is empty. Refill the sheath fluid tank following the instructions in the operator's guide. (Error ID : 00 - 106)</p>	<p>The sheath tank is empty. Refill the sheath tank with sheath fluid. <i>See "Refilling the Sheath Tank" (page 164).</i></p>
<p>Waste Fluid Tank is full. Empty the waste tank following the instructions in the operator's guide. (Error ID : 00 - 107)</p>	<p>Empty the waste tank. <i>See "Emptying/Changing the Waste Tank" (page 166).</i></p>
<p>Compressed air is not supplied. Confirm if the compressed air hose is connected or the sheath tank is set up properly. (Error ID : 00 - 110)</p>	<p>Check that the compressed air supply is supplying rated pressure.</p>

Error message	Solution
Fluidics cart is not connected. Connect the fluidics cart correctly following the instructions in the operator's guide. (Error ID : 00 - 111)	Connect or reconnect the cart connection cable between the main unit and the fluidics cart.
The ambient temperature exceeds the recommended operating range. Shut down the cytometer. (Error ID : 00 - 112)	The main unit may be overheating. Shut down the main unit, and allow the temperature to fall within the rated range before restarting. If the problem persists at temperatures not exceeding 35 °C (95 °F), contact your Sony distributor for service.
The temperature of the sample area is too high. (Error ID : 00 - 113)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
The temperature of the collection area is too high. (Error ID : 00 - 114)	
Leakage is suspected in the drip tray. (Error ID : 00 - 115)	Open the fluidics maintenance door and check for leakages. When doing so, do not touch anything other than the drip tray. If the problem persists, contact your Sony distributor for service.
The sheath tank on the fluidics cart is not detected. Check that the sheath tank is attached. (Error ID : 00 - 116)	Check that the sheath tank is positioned correctly in the fluidics cart. If the problem persists, contact your Sony distributor for service.
The waste tank on the fluidics cart is not detected. Check that the waste tank is attached. (Error ID : 00 - 117)	Check that the waste tank is empty and is positioned correctly in the fluidics cart. If the problem persists, contact your Sony distributor for service.
The temperature control of the sample area failed. Shut down the cytometer. (Error ID : 00 - 118)	Check that there is no object shielding the base of the main unit. If the problem persists, contact your Sony distributor for service.
The temperature control of the collection area failed. Shut down the cytometer. (Error ID : 00 - 119)	
The laser initialization timed out. (Error ID : 00 - 120)	The main unit may be overheating. Shut down the main unit, and allow the temperature to fall within the rated range before restarting. If the problem persists at temperatures not exceeding 35 °C (95 °F), contact your Sony distributor for service.

Error message	Solution
System power supply error on the PCT board. (Error ID : 00 - 121)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
AC/DC power supply error on the PCT board. (Error ID : 00 - 122)	
250 V power supply error on the HVC board. (Error ID : 00 - 123)	
+5 V power supply error on the HVC board. (Error ID : 00 - 124)	
-5 V power supply error on the HVC board. (Error ID : 00 - 125)	
-12 V power supply error on the HVC board. (Error ID : 00 - 126)	
Piezo short error on the HVC board. (Error ID : 00 - 127)	
12 V power supply error on the MD1 board. (Error ID : 00 - 128)	
24 V power supply error on the MD1 board. (Error ID : 00 - 129)	
5 V power supply error on the MD1 board. (Error ID : 00 - 130)	
Temperature error from the X axis in the chip stage motor on the MD2 board. (Error ID : 00 - 131)	
Temperature error from the Z axis in the chip stage motor on the MD2 board. (Error ID : 00 - 132)	
Temperature error from the chip load motor on the MD2 board. (Error ID : 00 - 133)	
Temperature error from the collection load motor on the MD2 board. (Error ID : 00 - 134)	
Temperature error from the LDS motor on the MD2 board. (Error ID : 00 - 135)	
12 V power supply error on the MD2 board. (Error ID : 00 - 136)	
24 V power supply error on the MD2 board. (Error ID : 00 - 137)	
5 V power supply error on the MD2 board. (Error ID : 00 - 138)	
12 V power supply error on the MD3 board. (Error ID : 00 - 139)	
24 V power supply error on the MD3 board. (Error ID : 00 - 140)	
5 V power supply error on the MD3 board. (Error ID : 00 - 141)	
Temperature error from the agitation motor on the MD3 board. (Error ID : 00 - 142)	
Pump error on the MD1 board. (Error ID : 00 - 143)	
Temperature error from the droplet camera motor on the MD2 board. (Error ID : 00 - 144)	The main unit may be overheating. Shut down the main unit, and allow the temperature to fall within the rated range before restarting. If the problem persists at temperatures not exceeding 35 °C (95 °F), contact your Sony distributor for service.

Error message	Solution
Power supply error on the MEC board. (Error ID : 00 - 146)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Error from FAN1 on the MD1 board. (Error ID : 00 - 148)	
Error from FAN2 on the MD1 board. (Error ID : 00 - 149)	
Error from FAN3 on the MD1 board. (Error ID : 00 - 150)	
Error from FAN on the MD3 board. (Error ID : 00 - 151)	
Circuit error on the LD1 board. (Error ID : 00 - 152)	
Circuit error on the LD2 board. (Error ID : 00 - 153)	
Circuit error on the LD3 board. (Error ID : 00 - 154)	
Temperature control of the LD1 failed. (Error ID : 00 - 155)	
Temperature control of the LD2 failed. (Error ID : 00 - 156)	
Temperature control of the LD3 failed. (Error ID : 00 - 157)	
Temperature control of the prism failed. (Error ID : 00 - 158)	
Power supply error for LD. (Error ID : 00 - 159)	
Power supply error for peltier. (Error ID : 00 - 160)	
Thermistor error for LD1. (Error ID : 00 - 161)	Shut down the main unit, then try again. If the problem persists, contact your Sony distributor for service.
Thermistor error for LD2. (Error ID : 00 - 162)	
Thermistor error for LD3. (Error ID : 00 - 163)	
Thermistor error for prism. (Error ID : 00 - 164)	
Temperature error of the 561 nm laser. (Error ID : 00 - 165)	The main unit may be overheating. Shut down the main unit, and allow the temperature to fall within the rated range before restarting. If the problem persists at temperatures not exceeding 35 °C (95 °F), contact your Sony distributor for service.
Error on the eeprom. (Error ID : 00 - 166)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
The ambient temperature is too low. (Error ID : 00 - 167)	The main unit may be too cold. Shut down the main unit, and allow the temperature to rise within the rated range before restarting. If the problem persists at temperatures exceeding 10 °C (50 °F), contact your Sony distributor for service.

Error message	Solution
Communication error on the PCT board. (Error ID : 00 - 168)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Communication error on the MD1 board. (Error ID : 00 - 169)	
Communication error on the MD2 board. (Error ID : 00 - 170)	
Communication error on the MD3 board. (Error ID : 00 - 171)	
Communication error on the CLM board. (Error ID : 00 - 172)	
Communication error on the HVC board. (Error ID : 00 - 173)	
No Reply error on the PCT board. (Error ID : 00 - 174)	
No Reply error on the MD1 board. (Error ID : 00 - 175)	
No Reply error on the MD2 board. (Error ID : 00 - 176)	
No Reply error on the MD3 board. (Error ID : 00 - 177)	
No Reply error on the CLM board. (Error ID : 00 - 178)	
No Reply error on the HVC board. (Error ID : 00 - 179)	
Load Chip Error. Wrong side out. (Error ID : 00 - 180)	Insert a sorting chip in the correct orientation, and try to load the chip again. If the problem persists, contact your Sony distributor for service.
Load Chip Error. Up side down. (Error ID : 00 - 181)	
Load Chip Error. Failed to hold the chip. Please reinsert the chip. (Error ID : 00 - 500)	
Load Chip Error. Failed to detect the chip. Please reinsert the chip. (Error ID : 00 - 182)	
Leakage is suspected in the collection area. (Error ID : 00 - 183)	Search the collection area and check to make sure there is no fluid leakage. If there is fluid leakage, wipe clean and wait a while before restarting. If the problem persists, contact your Sony distributor for service.
Leakage is suspected in the drip tray. Check if there is leakage in the drip tray and clean it off. If the detection will be continued, the cytometer will be disabled in two minutes. (Warning ID : 00 - 245)	Inspect the drip tray and check for any fluid leakages. When doing so, do not touch anything other than the drip tray. If the problem persists, contact your Sony distributor for service.
Leakage is suspected in the collection area. Check if there is leakage in the collection area and clean it off. If the detection will be continued, the cytometer will be disabled in two minutes. (Warning ID : 00 - 246)	Search the collection area and check to make sure there is no fluid leakage. If there is fluid leakage, wipe clean and wait a while before restarting. If the problem persists, contact your Sony distributor for service.
Failed to start feeding sheath fluid. (Error ID: 00-184) Click [Troubleshooting] to begin check the line connection and compressed air supply.	Click [Troubleshooting] to launch the troubleshooting wizard, and follow the on-screen instructions. <i>See "If fluid flow does not start" (page 48).</i>
Failed to fit the chip. (Error ID : 00 - 185)	Insert a sorting chip in the correct orientation, and try to load the chip again. If the problem persists, contact your Sony distributor for service.

Error message	Solution
Failed to release the chip. (Error ID : 00 - 186)	Shut down the main unit and turn off the compressed air supply. Pull the sample line out from the chip connector, and remove the sorting chip manually.
Unload Chip Error. Failed to hold the chip. (Error ID : 00 - 187)	
Unload Chip Error. Failed to release the chip. (Error ID : 00 - 188)	
The model number is invalid. (Error ID : 00 - 189)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
The laser configuration is invalid. (Error ID : 00 - 190)	
The ADC configuration is invalid. (Error ID : 00 - 191)	
Power supply error on the ADC board. (Error ID : 00 - 192)	
Power supply error on the ADC2 board. (Error ID : 00 - 193)	
Timed out to close the sample door. (Error ID : 00 - 194)	
Timed out to control sheath pressure. (Error ID : 00 - 195)	
Flash write error. (Error ID : 00196)	
Temperature error from the collection load motor on the MD2 board. (Error ID : 00 - 197)	
Timed out to control sheath pressure. (Error ID : 00 - 198)	
Timed out to control sample pressure. (Error ID : 00 - 199)	
Collection load motor temperature error on the MD2 board. (Error ID : 00-197)	
Sheath pressure control error. (Error ID : 00-198)	
Sample pressure control error. (Error ID : 00-199)	
Sheath Fluid Tank is almost empty. Refill the sheath fluid tank following the instructions in the operator's guide. (Warning ID : 00 - 200)	
Waste Fluid Tank is almost full. Empty the waste tank following the instructions in the operator's guide. (Warning ID : 00 - 201)	Empty the waste tank. <i>See "Emptying/Changing the Waste Tank" (page 166).</i>
Collection area door is open. (Warning ID : 00 - 202)	Close the collection area door.
Chip loader flip-up door is open. Close the chip loader flip-up door. (Warning ID : 00 - 203)	Close the flip-up door.
Warning for the 561 nm laser. Chip loader flip-up door is open. Close the chip loader flip-up door. (Warning ID : 00 - 210)	Laser has reached the end of its service life. The laser output does not drop immediately. The laser is still operational, but should be replaced as soon as possible.
Warning for the unsuccessful communication on MD2 board. (Warning ID : 00 - 214)	This warning does not affect automatic calibration operation, although it may appear during automatic calibration. In this case, continue the automatic calibration process.

Error message	Solution
Warning for the unsuccessful communication on PCT board. (Warning ID : 00 - 212)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Warning for the unsuccessful communication on MD1 board. (Warning ID : 00 - 213)	
Warning for the unsuccessful communication on MD3 board. (Warning ID : 00 - 215)	
Warning for the unsuccessful communication on CLM board. (Warning ID : 00 - 216)	
Warning for the unsuccessful communication on HVC board. (Warning ID : 00 - 217)	
LD1 laser power warning. (Warning ID : 00 - 229)	A reduction in laser output power has been detected. Contact your Sony distributor for service.
LD2 laser power warning (Warning ID : 00 - 230)	
LD3 laser power warning (Warning ID : 00 - 231)	
LD4 laser power warning (Warning ID : 00 - 232)	
LD1 laser current warning. (Warning ID : 00 - 233)	A laser error has been detected. Contact your Sony distributor for service.
LD2 laser current warning. (Warning ID : 00 - 234)	
LD3 laser current warning. (Warning ID : 00 - 235)	
The LD1 laser is approaching its end of life expiration. (Warning ID : 00 - 236)	Laser is approaching the end of its service life. The laser output does not drop immediately. The laser is still operational, but should be replaced as soon as possible.
The LD2 laser is approaching its end of life expiration. (Warning ID : 00 - 237)	
The LD3 laser is approaching its end of life expiration. (Warning ID : 00 - 238)	
The LD4 laser is approaching its end of life expiration. (Warning ID : 00 - 239)	
Failed to record LD1 laser total hours of use. (Warning ID : 00 - 240)	Total irradiation time was not recorded correctly. The laser is still operational, but you should contact your Sony distributor for service.
Failed to record LD2 laser total hours of use. (Warning ID : 00 - 241)	
Failed to record LD3 laser total hours of use. (Warning ID : 00 - 242)	
Failed to record LD4 laser total hours of use. (Warning ID : 00 - 243)	
DI Water Tank is almost empty. Refill the DI water tank following the instructions in the operator's guide. (Warning ID : 00 - 244)	Refill the DI water tank. <i>See "Refilling the DI Water Tank" (page 167)</i>
Waste catcher position warning. (Warning ID : 00-247)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
The temperature control of the collection area has failed and has turned off. Wait a while before turning on it again. (Warning ID : 00-260)	
Motor position has reached to the limit. (Motor limit over) (Warning ID : 00-261)	
Calculated well position has reached to the limit. These position cannot be updated. (Warning ID : 00-262)	
Power supply error on the LCD board. (Warning ID : 00-263)	An error occurred on the LCD board. The laser is still operational, but you should contact your Sony distributor for service.
The cytometer has been disconnected. Acquiring and sorting has stopped.	Check the connections to the main unit.

Error message	Solution
Communication to the cytometer has been lost. Exit and relaunch the software.	Check the connections to the main unit and relaunch Cell Sorter Software.
The initial setup did not complete successfully. Exit and relaunch the software.	Exit Cell Sorter Software and try starting again.
Acquiring and sorting has stopped because available free size of Hard Disk Drive is not enough.	Insufficient free space on the hard disk. Delete unnecessary files to increase the free capacity.
The application is already running. This instance will now exit.	Shut down the Cell Sorter Software instance that is already running.
Failed to save settings.	Try saving the settings again.
Parameter count in a selected file is different from current settings.	Select a file with the same number of parameters as the current settings.
A selected file has the wrong format.	Select a file in the correct format.
Unable to maintain stable droplet breakoff.	An error occurred when the control breakoff function was turned on. If a droplet formation problem is detected during measurement, acquisition stops. Try running [Sort Calibration]. If the problem persists after running [Sort Calibration] again, contact your Sony distributor for service.
The selected drive is not external storage device. Please select an external storage device.	Specify an external storage device.
The drive is not formatted as NTFS. Please use an NTFS formatted drive.	Use only NTFS-formatted drives.
You do not have write permission for the selected folder. Please add write permission to the folder or select another folder with write permission.	You do not have write access privileges for the selected folder. Change the folder access privileges or specify a folder with write access privileges.
The available disk free space is low.	Insufficient free space on the hard disk. Delete unnecessary files to increase the free capacity.
This chip has been used over the duration limit (12 hours) with the 405 nm laser. Please exchange for a new chip.	Replace the sorting chip.
This chip has been used over the duration limit (12 hours) with the 405 nm laser. You cannot use this chip any longer with the 405 nm laser. Please exchange for a new chip.	
System resources are scarce and may cause the application performance to be worse. Would you like to close unnecessary worksheets?	Close any unnecessary worksheets.
A number of plots are opened and may cause the application performance to be worse. Please close unnecessary plots or worksheets.	
System resources are scarce and may cause the application to shutdown. Would you like to start the recovery process?	Stop data acquisition and measurement, and shut down Cell Sorter Software.
System resources are scarce because of multiple applications/processes running in the background. This may impact the performance of the Cell Sorter Software. Please close unnecessary applications/processes.	Close any unnecessary running applications and processes.
Invalid biexponential parameters were selected.	Specify valid biexponential parameters.
Cannot assign this tube because available free space on the HDD is insufficient. Please increase the available free space on the HDD.	Cannot assign due to insufficient free space on HDD. Delete unnecessary files to increase the free capacity.
Detected UNKNOWN state. Please check the Cell Sorter.	An unknown problem has been detected. Check the SH800.
Unable to maintain stable droplet breakoff. The measurement will be stopped and Control Breakoff will be disabled.	Data acquisition stops. Check the instrument settings, as required.
This chip has been used over 9 hours with the 405 nm laser; the limit with a 405 nm laser is 12 hours. Please exchange the chip before the 12 hour limit.	This sorting chip is approaching the maximum recommended use period. Replace the sorting chip.
Cannot receive droplet camera image continuously. If you have some trouble, please stop acquisition.	The droplet camera image cannot be acquired. If a problem occurs, stop data acquisition.
Cannot receive side stream camera image continuously. If you have some trouble, please stop acquisition.	The side stream monitor image cannot be acquired. If a problem occurs, stop data acquisition.

Error message	Solution
<p>Cannot make appropriate droplet for sorting.</p>	<p>It is impossible to maintain the droplet shape determined during auto alignment. De-bubble the sorting chip to remove any trapped air. If the problem persists, run [Sort Calibration] again. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration].</p>
<p>Unable to perform ethanol cleaning with full waste tank. Please empty the waste tank and retry.</p>	<p>Cannot start ethanol cleaning due to insufficient free capacity in the waste tank. Empty the waste tank.</p>
<p>Droplet condition has changed. Please re-perform Sort Calibration for stable sorting. (Warning ID : W1001) (Warning ID : W1011) (Warning ID : W1022)</p>	<p>The droplet shape has changed. Run [Sort Calibration] again to restore stable sorting. Use the unit within the specified ambient temperature range. Use only in environments where the ambient temperature does not change suddenly. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration]. A message indicating that the chip is being used a second time appears, but the chip can be used.</p>
<p>Sample flow is paused due to change in droplet condition. Click [Stop] and re-perform Sort Calibration for stable sorting. (Warning ID : W1002)</p>	<p>The sample flow was paused due to changes in the droplet shape. Click [Stop] and run [Sort Calibration] again to restore stable sorting. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration]. A message indicating that the chip is being used a second time appears, but the chip can be used.</p>
<p>Sample flow is paused due to change in droplet condition. Click [Stop] and re-perform Sort Calibration for stable sorting. To continue sorting without re-calibration, deselect [Detect Satellite Droplets] under [Settings] - [Advanced Settings]. (Warning ID : W1012)</p>	<p>The sample flow was paused due to changes in the droplet shape. Click [Stop] and run [Sort Calibration] again to restore stable sorting. To continue sorting without running the calibration again, click [Settings] on the [Cytometer] tab of the ribbon to display the [Advanced Settings] dialog, and disable [Detect Satellite Droplets]. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration]. A message indicating that the chip is being used a second time appears, but the chip can be used. <i>See "[Detect Satellite Droplets]" (page 140).</i></p>
<p>Control Breakoff initialization timed out. Please re-perform Sort Calibration for stable sorting. (Warning ID : W1021)</p>	<p>Control breakoff could not be enabled. Run [Sort Calibration] again. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration]. A message indicating that the chip is being used a second time appears, but the chip can be used.</p>
<p>Control Breakoff initialization timed out. Please re-perform Sort Calibration for stable sorting. Click [Stop] and re-perform Sort Calibration for stable sorting. (Warning ID : W1023)</p>	<p>Control breakoff could not be enabled. Click [Stop] and run [Sort Calibration] again to restore stable sorting. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration]. A message indicating that the chip is being used a second time appears, but the chip can be used.</p>
<p>Sample flow is paused due to change in droplet condition. (Warning ID : W1024)</p>	<p>The sample flow was paused due to changes in the droplet shape. Run [Sort Calibration] again to restore stable sorting. If the error message is displayed repeatedly, remove the chip using [Chip Exchange], then re-insert the same chip and run [Sort Calibration]. A message indicating that the chip is being used a second time appears, but the chip can be used.</p>
<p>Failed to initialize the sheath pressure regulator. Sheath pressure is too high. Please restart the cytometer and software. If the error reoccurs, please contact your service representative. (Error ID : 00 - 510)</p>	<p>Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.</p>
<p>Failed to initialize the sample pressure regulator. Sample pressure is too high. Please restart the cytometer and software. If the error reoccurs, please contact your service representative. (Error ID : 00 - 511)</p>	

Error message	Solution
High sheath pressure is detected while initializing the regulator. Please contact your service representative. (Warning ID : 00 - 264)	A pressure value outside of rated values has been detected. The laser is still operational, but you should contact your Sony distributor for service.
High sample pressure is detected while initializing the regulator. Please contact your service representative. (Warning ID : 00 - 265)	
Low sheath pressure is detected while initializing the regulator. Please contact your service representative. (Warning ID : 00 - 268)	
Low sample pressure is detected while initializing the regulator. Please contact your service representative. (Warning ID : 00 - 269)	
Failed to move the waste tray. (Warning ID : 00 - 266)	The waste fluid tray is not operating correctly. Contact your Sony distributor for service.
Failed to agitate the sample tube. (Warning ID : 00 - 267)	The agitation unit of the sample loader is not operating correctly. Contact your Sony distributor for service.
Low chip chuck pressure is detected. Please contact your Sony distributor for service. (Warning ID : 00 - 274)	A pressure value outside of rated values has been detected. The laser is still operational, but you should contact your Sony distributor for service.
High chip chuck pressure is detected. Please contact your Sony distributor for service. (Warning ID : 00 - 275)	
Failed to lock the sample area door. (Warning ID : 276)	The sample area door could not be locked. Check the condition of the sample area door.
Invalid deflection voltage is detected. Check if deflection plate is wet or dirty and clean it off. (Warning ID : 00 - 277)	A voltage error was detected with the deflection plates. Check the deflection plates for any dust or moisture, and wipe clean as necessary. If the problem persists, contact your Sony distributor for service.
Sheath tank will be empty if you use automatic bleach/DI cleaning at shutdown. If you want to continue a current sorting, manual bleach/DI cleaning should be execute in the Instrument Cleaning wizard at shutdown. (Warning ID: 278)	The remaining volume of sheath fluid is insufficient if automatic cleaning at shutdown is performed. To continue measurement, perform cleaning manually using the cleaning wizard at shutdown.
Sorting error has been detected. Acquisition was stopped.	A sorting error has occurred. Check that there are no abnormal objects in the collection area. Wait for a short while before restarting sorting. If the problem persists, contact your Sony distributor for service.
Failed to collection load!!	Loading of the collection stage failed. Check that there are no abnormal objects in the collection area. Wait for a short while before loading again. If the problem persists, contact your Sony distributor for service.
Laser output error of the 561nm laser. (Error ID : 00 - 517)	A laser output error has been detected. Contact your Sony distributor for service.
The MD1 board configuration is invalid. (Error ID : 00 - 518)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Emergency stop was activated. Shut down the cytometer. (Error ID : 00 - 519)	The [Emergency Stop] button was pressed. Shut down the main unit.
Chip chuck pressure error. (Error ID : 00 - 520)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Sample door error. (Error ID : 00 - 521)	
Sample door error. (Error ID : 00 - 522)	
Failed to initialize instrument. (Error ID: E1001)	

Error message	Solution
No sample events detected in Coarse mode. (Error ID: E1010) Click [Troubleshooting] to clean sample line.	Click [Troubleshooting] to launch the troubleshooting wizard, and follow the on-screen instructions. See <i>"If an error occurs during auto calibration"</i> (page 51).
No sample events detected in Fine mode. (Error ID: E1011) Click [Troubleshooting] to clean sample line.	
No sample events detected in SuperFine mode. (Error ID: E1012) Click [Troubleshooting] to clean sample line.	
The statistics on the found point is not proper value. (Error ID: E1013)	Stable measurement could not be obtained of the automatic setup beads signal. Check the state of the fluidics lines and automatic setup beads, then run auto calibration again. If the problem persists, contact your Sony distributor for service.
The statistics on the found point is not proper value. (Error ID: E1014)	
Failed to initialize instrument. (Error ID: E2000)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Failed to initialize instrument. (Error ID: E2001)	An error occurred when moving the collection stage. Check that there are no abnormal objects in the collection area, then run auto calibration again. If the problem persists, contact your Sony distributor for service.
Failed to initialize instrument. (Error ID: E2003)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Failed to form appropriate droplet pattern. (Error ID: E3000)	A droplet shape suitable for sorting could not be obtained. Check the state of the fluidics lines, then de-bubble the sheath filter and sorting chip. See <i>"Releasing Air in the Sheath Filter and Sorting Chip"</i> (page 202).
Failed to calibrate the side stream position. (Error ID: E4000)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Failed to calibrate the side stream position. (Error ID: E4010)	Click [Troubleshooting] to launch the troubleshooting wizard, and follow the on-screen instructions. See <i>"If an error occurs during auto calibration"</i> (page 51).
Failed to calibrate the side stream position. (Error ID: E4020)	
Failed to calibrate the side stream angle. (Error ID: E4050)	
Failed to calibrate the side stream position. Droplet is unstable. (Error ID: E4060)	A droplet shape suitable for sorting could not be obtained. Check the state of the fluidics lines, then de-bubble the sheath filter and sorting chip. See <i>"Releasing Air in the Sheath Filter and Sorting Chip"</i> (page 202).
Failed to calibrate the sort delay. (Error ID: 5000)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Failed to calibrate the sort delay. Event rate is unstable. (Error ID: 5010)	An error occurred during sort delay calibration. Check the state of the fluidics lines and automatic setup beads, then run auto calibration again. If the problem persists, contact your Sony distributor for service.
Failed to calibrate the sort delay. Captured beads image is too bright. (Error ID: 5020)	
Failed to calibrate the sort delay. Captured beads image is too dim. (Error ID: 5030)	
Failed to calibrate the sort delay. Captured beads image is too bright. (Error ID: 5040)	
Failed to calibrate the sort delay. Captured beads image is too dim. (Error ID: 5050)	
Failed to calibrate the sort delay. Captured beads image is unstable. (Error ID: 5060)	
Failed to calibrate the sort delay. No sample event is detected. (Error ID: 5070)	

Error message	Solution
Failed to adjust droplet image. (Error ID: E6000)	Shut down the main unit, and wait a while before trying again. If an error with a "00-" prefix is displayed, try the solution provided for the corresponding error ID. If the problem persists, contact your Sony distributor for service.
Failed to adjust droplet image. (Error ID: E6100)	A stable droplet rate suitable for sorting could not be obtained. Check the state of the fluidics lines, then de-bubble the sorting chip. <i>See "Releasing Air in the Sheath Filter and Sorting Chip" (page 202).</i>
Automatic calibration has failed. (Error ID : E0001)	Shut down the main unit, and wait a while before trying again. If the problem persists, contact your Sony distributor for service.
Automatic calibration has failed. (Error ID : E1000)	
Failed to form appropriate droplet pattern. (Error ID : E3100)	A droplet shape suitable for sorting could not be obtained. Check the state of the fluidics lines, then de-bubble the sorting chip. <i>See "Releasing Air in the Sheath Filter and Sorting Chip" (page 202).</i>
Failed to form appropriate droplet pattern. (Error ID : E4100)	



Optional Accessories

- 2) Use with the LE-C3110 100 µm sorting chip is also supported.
 3) Use with the LE-C3113 130 µm sorting chip is also supported.
 4) Use with the LE-C3100 cleaning chip is also supported.
 5) The sheath filter and DI water filter are autoclavable.

For details about purchasing optional accessories, contact your Sony distributor.

Item	Product Number
Sorting chip (70 µm) ¹⁾	LE-C3207 ¹⁾
Sorting chip (100 µm) ²⁾	LE-C3210 ²⁾
Sorting chip (130 µm) ³⁾	LE-C3213 ³⁾
Cleaning chip ⁴⁾	LE-C3200 ⁴⁾
Automatic setup beads	LE-B3001
Sample line	LE-U3F001A
Sample probe	LE-U3F002
PEEK sample line	LE-U3F021
Probe adapter	LE-U3F023
Sheath filter ⁵⁾	LE-U3F003A
Sheath filter bypass line (yellow)	LE-U3F004A
DI water filter bypass line (green)	LE-U3F020
DI water tank kit (with DI water tank filter)	LE-U3F005
Sheath tank	LE-U3F007A
Sintered sheath filter	LE-U3F008
Waste tank	LE-U3F009
Waste tank air filter	LE-U3F011A
Ethanol tank kit (with ethanol tank air filter)	LE-U3F012
Air filter for ethanol and DI water tanks	LE-U3F014
DI water tank filter ⁵⁾	LE-U3F019A
Pressurized air line connection	LE-U3F016
Waste line connection	LE-U3F017
Ethanol line connection	LE-U3F018
Sheath line connection	LE-U3F022
Sample loader O-ring	LE-U3M001
Deflection plate kit	LE-U3M006
Waste catcher	LE-U3M008A
Waste catcher O-ring	LE-U3M009
Rubber packing for collection area door	LE-U3M012
Fluorescent protein filters	LE-U3L001
Splash guard	LE-U3M013
Two-way collection tube holder with embedded cooling agent (5 ml)	LE-A3109
96-well plate holder	LE-A3103
96-well PCR plate holder	LE-A3104
384-well plate holder kit	LE-K3103
384-well plate holder	LE-A3107
384-well PCR adapter	LE-A3108

1) Use with the LE-C3107 70 µm sorting chip is also supported.

Software License

LGPL

This product uses the following LGPL software. The customer is entitled to obtain, modify, and redistribute the source code for the software.

DirectShow.NET

Source codes can be obtained from the “\Others\SourceCode” folder on the supplied installation disc.

However, Sony cannot accept inquiries regarding the content of these source codes.

For license information, refer to the “\Others\License” folder on the supplied installation disc.

Apache License

This product uses the Apache license software, ZXing and Unity.

For license information, refer to the “\Others\License” folder on the supplied installation disc.

Microsoft Public License (Ms-PL)

This product uses the Microsoft Public License (Ms-PL) software, WPF Toolkit, DotNetZip, and Fluent Ribbon Control Suite.

For license information, refer to the “\Others\License” folder on the supplied installation disc.

Third-party Software

This product uses third-party software Entity Framework, .NET Framework, Visual C++ Redistributable Packages, and Microsoft SQL Server. For license information, refer to the “\Others\License” folder on the supplied installation disc.

Environmental Notices

Disposal of Old Electrical & Electronic Equipment (Applicable in the European Union and other European countries with separate collection systems)



This symbol on the product or on its packaging indicates that this product shall not be treated as household waste. Instead it shall be handed over to the applicable collection point for the recycling of electrical and electronic equipment. By ensuring this product is disposed of correctly, you will help prevent potential negative consequences for the environment and human health, which could otherwise be caused by inappropriate waste handling of this product. The recycling of materials will help to conserve natural resources. For more detailed information about recycling of this product, please contact your local Civic Office, your household waste disposal service or the shop where you purchased the product.

Index

Numerics

8-well strips, collection 21, 35

A

AC power connector 24
Air filter, tank 186
Analyzing data 91
Automatic calibration 50

B

Back scatter (BSC) 224
Backup 76
Bandpass optical filter (BPF) 225
Bleach cleaning 170

C

Cart connection cable 25, 26
Chip insertion slot 19
Chip loader 19
Cleaning 172
 Internal sheath line and DI water
 line using ethanol 209
 Optical filter 207
 Reminder 180
 Sample fluidics system using
 bleach 170
Collection area
 Door 18
 Door button 18
Compensation
 introduction 227
Compressed air supply input
 connector 25
Custom sorting mode 117

D

Data analysis 56, 91
Deflection plates
 cleaning 174
Detection matrix, fluorochrome 232
Detection module 20, 225
DI water filter
 autoclaving 204
 De-bubble 203
 replacing 198
DI water tank 23
 cleaning 187
 refilling 167
DISPLAY MODE button 19
Drip tray 23
Droplet
 sorting 222
Droplet camera
 Image 51

Droplet stream calibration 50

E

Environmental notices 258
Error messages 244
Ethanol cleaning 209
Ethanol tank 26
 refilling 169
Experiment
 Export 82
 Import 83
 Searching 84
 Sharing 83

F

Flip-up door 18
Fluidics cart 25
 Cart connection cable 26
 Ethanol tank 26
 Fluidics cart air line connector
 26
 Sheath tank 26
 Waste tank 25
Fluidics cart air line connector 25, 26
Fluidics cart connectors 25
Fluidics maintenance door 23
Fluorescence compensation, see
 compensation
Fluorescent light (FL) 224
Fluorochrome
 detection matrix 232
Forward scatter (FSC) 224

H

Heat exhaust vent 24
Hydrodynamic focusing 221

I

[Initial Instrument Setup] screen 46
Installation disc 11

L

Laser
 configuration 47
 light collection 225
 source 225
 wavelength 229
LCD monitor 19, 30
Login 45
Longpass optical filter (LPF) 225

M

Main power switch 24
Maintenance 161
Maintenance mode 177

O

Operation
 principles 219
Operation flow 43
Optical filter
 cleaning 207
 pattern 230
Optional accessories 257
Overlay display 101

P

Parts
 replacement 257
Password, changing 41
PC connection cable connector 24
Pinch valve 19
Power, turning on 44
POWER/STANDBY button 18
Pulse parameter
 description 226

Q

QR code 46

R

Replacement parts 257

S

Sample fluid
 flow 220
Sample line 19
 changing 192
Sample loader
 Door 18
 Door button 18
 O-ring 181
Sample probe
 changing 192
Scatter light 224
Sheath filter 23
 autoclaving 204
 de-bubble 202
 replacing 196
Sheath fluid
 flow 220
Sheath line
 cleaning 183
Sheath line filter 200
Sheath tank 26
 cleaning 183
 refilling 164
Side stream calibration 50
Side stream monitor
 Image 51
Slide glass, collection 21, 35
Software license 258
Sort delay calibration 50

- Sort parameters 117
 - Manual adjustment 117
- Sorting 58
 - Adjusting 113
 - modeSingle Cell 108
- Sorting chip 14, 27
 - Alignment 50
 - Exchanging 73
 - Loading 46
 - QR code 46
 - thumbwheel 206
- Sorting chip thumbwheel 19
- Sorting mode 106
 - Normal 109
 - Purity 110
 - Semi-Purity 110
 - Semi-Yield 111
 - Ultra Purity 111
 - Ultra Yield 112
 - Yield 112
- Starting 44
- System components 13

T

- Troubleshooting 241

U

- User preferences 72
- Users, adding 40

W

- Waste catcher
 - cleaning 176
 - O-ring 177
- Waste tank 25
 - emptying or replacing 166
- Wavelength, laser 229
- Window extension 227
- Workflow 43



The material contained in this manual consists of information that is the property of Sony Corporation and is intended solely for use by the purchasers of the equipment described in this manual.

Sony Corporation expressly prohibits the duplication of any portion of this manual or the use thereof for any purpose other than the operation or maintenance of the equipment described in this manual without the express written permission of Sony Corporation.

Trademarks

- Microsoft and Windows are registered trademarks of Microsoft Corporation in the United States and/or other countries.
- Intel and Intel Core are trademarks or registered trademarks of Intel Corporation or its subsidiaries in the United States and other countries.
- QR Code is a registered trademark of DENSO WAVE INCORPORATED.
URL: <http://www.qrcode.com/faqpatent.html>
- Alexa Fluor, Qdot, and PE-Texas Red are registered trademarks of Life Technologies Corporation.
- eFluor is a registered trademark of eBioscience.
- Cy and CyDye are trademarks of GE Healthcare, Ltd.
- Brilliant Violet is a trademark of Sirigen, Ltd.
- Pacific Blue and Pacific Orange are trademarks of Life Technologies Corporation.
- S.A.F.E. System is a trademark of Edge Systems LLC.

The products or system names appearing in this document are trademarks or registered trademarks of their respective owners.

Further, the © or ™ symbols are not used in the text.

