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VWR International bvba / spr
B-3001 Leuven

ENGLISH

You will find the Operating Instructions in English language beginning at page 3.

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Dear Customer

Thank you for purchasing a USC ultrasonic cleaning unit from VWR International. The product you have purchased is one of the most innovative in ultrasonics and provides a maximum of performance and comfort.

This documentation was created in order to make all information concerning the installation, handling and operation of ultrasonic cleaning unit available to the operator. This documentation corresponds to the latest release of the date of publication and is not included in revision service. We reserve the right to make technical changes and improvements at any time.

All devices of the USC-series comply with the regulatory affairs and the normative standards. This is indicated by the -symbol on the label.

Any enclosed documentation of other vendors serves only for information. We assume no liability for its completeness or correctness.

If you require information extending beyond this documentation, we will be pleased to offer you technical advice at any time.

Technical service

Web Resources

Visit the VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

Contact us for information or technical assistance contact your local VWR representative or visit www.vwr.com.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you.

VWR International


1. SAFETY INFORMATION	5
2. GENERAL	8
2.1. Intended purpose / lifetime	8
2.2. Cleaning liquid	8
3. START-UP	8
4. OPERATION	9
4.1. Operating elements	9
4.2. Ultrasonic Cleaning	9
4.3. Emptying the tank	9
4.4. Accessories	10
5. MAINTENANCE	10
6. APPENDIX	11
6.1. Technical Data	11
6.2. Operating elements / front view	12
6.3. Connections	12
6.4. Dimensions and weights	12
6.5. Persistency of the front overlay to solvents	13
7. NOTES	38

1. SAFETY INFORMATION

Explanation of the used symbols:

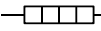Symbol	Explanation
 or 	<p>This symbol points to some danger for persons or the unit, maximum attention is needed in this case. Read the referring chapters especial carefully and follow the requirements strictly.</p> <p>Globally this symbol is also used as reference to follow the operating instructions resp. special advices.</p>
	<p>This symbols shows, that the unit can have hot surfaces.</p>
 or 	<p>Information on the product, that the operating instructions are a part of the unit and must be read in all details.</p>
	<p>CE-sign (Communauté Européenne). The product which carries this symbol meets the requirements of the referring EU-directives (European standards).</p>
	<p>This symbols shows, that the unit must not be disposed with the domestic waste, but collected at special collection points for disposal.</p>
	<p>This symbol marks the unit as unit protection class 1. All electrical conductive parts of the housing are connected to the protective earthing system of the installation, which has earth potential.</p>
	<p>This symbol indicates that the unit has a heating system to warm up the cleaning liquid.</p>

E

Safety Information

Following safety information must be read and respected before start-up and operation of the ultrasonic cleaning unit:

- Any person using this product must have read and understood this manual. Keep this manual close to the device; it must be permanently accessible to the operator. This will prevent operator error and ensure operation free of trouble.
- Check your device for any transport damage. If you find shipping damage, the device must not be connected to the mains.
- The power cord must be removed from the socket only by pulling the plug, do not pull at the cable. The plug must not be touched with wet / damp hands.
- Power cords must be routed without loops and must not be bent or wrapped around sharp edges.
- The intended use of the ultrasonic cleaning unit is described in this documentation and must be followed.
- Insufficient information about handling or operation can cause irreparable damage and injuries. As a result, the unit must be operated by trained personnel only.
- The ultrasonic cleaning unit must be located on a dry, hard and flat horizontal surface.
- The ultrasonic cleaning unit must be connected to the supply voltage specified on the rating label only. The mains connection must be designed for the rated power consumption.

Always use the correct power cord or –adapter for the available socket outlet.

- Only use cleaning liquids which are approved for use with ultrasonic energy.
- Cleaning liquid in use becomes hot. When cleaning temperature sensitive parts check temperature of cleaning liquid frequently.
- Flammable or explosive liquids and liquids which will corrode the tank material (CrNi steel 1.4301) must not be used. Liquids which contain halogens (chlorine, bromine or iodine ions) can cause irreparable damages.
- Never switch on the ultrasonic cleaning unit without cleaning liquid. The cleaning tank must be filled at least to 7/8 of volume (approx. up to the seam). Also ensure that the tank is filled enough during operation.
- The temperature rise of the cleaning liquid caused by heaters or ultrasonic energy can result in steaming. Ensure adequate ventilation at the environment of installation.
- Fluid must not get inside the unit. Wipe off any fluid which has slopped over and is leaking down the side walls as soon as possible
- The items to be cleaned must not be placed directly on the bottom of the tank. Use a suitable cleaning basket.
- Never leave the ultrasonic cleaning unit unattended during operation.
- If the unit does not have to be ready for operation, disconnect it from the power supply (unplug it from mains).
- Never operate the unit if it has been damaged. Inspect the device and the power cord for outer damages frequently. Repairs must only be carried out by the manufacturer.
- Observe the statutory regulations when disposing of the cleaning liquid and disposing the unit.

WARNING:

To avoid risk of **electric shock**, this equipment must only be connected to supply mains with **protective earth (PE)**.

Take special note on using the **proper power cable and/or adapter** for the available power outlet.

WARNING:

There is a danger of **injury and burns** through:

- Ultrasonic energy!
Never grab in the tank with your hands when the ultrasonic energy is switched on. Always switch off the ultrasonic when inserting or removing the items to be cleaned!
- Hot liquid!
- Hot surfaces!
- Toxic vapor!

During operation it is permitted to touch or to handle the unit with the handles or at the control panel only.

After use the unit must cool down min. 30min prior to touching the housing.

WARNING:

For usage of the cleaning unit and handling with cleaning agent you generally have to wear adequate personal protective equipment (PPE) corresponding to the standards:

- Safety gloves
- Safety glasses
- Ear protection
- Breathing protection
- Protective clothing
- Protective boots

WARNING:

No modification of this equipment is allowed.

Do not carry out any repair. In case of service, send the unit back to the manufacturer.

General / Start-up

2. GENERAL

2.1. Intended purpose / lifetime

The USC ultrasonic cleaning units were especially developed for cleaning tasks in laboratory facilities of commercial use.

The expected life time of the USC ultrasonic cleaning units depends basically on the quality of maintenance and cleaning. With proper use, maintenance, cleaning and any necessary repairs a lifetime of five years can be expected. This does not include consumables such as cable / connectors, cleaning baskets and similar.

Only accessories which are approved by or can be obtained from the manufacturer have to be used.

2.2. Cleaning liquid

The cleaning liquid which has to be used depends on the material of the items to be cleaned and the type of soiling. Please contact the supplier of the cleaning agent.

3. START-UP

- Check if the mains connection, which must have a protective earth conductor, confirms with the specification on the rating label of your device.
- Connect the power socket connector at the rear of the unit. If the power cord is already connected, check that this is done properly.
- The ultrasonic cleaning unit must be located on a dry, hard and flat horizontal surface.
- Fill the cleaning tank with a suitable cleaning liquid depending on the application. The level must be at least 7/8 of volume (approx. up to the seam).

Take note of the respective safety regulations for handling chemicals.

- Connect the ultrasonic cleaning unit to the mains, the unit is ready for operation.

WARNING:

To avoid risk of **electric shock**, this equipment must only be connected to supply mains with **protective earth**.

Take special note on using the **proper power cable and/or adapter** for the available power outlet. As a result of the design, the EU connector (CEE 7/7) fits mechanically into the DK socket outlet (Denmark) but does not connect the protective earth conductor (PE).

**Never use the EU connector in combination with DK socket outlets
DANGER TO LIFE AND LIMB!**

Adapters are one-time mountable only and cannot be removed after applying to the connector. Therefore be sure to mount the proper adapter.

WARNING:

The power plug is used to disconnect the unit from mains.

The used **power outlet** must be **permanently accessible** and must not be covered by the unit or any other objects, so the mains plug can be unplugged at any time.

4. OPERATION

Following configurations were available:

Configuration	Series	
	T	TH
Table top cleaning unit, ultrasonic frequency approx. 45kHz	X	X
Timer for ultrasonic operation adjustable up to 30 minutes or continuous	X	X
Integrated heater		X

4.1. Operating elements

The devices have a mechanical timer which allows an adjustable time set of ultrasonic operating up to 30 minutes or continuous operation. For on-times below 2 minutes, first of all turn the knob to a longer time and then turn it back to the desired value. The continuous ultrasonic operation is obtained by turning the rotary knob anti-clockwise from the '0' position over the latching point.

The ultrasonic energy applied to the fluid is visible in the form of undulations on the surface of the fluid.

USC100TH: An orange indicator in the heater switch lights up when the heater is enabled. The temperature is preset fixed to 65degC.

USC200T, TH: During ultrasonic operation a green indicator lights up. The USC200TH has additionally a green indicator when the heater is enabled and a yellow indicator if the set temperature is not reached. To adjust the desired set temperature turn the knob to the desired value (in degC).

4.2. Ultrasonic cleaning

- Immerse the items to be cleaned in the cleaning bath, the distance of the items from the bottom must be at least 20 mm (0.8 inches). If required, use the cleaning basket which is available as an accessory.
- Set the required cleaning time and temperature for the cleaning process (if applicable). The value excessively depends on the cleaning fluid, temperature, type and degree of soiling of the items requiring cleaning. The time can vary from a few seconds to a few minutes. The best combination of the parameters may have to be experimentally determined.
- Start the cleaning process by switching on the ultrasonic operation.
- After the cleaning process remove the items out of the liquid with suitable care. We recommend rinsing the parts in clean water.
- Disconnect the ultrasonic cleaning unit from mains if the unit is not needed.

WARNING:

Consider the **temperature tolerance** of the material to be cleaned!

Please note that the bath temperature rises considerably with duration of the ultrasonic operation and can exceed the set temperature.

Never put the items to be cleaned directly on the bottom of the tank.

Never clean too many items at once, the cleaning result can deteriorate.

4.3. Emptying the tank

If the cleaning result becomes worse or if the cleaning liquid has got contaminated, the cleaning liquid must be replaced.

Prior to any emptying or cleaning the unit must be disconnected from mains.

4.4. Accessories

- **Tank cover (stainless steel)**
To prevent evaporation of the tank contents and odor nuisance.
- **Cleaning basket (stainless steel)**
For carrying the items to be cleaned. The basket can be suspended in the cleaning tank and it also can be placed on the edge of the tank with the handles for draining.
- **Cover with openings for beakers (different sizes)**
 - easy cleaning of small parts
 - less needs of cleaning solvent
 - cleaning in different cleaning solvents

The ultrasonic energy is coupled inside the filled beaker by the liquid around through the glass. Fill the tank with water and add a surface-active agent (e.g. washing-up liquid).

5. MAINTENANCE

WARNING:

The ultrasonic cleaning unit needs low maintenance.
Regard following notes:

- Prior to any emptying or cleaning the unit must be disconnected from mains.
- Keep the device clean. Wipe the housing frequently with a damp cloth. Remove cleaning agent residues immediately.
- Ensure that the cleaning liquid does not soil the housing during cleaning.
Clean, replace and dispose the cleaning liquid in accordance with the supplier's recommendations.
- Avoid scratching the inner tank surface, this accelerates cavitation (erosion). Ensure the items to be cleaned have no direct contact with the bottom or inner sidewalls of the tank and try to keep none or only small amounts of pollution on the bottom of the cleaning tank. If necessary, the cleaning liquid should be drained off prematurely, so the tank can be cleaned.
- Check the unit and the power cord for damages frequently and prior to any use.
A visibly damaged unit or cable must not be put in operation and must be disconnected from the mains.
Do not carry out repairs yourself. Send the unit to the manufacturer for servicing or repair.
- There are any parts to maintain or which have to be changed by the user inside the unit.

If the power cord is damaged, it must be replaced with the manufacturer's original part only.

6. APPENDIX

6.1. Technical Data

Intended Use Cleaning applications in laboratory facilities of commercial use.

Executions Table top cleaning unit

Unit series	Execution
USC - T	<ul style="list-style-type: none"> ultrasonic energy with approx. 45kHz
USC - TH	<ul style="list-style-type: none"> ultrasonic energy with approx. 45kHz built-in liquid heater

Output ultrasonic energy with frequency of approx. 45kHz
switch-on duration: 100% (continuous operation)

Environment conditions

Operating ambient temperature: 0degC ... +40degC (32F ... 104F)
 Transportation & storage temperature: - 20degC ... +65degC (- 4F ... 149F)
 Humidity: 20% ... 70%, non-condensing
 Altitude (sea level): max. 2000m (a.s.l.)

Classification:

Protection class: class I equipment (earthed housing)
 Protection category according IEC 529: IP 21
 Moisture degree 2
 Over voltage category II


Mains connection 230V-versions: AC230V ±10%, 47...53Hz, 1~, N, PE / 16A

Unit type (230V)		HF-power		connected load		heater power
		peak	effective	'T'	'TH'	'TH'
USC 100	T	60W	30W	60VA	120VA	60W
USC 200	TH	120W	60W	100VA	180VA	90W


6.2. Operating elements / front view


USC100T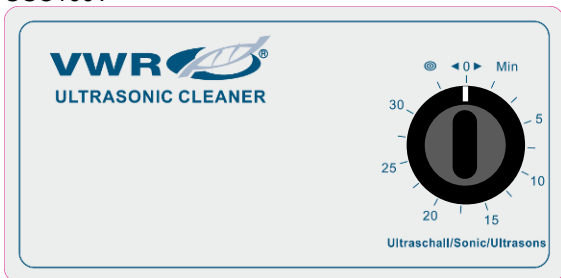


USC100TH

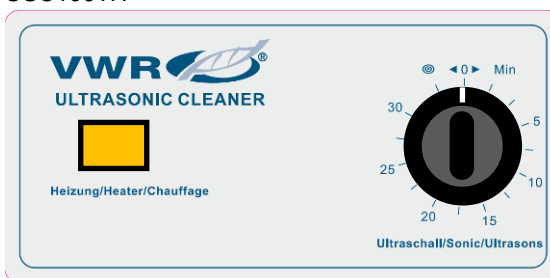USC200T

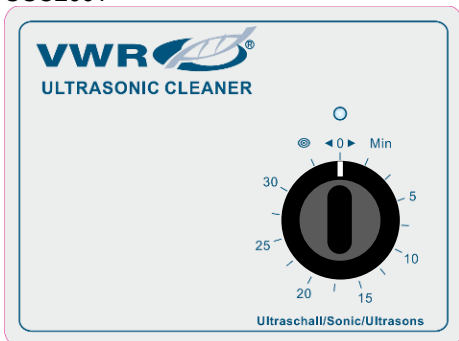USC200TH

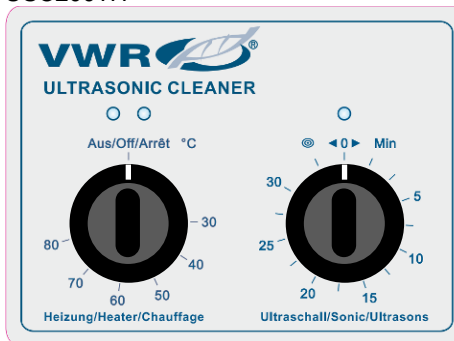6.3. Connections

Mains connection, single phase 230V:
 power cord detachable with power plug CEE 7/7
 cord length approx. 2.0m (6.5ft)

- Adapters available for
- United Kingdom – BS-1363
 - Switzerland – SEV1011
 - Denmark – DS 60884-2-D1
 - China CPCS-CCC / Australia – AS 3112

Attention:

Adapters are one-time mountable only and cannot be removed after applying to the connector. Therefore be sure to mount the proper adapter.

6.4. Dimensions and weights

Unit type	Tank capacity (liters)	inner dimensions (LxWxD / mm)	outer dimensions (LxWxD / mm)	weight (kg)
USC 100	0.8	190 x 85 x 60	205 x 100 x 155	1.8
USC 200	1.8	150 x 140 x 100	175 x 165 x 270	2.6

6.5. Persistency of the front overlay to solvents

The front overlay is based on a polyester foil with a biaxial alignment and has therefore a higher persistence against solvents compared with e.g. polycarbonate or PVC. It is stronger and more durable than other foils generally used for plastic foil keyboards and front panels.

The material is resistant against following chemicals over a period of more than 24 hours without any perceptible changes:

Ethanol Cyclohexane Diaceton alcohol Glycol Isopropanol Glycerin Methanol Triacetine Dowandol DRM/PM	Formaldehyde 37 – 42 % Acetaldehyde Aliphatic Toluene Xylene Thinning agent (white spirit)	Trichloroethylene Ethyl acetate Diethyl ether N-butyl acetate Amyl acetate Butylcellosolve Ether
Acetone 2-butanone Dioxane Cyclohexane MIBK Isophorone	Formic acid < 50 % Acetic acid < 50 % Phosphoric acid < 30 % Hydrochloric acid < 36 % Nitric acid < 10 % Trichloroacetic acid < 50 % Sulphuric acid < 10 %	Soda chlorine < 20 % Hydrogen peroxide < 25 % Potash soft soap Detergents Tensides Softening agents Iron (II) chloride (FeCl ₂) Iron (III) chloride (FeCl ₃) Dibutyl phthalate Dioctyl phthalate Sodium carbonate
Ammonia < 40 % Sodium hydroxide solution < 40% Potassium hydroxide Alkali carbonate Dichromate Potassium hexacyanoferrate Acetonitrile Sodium bisulphate	Cutting emulsions Diesel oil Varnish Paraffin oil Castor oil Silicone oil Turpentine substitute Brake fluid Decon Aviation fuel Petrol Water Salt water	

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India

VWR Lab Products Private Limited
No.139. BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India – 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR
International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02-3320311/02-487791
Email: info.it@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 020 4808 400
Email: info.nl@vwr.com

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: +47 22 90 00 00
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 058 32 38 210
Email: info.pl@vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Alfragide
Tel.: +351 21 3600 770
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnologia 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 08 621 34 00
Email: info.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 044 745 13 13
Email: info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
Email: uksales@vwr.com

**GO TO VWR.COM FOR THE
LATEST NEWS, SPECIAL OFFERS
AND DETAILS OF YOUR LOCAL
VWR DISTRIBUTOR**